

PrimePCR SARS-CoV-2 Multiple Mutation RT-qPCR Assay Using Reliance One-Step Multiplex RT-qPCR Supermix Protocol

Description

The PrimePCR SARS-CoV-2 Multiple Mutation RT-qPCR Assay* is a single-well multiplexed panel which can be validated by researchers to detect and differentiate four key mutations identified in the severe acute respiratory syndrome coronavirus 2 (SARS-CoV-2) genome. The mutations targeted by the assay are believed to confer evolutionary advantages to the virus, such as increased transmissibility, disease severity, and ability to evade vaccinations and treatments. The assay's high sensitivity can be validated by researchers to assess low-level samples with SARS-CoV-2 quantification cycle (Cq) values up to 40. The single tube assay is 20x concentrated and contains a primer set plus up to six probes, including one probe that is the multiplexed control (MPC) (Tables and Figures 1–2). This probe has homology to a conserved region between the primers and serves as an internal RNA integrity and experimental setup control. The other probes are perfect matches to the four mutations of interest in the spike protein gene. Specific mutation(s) in the sample can be evaluated by researchers to assess the likelihood of a specific variant in the sample. The assay format is optimal for use with the Reliance One-Step Multiplex RT-qPCR Supermix and should be run on a real-time PCR detection system equipped with five fluorescence channels, such as the CFX96 Touch or CFX Opus 96 Real-Time PCR System.

* For research use only (RUO). Not for use in diagnostic procedures. ** The multiplexed control (MPC) is designed in a conserved region of the S gene. MPC functionality may be compromised if the virus acquires mutations in this region.

Table 1. Detectable mutations for assay 1.

SARS-CoV-2 Variant	K417N	L452R	MPC	E484K	E484Q	N501Y
Alpha (B.1.1.7)			•			•
Beta (B.1.351)	•		•	•		•
Delta (B.1.617.2)		•	•			
Epsilon (B.1.427/B.1.429)		•	•	•		
Gamma (P.1)			•	•		•
Kappa (B.1.617.1/ B.1.617.3)		•	•		•	

MPC, multiplexed control.

Assay 1

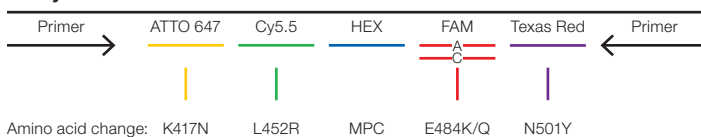


Fig. 1. Schematic depiction of the PrimePCR SARS-CoV-2 Multiple Mutation Assay (assay 1). MPC, multiplexed control; RT-qPCR, reverse transcription quantitative PCR.

Table 2. Detectable mutations for assay 2.

SARS-CoV-2 Variant	K417T	L452R	MPC	T478K	N501Y
Alpha (B.1.1.7)			•		•
Beta (B.1.351)			•		•
Delta (B.1.617.2)		•	•	•	
Epsilon (B.1.427/B.1.429)		•	•		
Gamma (P.1)	•		•		•
Kappa (B.1.617.1/ B.1.617.3)		•	•		

MPC, multiplexed control.

Assay 2

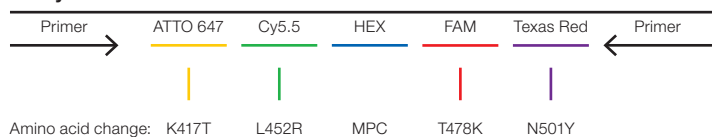


Fig. 2. Schematic depiction of the PrimePCR SARS-CoV-2 Multiple Mutation Assay (assay 2). MPC, multiplexed control; RT-qPCR, reverse transcription quantitative PCR.

Additional assays will become available as the virus continues to evolve and the need to identify novel variants arises. Visit [bio-rad.com/SC2Variants](https://www.bio-rad.com/SC2Variants) for a complete list of assays available.

Storage and Stability

The assays are stable for 12 months when stored at 4°C, protected from light. The 20x assay mix can be kept at -20°C for long-term storage.

Product Contents

This product is a 20x concentrated, ready-to-use primer/probe mix, which can be validated for use with the Reliance One-Step Multiplex RT-qPCR Supermix for detection of four mutations in a single well. It is provided as 200, 1,000, or 2,500 µl volumes, sufficient for 200, 1,000, or 2,500 x 20 µl reactions.

Suggested Reagents and Equipment*

- Reliance One-Step Multiplex RT-qPCR Supermix (catalog #12010220, 12010221)
- Reference control: SARS-CoV-2 Standard (#COV019)
- SARS-CoV-2 Variant Controls:
 - SARS-CoV-2 S Gene Alpha Variant Control (B.1.1.7) (#COVA)
 - SARS-CoV-2 S Gene Beta Variant Control (B.1.351) (#COVB)
 - SARS-CoV-2 S Gene Epsilon Variant Control (B.1.427 and B.1.429) (#COVE)
 - SARS-CoV-2 S Gene Gamma Variant Control (P.1) (#COVG)
- Negative control: SARS-CoV-2 Negative (#COV000)
- CFX Opus 96 or CFX96 Touch Real-Time PCR System (#12011319 or 1855195, respectively) or equivalent 5-color instrument
- Hard-Shell 96-Well PCR Plates (#HSP9655, HSP9955) or equivalent
- Microseal 'B' PCR Plate Sealing Film (#MSB1001)

* Bio-Rad reagents and equipment, unless otherwise noted. Reference, variant, and negative controls are from Exact Diagnostics, a Bio-Rad company.

Sample Preparation

Sample Type

The optimal sample type is total RNA extracted from human specimen samples. Alternative sample types may be compatible but should be optimized and validated before use by the researcher. Assay performance depends on the amount and quality of template RNA validated. The PrimePCR SARS-CoV-2 Multiple Mutation Assay is for research use only and is not intended for patient diagnosis from clinical samples.

Reference Control

The SARS-CoV-2 Standard (#COV019) from Exact Diagnostics is manufactured with synthetic RNA transcripts containing five gene targets (the *E*, *N*, *ORF1ab*, *RdRp*, and *S* genes of SARS-CoV-2). The product is optimal to validate reverse transcription, amplification, and detection of SARS-CoV-2 RNA. This product also contains human genomic DNA, which can serve as an amplification control (*RPP40*). Add 10 µl per reference control reaction; extraction or dilution should not be required.

Positive Variant Controls

The SARS-CoV-2 S Gene Variant Controls from Exact Diagnostics are individually manufactured with synthetic RNA transcripts of the S gene from SARS-CoV-2 variants of interest or concern.

These controls contain human genomic DNA, which serves as an amplification control. These variant controls can be validated for use to monitor the reverse transcription, amplification, and detection steps in assays that target S gene mutations. Add 10 µl per variant control reaction; extraction or dilution should not be required.

gBlocks Gene Fragments from Integrated DNA Technologies, Inc. are DNA oligos, which are optimal as positive controls in assays that target mutations other than the S gene. Ordering information can be found in the Assay Control tab of each PrimePCR SARS-CoV-2 Multiple Mutation Assay product, or the PrimePCR SARS-CoV-2 Multiple Mutation Assay product page. Dilute gBlocks Controls in 1x Tris-EDTA (TE) buffer to 10⁵–10⁶ copies/µl (50–500 fM). Use 5 µl per variant control reaction.

Negative Control

The SARS-CoV-2 Negative (#COV000) from Exact Diagnostics is formulated in a synthetic matrix and contains genomic DNA. It should be used as negative reference material when testing for SARS-CoV-2. Add 10 µl per negative control reaction; extraction or dilution should not be required.

No Template Control (NTC)

This control can be validated by researchers to monitor for reagent and environmental contamination of the reactions. Use 10 µl of nuclease-free water per NTC reaction. A minimum of two NTC wells per plate is required to analyze data using the allelic discrimination module in CFX Maestro Software.

Bio-Rad Instrument Setup Example

The following instructions are suggested for researchers to develop a validation protocol in their laboratory. The example shown here is for the CFX Opus 96 and CFX96 Touch Real-Time PCR Systems, which are research instruments equipped with five fluorescence channels. For more detailed information, refer to the respective instrument manual.

Cycling Protocol Setup

1. Click **File > New > Protocol** in the menu bar to open the Protocol Editor.
2. Change Sample Volume to 20 µl.
3. Modify the thermal cycling protocol according to the guidelines in Table 3.

Note: Up to 40 cycles is optimal as less than 40 cycles can be a positive result.

Table 3. Example of thermal cycling protocol.*

Step	Description	Temperature, °C	Time	Number of Cycles	
1	Reverse transcription	50	10 min	1	
2	DNA polymerase activation and template denaturation	95	10 min	1	
3		Template denaturation	95	10 sec	45
4	Amplification	Annealing/extension and plate read	58	30 sec	

* Set the lid temperature to 105°C.

- Confirm the annealing/extension step includes a plate read, as indicated by a camera symbol in the step.
- To add a plate read, click on the step to highlight, then click **Add Plate Read to Step**.
- Save the protocol by clicking **File > Save As**, assign a name, and click **Save**.
- Click **OK** to exit the Protocol Editor.

Plate Setup

- Click **File > New > Plate** in the menu bar to open the Plate Editor.
- Select **Settings > Plate Size > choose 96 Wells**.
- Select **Settings > Plate Type > choose BR White**.
- Expand the dropdown menu to the right of Scan Mode and select **All Channels**.
- Highlight the wells where samples and controls will be on the plate. To highlight all the wells, click the upper left corner of the plate graphic.
- Click **Select Fluorophores** and choose FAM, HEX, Texas Red, Cy5,* and Cy5.5 by checking the box to the right of the selected fluorophore (uncheck **SYBR**). Click **OK** to apply changes.

- Define the sample type for each well by highlighting the wells, then choosing the appropriate identifier from the Sample Type dropdown menu.

Note: If samples contain very low amounts of viral RNA, consider adding a second set of control samples that contain SARS-CoV-2 Variant Controls and the SARS-CoV-2 Standard diluted in TE buffer at 1:100, and undiluted.

- Apply target names and fluorophores to all wells by highlighting the wells, then checking the Load box to the left of each of the fluorophores listed in the Target Name section. To include the target names, replace <none> in the open text box to the right of each fluorophore. For example, if using PrimePCR SARS-CoV-2 Multiple Mutation Assay 1:
 FAM — 484K/Q
 HEX — MPC
 Texas Red — 501Y
 Cy5 — 417N
 Cy5.5 — 452R
- Save the file by clicking **File > Save As**.
- Name the plate file and click **Save**.
- Click **OK** to exit the Plate Editor (Figure 3).

* Assays utilize an ATTO 647 fluorophore, which has a similar excitation and emission spectra to Cy5.

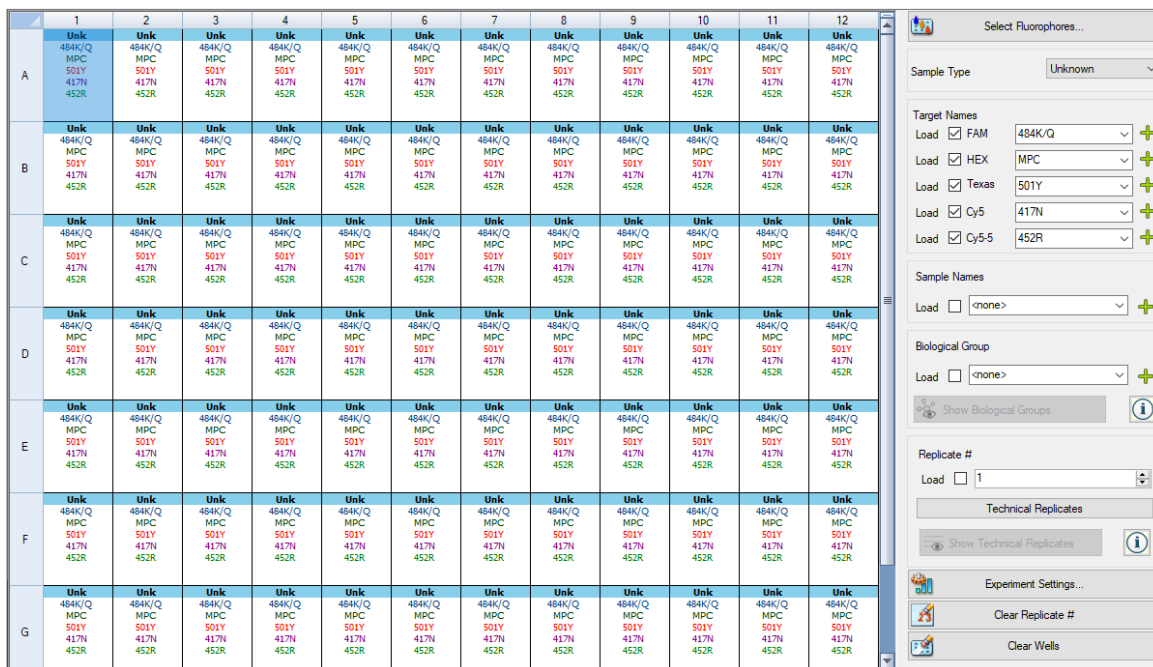


Fig. 3. Plate layout in CFX Maestro Software Plate Editor.

Suggested Reaction Setup

1. Thaw the Reliance One-Step Multiplex RT-qPCR Supermix on ice. Mix thoroughly by vortexing the tube to ensure homogeneity and resuspend any precipitates that may have formed during storage. Centrifuge briefly to collect contents at the bottom of the tube. Store on ice protected from light until ready to use.

Note: Reliance One-Step Multiplex RT-qPCR Supermix is a 4x ready-to-use solution. The viscosity of this solution is greater than a 2x mix. If using an automated liquid handling system, allow the supermix to equilibrate at room temperature for 10 min to reduce its viscosity.

2. Thaw the PrimePCR SARS-CoV-2 Multiple Mutation Assay at room temperature. Mix thoroughly by vortexing the tube to ensure homogeneity. Centrifuge briefly to ensure that all components are at the bottom of the reaction tube. Store protected from light.
3. Thaw samples on ice. Place the SARS-CoV-2 Standard, Negative, and Variant Controls (stored at 4°C) on ice. Mix thoroughly by vortexing each tube to ensure homogeneity. Centrifuge briefly to ensure that all components are at the bottom of each reaction tube.

Note: The controls are supplied at 200,000 copies/ml and can be validated for use in these assays without dilution and without extraction. However, if samples contain very low amounts of viral RNA (<300 copies/ml or have Cq values ≥30), it is recommended to prepare diluted SARS-CoV-2 Variant Controls (1:100) so that the control wells cross the threshold at approximately the same cycle as the samples.

4. Prepare a master mix volume sufficient for the number of samples and controls to be tested, plus 10% more to account for pipetting variability (Table 4). The volumes in Table 4 are based on 10 µl of sample being added to a final volume of 20 µl.

Table 4. Example of preparation of master mix.

Component	Volume per Reaction, µl	Volume for n Samples, µl
Reliance One-Step Multiplex RT-qPCR Supermix	5 µl	= (5 x n) x 1.1
20x PrimePCR SARS-CoV-2 Multiple Mutation Assay	1 µl	= (1 x n) x 1.1
Nuclease-free water*	4 µl	= (volume x n) x 1.1

* Adjust volume to accommodate larger sample volume. We recommend using 10 µl of sample. In each 20 µl reaction, the volume of sample plus water should equal 14 µl.

5. To obtain a final volume of 20 µl, pipet 10 µl of master mix into each well of a Hard-Shell 96-Well PCR Plate.
6. Add 10 µl of RNA sample, reference control, variant control, negative control, or nuclease-free water (no template control) to the appropriate well of the reaction plate for a final volume of 20 µl. Adjust volume, if necessary.

7. Seal the plate with Microseal 'B' PCR Plate Sealing Film and vortex to mix reagents. Centrifuge the plate briefly to ensure that all components are at the bottom of the reaction tube.
8. Load the PCR plate into the CFX96 Touch Real-Time PCR System, select the plate and protocol files defined in earlier steps, and start the run.


Suggested Data Analysis

Analysis should be performed in a three-step process. First, adjust the baseline and threshold setting for each fluorophore. Second, confirm that reactions were set up properly and that all samples contain amplifiable RNA by assessing the MPC amplification curve in the HEX channel for each sample. Third, assess the presence or absence of each mutation by determining whether the amplification plot indicates exponential growth and crosses the threshold for each fluorophore.

Baseline-Threshold Setting

The baseline and threshold settings should be confirmed and manually defined, if necessary, in the Quantification tab of CFX Maestro Software.

1. All detection channels should be selected by default during plate setup; confirm before proceeding to step 2.
2. Select **Settings > Cycles to Analyze**.
3. Change lower cycles to 5 and upper cycles to 45.
4. Click **OK** to apply changes and exit.

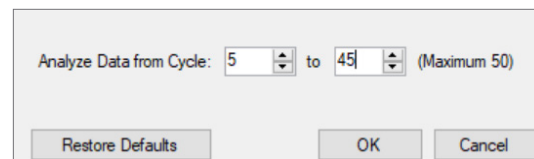

5. Select **Settings > Baseline Setting > Apply Fluorescence Drift Correction**.
6. Deselect the check boxes for HEX, Texas Red, Cy5, and Cy5.5 below the amplification plot (only **FAM** should be selected).
7. Select **Log Scale** by checking the box in the lower right corner of the amplification plot. Change the log scale to a minimum value, such as 0.5, so that the baseline signals are visible.
8. Click and drag the bold horizontal threshold line until it lies in the exponential phase of the fluorescence curve and above any baseline signal.
9. Repeat steps 4–8, for each of the remaining fluorophores (HEX, Texas Red, Cy5, and Cy5.5).

Confirm Expression for Multiplexed Control

MPC (HEX) can be an internal sample control. The probe has homology to a highly conserved region between the primers. Failure to amplify MPC may suggest incorrect reaction setup, inadequate reagent quality, or insufficient intact RNA by the researcher's method. Alternatively, the sample may contain a new mutation that impacts the ability of the probe to bind.

1. Select **HEX** by unchecking all other fluorophores in the boxes under the amplification plot.
2. Confirm that all wells identified by researchers as unknowns cross the threshold before cycle 45 in the HEX channel.
3. Exclude from analysis any well containing an unknown that fails to cross the threshold before cycle 45 in the HEX channel.

Assess Presence or Absence of Mutations

1. Select **FAM** by unchecking all other fluorophores in the boxes under the amplification plot.
2. Identify unknowns that may contain the mutation associated with the FAM fluorophore by determining whether the amplification curves in the FAM channel display exponential growth and cross the threshold before cycle 40.
3. Repeat step 1 and 2 for fluorophores Texas Red, Cy5 (ATTO 647), and Cy5.5 sequentially to identify unknowns that contain mutations that encode the other targeted mutations.

Data Interpretation

Researchers can determine the likelihood that a sample contains a variant of interest by considering which of the mutations are assessed as positive in the sample. The currently reported most common SARS-CoV-2 variants of interest or concern that can be identified using the PrimePCR SARS-CoV-2 Multiple Mutation Assay are listed in Table 5. To determine the likelihood that a sample contains a specific variant, researchers should identify which fluorophore(s) cross the threshold before cycle 45 in a sample, and then confirm the sample is positive for MPC. Data can be compared to the information about currently reported variants of interest or concern in Table 5. For example, using PrimePCR SARS-CoV-2 Multiple Mutation Assay 1, if a sample is determined to be positive for FAM (E484K/Q), Texas Red (N501Y), ATTO 647 (K417N), and HEX (MPC), then the sample mutation profile may be consistent with the Beta variant (B.1.351).

Table 5. Example of mutation fingerprint of SARS-CoV-2 variants of interest or concern for assay 1.

WHO Label	Pango Lineage	E484K/Q (FAM)	N501Y (Texas Red)	K417N (ATTO 647)	L452R (Cy5.5)	MPC (HEX)
Alpha*	B.1.1.7	—	•	—	—	•
Beta	B.1.351	•	•	•	—	•
Gamma	P.1	•	•	—	—	•
Delta**	B.1.617.2	—	—	—	•	•
Epsilon**	B.1.427/B.1.429	—	—	—	•	•
Kappa	B.1.617.1	•	—	—	•	•
No designation***	B.1.617.3	•	—	—	•	•

WHO, World Health Organization.

* Some Alpha strains contain the E484K mutation, but the majority do not. Those that do contain the E484K mutation will have a mutation pattern similar to the Beta variant. In such cases, the PrimePCR SARS-CoV-2 Single Mutation Assay targeting the D80A or A701V assay for Beta or the HV69-70 deletion assay for Alpha can be used by researchers to further distinguish the viral strain present in the sample.

** In this panel, it is not possible to distinguish the Delta variant from the Epsilon variant. To differentiate between these strains, the singleplex T478K assay can be used by researchers.

*** Additional PrimePCR SARS-CoV-2 Multiple Mutation Assays will become available as the virus undergoes additional mutations and a need to survey for novel mutations arises. Visit bio-rad.com/SC2Variants for a complete list of assays available and evaluation of results.

Bio-Rad PCR Plastics, Reagents, and Equipment

Catalog #	Description
MSB1001	Microseal 'B' PCR Plate Sealing Film , adhesive, optical, pkg of 100
HSP9655	Hard-Shell 96-Well PCR Plates , low profile, thin wall, skirted, white shell/white well
HSP9955	Hard-Shell 96-Well PCR Plates , low profile, thin wall, skirted, white shell/white well, barcoded
12010220	Reliance One-Step Multiplex RT-qPCR Supermix , 1,000 x 20 µl reactions
12010221	Reliance One-Step Multiplex RT-qPCR Supermix , 2,000 x 20 µl reactions

Instruments

12011319	CFX Opus 96 Real-Time PCR System
1855195	CFX96 Touch Real-Time PCR System

Exact Diagnostics SARS-CoV-2 Reference Control (RUO)

COV019	SARS-CoV-2 Standard
--------	---------------------

Exact Diagnostics SARS-CoV-2 Variant Mutation Controls (RUO)

COVA	SARS-CoV-2 S Gene Alpha Variant (B.1.1.7)
COVB	SARS-CoV-2 S Gene Beta Variant (B.1.351)
COVE	SARS-CoV-2 S Gene Epsilon Variant (B.1.427 and B.1.429)
COVZ	SARS-CoV-2 S Gene Gamma Variant (P.1)

Exact Diagnostics SARS-CoV-2 Negative Control (RUO)

COV000	SARS-CoV-2 Negative
--------	---------------------

PrimePCR SARS-CoV-2 Multiple Mutation Assays

Refer to [bio-rad.com/SC2Variants](https://www.bio-rad.com/SC2Variants) for assays and ordering information.

Visit [bio-rad.com/SC2Variants](https://www.bio-rad.com/SC2Variants) for a complete list of PrimePCR SARS-CoV-2 Mutation Assays and ordering information.

Visit [bio-rad.com/PCRPlastics](https://www.bio-rad.com/PCRPlastics) for a full range of PCR plastics and seals.

Visit [bio-rad.com/RelianceOneStep](https://www.bio-rad.com/RelianceOneStep) for more information.

Visit [ExactDiagnostics.com](https://www.ExactDiagnostics.com) for more information about these controls. [Q1678 COVID-19 Quality Control Solutions](#) contains a complete list of SARS-CoV-2 controls.

Bio-Rad Technical Support

The Bio-Rad Technical Support department in the U.S. is open Monday through Friday, 5:00 AM to 5:00 PM, Pacific time.

Phone: 1-800-424-6723, option 2

Email: Support@Bio-Rad.com (U.S./Canada only)

For technical assistance outside the U.S. and Canada, contact your local technical support office or click the Contact us link at [bio-rad.com](https://www.bio-rad.com).

BIO-RAD, HARD-SHELL, and MICROSEAL are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions. SYBR is a trademark of Thermo Fisher Scientific Inc. All trademarks used herein are the property of their respective owner.

For research use only. Not for use in diagnostic procedures.



**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Website [bio-rad.com](https://www.bio-rad.com) **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399
Canada 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23
France 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300 **Israel** 0 3 9636050
Italy 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 2 3473 4460 **Luxembourg** 00 800 00 24 67 23 **Mexico** 52 555 488 7670
The Netherlands 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23 **Portugal** 00 800 00 24 67 23
Russian Federation 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23 **Spain** 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23
Switzerland 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311 **United Arab Emirates** 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23

