

Plate Lifter for 384-Well Reaction Module



The Plate Lifter for 384-Well Reaction Module is designed for use with 384-well plates in the following instruments:

- PTC Tempo 384 Thermal Cycler
- CFX Opus 384 Real-Time PCR System
- C1000 Touch Thermal Cycler with 384-Well Reaction Module
- S1000 Thermal Cycler with 384-Well Reaction Module

The Plate Lifter for 384-Well Reaction Module aids in the removal of a 384-well plate by gently lifting it slightly out of the reaction block when the heated lid is opened. Install the Plate Lifter for 384-Well Reaction Module as necessary. Save the 384-Well Plate Lifter for all potential future uses.

Instructions

1. Adjust orientation.


Place the plate lifter around the outside of the 384-well block, with the lifters facing up.

2. Install the plate lifter.

Gently press down on the long edges of the plate lifter until the four notches on the short side of the plate lifter push against the 384-well block.

3. Ensure proper installation.

Examine the plate lifter to ensure that it is properly installed.

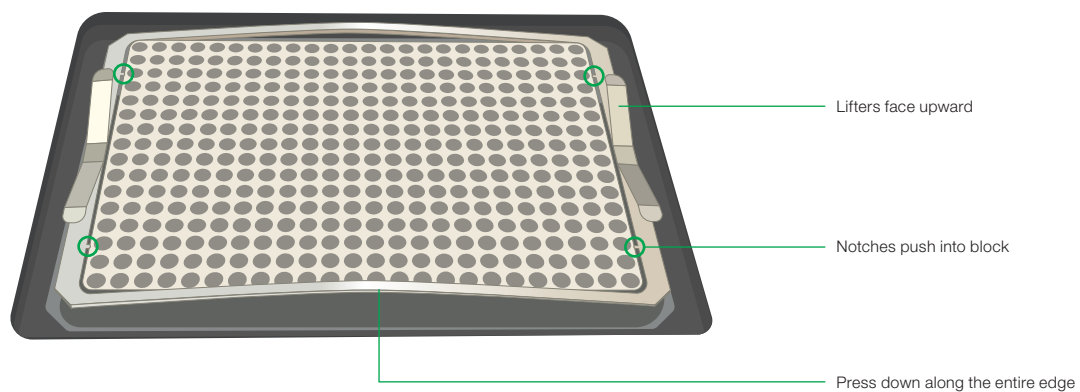



Bio-Rad PCR reagents and analytical instruments are manufactured under an ISO 13485:2016 certified Quality Management System and are quality control tested to ensure consistent product performance you can trust.

BIO-RAD is a trademark of Bio-Rad Laboratories, Inc. All trademarks used herein are the property of their respective owner.
© 2024 Bio-Rad Laboratories, Inc.



**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Website bio-rad.com **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399
Canada 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23
France 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300
Israel 0 3 9636050 **Italy** 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 080 007 7373 **Luxembourg** 00 800 00 24 67 23
Mexico 52 555 488 7670 **The Netherlands** 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23
Portugal 00 800 00 24 67 23 **Russian Federation** 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23
Spain 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23 **Switzerland** 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311
United Arab Emirates 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23

