



Getting Started with ChemiDoc™ Go and GelDoc™ Go Sample Trays

Quick Guide

The following trays, accessories, and applications are supported for use with the ChemiDoc Go and GelDoc Go Imaging Systems. The UV/Stain-Free Tray ships with the instrument; all other trays are optional accessories.

Tray Type or Accessory	Description
UV/Stain-Free Tray for ChemiDoc Go and GelDoc Go Imaging Systems (catalog #12012189)	<p>Use with the following applications:</p> <ul style="list-style-type: none"> ▪ Chemiluminescence ▪ Colorimetric blots ▪ Coomassie Fluor Orange Stain ▪ Diamond Stain ▪ Ethidium bromide (EtBr) ▪ Flamingo Stain ▪ Gel excision ▪ GelGreen Stain ▪ GelRed Stain ▪ Oriole Gel Stain ▪ Ponceau blots ▪ Stain-Free gels and blots ▪ SYBR® Stains ▪ SYPRO Ruby Stain ▪ UView Stain <p>Caution: You must use the accessory UV Shield for ChemiDoc Go and GelDoc Go Imaging Systems with the UV/Stain-Free Tray during any band excision procedure. This ultraviolet (UV) shield reduces exposure to UV rays.</p> <p>Caution: Before you perform band excision, you must protect yourself from UV radiation by wearing protective gear, including UV-protective eyewear, lab coat, and gloves. Follow instructions in the Band Excision Using the ChemiDoc Go and GelDoc Go Imaging Systems Quick Guide (10000122134).</p>
White Tray for ChemiDoc Go Imaging System (#12019873) White Tray for GelDoc Go Gel Imaging System (#12012165)	<p>Use with protein applications that call for colorimetric stains:</p> <ul style="list-style-type: none"> ▪ Coomassie Blue Stain ▪ Fast Blast DNA Stain ▪ Silver stain
Blue Tray for ChemiDoc Go Imaging System (#12019748) Blue Tray for GelDoc Go Gel Imaging System (#12012160)	<p>Use with all nucleic acid applications that use blue-excitable stains:</p> <ul style="list-style-type: none"> ▪ Diamond Stain ▪ Gel excision ▪ GelGreen Stain ▪ GelStar Stain ▪ StarBright™ Blue 520 Dye (only for ChemiDoc Go Imaging System) ▪ StarBright Blue 700 Dye (only for ChemiDoc Go Imaging System) ▪ SYBR® Gold Stain ▪ SYBR® Green Stain ▪ SYBR® Safe Stain <p>Note: Because the Blue Tray does not use UV, it avoids damage to nucleic acids during imaging and is safer for the user during band excision. XcitaBlue Viewing Goggles (#1708185) are available.</p>
UV Shield for ChemiDoc Go and GelDoc Go Imaging Systems (#12012164)	To be used with the UV/Stain-Free Tray when performing band excision.

Using the Trays

1 Clean the Tray

Use water and standard laboratory detergents or mild solvents such as ethanol or methanol.

2 Dry the Tray

Use lint-free wipe as dust particles or lint on a sample tray can glow under UV illumination.

3 Turn On the Instrument and Use Tray for Imaging

Center the sample (gel/blot) on the tray glass. Turn on the instrument. Pull the drawer door to open. Place tray on the transilluminator and close door fully.

Caution: Clean tray with water and standard laboratory detergents or mild solvents after use. Do not completely submerge tray and do not wash it in a dishwasher.

BIO-RAD, CHEMIDOC, GELDOC, and STARBRIGHT are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions. All trademarks used herein are the property of their respective owner. © 2024 Bio-Rad Laboratories, Inc.

SYBR is a trademark of Thermo Fisher Scientific Inc.

The GelDoc Go Imaging System uses the Linux operating system, which contains software licensed under the following licenses as well as others:

Lesser Gnu Public License v 2.0, 2.1, 3.0

Apache License 2.0

BSD License



**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Website bio-rad.com **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399
Canada 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23
France 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300 **Israel** 0 3 9636050
Italy 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 080 007 7373 **Luxembourg** 00 800 00 24 67 23 **Mexico** 52 555 488 7670
The Netherlands 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23 **Portugal** 00 800 00 24 67 23
Russian Federation 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23 **Spain** 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23
Switzerland 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311 **United Arab Emirates** 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23

