


# Gel Alignment Templates for ChemiDoc™ Go and GelDoc™ Go Imaging Systems

## Quick Guide


Bio-Rad Gel/Blot Alignment Templates enable you to quickly and consistently place gels and blots on ChemiDoc Go and GelDoc Go Sample Trays.

The ChemiDoc Go and GelDoc Go Gel Alignment Template Kit (catalog #12012190) contains the following:

- Alignment template for small gels, 10 x 7 cm

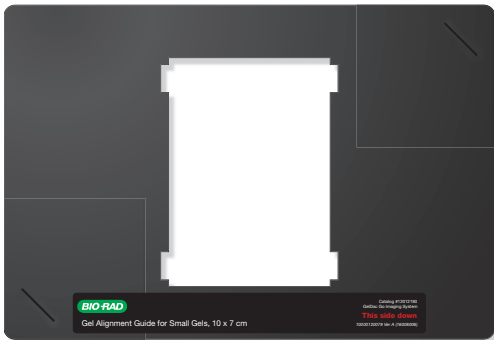

- Alignment template for large gels, 15.5 x 10 cm

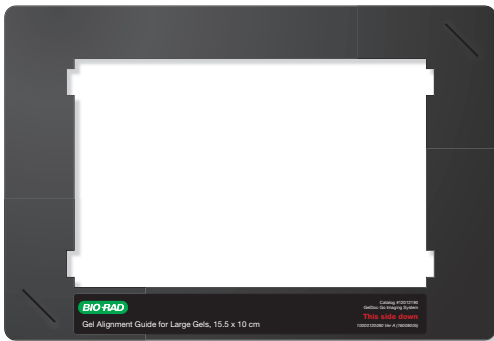

You can use the gel/blot alignment templates with all supported tray types. For more information about supported tray types, see the Getting Started with ChemiDoc Go and GelDoc Go Sample Trays Quick Guide (10000122130).

### Prepare the Gel/Blot Alignment Template

- 1 Select the gel/blot alignment template that matches the size of your sample.
- 2 To avoid the alignment template label getting in your image, turn the label side to face down.
- 3 Bend a corner at the precut slit to make a tab.  
The tab will be used to remove the template from the tray after use.
- 4 Place the gel/blot alignment template on the sample tray.
- 5 Place the gel or gel tray in the open area of the template.



### Clean and Store Templates

It is important to keep the templates clean and to store them flat.

- 1 After using a gel/blot alignment template, rinse it with water and allow it to air dry.
- 2 Store gel/blot alignment templates flat to prevent warping.

BIO-RAD, CHEMIDOC, and GELDOC are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions. All trademarks used herein are the property of their respective owner. © 2024 Bio-Rad Laboratories, Inc.

The GelDoc Go Imaging System uses the Linux operating system, which contains software licensed under the following licenses as well as others:

Lesser Gnu Public License v 2.0, 2.1, 3.0

Apache License 2.0

BSD License



**Bio-Rad  
Laboratories, Inc.**

Life Science  
Group

**Website** [bio-rad.com](http://bio-rad.com) **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399  
**Canada** 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23  
**France** 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300 **Israel** 0 3 9636050  
**Italy** 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 080 007 7373 **Luxembourg** 00 800 00 24 67 23 **Mexico** 52 555 488 7670  
**The Netherlands** 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23 **Portugal** 00 800 00 24 67 23  
**Russian Federation** 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23 **Spain** 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23  
**Switzerland** 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311 **United Arab Emirates** 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23

