

	<b>TYPE :</b>	<i>Standard working instructions (SWI)</i>		
	<b>TITLE</b>	<i>Decontamination procedure for 6-color naica® system (N20001.6)</i>		
	<b>DOC NO. :</b>	SWI- 00102	<b>DATE :</b>	20220818
	<b>REVISION :</b>	B		

#### DOCUMENT HISTORY TABLE

Rev.	Reason for the change	Author	Change Date
A	Creation	ABI	20210329
B	Updates document : - Note about disinfection products	ABI	20220818



 	<b>TYPE :</b>	<i>Standard working instructions (SWI)</i>		
	<b>TITLE</b>	<i>Decontamination procedure for 6-color naica® system (N20001.6)</i>		
	<b>DOC NO. :</b>	SWI- 00102	<b>DATE :</b>	20220818
	<b>REVISION :</b>	B		

## Decontamination procedure for 6-color naica® system (N20001.6)

### 1. Overview and context of the document

- This document provides the instructions to perform the decontamination of all the components of the naica® system in case of return at Stilla Technologies.
- The actions described in this protocol can be performed by a Stilla Technologies’ service specialist, by a distributors’ service specialist or by a client under the supervision of a Stilla Technologies’ or distributors’ service specialist.
- In case of a return of one or more instruments of the naica® system to Stilla Technologies, the following actions are required:
  - Clean and disinfect the instrument(s) that are supposed to return at Stilla Technologies. Be sure that all the procedures are carried out correctly, following the instruction present in the second section of this document “Decontamination Procedure – Instructions”,
  - Fill out and sign the “*Decontamination certificate for naica® system (Prism6)*”.
  - Pack the instrument to be returned in an adequate and protective packaging. Use the original boxes the naica® system was delivered in. Stilla Technologies will not be responsible for any damages incurred during transport of the instrument.
  - Stick the “*Decontamination Certificate*”, firmly on the external surface of the box along with any other documentation required for shipment.
  - Depending on the location of the naica® system, return the instrument at the following address:

<input type="checkbox"/> From Europe, China and Africa:  <b>Stilla Technologies</b> Villejuif Biopark 1, Mail du Professeur Georges Mathé 94800 Villejuif, FRANCE  Phone : (+33) 09 82 27 47 47 Email : <a href="mailto:support@stilla.fr">support@stilla.fr</a>	<input type="checkbox"/> From America, Asia (without China), Oceania  <b>Stilla Technologies Inc.</b> 66 Cherry Hill Drive, Suite 110 Beverly MA 01915 USA  Tel : 1-833-888-0150 Email : <a href="mailto:support@stilla-inc.com">support@stilla-inc.com</a>
---	--

 	<b>TYPE :</b>	<i>Standard working instructions (SWI)</i>		
	<b>TITLE</b>	<i>Decontamination procedure for 6-color naica® system (N20001.6)</i>		
	<b>DOC NO. :</b>	SWI- 00102	<b>DATE :</b>	20220818
	<b>REVISION :</b>	B		

## Reference Documents & standards

Reference	Document title	Date
SWI-00108	How to re-package a pump used for the naica® system	20210831
SWI-00109	How to pack a Geode instrument	20210916
SWI-00132	Unboxing instructions of the Prism6 instrument	20210125
QFD-00122	Decontamination certificate for naica® system (Prism6)	20210329

## Decontamination procedure – Instructions

### Material necessary for the decontamination procedure:

- Gloves;
- Glasses;
- Mask;
- Laboratory coat;
- Hydroalcoholic disinfectant solution for device’s surfaces commonly used for biological and medical devices (Ex. Phagospray).  
 Note: Certain disinfection products have been observed to leave white stains on the Prism6 lid surface.  
 In case of anticipation of using a similar disinfection product due to non-availability of the above specified Phagospray reference, Stilla Technologies recommends to first test another alcohol-based product on a small, less visible surface of the top cover of the Prism6 instrument.  
 Stilla Technologies does not recommend using a disinfection agent based on isopropanol and glutarate e.g., Bacillo® Plus.
- Decontaminant solution for device’s surfaces used for biological and medical devices, specifically targeting nucleases and DNA contaminations (Ex. RNase away).
- Standard laboratory paper towels.

Select the instrument (or the instruments) that will undergo the decontamination procedure and follow the instructions.



#### Instrument : Pump

##### Prerequisites:

- Switch off and unplug the pump from the power socket.  
(For more information check the procedure “How to pack a pump of the naica® system”).

##### Removal and cleaning/sterilization of the detachable parts:

- Moisten a paper towel with the disinfectant/decontaminant cleaner and pass it through the external surfaces of all the detachable parts (they could be different based on the model of the pump).

 	<b>TYPE :</b>	<i>Standard working instructions (SWI)</i>		
	<b>TITLE</b>	<i>Decontamination procedure for 6-color naica® system (N20001.6)</i>		
	<b>DOC NO. :</b>	SWI- 00102	<b>DATE :</b>	20220818
	<b>REVISION :</b>	B		

**Cleaning/sterilization of the instrument's surface:**

- Moisten a paper towel with the disinfectant/decontaminant cleaner and pass it through all the surface of the pump, paying attention to avoid the power connector.

**Instrument : Geode**

**Prerequisites:**

- Shut off and unplug the Geode from the power socket.
- Remove the tube coming from the pump, from the rear side of the Geode.  
(For more information check the procedure "*How to pack a Geode instrument*").

**Removal and cleaning/sterilization of the detachable parts:**

- Open the lid of the Geode and take away eventual residual chips.
- Remove the rubber gasket laying on the edges of the thermo-block and trash it to an appropriate bin.
- Remove the magnetic frame laying on the thermo-block surface.
- On a bench, put several layers of paper towel and place the magnetic frame on top.
- Spray the disinfectant/decontaminant cleaner on the whole surface of the magnetic frame, on both sides and leave it for 5 minutes.
- With a clean and dry paper towel absorb the liquid in excess from the magnetic frame and rinse it if necessary.
- Please take notice that a Geode instrument may include several magnetic frames for the different naica® system chip consumable types. Ensure to perform the above-described procedure for all magnetic frames and for all magnetic adaptor(s) fitting in the internal part of the lid.
- At the end of the procedure, put all the disinfected/decontaminated tools in clean zipper bag (not provided).



**Cleaning/sterilization of the instrument's surface:**

- Moisten a paper towel with the disinfectant/decontaminant cleaner and pass it through all the surface of the Geode, paying attention to avoid the power connector.
- Clean the surface of the thermo-block and the internal surface of the lid in the same way.

**Instrument : Prism6**

**Prerequisites:**

- Switch on the Prism6, launch Crystal Reader software to open the Prism6's flap. Take away eventual residual chips and remove the Prism6 tray holder. Close the flap and shut off the instrument.
- Unplug the Prism6 from the power socket.
- Remove all the cables coming plugged on the rear side of the Prism6.  
(For more information check the procedure "*Unboxing instructions of the Prism6 instrument*").

 	<b>TYPE :</b>	<i>Standard working instructions (SWI)</i>		
	<b>TITLE</b>	<i>Decontamination procedure for 6-color naica® system (N20001.6)</i>		
	<b>DOC NO. :</b>	SWI- 00102	<b>DATE :</b>	20220818
	<b>REVISION :</b>	B		

**Removal and cleaning/sterilization of the detachable parts:**

- On a bench, put several layers of paper towel and place the Prism6 tray holder on top.
- Spray the disinfectant/decontaminant cleaner on the whole surface of the Prism6 tray holder, on both sides and leave it for 5 minutes.
- With a clean and dry paper towel absorb the liquid in excess from the frame, rinse it if necessary.
- Please take notice that a Prism6 instrument may include several Prism6 tray holders for the different naica® system chip consumable types. Ensure to perform the above-described procedure for all Prism6 tray holders.
- At the end of the procedure, put all the disinfected/decontaminated tools in a clean zipper bag (not provided).
- Unplug the mouse, the keyboard, and the monitor connected to the Prism6. Moisten a paper towel with the disinfectant/decontaminant cleaner and pass it through all the surface of these instrument accessories.  
(For more information check the procedure “How to pack the Prism6”).

**Cleaning/sterilization of the instrument’s surface:**

- Carefully read the procedure “Unboxing instructions of the Prism6”
- It is important to lock the optical unit of the Prism6 instrument before moving it. Stilla Technologies will not be responsible for any damages incurred during packaging and / or transport of the instrument, if the transportation locks have not been put in place correctly.
- Moisten a paper towel with the disinfectant/decontaminant cleaner and pass it through all the surface of the Prism6, paying attention to avoid the power connector.

Following the decontamination procedure continue to package the instruments as instructed in the documents listed in the section “Documents reference & standards”.

Fill the “Decontamination certificate for naica® system (Prism6)” and add it on the instrument boxes as described on page 2 of this document.