



**Prism6 & Crystal Reader
software
for the
6-color naica® system**



Prism6 H24000.6

Crystal Reader software v4.0

Sapphire Chip & Ruby Chip

TECHNICAL SUPPORT

For help and technical advice, please contact the Technical Support Department at Stilla Technologies. The Technical Support Department:

For European Customers:

Monday to Friday, 9:30 AM - 6:30 PM Central European Time (CET).

Closed on French bank holidays.

Phone: (+33) 9 82 27 47 47

Email: support@stilla.fr.

For North American Customers:

Monday to Friday, 8:00 AM – 6:00 PM, EST.

Closed on American bank holidays.

Phone: 1-833-888-0150 ext. 1

Email: support@stilla-inc.com

Online Technical Support is also available at: www.stillatechnologies.com/technical-support/

LEGAL NOTICES

No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopy, recording, or any information storage or retrieval system, without permission in writing from Stilla Technologies.

Stilla Technologies reserves the right to modify its products and services at any time. Information contained in this user manual is subject to change without notice. Stilla Technologies assumes no liability for errors or any damage resulting from the use of this information.

The naica® system and/or its uses are covered by US and corresponding non-US patents owned by or under license to Stilla Technologies.

The naica® system and/or its use is licensed under U.S. and/or non-U.S. patents, and/or pending U.S. and non-U.S. patent applications owned by or under license to Bio-Rad Laboratories, Inc., including, but not limited to, U.S. Patent Nos. 9,968,933; 9,127,310; RE41,780; and 8,871,444.

Details available at: www.stillatechnologies.com/patents.

Purchase of the naica® system includes a limited, non-transferable right under such intellectual property for use of the products for research and applied applications, including commercial services, but excluding any commercial NIPT services. The exclusion of commercial NIPT services applies only in the United States and the European Union.

No rights are granted for diagnostic uses.

Further information on purchasing licenses may be obtained by contacting Stilla Technologies, 1 Mail du Professeur Georges Mathé, 94800 Villejuif, France. It is the responsibility of the purchaser/end user to acquire any additional intellectual property rights that may be required from Bio-Rad Laboratories, Inc. It is the responsibility of the purchaser/end-user to acquire any additional intellectual property rights that may be required.

TRADEMARKS

Crystal Digital PCR™, naica® and Stilla® are trademarks of Stilla Technologies. All other trademarks are the sole property of their respective owners.

COPYRIGHT & INTENDED USE

Copyright 2018-2022, Stilla Technologies. All rights reserved.

For Research Use Only. Not for use in diagnostic procedures.

Table of Contents

1. INTRODUCTION TO CRYSTAL DIGITAL PCR™	3
1.1. Purpose of the document	3
1.2. Intended use of the naica® system	5
1.3. Citing the naica® system in scientific publications, presentations, seminars, etc.,	6
2. MATERIALS AND EQUIPMENT	7
2.1. Prism6 packaging	7
2.2. Safety instructions for the naica® system	8
2.3. Prism6 instrument safety information	8
2.3.1. Prism6 instrument labeling	11
2.3.2. Prism6 declaration of conformity	13
2.3.3. Prism6 instrument specifications	13
2.3.4. Prism6 installation instructions	17
3. OPERATING THE PRISM6 INSTRUMENT	18
3.1. General considerations and prerequisites for using Crystal Reader software	18
3.2. Prism6 instrument initialization	18
3.3. Prism6 instrument availability	19
3.4. Instructions for chip positioning	20
3.5. Definition of a Crystal Digital PCR™ experiment	27
3.6. Experimental set-up with Crystal Reader software	28
3.6.1. Step 1 - Create a new experiment by selecting the scanning template file	28
3.6.2. Step 2 - Confirm or modify the experiment context	31
3.6.3. Step 3 - Select the chambers to scan and define chamber details	36
3.6.4. Step 4 - Launch the scan	39
3.6.5. Step 5 - Check image quality	43
3.6.6. Perform quality control:	45
4. ADVANCED FUNCTIONALITIES	47
4.1. How to rescan existing chambers and how to scan new ones	47
4.1.1. How to rescan all the chambers of the experiment	48
4.1.2. How to rescan only a subset of chambers	48
4.2. How to edit the Prism6 instrument settings	48
4.3. How to perform focus calibration	49
4.4. How to Pool/Unpool Chambers	51
5. MAINTENANCE AND TECHNICAL SUPPORT	53
5.1. Cleaning and Decontamination	54
5.2. Disposal	55
5.3. To ship the instrument back to Stilla®	55
6. TROUBLESHOOTING	56
6.1. Prism6 and Crystal Reader software	56
6.2. Sapphire Chip	57
6.3. Ruby Chip	63
7. SOFTWARE LICENSE INFORMATION	70
7.1. How to view the software version	70
7.2. Third-party licenses	71

1. Introduction to Crystal Digital PCR™

1.1. Purpose of the document

This document provides general information for the use of the Prism6 instrument and Crystal Reader software for Crystal Digital PCR™. The Prism6 workflow, using the Crystal Reader software and the different hardware components, is described in detail. It is essential to read the User Manual carefully and pay attention to the safety information provided. The instructions and safety information in the User Manual must be followed to ensure the safe operation of the instrument and to maintain the instrument in a safe condition.

All documents referenced in this User Manual can be accessed here:

<https://www.stillatechnologies.com/digital-pcr/naica-system-support/technical-resources/>

Overview of the Crystal Digital PCR™ Workflow

Crystal Digital PCR™ is Stilla Technologies' next-generation technology for the absolute quantification of nucleic acids.

Using cutting-edge microfluidic innovations, this technology integrates the digital PCR process in a single consumable. The sample is first flowed through a network of microchannels and partitioned into a large 2D array of 12,900 to 28,500 individual droplets (Sapphire Chip)¹ and 10,000 to 17,000 droplets (Ruby Chip)², also called a Droplet Crystal. PCR is then performed within the chips and the Droplet Crystal is imaged to reveal the droplets that contain amplified targets. The last step consists of counting the number of these positive droplets to precisely calculate the absolute quantity of nucleic acids.

With Crystal Digital PCR™, the combination of powerful image analysis and intuitive visual inspection offers an unmatched level of confidence in the digital PCR measurement, yielding data you can trust.

¹ 12,900 to 25,800 for naica® multiplex PCR MIX and naica® PCR MIX

15,000 to 30,000 for qScript™ XLT One-Step RT-qPCR ToughMix®

² naica® multiplex PCR MIX and naica® PCR MIX & qScript™ XLT One-Step RT-qPCR ToughMix®

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip




SAMPLE LOADING	GENERATING DROPLET CRYSTAL AND PCR	READING AND ANALYSIS
		
<p>Prepare the sample for the reaction mix. Stilla Technologies recommends naica® PCR MIX reagents, specifically developed for Crystal Digital PCR™. Load the reaction mix into the wells of the selected chip, and seal with the provided caps.</p>	<p>Place the prepared chips into the Geode. Launch the combined partitioning and amplification program; Droplets are generated by partitioning of each sample and PCR amplification is performed immediately after Droplet Crystal generation.</p>	<p>After PCR, transfer the chips to the Prism6 instrument. Set up the read-out using Crystal Reader software for data acquisition of the Droplet Crystal using up to 6 fluorescent channels (Blue, Teal, Green, Yellow, Red, and Infra-Red). Image analysis and data extraction are performed using the Crystal Miner software.</p>

Figure 1: Overview of Crystal Digital PCR™ Workflow.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

1.2. Intended use of the naica® system

The naica® system for Crystal Digital PCR™ is composed of two instruments: the Geode, which performs droplet generation and amplification, and the Prism6, which enables imaging of the Droplet Crystal in up to six detection channels. The Crystal Reader software is used to control and set-up Crystal Digital PCR™ experiments on the Prism6 instrument.

The Crystal Reader software functions as the user interface to prepare the experiment for the Prism6 and it is dedicated to the image acquisition in the Crystal Digital PCR™ workflow.

The Crystal Reader software:

- Allows defining the analytical context of the experiments. Experiments can be set up on-demand or dedicated experimental templates can be created for recurring experimental setups.
- Controls the Prism6 instrument for the acquisition of the fluorescence images of Sapphire Chip or Ruby Chip.
- Applies pre-analysis treatments to the acquired images and provides the initial quality control in preparation for the detailed downstream experiment analysis performed by the Crystal Miner software.

The Crystal Miner software is used to extract data from the images acquired using the Prism6 instrument and to calculate the absolute concentrations of the targeted nucleic acids.

Crystal Reader and Crystal Miner software are pre-installed on the Prism6 instrument and delivered as part of the naica® system.

The naica® system performs Crystal Digital PCR™ within microfluidic chips (Sapphire Chip & Ruby Chip). naica® PCR MIX reagents are recommended to achieve optimal Crystal Digital PCR™ performance on the naica® system.

For detailed instructions for the Geode and for the Crystal Miner software, please refer to the respective User Manuals. For detailed instructions for Sapphire Chip, Ruby Chip, naica® PCR MIX, and naica® multiplex PCR MIX, refer to the respective Instructions for Use (IFU).

The naica® system is a laboratory instrument to be used by qualified personnel in a controlled environment. Before using the naica® system, the user must be trained by a Stilla Technologies representative.

In general, Crystal Digital PCR™ can be performed on the naica® system with all types of DNA sample types. However, individual sample-type compatibility for digital PCR applications may require dedicated assay validation by the end-user. Please note that the extraction method used, and sample purity, can also influence sample compatibility for digital PCR applications.

The Prism6 instrument and Crystal Reader software is intended for use by professional users trained in molecular biology techniques and in the operation of the Prism6 instrument and Crystal Reader software.

The Prism6 instrument and Crystal Reader software are part of the naica® system.

The naica® system is intended for Research Use Only. Not for use in diagnostic procedures.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

1.3. Citing the naica® system in scientific publications, presentations, seminars, etc.,

To cite the use of the naica® system use:

Crystal Digital PCR™ (Stilla Technologies, France)

naica® system (Stilla Technologies, France)

naica® system component names:

- Geode
- Prism6
- Sapphire Chip
- Ruby Chip
- Crystal Reader software
- Crystal Miner software
- naica® PCR MIX reagents:
 - naica® PCR MIX (Stilla Technologies, France)
 - naica® multiplex PCR MIX (Stilla Technologies, France).

2. Materials and Equipment

2.1. Prism6 packaging

- Prism6 instrument:
 - Prism6 instrument (H24000.6)
 - 1 Prism6 tray holder Sapphire Chip – H24002
 - 1 Prism6 tray holder Ruby Chip - H24001
 - Certificate of Conformity
- For the peripheral box :
 - Mini display cable
 - Monitor
 - Keyboard
 - Mouse
 - Mousepad
 - Power cord for Prism6
 - Cable for Monitor
 - External USB Barcode Scanner
 - Precision Wipes (Kimtech™ Science, 7552, 1 ply, 213x114 mm)
 - Antistatic wetted wipes (ACL Staticide®, Reference: SW12)
 - USB key

Note- Replacement Precision Wipes must be purchased using the reference - Precision Wipes (Kimtech™ Science, 7552, 1 ply, 213x114 mm) from global standard laboratory suppliers

Note - Replacement Antistatic wetted wipes (ACL Staticide®, Reference: SW12) can be ordered through Stilla Technologies using product reference H10000.472 or directly from the supplier Digi-Key using the reference ST1059-ND. [SW12 ACL Staticide Inc | Static Control, ESD, Clean Room Products | DigiKey](#)





2.2. Safety instructions for the naica® system

Before the first use of the naica® system, please read the User Manual entirely.


Special notice should be taken of the following.

2.3. Prism6 instrument safety information

The following warning labels are located on the Prism6 instrument:

SYMBOL	MEANING
	<p>Caution - the possibility of pinch risk</p> <p>Please ensure to correctly insert the tray holders onto the movable tray of the Prism6 instrument to avoid pinch risk.</p> <p>Avoid any manual operation of the Prism6 instrument flap to access the movable tray to avoid pinch risk.</p>
	<p>Caution - optical radiation</p> <p>High power light sources used for fluorescence imaging. Do not stare directly at the beam.</p> <p>Do not open the Prism6 instrument flap manually while the Prism6 instrument is switched on.</p> <p>Do not open the Prism6 instrument flap manually while the Prism6 instrument is in operation.</p> <p>Do not stare into the light beam when using a handheld bar code scanner.</p>
	<p>Grounding point</p> <p>This point is connected to the earth. Use this connection if grounding is required.</p>
	<p>Caution – attention required</p> <ul style="list-style-type: none"> ○ The Prism6 instrument must be placed on a stable surface that is clean and vibration-free. ○ Do not block the area in front of the Prism6 instrument and ensure that the movable tray has sufficient space to move while the Prism6 instrument flap is opening and closing. ○ A sufficient clearance area around the Prism6 instrument is required to allow access to the main power switch. ○ A 10cm (4 inches) clearance must be left at the back of the device to ensure proper ventilation. ○ Before transportation, the respective transportation locks must be put in place. ○ Before transportation, the respective tray holder (Sapphire Chip & Ruby Chip) must be removed.

Additional warning for the Prism6 instrument:

SYMBOL	MEANING
	<p>Warning- heavy object</p> <p>Do not try to lift the Prism6 instrument alone.</p> <p>Two persons are required to lift the Prism6 instrument.</p>

General safety instructions for the Prism6 instrument:









- Please ensure that the Prism6 instrument is operated as instructed in the provided User Manuals. User Manuals are subjected to changes. The latest version can be accessed at the Technical Resources webpage.
- Stilla Technologies cannot be held responsible for any damage or injuries arising from improper operation.
- Do not use the Prism6 instrument if any parts are broken, chipped, rusty, or if the power cables are damaged in any way.
- Do not open the Prism6 instrument flap manually while the Prism6 instrument is switched on.
- Do not open the housing of the Prism6 instrument unless explicitly instructed by a Stilla Technologies representative.
- Opening the Prism6 instrument housing may breach any warranty for the Prism6 instrument.
- Only operate the Prism6 instrument with the provided and specified detachable main supply cords. Do not replace the components with non-specified cords.
- Do not attempt any repairs or alterations except as expressly described in this User Manual or as instructed by a Technical Support Representative.
- Always disconnect all instruments from their power source before cleaning or moving the Prism6 instrument.
- For the Prism6 instrument, the mains supply source must meet the national regulatory requirements.
- The power cord of the Prism6 instrument must be connected to a wall outlet with a grounded conductor.
- For the Prism6 instrument, the mains voltage must correspond to the range given in the product specification.
- Keep liquids away from the Prism6 instrument; avoid percolation of liquids inside the Prism6 instrument.
- For optimal use of the Prism6 instruments, the room temperature must be between 15°C and 30°C.
- For optimal performance, ensure to operate the Prism6 instrument in a dust-free environment.
- Do not block the area in front of the Prism6 instrument and ensure that the tray has sufficient space to move while the Prism6 instrument flap is opening and closing.
- A sufficient clearance area around the Prism6 instrument is required to allow access to the main power switch.
- Samples can be infectious or cause other damage to health. Safety regulations issued for the handling of sample material in the laboratory must be followed by wearing the proper Personal Protective Equipment (e.g., gloves and protective clothing).

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip






- For instrument cleaning and decontamination, follow the provided instructions. Do not use any cleaning agents other than those specified.
- For instrument shipment, follow only the detailed packaging and shipment procedures provided by Stilla Technologies. Only package and ship the instrument when instructed by a Stilla Technologies representative.
- Regular instrument service maintenance is recommended to ensure optimal system performance at all times. Please contact Stilla Technologies for the respective service offers.

2.3.1. Prism6 instrument labeling

Labeling symbols present on the Prism6 instrument:

SYMBOL	MEANING
	<p>Manufacturer</p>
	<p>Product reference (part number)</p>
	<p>Product serial number</p>
	<p>Read the User Manual before using the product</p>
	<p>Caution: documentation must be consulted in all cases where this symbol is marked. Using the product without applying the instructions explained in the User Manual may result in personal injury or damage the equipment and facilities.</p>
	<p>Alternating current.</p>
	<p>This device has been confirmed by the global safety certification company UL that the product is safe for use.</p>
	<p>Restriction of Hazardous Substances (Directive 2002/95/EC on the restriction of the use of certain hazardous substances in electrical and electronic equipment).</p>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

	<p>The product should be disposed of in an appropriate recovery and recycling structure.</p>
	<p>CE marking (manufacturer's declaration that the product meets the requirements of the applicable EC directives).</p>
	<p>Power on.</p>
	<p>Power off.</p>
	<p>Protective conductor terminal.</p>
<p>UDI</p>	<p>Unique device identifier</p>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

2.3.2. Prism6 declaration of conformity

To view the declaration of conformity certificate, please visit the Technical Resources webpage <https://www.stillatechnologies.com/digital-pcr/naica-system-support/technical-resources/>.

2.3.3. Prism6 instrument specifications

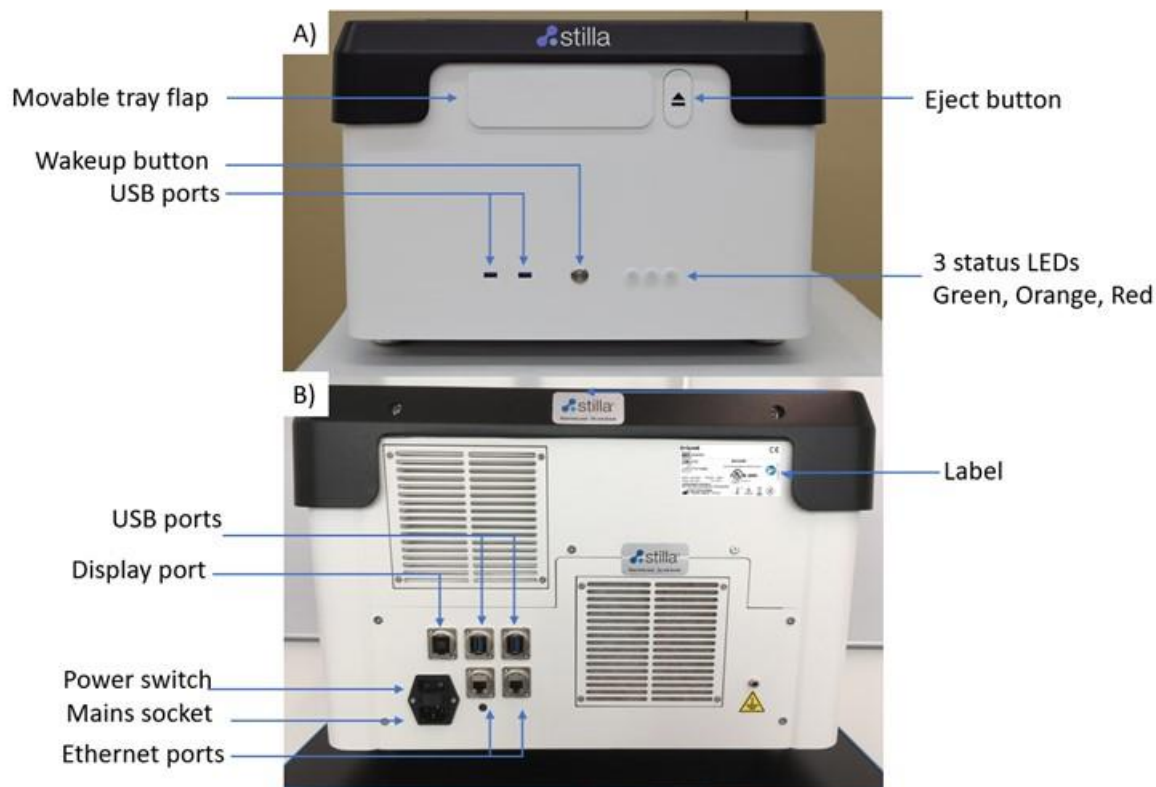


Figure 2: A) Front panel of the Prism6 instrument. B) Back panel of the Prism6 instrument.

Further details about Fig. 2 A

The front panel of the Prism6 instrument contains:

- a covering flap where chips are loaded
- an eject button to eject the movable tray from the Prism6
- Three status LEDs (Green Orange, Red)
 - Green: Prism6 instrument ready for operation
 - Orange: Prism6 instrument busy (instrument initialization, instrument operation)
 - Red: Prism6 instrument in failure mode
- Two USB 3.0 ports, for data transfer
- One soft power button, to revive the instrument if it has been powered up beforehand

Note: There can be differences in USB key compatibility with the Prism6 front panel USB ports. Stilla Technologies recommends using a USB key with a maximum capacity of 8Gb. The compatibility of individual USB keys needs to be validated by the end user for functional data transfer.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Further details about Fig. 2 B

The back panel of the Prism6 instrument has:

- 2 USB 3.0 ports
- 1 display port to connect the Prism6 monitor
- 2 Ethernet ports
- 1 label
- 1 mains socket
- 1 power switch

Technical Specifications

TECHNICAL SPECIFICATIONS	
INSTRUMENT DIMENSIONS AND WEIGHT	
Instrument dimensions (W x D x H in mm)	514 x 564 x 365
SAMPLE THROUGHPUT CAPACITY AND SCANNING TIMES	
Sample throughput capacity per run	Up to 3 chips/run: Sapphire Chip: up to 12 samples/run Ruby Chip: up to 48 samples / run
Scan time per sample	~3-4 min/sample for Sapphire Chip <2 min/sample for Ruby Chip
Scan time per run	Depending on the number of LEDs selected: <u>3-color scanning:</u> 30 min/run for 3 Sapphire Chip 60 min/run for 3 Ruby Chip <u>6-color scanning:</u> 45 min/run for 3 Sapphire Chip 80 min/run for 3 Ruby Chip
Specifications for chip re-scanning	Up to 3 scans of the same chip within 48h

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

IMAGING SYSTEM		
LED	Excitation wavelengths [nm]	Emission wavelengths [nm]
Blue	450-490	505-535
Teal	509-519	530-551
Green	533-557	568-593
Yellow	564-586	600-640
Red	625-643	655-685
Infrared	645-695	708-753
Recommended fluorophores*	FAM, Yakima Yellow®, Atto®550, ROX, Cy®5, Atto®700/Cy®5.5	
*Additional fluorophores within the specified wavelengths can be used, fluorophore selection requires validation depending on the digital PCR assay design.		
Result data format	16bit TIFF grayscale, CSV spreadsheet, XML files	
EMBEDDED INSTRUMENT PC		
Motherboard	Advantech AIMB 276	
CPU	intel i5-8500	
RAM	2x16GB	
System drive (SSD)	512GB	
Data drive (SSD)	1TB	
AMBIENT CONDITION		
Pollution degree	2	

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

OPERATION	
Environmental conditions	Clean indoor laboratory
Temperature	+15°C to +30°C
Relative humidity	20% to 90% non-condensing
Altitude	Operating at max. 1000m above sea level
Barometric pressure	795 hPa to 1060 hPa
Noise	<45 dB at a distance of 1 meter
ELECTRICAL	
AC input	100-240V~, 50/60Hz
Installation category	II
Input mains fluctuation	Max. 10%
Input power	Max. 400W

2.3.4. Prism6 installation instructions

Operating requirements

Proper infrastructure requirements for Prism6 installation :

- A clean laboratory environment
- A sturdy surface, for supporting up to 40 kg
- A minimum distance of 10 cm to neighboring objects around the Prism6 instrument
- Room temperature between 15°C and 30°C

Installation and performance validation

The initial naica® system installation, including the Prism6 instrument, must always be executed by a Stilla Technologies representative.

The naica® system installation and operational qualification (IQOQ), according to Stilla Technologies specifications, will be documented with a certificate.

The certification is required to release the naica® system for customer operation.

3. Operating the Prism6 instrument

3.1. General considerations and prerequisites for using Crystal Reader software

The Crystal Reader software is pre-installed on the Prism6 instrument. Future software version updates for Crystal Reader software are possible. Stilla Technologies will communicate about any required and available software update.

To download the latest version of the Crystal Reader software, visit the Technical Resources webpage: <https://www.stillatechnologies.com/digital-pcr/naica-system-support/technical-resources/>

Note: 2 GB of storage is required for software installation.

Note: Only install the required Windows 10 Operating System security updates for the naica® system. DO NOT update full Windows 10 Operating System updates, as they may have a negative impact on Crystal Reader & Crystal Miner software robustness and stability. Contact Technical Support with any questions.

The officially supported scanning templates for the Crystal Reader software are located on the Prism6 instrument by default at, “C:\Program Files\Stilla\CrystalReader\templates”.

Stilla Technologies does not recommend modifying or deleting any of the default scanning templates provided. Instructions to create a new scanning template from the default scanning templates provided are detailed below in section 3.6.1. Create a new experiment by selecting the scanning template file

All scanning templates are named as described below:

FileType_InstrumentName_ChipType_PCR-Mix_Name_AssayIndicator_version.ncx

3.2. Prism6 instrument initialization

The Prism6 instrument is operated through the monitor, keyboard, and mouse connected to the back panel of the Prism6 instrument.

To turn on and initialize the Prism6 instrument, and log into the Crystal Reader software, follow these steps:

1. Switch on the Prism6 instrument.

First, switch on the power button located at the back of the Prism6, then press the wake-up power button located on the front panel of the Prism6 instrument. Wait for instrument initialization. The left-most light on the front LED panel of the Prism6 instrument lights up green.

2. Turn on the Monitor of the Prism6 instrument. With instrument initialization, the Microsoft Windows Login window appears. The default password is “admin” for “admin” user. Refer to Microsoft Windows User Manual regarding account features.
3. The Crystal Reader software application is automatically launched at start-up. When the software is launched, the Crystal Reader software homepage is displayed.
4. To download the latest version of the Crystal Reader software, visit the Technical Resources webpage: <https://www.stillatechnologies.com/digital-pcr/naica-system-support/technical-resources/>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

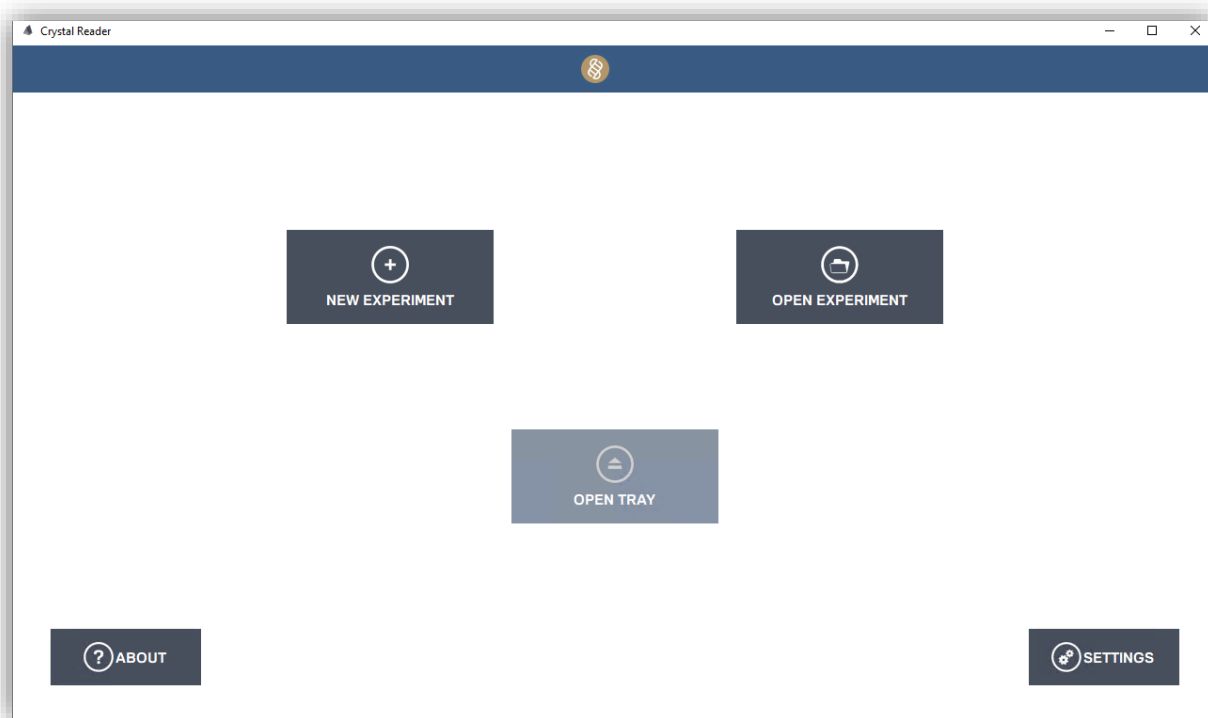


Figure 3: Crystal Reader software homepage with disabled “OPEN TRAY” button during Prism6 initialization or operation mode.

The Crystal Reader software homepage displays the following buttons:

- LOGIN: log in to apply user preference.
- NEW EXPERIMENT: create a new experiment
 - Embedded drop-down menu for chip-specific experimental templates
 - New Sapphire Chip experiment
 - New Ruby Chip experiment
- OPEN EXPERIMENT: open a previously saved experiment file to use identical experimental settings.
- OPEN TRAY/CLOSE TRAY: engagement button to open/close the Prism6 instrument flap to eject/retract the movable tray.
 - If the movable tray is fully retracted, this button displays “OPEN TRAY”.
 - If the movable tray is fully ejected, this button displays “CLOSE TRAY”.
 - If the moveable tray is in initialization mode, or in operation, the button is greyed out and inactive for engagements.
- ABOUT: detailed information about the Crystal Reader software.
- SETTINGS: edit user preferences, and the settings of the Crystal Reader software and the Prism6 instrument. For detailed instructions please refer to section 4.2 How to edit the settings.

3.3. Prism6 instrument availability

Only one instance of Crystal Reader can operate the Prism6 instrument at a time. When Crystal Reader is launched, the application takes control of the instrument and locks the access for other Crystal Reader instances. To unlock the access of the instrument, the Crystal Reader instance using the instrument must be closed.

Note: Always close Crystal Reader after scanning your experiment to make the Prism6 instrument available for the next user.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

If another Crystal Reader instance is running when Crystal Reader is started, the software will detect that the Prism6 instrument is unavailable. Crystal Reader will display a warning message while opening and the scan button will be disabled. Until the instrument is made available, it will only be possible to view and edit parameters to prepare the next experiment.

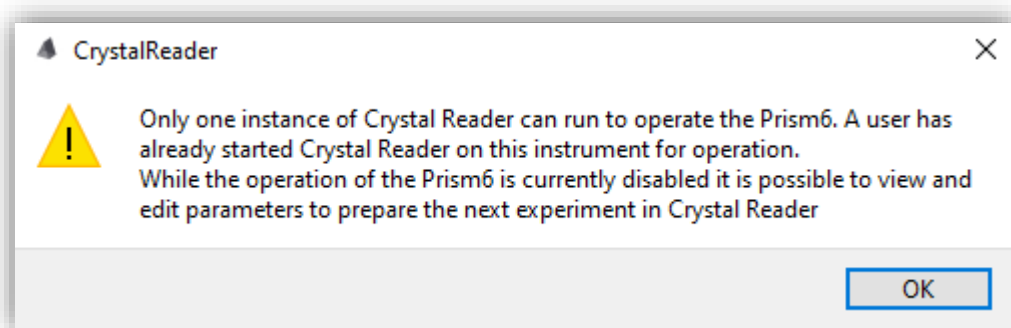


Figure 4 Warning message at Crystal Reader launch when the Prism6 instrument is unavailable

To take control of an instrument that is locked, first make sure that the Prism6 instrument is not in use by another user. The Crystal Reader software instance holding the Prism6 instrument resource can be running on the current Windows user session or in another user session. Connect to the user session running Crystal Reader and close the software. Restart Crystal Reader and check that the instrument has connected successfully.

3.4. Instructions for chip positioning

Note: Prior to chip positioning for scanning, read the respective Sapphire Chip and Ruby Chip Instructions for Use.

Eject the moveable tray for chip positioning

- Press the "OPEN TRAY" button on the Crystal Reader software homepage.
- Alternatively, press the "Eject" button located at the front of the Prism6 instrument for at least 3 seconds (Figure 2).

Wait until the moveable tray is fully ejected before placing the chips for scanning.



Figure 4: Fully ejected position of the movable tray. In this position, the moveable tray is ready to be loaded with the tray holder for chip scanning.

CAUTION!

Do not manually open the Prism6 instrument flap.

Do not manually engage with the movable tray.

Do not manually pull the moveable tray out any further.

Make sure that the moveable tray is clean and free from dust particles (clean it occasionally with ethanol or water)

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip



Figure 5: Prism6 tray holder Sapphire Chip H24002 (left) & Prism6 tray holder Ruby Chip H24001 (right)

Up to three chips can be fitted into one tray holder for simultaneous scanning. Two kinds of tray holders are provided with each Prism6 instrument:

- Prism6 tray holder Sapphire Chip – H24002, to be used with Sapphire Chip only
- Prism6 tray holder Ruby Chip– H24001, to be used with Ruby Chip only.

Note: The Prism6 tray holder Ruby Chip may display the following label: Prism6 tray holder Opal Chip – H24001 for the previous generation Opal Chip. The Opal Chip labeled tray holder is compatible for Ruby Chip positioning and operation for Crystal Reader software version v4.0.

The direction, for insertion of the tray holders onto the moveable tray, is indicated on the individual tray holder labels. Stilla Technologies recommends cleaning all tray holders occasionally with 70% ethanol solution, then rinsing with water using a Precision Wipe.

Position one of the chips (Sapphire Chip or Ruby Chip) in the individual slots of the corresponding tray holder (Prism6 tray holder Sapphire Chip or Prism6 tray holder Ruby Chip). Check for correct positioning. The respective chips should fit perfectly into the individual slots.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Instructions to position the chips in the tray holder.

Example provided for Sapphire Chip and 'Prism6 tray holder Sapphire Chip' only.


Similar instructions apply for Ruby Chip placement in the 'Prism6 tray holder Ruby Chip'. Differences are indicated in the provided text description with the respective image instructions.



Place the tray holder on a clean surface.



Release the chip clamp, to allow the chip to pass, by rotating it away from the slot.

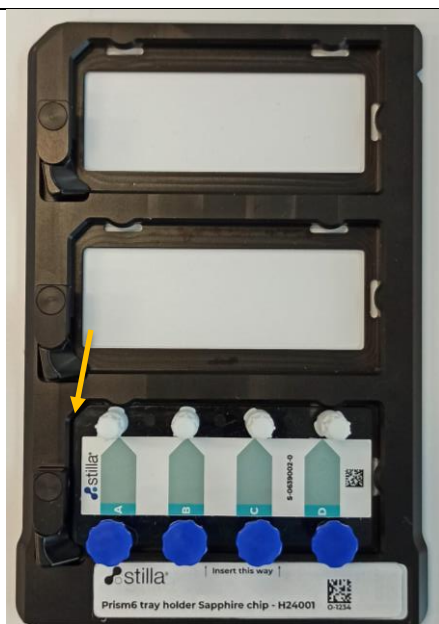



Insert the chip in the tray holder.

Each type of chip has a poka-yoke (cut-off) corner matching the respective tray holder.

On the Prism6 tray holder Sapphire Chip, the cut corner is at the top-right of the chip. Here the arrow points to the top right corner in the orientation of the chip (rotated 90 degrees in the photo, which is the orientation of the tray holder).

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

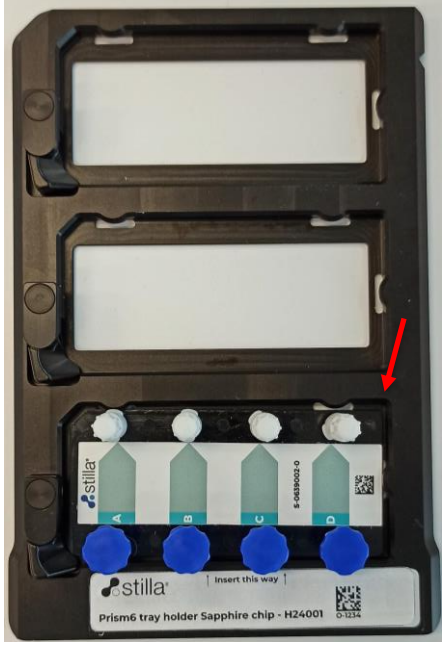	<p>Caution: this image shows <u>incorrect positioning</u> of a chip in the Prism6 tray holder Sapphire Chip. The chip is lying at an angle in the tray holder, i.e. it is not well inserted and is not flush at the corners (arrow).</p>
	<p>Correctly loaded Prism6 tray holder Ruby Chip.</p> <p>For the Prism6 tray holder Ruby Chip, the cut corner is at the bottom-left of the chip (arrow).</p>
<p>Repeat the described chip positioning operation for all the tray holder slots to be used.</p>	

Note: Make sure to only use the specific Prism6 tray holder Sapphire Chip or Prism6 tray holder Ruby Chip with the respective chip types:




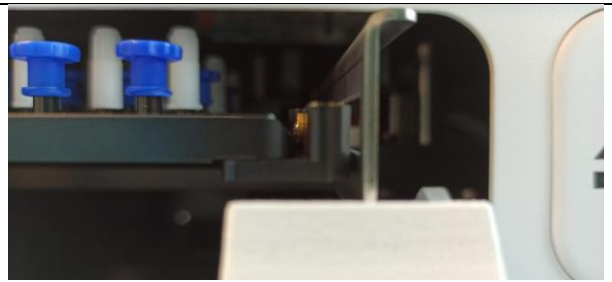
- Do not operate Sapphire Chip on Prism6 tray holder Ruby Chip.
- Do not operate Ruby Chip on Prism6 tray holder Sapphire Chip.
- Do not mix chip types on one tray holder.
- The tray holder is clearly identifiable by the tray holder label.

Note! The Prism6 tray holder may display the following label: Prism6 tray holder Opal Chip – H24001 for the previous generation Opal Chip. The Opal Chip labeled tray holder is compatible for Ruby Chip positioning and operation for Crystal Reader software version v4.0.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Instructions for loading the tray holder in the movable tray of the Prism6 instrument:

Example provided for Sapphire Chip, and Prism6 tray holder Sapphire Chip, only. Identical instructions apply for Prism6 tray holder Ruby Chip placement.

	<p>Place the tray holder on the movable tray</p>
	<p>Gently push the tray holder towards the back of the moveable tray.</p>
 <p>ball bearing in contact with tray holder</p>	<p>The tray holder is correctly positioned after the spring-loaded ball bearing (blue circle) is set in position with a distinct click.</p>
	<p>Ensure the tray holder is lying flat on the movable tray.</p> <p>Correct placement of the tray holder on the moveable tray.</p>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip


	<p>Incorrect placement of the tray holder on the moveable tray. Tray holder not sitting cleanly.</p>
<p>Retract the moveable tray to allow for chip scanning.</p> <ul style="list-style-type: none">a) Press the “CLOSE TRAY” button on the Crystal Reader software homepage.b) Alternatively, press the “Eject” button located at the front of the Prism6 instrument (Fig 1). <p>Ensure that the moveable tray is fully retracted and the Prism6 instrument flap is fully closed before proceeding to chip scanning.</p>	

Figure 6 - Instructions for correct position of chips in chip trays and of chip trays in the Prism6.

Note: In the event that the tray holder, a chip or multiple chips or parts thereof, fall into the Prism6 instrument interior, turn off the Prism6 instrument immediately and contact Technical Support.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

3.5. Definition of a Crystal Digital PCR™ experiment

An experiment is defined as being a set of samples processed during an individual Crystal Digital PCR™ workflow on the naica® system:

- The samples of an individual experiment were all processed on the same respective chip type (Sapphire Chip or Ruby Chip). Chip types cannot be mixed in a single experiment, for image acquisition with the Prism6 instrument.
- The samples of the individual experiment were prepared using the same PCR mastermix and the same PCR program for partitioning and amplification in the Geode.
- The same scanning parameters and the experimental details will be applied to all samples for an individual experiment during a Prism6 run, i.e.
 - the same exposure times (in ms) for each of the selected LEDs;
 - the same focus value (in mm).

Once an experiment is scanned by the Prism6, the Crystal Reader software automatically records the data in a ZIP format in a “.ncx” file (naica® crystal experiment). This generated “.ncx” file contains all the chamber images as well as all the contextual and pre-analysis data.

The indicative sizes of an individual “.ncx” file is:

Chip type (6-plex data)	.ncx file size / chamber
Sapphire Chip	237 +/- 12 MB
Ruby Chip	111 +/- 2 MB

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

3.6. Experimental set-up with Crystal Reader software

The 5 main steps of the experimental set up, for chip scanning, using Crystal Reader software include:

1. Create a new experiment by selecting the scanning template file
2. Confirm or modify the experiment context
3. Select the chambers to scan and define chamber details
4. Launch the experiment scan
5. Check chamber image quality

3.6.1. Step 1 - Create a new experiment by selecting the scanning template file

Each experiment starts with the selection of the scanning template (".ncx" format). The scanning template contains all the information and components required to set up an experiment consistently and quickly.

On the Crystal Reader software homepage click on "NEW EXPERIMENT". A pop-up window appears for scanning template selection (Figure 7).

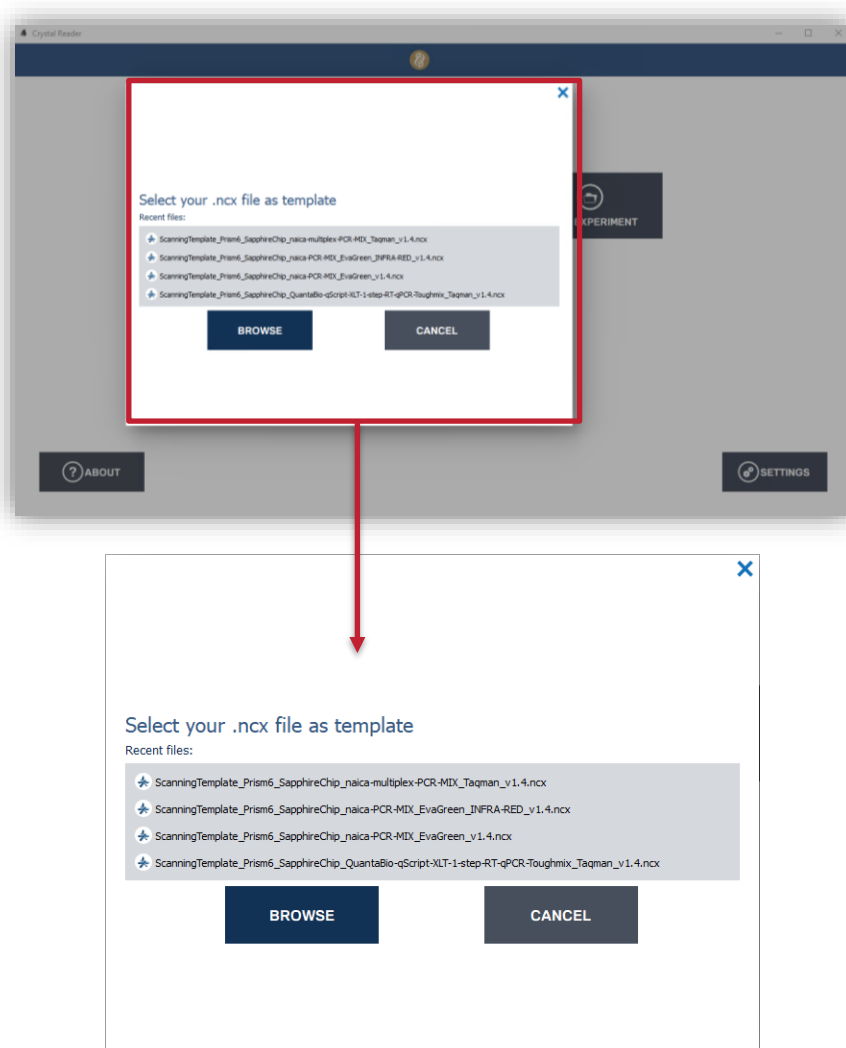


Figure 7: Crystal Reader software prompt to select the scanning template (".ncx" file) for the experiment.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Recently used scanning templates are provided for direct selection from the displayed menu “Recent files”.

To select a different scanning template, proceed to “BROWSE” and select the respective scanning template from the set of default scanning templates provided at the location: **C:\Program Files\Stilla\CrystalReader\templates**.

Different default scanning templates are available based on the parameters of the experiment:

- Chip type
- PCR mastermix

All scanning templates and analysis configuration files are named as described below:

FileType_InstrumentName_ChipType_PCR-Mix-Name_AssayIndicator_version.ncx

Examples:

- For an experiment, using the naica® multiplex PCR MIX, compatible with TaqMan® probe and the Sapphire Chip select:
ScanningTemplate_Prism6_SapphireChip_naica-multiplex-PCR-MIX_Taqman_v1.4.ncx
- For an experiment using the naica® PCR MIX compatible with EvaGreen® fluorescent intercalating dye (with droplet recognition in the blue channel) and the Ruby Chip select:
ScanningTemplate_Prism6_RubyChip_naica-PCR-MIX_Evagreen_v1.0.ncx.

Note: Stilla Technologies does not recommend modifying the provided original template files. Always use the “save as” function to ensure to keep the original template files. If a modification of any template file is required, please contact Technical Support.

After selecting the scanning template, the user is prompted to choose to use the default sample details (i.e. samples are simply named “Sample 1” to “Sample 12” for Sapphire Chip and “Sample 1” to “Sample 48” for Ruby Chip) or load the chamber details that are embedded in the scanning template (Figure 8). This feature provides the user a convenient way to set up chamber settings that will be run routinely.

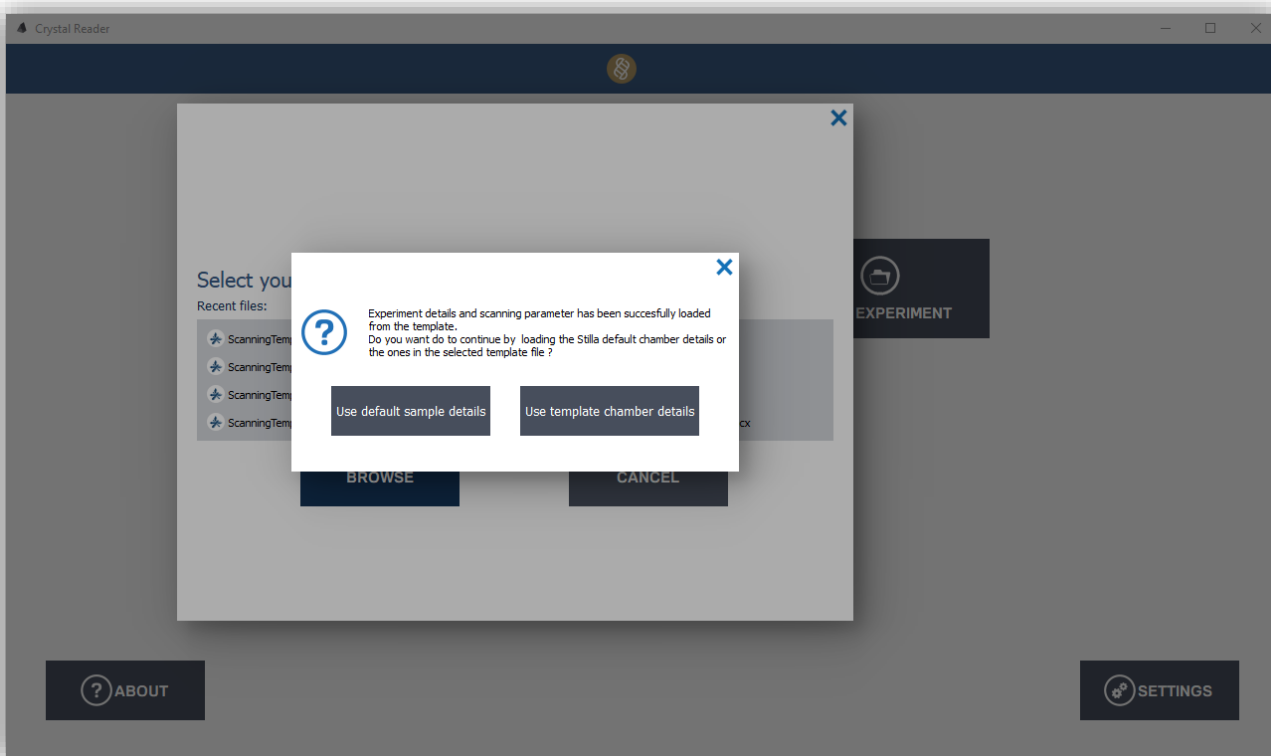


Figure 8: After the template has been loaded a pop-up window proposes the chamber details configurations that the user can load.

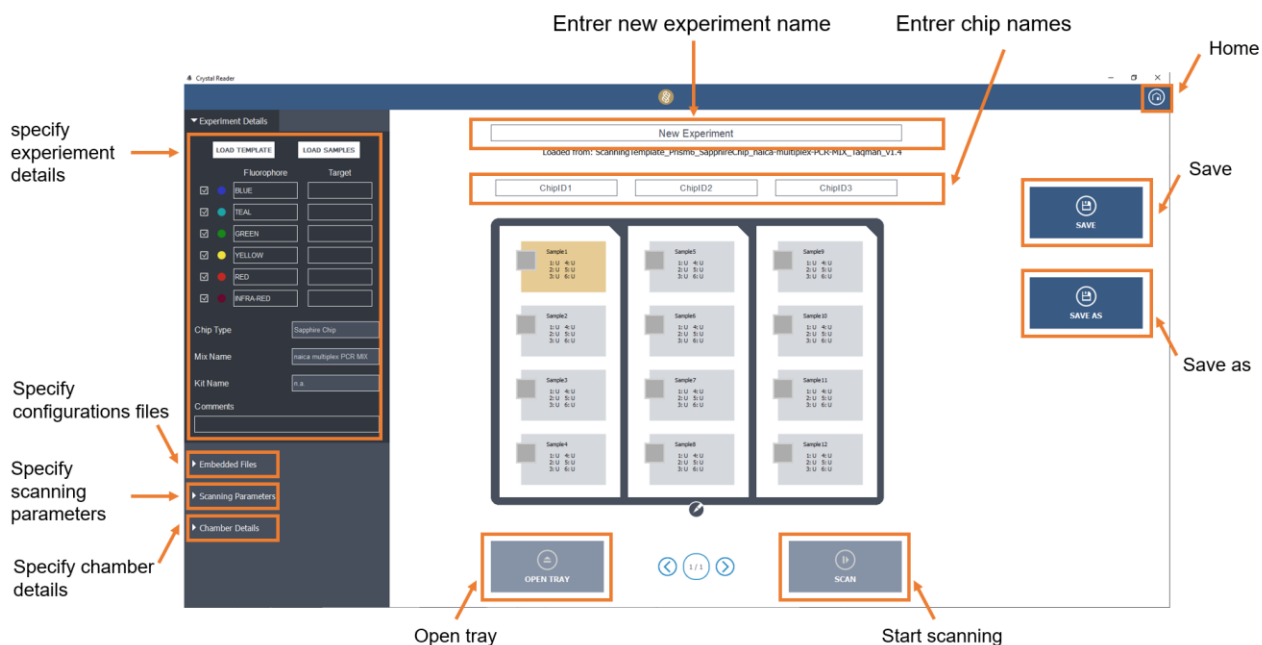


Figure 9: The Crystal Reader software “New Experiment” page for Sapphire Chip.

The New Experiment page display allows the user to check and (if necessary) modify the experiment context (Figure 9 for Sapphire Chip and Figure 10 for Ruby Chip).

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

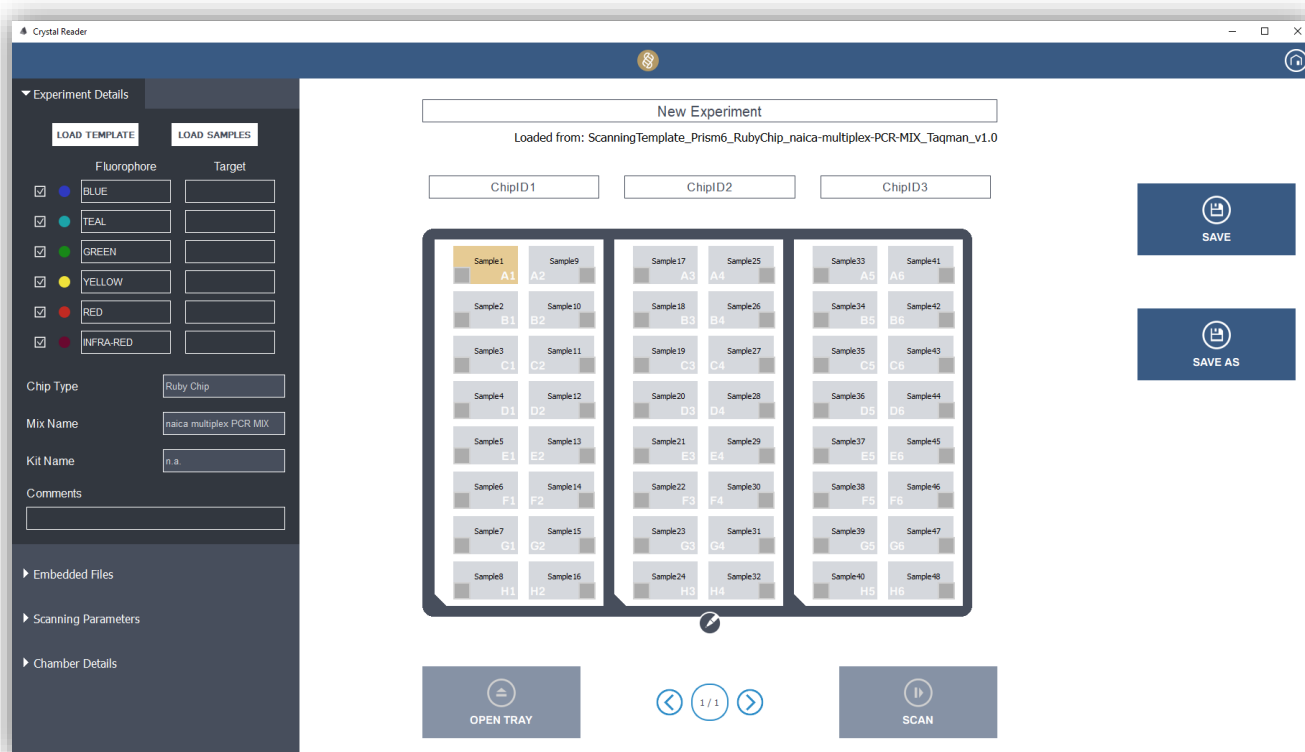


Figure 10: The Crystal Reader software “New Experiment” page for Ruby Chip.

The Experiment Details, Embedded Files, Scanning Parameters and Chamber Details are pre-filled according to the selected scanning template; some settings can be checked and modified if needed at the next step.

3.6.2. Step 2 - Confirm or modify the experiment context

Note: All subsequent sections provide instructions for Sapphire Chip that can also be applied to Ruby Chip.

Name the experiment

Enter the experiment name in the “New Experiment” field.

Note: Stilla Technologies recommends establishing a traceability rule for experiment naming (e.g., date prefix, chip IDs, user initials, objective of the experiment; 20210507_04527983-04568924-04698435_ABC_valid_6plex).

Enter the Chip IDs

The chip ID is a unique identifier that contains traceability information. It is located on the chip box label in both text and QR code formats. Stilla Technologies strongly recommends entering the chip IDs for each experiment. To do so, select Chip ID1, Chip ID2 or Chip ID3 according to the position of the chip in the tray holder and scan the QR code of the corresponding chip with the external barcode reader provided with the naica® system.

The Chip ID can also be manually written, although this is more error prone.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Edit experiment details

In the left panel, drop down the “Experiment Details” menu (Figure 11).

Use the “LOAD TEMPLATE” button to load a new scanning template.

Use the “LOAD SAMPLES” button to load new sample information (see section “ 3 Step 3” for more details about sample information).

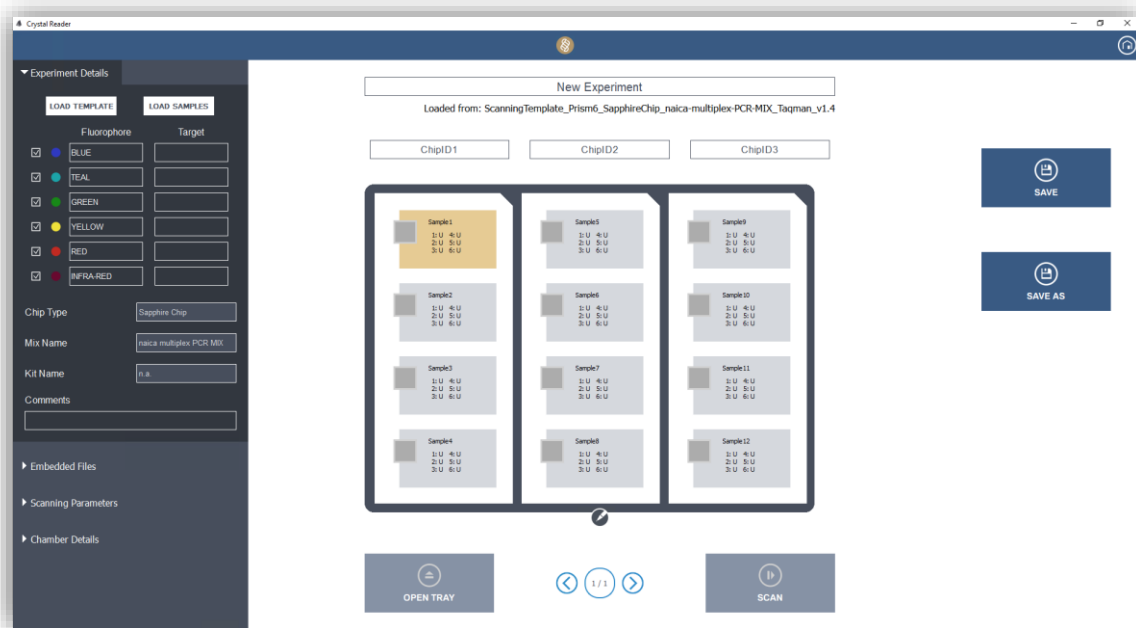


Figure 11: Enter Experimental Details in the “New Experiment” interface.

For each detection channel (“Blue”, “Teal”, Green”, “Yellow” “Red” and “Infra-Red” LED) enter the following (Figure 12):

- The name of the fluorophore excited by the LED.
- The name of the target associated with this fluorophore.
- Select/unselect the checkbox to include/exclude a detection channel for/from scanning.

It is mandatory to select a minimum of three detection channels and to include the blue channel, as it is the main detection channel.

Note: Ensure to select the appropriate detection channel for the experiment to avoid any false-positive results. If the wrong detection channel was selected, re-scan the chips.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

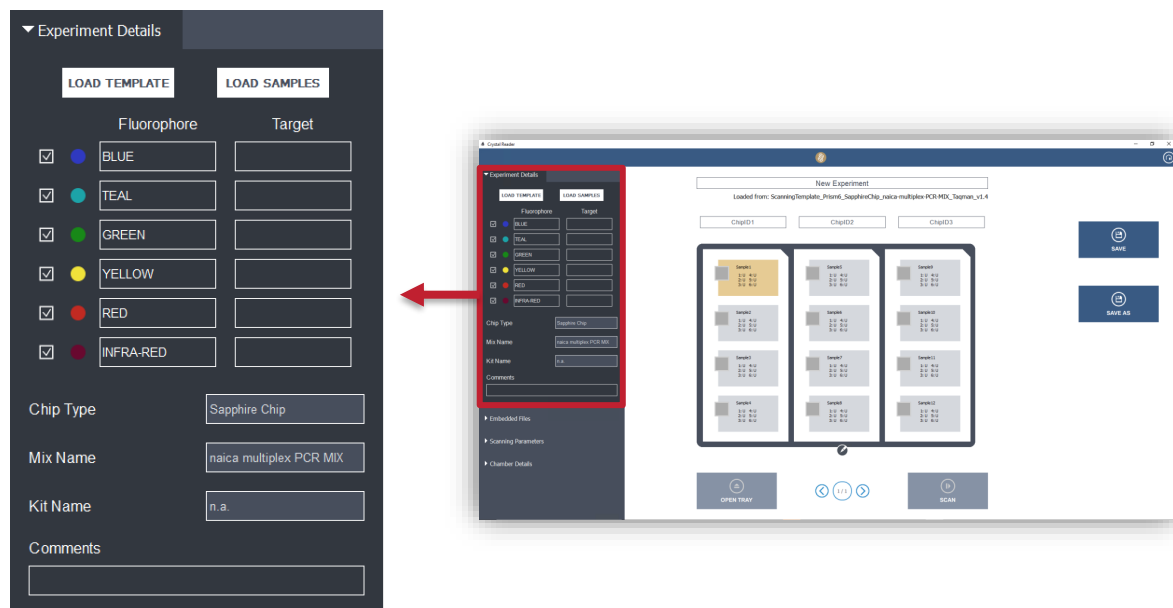


Figure 12: Enter fluorophore and target names in the “Experiment Details” interface.

The “Comments” field can be used to add any information pertaining to the experimental record.

Check the Embedded Files

The embedded file selection menu concerns the supported files that can be pre-loaded in the Crystal Reader software user interface. The only mandatory file is the Image Analysis Configuration File (.yaml), which defines the set of chip type, scanner type and mix type.

The three other files are optional:

- The Spillover Compensation Matrix (.ncm)
- The Plots configurations (.ncp) containing favorite views and the settings of dot-plots.
- The Analysis Configuration (.nca) containing the zoning (thresholds or polygons) and populations settings.

When these files are loaded in Crystal Reader software, they will be saved in the .ncx file together with the raw data. When opening the .ncx file in Crystal Miner software, the display settings of the histograms plots, 1D and 2D dot-plots will be pre-set from the .ncp file and the thresholds and populations will be pre-defined from the .nca file. This is especially useful for routine use and allows the user to directly access the quantification without having to reset the display and population settings for every analysis.

Please refer to the Crystal Miner software User Manual for additional information.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

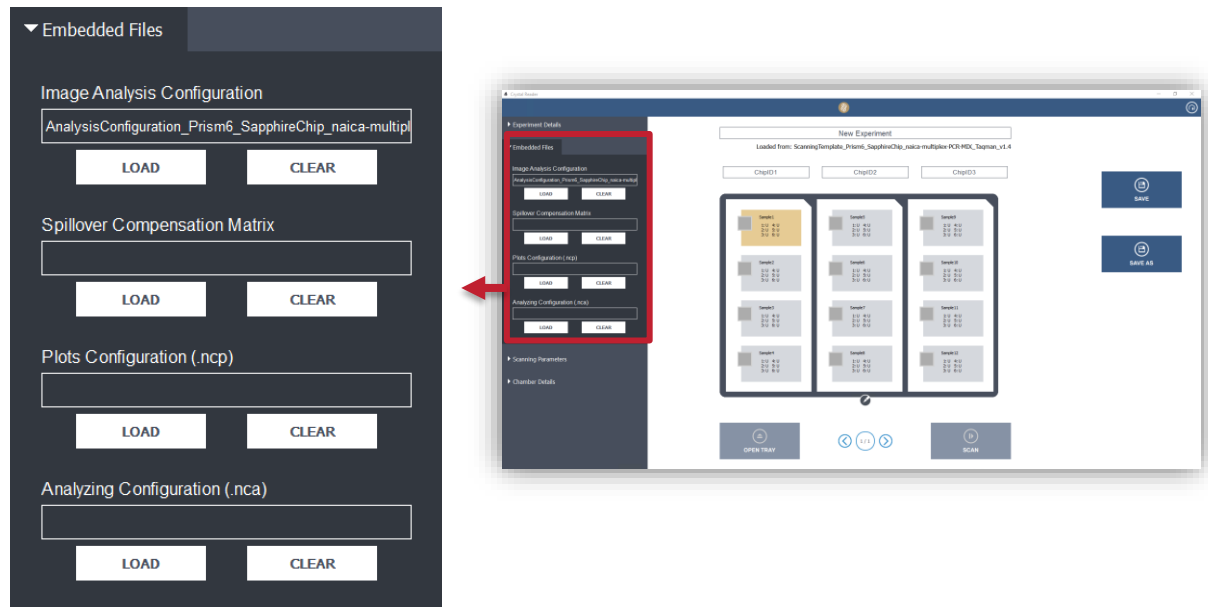


Figure 13: Select the relevant “Embedded Files” for the experiment.

Use the LOAD button to add a file (e.g. to add a spillover compensation matrix if the scanning template does not contain one).

Use the CLEAR button to remove a file if needed.

If the “Spillover Compensation Matrix” field is empty, it is possible to load a spillover compensation matrix that has been created for a previous experiment with the Crystal Miner software matrix computation tool. To load a pre-existing matrix, just click on the “Load” button in the “Spillover Compensation Matrix” widget and select the desired “.ncm” file.

If a spillover compensation matrix is loaded, the fluorescence spillover will be automatically applied, when opening the data later, in the Crystal Miner software.

CAUTION!

The spillover compensation matrix (.ncm), is specific to the set of fluorophores used. However, the matrix may also depend on the biological setup (mastermix used and PCR program) and the scanning parameters (exposure times of the LEDs). When changing an assay parameter or a scanning parameter, Stilla Technologies recommends to either redefine the compensation matrix or to verify that any compensation matrix, defined for the previous parameters, still provides optimal spillover compensation, before validating its use on the assay with the new parameters.

Please refer to the Crystal Miner software User Manual for additional information about spillover compensation.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Check the scanning parameters

Click on the “Scanning Parameters” drop-down to edit the focus value and exposure time of each excitation LED (in ms):

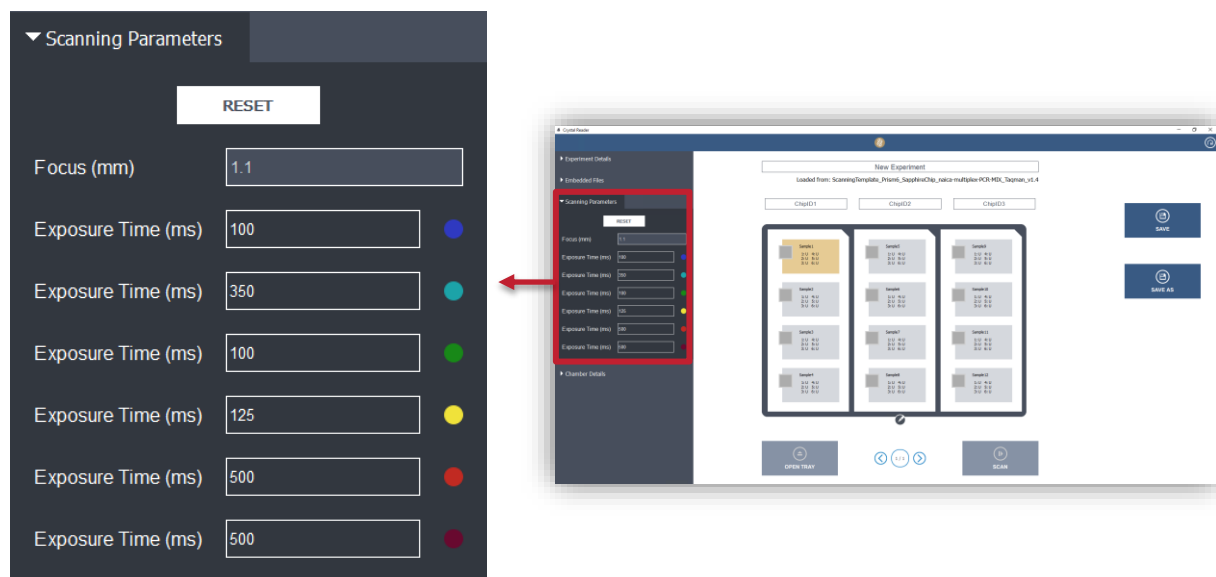


Figure 14: The “Scanning Parameters” menu.

Focus value




- The focus value is a set value based on the original instrument calibration and it does not need to be modified.
- The focus value represents the optimal z-distance of the imaging plane in mm.

Exposure times

- When a scanning template is selected, the default exposure times from the template will be displayed
- Exposure times can be modified to optimize experimental settings. Only positive integer values between 5 ms and 999 ms must be used for exposure times.
- Increasing or decreasing the exposure time of an LED will correspondingly increase or decrease the fluorescence intensity value of the droplets in the image in that detection channel.
- The exposure time of each LED should be high enough to discriminate between negative and positive droplets in the detection channel, but not so high that it saturates the positive droplets in the detection channel (intensity saturation incurs the loss of quantitative information).
- To find the optimal LED exposure times for new experimental conditions, the user may create a control experiment in which only one chamber is activated. In this case a few scans should be run until an optimal set of exposure times is found.
- When at least one chamber of the experiment has been scanned in the experiment, the LED exposure time should no longer be edited. This will ensure that all the chambers of the experiment share the same scanning parameters. If the user tries to change an exposure value, the software will give a prompt, asking if the user wishes to proceed to erase the images of the already scanned chambers.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

3.6.3. Step 3 - Select the chambers to scan and define chamber details

- Click on the “pencil icon”  (located below the chip holder display) to modify the selection of the chips & chambers to be scanned (by default the 12 chambers for the Sapphire Chip and 48 chambers for Ruby Chip are selected).
- Clicking on the “cross icon”  will deactivate the object (chamber, chip, chip holder).
- Clicking on the “plus icon”  will activate the object (chamber, chip).

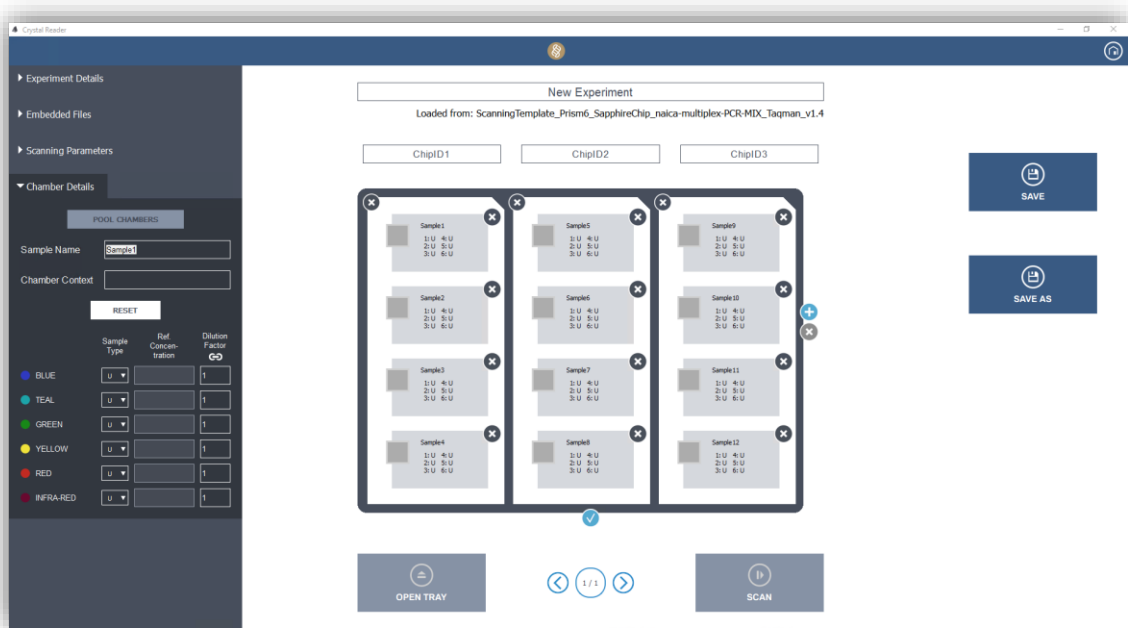




Figure 15: Click on the cross icons to deactivate objects; click on the “plus icon”  to activate objects.

- If an additional chip holder is required, click on the “plus icon”  on the right side of the chip holder display; another chip holder is added to the experiment.
- To navigate from one chip holder to another, click on the left or right arrows around the chip holder counter at the bottom (Figure).

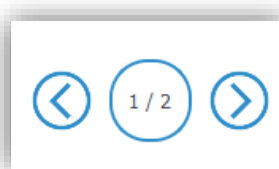


Figure 16: Chip holder counter.

- Click on the “validation icon”  (located below the chip holder layout) to validate the selected chambers.

The Chamber Details panel

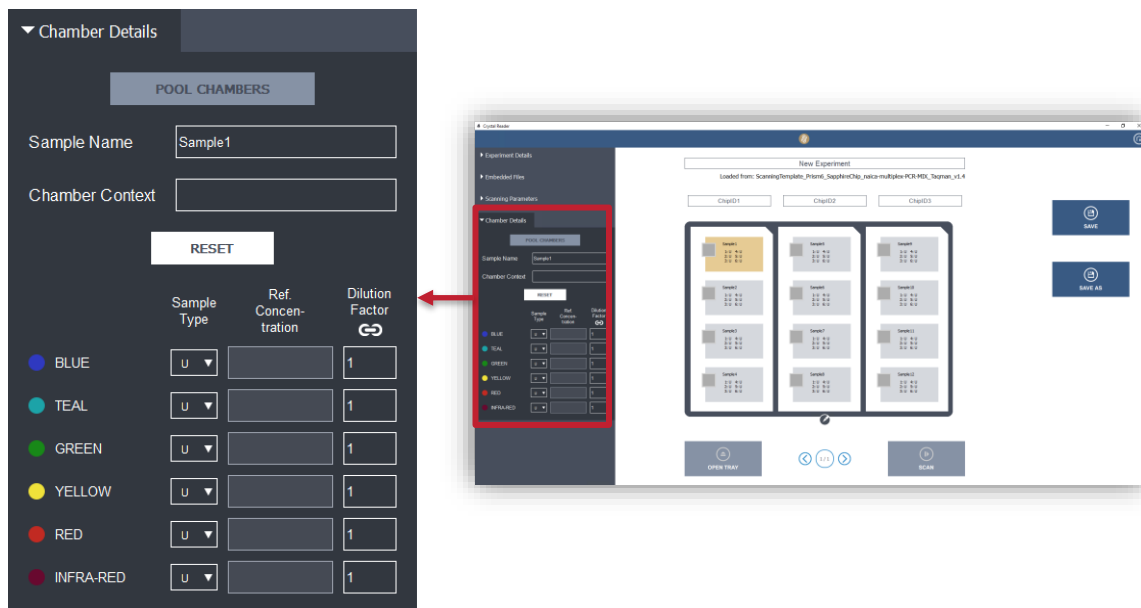



Figure 17: The “Chamber Details” form.

To enter the details of a given chamber; first click on the chamber in the rack layout (chamber rectangle becomes golden), then proceed to edit the “Chamber Details”:

- “Sample Name”: to identify sample loaded in the chamber.
- The USB barcode reader can be used to autofill the chip code to gain time.
- “Chamber Context” (optional): to enter additional information to make the scan data unique (useful if scanning the chamber multiple times).
- For each detection channel (“Blue”, “Teal”, Green”, “Yellow” “Red” and “Infra-Red” LED) enter:
 - “Sample Type” (see details below).
 - “Reference Concentration”. The Sample Type “S” stands for Standard. It is used when a sample of known concentration is processed. In that case, it is possible to edit the expected concentration, in cp/μL. By doing so, both the expected concentration and the measured concentration will be displayed close to each other in the Results page in Crystal Miner software.
 - “Dilution Factor” (e.g., 10 for a 10-fold dilution from the stock sample).
 - By default, the same dilution factor is applied to every channel. To make the dilution values independent from one channel to another, click on the “unlink icon” . Clicking again on this icon will reset the default configuration.
 - If no value is specifically added, the final concentration in copies/μL will correspond to the number of DNA copies for each μL in the whole 25 μL of reaction mix for Sapphire Chip or 7 μL of reaction mix for Ruby Chip.
 - If, for example, the volume of DNA solution is 4 μL out of 25 μL of reaction mix, it could be useful to insert a dilution factor of 6.25 (25 μL/4 μL) to obtain the effective concentration (in copies/μL) in the 4 μL of the initial sample.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

- Then, edit the type of the sample for each given channel, the possible values are:
 - U: Unknown
 - S: Standard. This type is used when a sample with a known concentration is run. In that case, it is possible to edit the expected concentration in the “reference Concentration” field (in copies/μL). This is not mandatory but by doing so, the expected concentration will display next to the measured concentration in the results page in Crystal Miner software.
 - N: Negative control (i.e., there are only negative droplets for the channel of interest)
 - P: Positive control (i.e., there are positive droplets for the channel of interest)

Note: It is possible to navigate between the chambers using the arrows of the keyboard by pressing the “ALT” key. This method may help to save some time editing the chamber context.

- The chamber details are displayed on the chamber button of the chips’ layout:
- The sample ID and sample types per channel are displayed on the chamber in the chips’ layout.

Note: Before the scan is completed, the flag button is grayed and deactivated. The user will be able to click on this button once the chamber has been scanned and analyzed.

The same “Experimental Details”, “Embedded Files” and “Scanning Parameters” will be applied to all samples for an individual experiment during a Prism6 run (see the previous three sections for more details).

If different “Experimental Details” or “Scanning Parameters” are to be applied to the same set of samples for an experiment, chips can be scanned up to three times within 48 hours following the first scan.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

3.6.4. Step 4 - Launch the scan

Click on the “SCAN” button, then validate the path to save scanned data (as a “.ncx” file), and confirm the loading of the first chip holder to launch the scan (Figure 18).

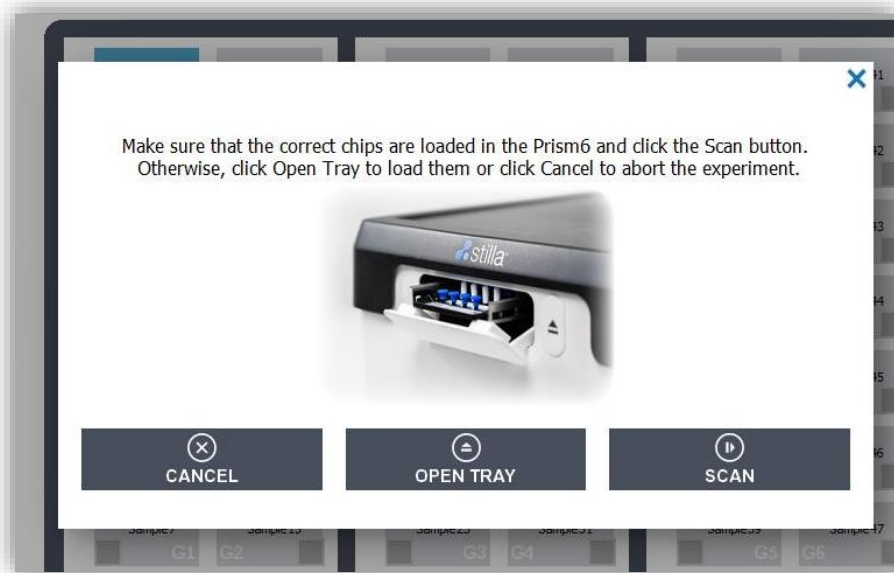


Figure 18: Pop-up confirmation before launching the scan.

When saving an “.ncx” file, the available disk space is checked and the user is warned if the available disk space is less than 10 GB (Figure 19).

If the chips to be scanned are loaded in the scanner, click on “OK”. Otherwise click on “CANCEL” and go back to the previous steps.

Once the scan is processing, a progress bar with the estimated remaining time is displayed.

CAUTION!

Do not push the power button on the front panel of the Prism6 instrument while it is still in operation. This would cause the instrument to shut down with a risk of data being lost, necessitating re-scanning of the chips.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

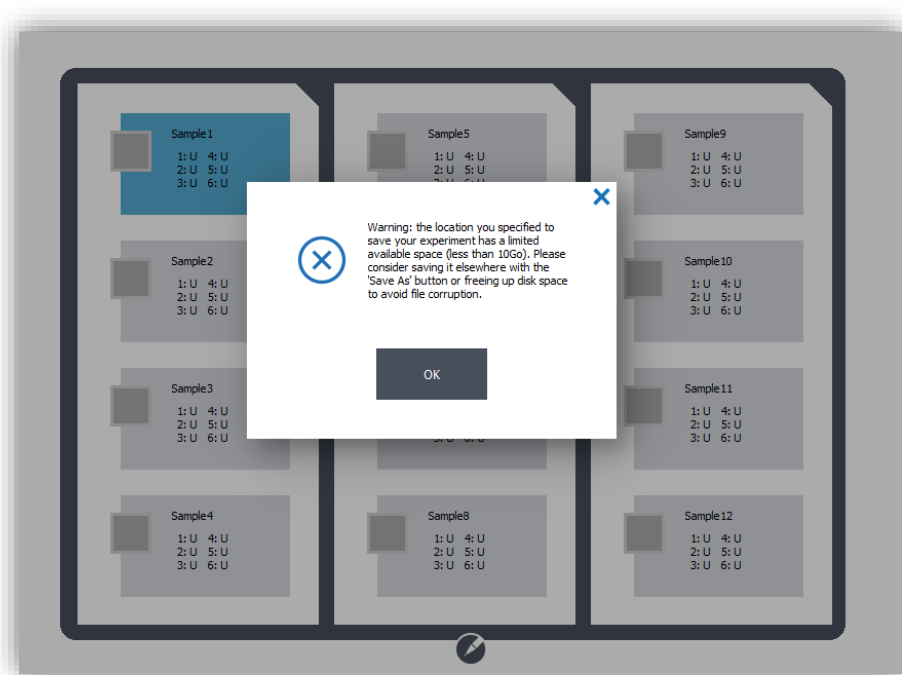


Figure19: Pop-up warning about low disk storage space

The scan can be canceled at any time, but you might lose some scanned data. Editing the experiment context is deactivated during the scan process.

- If several tray holders are to be scanned sequentially as part of a larger experiment, the scanning process is split into several steps:
 - Once the current chip holder is completely scanned, a pop-up window is displayed that instructs to load the next chips.
 - Click on “OPEN TRAY”, place the next chips to be scanned, then click on “CLOSE TRAY”.
 - Click on “Continue Scanning” to confirm the loading of the next chip holder to launch the next scan.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Check chamber image quality

As the individual chambers are scanned, the quality flags are displayed on the Crystal Reader software scan page.

A quality flag can be green or yellow.

Chamber Quality Flag

A chamber quality flag is displayed next to each chamber name on the top left side of the application window. This chamber quality flag provides visual feedback for chamber quality control. Chamber level: the chamber flag corresponds to the lowest quality flag, among the criteria computed for the chamber.

There are 2 types of chamber quality flags:



- “Green” quality flag indicates that the chamber and its picture are good quality.



- “Yellow” quality flag indicates that the software detected major deviations in the analyzed chamber quality parameters. Visual inspection of the Droplet Crystal is required to determine if the experiment may have to be repeated.
- For a visual pre-analysis, while the scanning is still in progress, clicking on one of these quality flags will display the corresponding chamber page.
- In case of non-green flags, a visual inspection of the chamber must be performed in every channel, directly in Crystal Reader software. This is because re-scanning with adjusted Scanning Parameters may suffice to improve the quality of the picture and get a green flag.
- For further details on quality flags, refer to the section 3.6.3 ‘Check image quality’.
- When the image of the last chamber of the experiment is acquired, the following message is displayed:

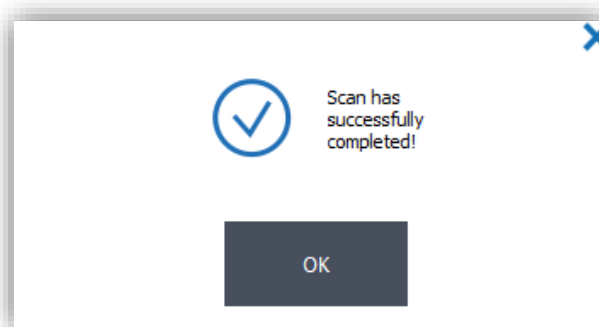


Figure 20: Pop-up signaling a completed scan.

- Click “OK” to display all results of the scanned experiment.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

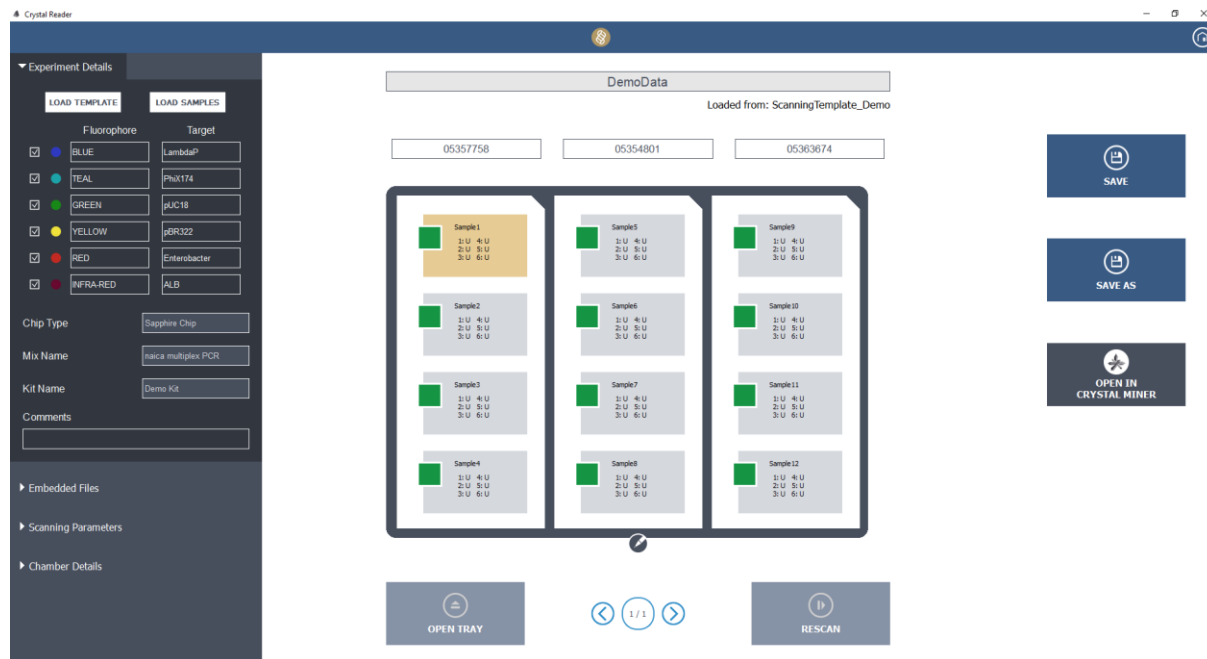


Figure 21: Example of a Sapphire Chip experiment layout with all green chamber quality flags at the end of the scan.

- Note: At the end of the scanning process, the experiment is automatically saved in “.ncx” format in the user-defined output directory under the name: “<ExperimentName>.ncx”.
- The default output directory is: “%USERPROFILE%\Documents\Stilla\CrystalReader”

Stilla Technologies recommends initially checking the scanning performance by inspecting the image quality. Image quality for the scanned experiment can either be directly inspected in Crystal Reader software or later in the Crystal Miner software after clicking on the “OPEN IN ” button:

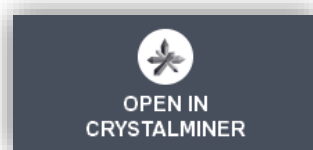


Figure 22: Button to open the scanned experiment with Crystal Miner software.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

3.6.5. Step 5 - Check image quality

Click on the quality flag of a chamber to explore the respective chamber image in detail.

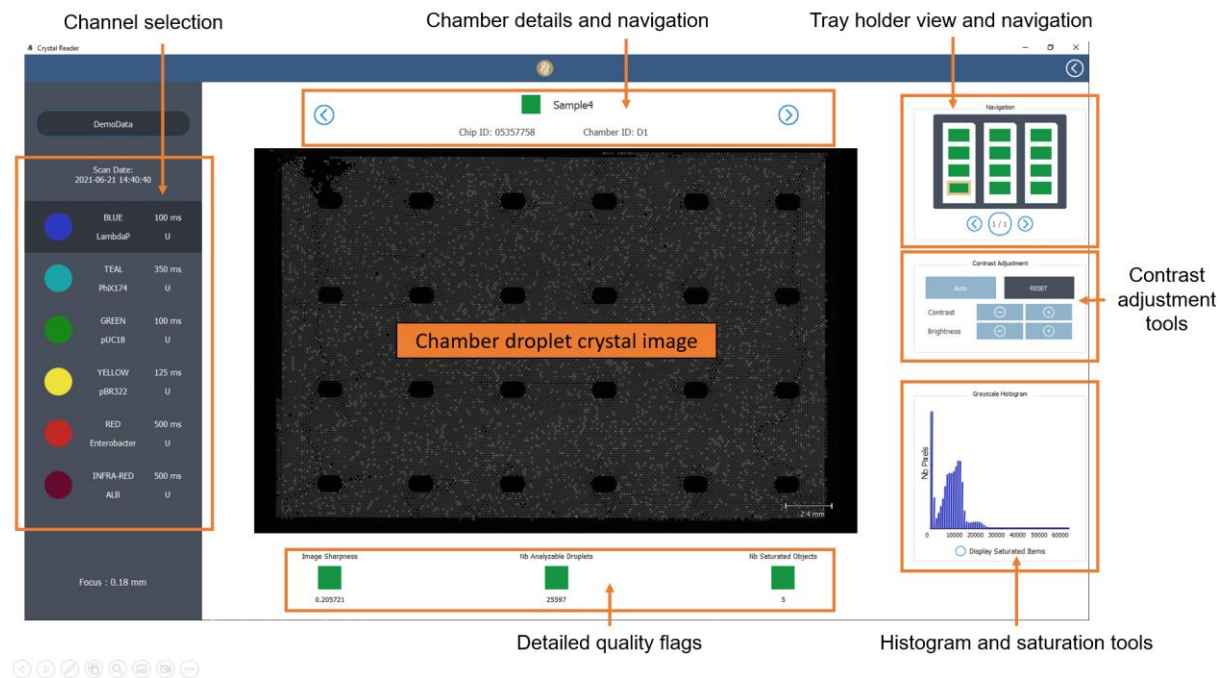


Figure 23: Overview of a Sapphire Chip chamber image in the “Blue Channel” detection channel in the Crystal Reader software.

To explore the images of a chamber:

- Click on the LED icons (on the left panel) to select the acquisition channel.
- Use the mouse scroll to zoom in or out in the image.
- Left mouse click and drag to span the image.
- Click on the “Auto” button to automatically adjust the image contrast.
- Click on the “Reset” button to reset the image contrast.
- It is possible to navigate across the chambers using:
 - the chip holder layout on the top right (click on the rectangle of a chamber to explore it)
 - The left/right arrow buttons, on the top, to explore the previous/next chamber.
 - The left/right/top/left arrows of the keyboard to explore the chamber in the specified direction.
- Check that the negative droplets are visible in the “Blue Channel” image via the reference fluorophore (e.g., fluorescein). If necessary, click on the “Auto” button to adjust the image contrast.

CAUTION!

For every chamber, the negative droplets should be visible in the Blue channel image for the experiment to be analyzable.

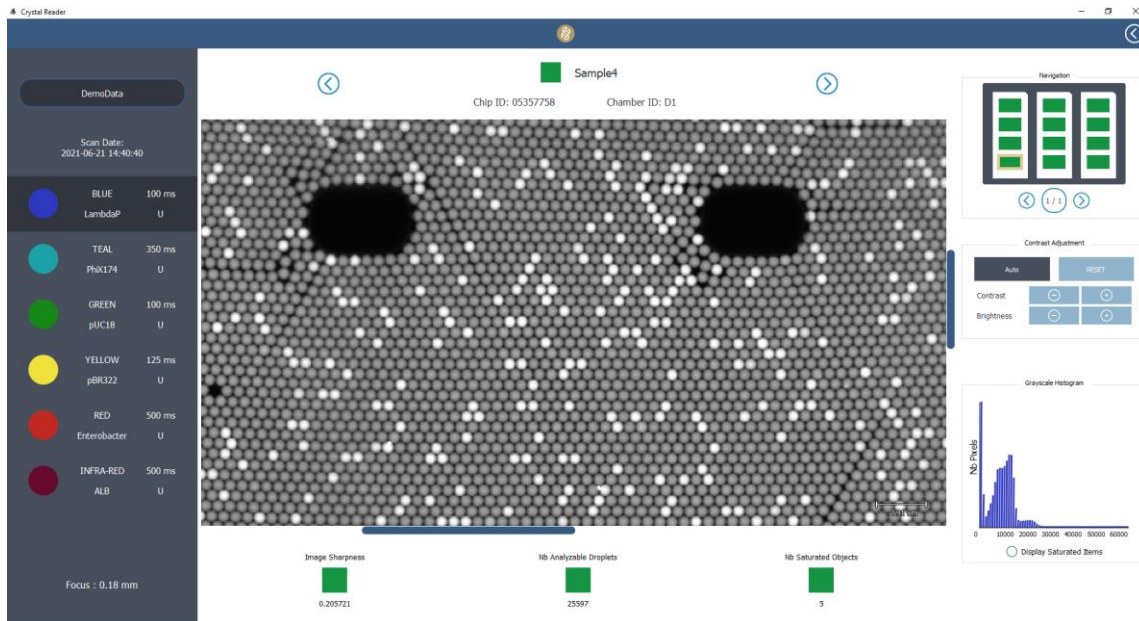


Figure 24: Zoomed chamber image in the Blue detection channel in the Crystal Reader software. All negative droplets should be visible in the Blue channel image.

- The Grayscale Histogram tool (bottom right) helps understand the distribution of pixel intensities expressed in RFUs (Relative Fluorescence Units), where zero RFUs means black and $216 - 1 = 65,535$ RFUs indicates saturation.
- By dragging & dropping in the histogram (with the left mouse button) you can set the minimum and maximum intensity cursors and draw an interval. The pixels with RFU values included in this interval will be highlighted in orange color in the chamber image.
- In the following example we select the interval 24576-37632 RFU. This interval is the selected orange band on the histogram that reveals the positive droplet population in the crystal image.

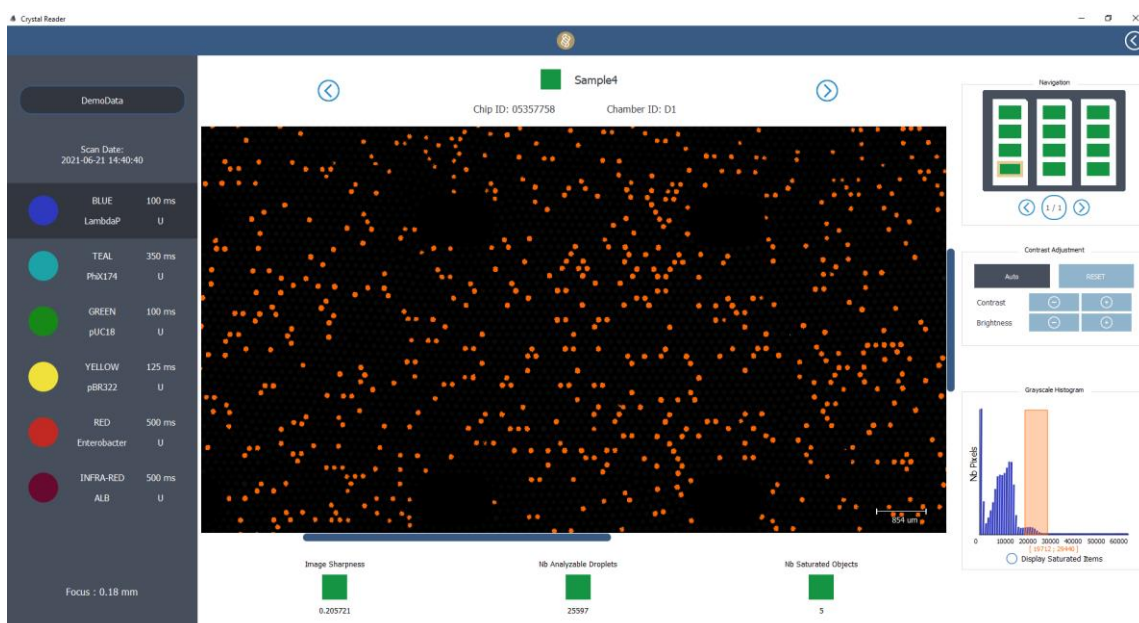


Figure 25: Use the histogram to place intensity cursors and reveal the pixels corresponding to the RFU interval in the Crystal Reader software.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Refer to the “Perform Quality Control” paragraph below for more details about the “Display Saturated Items” checkbox.

3.6.6. Perform quality control:

- Performing the quality control of a given chamber consists first in checking its quality flag. The quality flag of a given chamber is automatically computed based on three criteria, which are themselves also displayed in the form of quality flags. The three criteria are:
 - The “Image Sharpness” which reflects the sharpness of the image. A yellow flag may be caused by an out-of-focus issue. This can be corrected by checking that the chip holder is clean and the chip is flat in the tray holder. If the instrument is still out of focus after checking this, contact customer support for assistance, to run the focus calibration procedure.
 - “Number of analyzable droplets”, which corresponds to the total number of droplets included in the digital PCR computation, after image analysis of the chamber by Crystal Reader software. The higher the number the better the precision of the concentration results.
 - “Number of saturated items”, which are pixels whose fluorescence intensity is above 65,535 RFU. These pixels are automatically excluded from the analysis. By checking the “Display Saturated Items” box under the Grayscale Histogram tool, the saturated pixels are highlighted in orange color in the Droplet Crystal image (Figure 26). Counted saturated objects are either saturated droplets caused by a high exposure time or saturated artifacts located inside the chamber. They can be avoided by checking that exposure times are not too high.

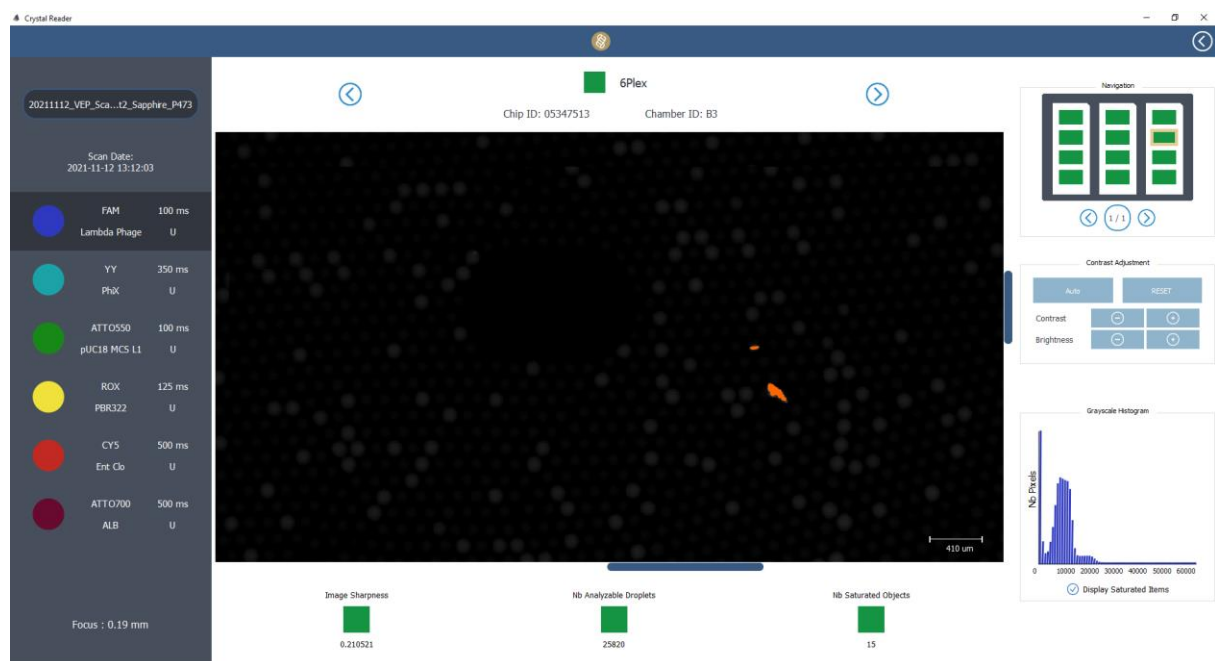


Figure 26: Overlay of the saturation map in the Blue detection channel in the Crystal Reader software (saturated pixels are displayed in orange color). In this case, they are dust particles.

CAUTION!

Ensure that none of the saturated pixels are located in positive droplets, in any channel. Otherwise, these positive droplets will be excluded from the analysis by the software. When positive droplets saturate in a given channel, it means that the chamber(s) has(have) been over-exposed. The exposure time of the corresponding channel must be reduced before rescanning the experiment.

- If exposure times need to be modified to improve the image quality, the chips should be rescanned (see the section “Advanced Functionalities” for more details about chamber rescanning).
- If the underside of a chip needs to be cleaned to improve the image quality of some chambers, the chip should be rescanned for these respective chambers (see the section “Advanced Functionalities” for more details about chamber rescan).
- In case of saturated artifacts with a yellow flag, clean the back surface of the chip by slowly sliding it once along the length of a Precision Wipe and rescan.

The saved experiment can be now analyzed with Crystal Miner software (please refer to Crystal Miner software User Manual available on the Technical Resources webpage for details).

- Before turning the Prism6 instrument off, click on “Open tray”/ “Close tray” to remove the chips. Then press the power button located at the front (the green light becomes orange) for a few seconds before extinguishing.

4. Advanced Functionalities

4.1. How to rescan existing chambers and how to scan new ones

Once a scan is finished, the SCAN button becomes a RESCAN button. It is then possible to, either:

- a) Rescan all the chambers of the experiment.
- b) Rescan only specific chambers of the experiment and keep the images of the other chambers.

To do so, Click on the RESCAN button

Note: If different scanning parameters or experimental details are to be applied to the same set of samples for an experiment, the chips can be scanned up to three times within 48 hours following the first scan.

Irrespective of the option chosen, if there is at least one already scanned chamber when the “SCAN” button is clicked, this window will be displayed:

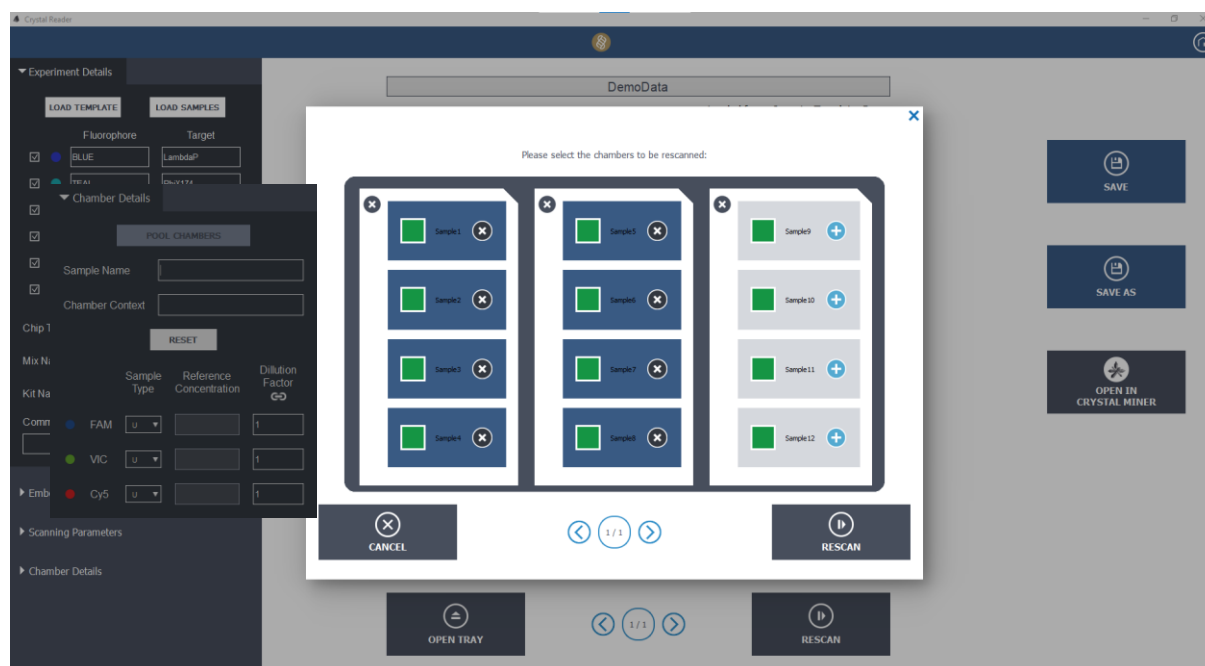


Figure 27: Default rescan window (chambers to be rescanned are highlighted in blue color) in the Crystal Reader software.

In the default configuration, all the previously scanned chambers are selected for rescanning (i.e., they are highlighted in blue color).

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

4.1.1. How to rescan all the chambers of the experiment

For this scenario: click on the Rescan button, and the scan will start.

4.1.2. How to rescan only a subset of chambers

For this scenario: unselect the previously scanned chambers to be kept and click on the new ones desired to be added to the experiment. Click on the Rescan button of the pop-up window and the scan will start.

4.2. How to edit the Prism6 instrument settings

To access the settings panel, click on the “SETTINGS” button from the HOME page in Crystal Reader software. The settings window will display different editable fields.

The following settings can be edited:

- Prism6 instrument settings:
 - Focus Value (mm): z-distance to the imaging plane used for image acquisition. If it is necessary to reset the focus settings, please contact Technical Support for guidance.
 - Prism6 Serial Number (SN): unique serial number of the Prism6 (set by Stilla Technologies).
 - “Focus Calibration” button: advanced functionality to automatically adjust the focus value for the image acquisition process (available only for the lab manager).
- System Settings:
 - Default paths: default output and input directory paths.

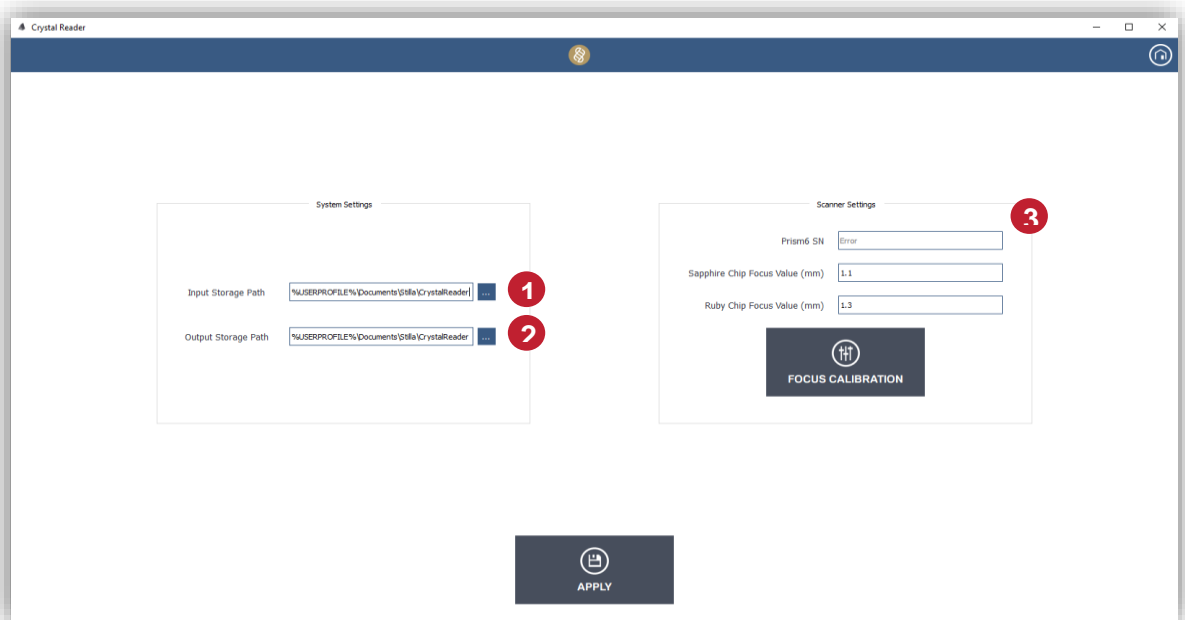


Figure 28: Settings page in the Crystal Reader software.

- A user can:
 - Modify the “Focus Value” used for experiment scanning in the “Scanner Settings”, either manually or via the “Focus Calibration” functionality (see next section), for each chip model (3).


Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

- Modify the default output and input directory paths. System preferences for the input/output storage directories (1 and 2) will be saved after quitting Crystal Reader
- In the path fields, “%USERPROFILE%” can be used; this is a Windows environment variable that redirects to the path of the current Windows session user.
- For example: if the user is logged into Windows with the account “John”, %USERPROFILE% has the following value: “C:\Users\John”
- For calibration purposes, it is possible to perform an automated estimation of the optimal acquisition focus (i.e., optimal z-value) by clicking on the “Focus Calibration” button (see next section).

4.3. How to perform focus calibration

The focus calibration is an advanced feature.

Caution!
It is strongly recommended not to alter focus settings.
If there is a requirement to reset the focus settings, please contact Technical Support .

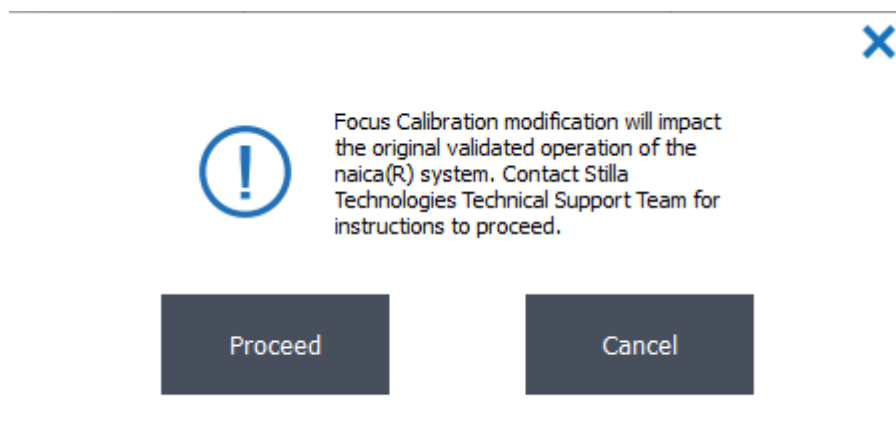

- Assuming that a chip with a chamber of interest is already inserted in the Prism6 instrument, access the SETTINGS menu, click on the “Focus calibration” button, confirm, then select the chip model of interest (Sapphire Chip or Ruby Chip).

Note: For optimal focus calibration on the Sapphire Chip, it is recommended to use at least one chip that includes a chamber with a good quality Droplet Crystal and to place this chip in the middle slot of the chip holder (i.e., slot 2).

Note: For optimal focus calibration on the Ruby Chip, it is recommended to use three chips including chambers with good quality Droplet Crystals in positions A, D, and H.

Select the chamber(s) of interest (i.e., the chamber for which the optimal acquisition focus needs to be calibrated) by clicking on the chamber rectangle in the chip holder layout (the rectangle becomes blue). It is also possible to modify the scanning times, as well as the min, max, and step values for the z-values to be tested.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

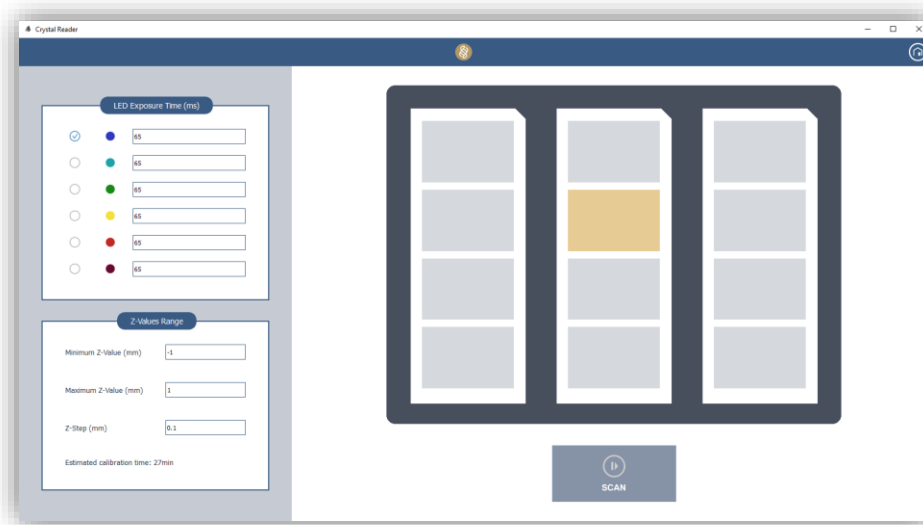


Figure 29: Example of selection of one chamber to calibrate the focus for the Sapphire Chip in the Crystal Reader software.

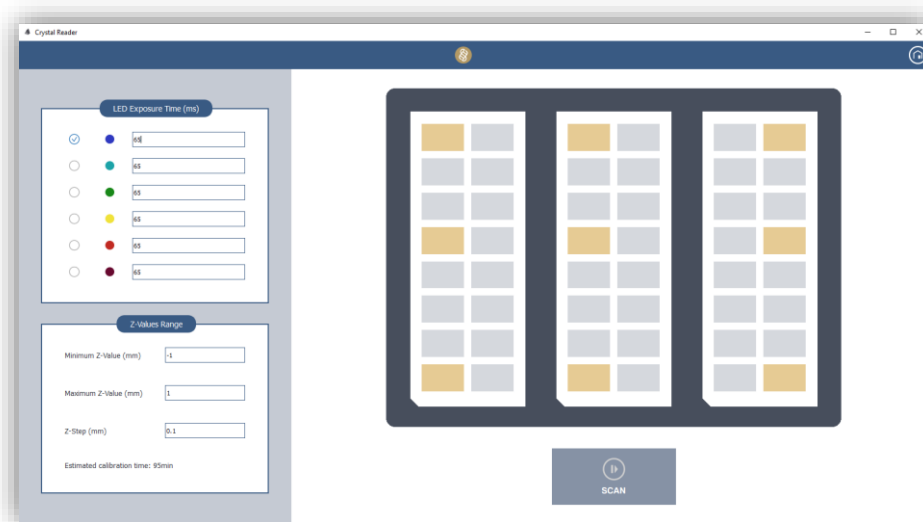


Figure 30: Example of selection of nine chambers to calibrate the focus for the Ruby Chip in the Crystal Reader software.

- Click on the “SCAN” button at the bottom to launch the focus calibration process.
- A progress bar is then displayed with the estimated remaining time.
- At the end of the focus calibration process, a result page is displayed, showing the value of the optimal focus (in mm). This value maximizes the average image sharpness score among all the selected chambers. The result page also shows the chamber image of the first selected chamber at this optimal focus.
- A table with all the computed sharpness scores for each selected chamber and each tested z-value is also provided, together with a graph showing the sharpness score for each chamber as a function of the tested z-value.
- The “manager” user may select the other tested focus values by clicking on them in the table, to check the sharpness of the chamber images acquired at different focus values.
- Once the optimal focus value has been obtained from the calibration experiment, please contact Technical Support for confirmation. The user may keep the optimal

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

focus value or select a different focus value, then click on “Apply this z-value” to apply this new focus value to all the next scanning processes for the chip model of interest.

- To cancel, click on the “Home” button on the top right.

4.4. How to Pool/Unpool Chambers

An additional feature, where users can pool multiple selected chambers containing the same sample, called “POOLING CHAMBERS”, is available in the Crystal Reader software.

Pooling a set of chambers, in which the same sample has been loaded, increases both detection sensitivity and quantification precision. Indeed, by considering each set of pooled chambers as one larger chamber, this pooling strategy allows the analyzed volume to be increased.



Figure 31: Pooling feature of the Crystal Reader software Chamber Details section.


CAUTION!

The chamber pooling functionality should only be used under the assumption that pure replicates have been loaded in the pooled chambers (e.g., the same sample has been loaded).

- If pooling (resp. unpooling) a set of chambers in which the same sample has been loaded is required, simply select these chambers in the chip layout using “Ctrl+Click” or “Shift+Click”, and then click on the “POOL CHAMBERS” (resp.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

“UNPOOL CHAMBERS”) button in the “Chamber Details” tab (

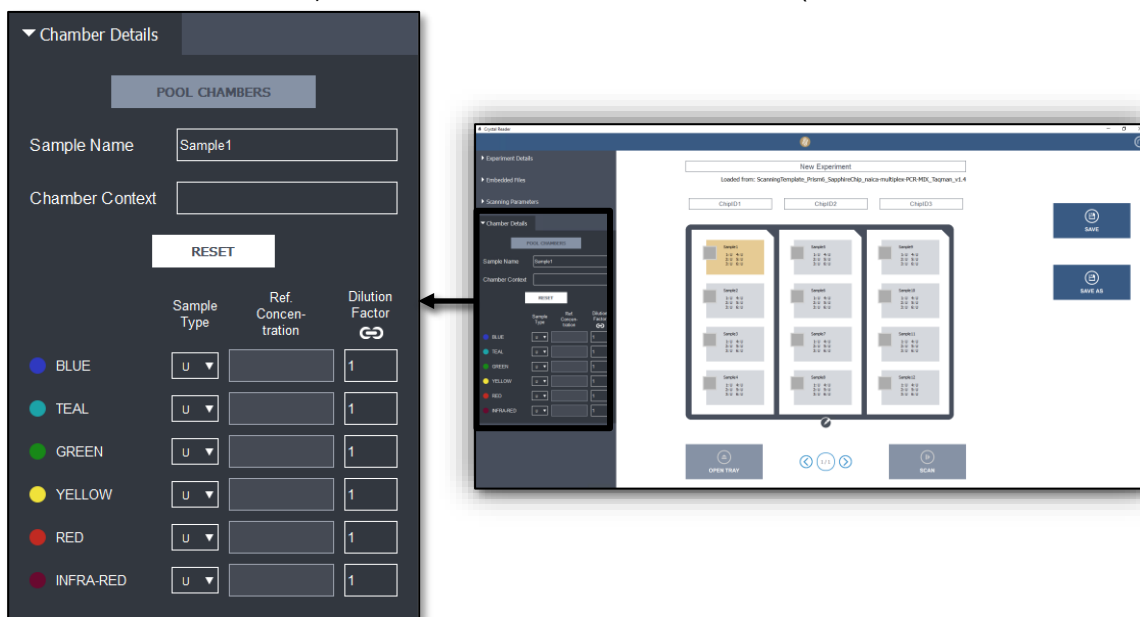

• Figure).

All chambers pooled together will automatically share the same “Pool ID” (displayed as an incremented number in the chip layout), as well as the same “sample name”, “chamber context” and, for each detection channel, the pooled samples will have the same: “sample type”, “reference concentration” and “dilution factor”.

Note: If pooled chambers have not been defined before scanning, it is still possible to pool the chambers after scanning, or during data analysis using the Crystal Miner software (see Crystal Miner software User Manual for more details).

5. Maintenance and Technical Support

Maintenance operations of the naica® system should be executed by a Stilla Technologies Technical Specialist during a visit on-site or by a return shipment of the device to Stilla Technologies premises. Stilla Technologies cannot be held responsible for any intervention or modification done by the user on devices of the naica® system. Before the intervention, we will request the user to decontaminate the instrument following instructions detailed in the Decontamination Protocol and to thereafter fill a Decontamination Certificate; both these documents are provided by a Stilla Technologies Service Specialist.

For technical questions or any issue regarding instrument or software malfunction contact us:


For customers from Europe, China and Africa:
Monday to Friday, 9:30 AM - 6:30 PM, Central European Time (CET).
Closed on french bank holidays.
Phone: (+33) 9 82 27 47 47
Email: support@stilla.fr.

For customers from America, Asia (excluding China) and Oceania:
Monday to Friday, 8:00 AM – 6:00 PM, Eastern Standard Time (EST).
Closed on American bank holidays.
Phone: 1-833-888-0150 ext. 1
Email: support@stilla-inc.com

Online Technical Support is also available at: www.stillatechnologies.com/technical-support/

We will try our best to answer as promptly as possible.

If applicable, Stilla Technologies offers to engage for remote troubleshooting by using the shared desktop application “Team Viewer”, which is installed on the naica® PC.

To enable remote software maintenance on the naica® PC provided by Stilla Technologies, check that the naica® PC is turned on and connected to the internet. Then, double-click on the “Team Viewer” desktop icon  and send both the user ID & password information by email to the Technical Support Team. Please have the following information ready for remote maintenance:

Please ensure to provide the following information when contacting the respective service & support team:

- The Crystal Reader software version number, which is available in “About” in the Home page
- The log files of the experiments which have been generated in the directory: “%USERPROFILE%\Stilla\CrystalReader”

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

5.1. Cleaning and Decontamination

Please see below for maintenance-related instructions for the Prism6 instrument.

For optimal performance, it is recommended to limit the contact of dust particles with the naica® system.

All naica® system devices should be switched off and unplugged before cleaning and decontamination operations. Make sure that there is no chips left in the instrument prior cleaning and decontamination.

External parts of the Prism6 instrument, can be cleaned using a dust-free cloth soaked with one of the following cleaning agents listed below and air-dried.

- In case of an accidental spill of liquid on the instrument, absorb the liquid with and then clean the instrument, using one of the following cleaning agents and air-dry:
Water
- DNase-away
- Ethanol (70%)
- RNase-away
- Bleach (final concentration 5%)
- Ethanol based disinfection solutions (e.g., Phagospray (reference: <https://www.nmmedical.fr/nettoy-desinfec-phago-spray-dm-750ml.html>))

Note: Certain disinfection products have been observed to leave white stains on the Prism6 lid surface.

If you wish to use a similar disinfection product, due to non-availability of the above specified Phagospray reference, Stilla® recommends to first test another alcohol-based product on a small, less visible surface of the top cover of the Prism6 instrument.

Stilla Technologies does not recommend using a disinfection agent based on isopropanol or glutarate e.g., Bacillol® Plus.

Ensure to regularly clean the Prism6 instrument to prevent any marks on the housing or the benchtop.

At least once a month, clean the tray holders of the Prism6 instrument of dust particles using a dust remover.

When decontamination for return shipments is required, please contact Technical Support for the respective protocol and instructions:

Material necessary for the decontamination procedure:

- Gloves
- Safety Glasses
- Mask
- Laboratory coat
- Hydroalcoholic disinfectant solution for device surfaces commonly used for biological and medical devices (e.g., Phagospray).

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

- Decontaminant solution for device surfaces used for biological and medical devices, specifically targeting nucleases and DNA contamination (e.g. RNase away).
- Standard laboratory paper towels.

5.2. Disposal

- Disposal of naica® system equipment

The disposal of the Prism6 at the end of the product's life should comply with the current legislation, in force in the country of use, regarding electrical and electronic waste.

5.3. To ship the instrument back to Stilla®

The Prism6 instrument should only be packed and shipped back to the manufacturer after a Stilla Technologies Service Specialist has deemed it necessary. If shipping the Prism6 is required, use all the original packaging and cables provided upon receipt of the naica® system. To prepare the Prism6 instrument for shipment, please contact Technical Support for detailed instructions.

Ensure that no chips (Sapphire Chip/Ruby Chip) or tray holders (Prism6 tray holder Sapphire Chip or Prism6 tray holder Ruby Chip) are left inside the Prism6 instrument for shipment.

6. Troubleshooting

6.1. Prism6 and Crystal Reader software

DISPLAY SCREEN ERROR MESSAGE	RECOMMENDATIONS
A red light appears in the third position at the front of the Prism6 instrument	Reboot the Prism6 instrument (press the power button on the front for a few seconds until the Prism6 instrument shuts down, then press again to start it). Restart Crystal Reader software.
The Crystal Reader software application stops responding	Go to “Control Panel” > “Task Manager” and stop the Crystal Reader process, then start it again.
Image acquisition issue (Black or non-existent flag)	Re-scan the concerned chamber.
Non-optimal image assembly in a chamber	Re-scan the concerned chamber.
Error message: Unable to scan chamber(s)	Re-scan the non-scanned chamber(s).
Error message: It is not possible to deactivate the blue channel, it is used for droplet detection.	Ensure to select at least 3 detection channels, including the Blue channel.
Error: it is necessary to scan three channels, including the Blue channel, to perform a scan.	
Error: the current Image Analysis Configuration File (.yaml) is not designed to be used on this scanner type. Please ensure the consistency between the “.yaml” file and the experiment settings.	<p>Please use the correct scanning templates provided by Stilla Technologies accessible at “C:\Program Files\Stilla\CrystalReader\templates”.</p> <p>Please ensure the consistency between the “.yaml” file and the experiment settings when editing “.yaml” files.</p>
Error: the selected Crystal Miner Plot configuration file (“.ncp”) uses different channels than the ones activated.	Please use the scanning templates provided by Stilla Technologies accessible at “C:\Program Files\Stilla\CrystalReader\templates”. Please ensure the consistency between the “.ncp” file and the experiment settings.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

<p>Error: the selected Spillover Compensation Matrix file (".ncm") uses a different number of channels than the "current experiment. Please check the consistency of these elements.</p>	<p>Please use the scanning templates provided by Stilla Technologies, accessible at "C:\Program Files\Stilla\CrystalReader\templates".</p>
<p>Error:The Prism6 scanner initialization failed. The instrument configuration may not support the selected channels. Please check the software license in place.</p>	<p>If the Prism6 license file is missing, the Prism6 software that communicates with Crystal Reader software will fail to initialize the instrument. An error message explaining the scenario will be displayed in Crystal Reader software and inform about the missing license. In this eventuality please contact Technical Support.</p>

If these observations persist, please contact Technical Support.



6.2. Sapphire Chip

OBSERVATION	POSSIBLE CAUSE	RECOMMENDATIONS
<p>No, or few, analyzable droplets</p> <p>No air bubble</p>	<p>The partitioning step did not occur</p>	<p>Check that the crucial partitioning step is included in the PCR program used (1st step).</p> <p>If the problem persists, please extract the live and system logs (refer to the section "How to use the Geode Software User Interface" in the Geode User Manual and contact Technical Support.</p>
	<p>Problem during the pressure release step</p>	<p>Check that the crucial pressure release step is included in the PCR program used (last step).</p> <p>If the problem persists, please extract the live and system logs (refer to the section "How to use the Geode Software User Interface" in the Geode User Manual and contact Technical Support.</p>
		<p>Make sure the experiment was not interrupted.</p>



Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

OBSERVATION	POSSIBLE CAUSE	RECOMMENDATIONS
Presence of a large droplet coalescence area in the chamber	Droplet Crystals may be sensitive to static shocks	Clean the foil of the chip with a Staticide anti-static wipe and a Precision Wipe before the loading step.
Droplets overlapping (typically in the bottom right corner of the chamber)	There are too many droplets within the chamber	Stacked droplets are considered artifacts and thus excluded from the analysis. To relax the crystal, bring back the chips to the Geode. In the menu, select Template > Protocols and launch the program named "Sapphire Protocol Unpacking Droplets". After the reading step, the neural network of Crystal Miner should recognize more droplets.
	Droplets are too soft (i.e., the surface tension is too low)	
	The chip has been tilted for a few minutes	
The droplets in the chamber have low background fluorescence	Droplets are not detected.	Check that the appropriate amount of reference fluorophore (e.g., fluorescein at 100 nM) was included in the reaction mix.
Presence of droplets of different sizes in the chamber.	Some microchannels are clogged.	Check the sample preparation conditions and/or the PCR mix for precipitates. Contact Technical Support for further information.
	There was a slight static shock on the chip	Clean the foil of the chip with an anti-static spray and a Precision Wipe before the loading step.
While using EvaGreen®, random exclusion of droplets.	The fluorophore EvaGreen® usually provides basic fluorescent background, necessary for droplet detection. In rare cases, this would not be sufficient.	Contact Technical Support.



Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

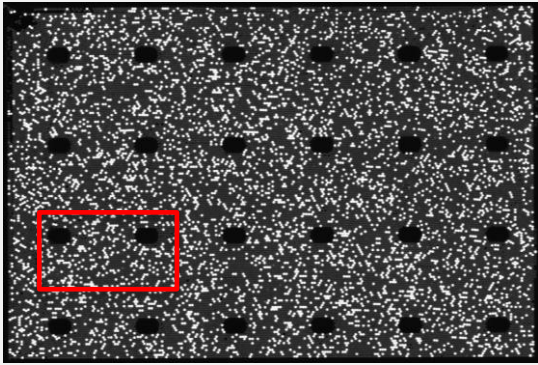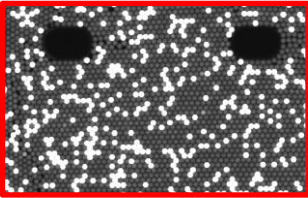OBSERVATION	POSSIBLE CAUSE	RECOMMENDATIONS
<p>Presence of an out of focus sub-image.</p>	<p>Autofocus failed during sub-image scanning</p>	<p>Rescan the concerned chamber.</p>
<p>Oversegmentation or undersegmentation of droplets.</p>	<p>There are two or more droplets (crosses) instead of one (oversegmentation) or one droplet instead of two or more (undersegmentation) which may lead to false positive or false negative results.</p>	<p>Manually inspect the droplets, in case of red crosses, the software automatically rejects the droplets, no action is required. However, if the crosses are green, then manual exclusion of the crosses is necessary.</p>
<p>Droplets with a saturated fluorescence level (saturation is indicated by the quality flags).</p>  	<p>The exposure time is too long for at least one channel</p>	<p>Saturated droplets have an intensity equal to $2^{16} - 1 = 65535$ RFU in at least one channel. They can be visible by clicking on "Display Saturation Map". Decrease the exposure time (usually by a factor 2) set for the detection channel where saturation is observed and re-scan the experiment.</p>



Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip
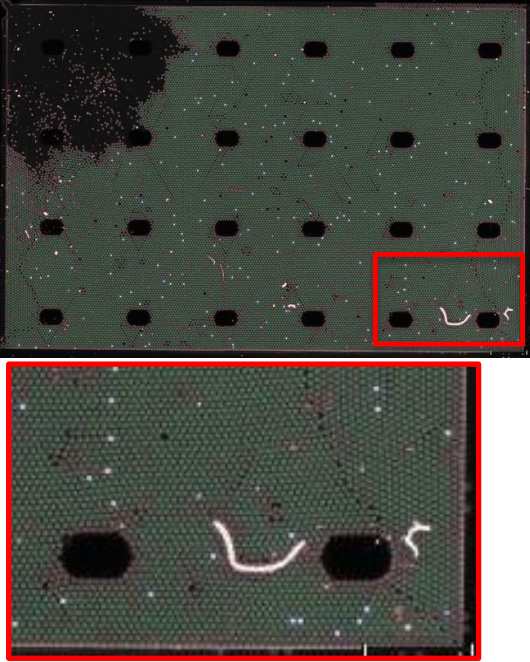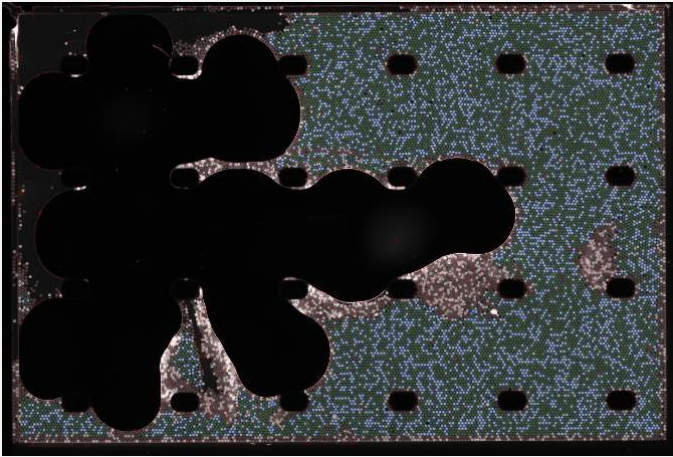
<p>Shiny dust particles on the external film of the chip.</p> 	<p>Presence of dust particles on the chip foil</p>	<p>Clean the foil with a Precision Wipe and re-scan the concerned chamber.</p>
<p>Large air bubbles in the chamber</p> 	<p>Air was injected into the chamber during partitioning</p>	<p>Ensure to pipet 25 μL of the reaction mix. The volume should reach the top of the inlet.</p> <p>Be sure to not introduce air bubbles at the top of the inlet port or between oil and PCR mix during the pipetting step.</p> <p>Check the expiry date of the used chips listed on the box.</p> <p>Make sure the white PCR caps are on the inlet port during PCR.</p>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

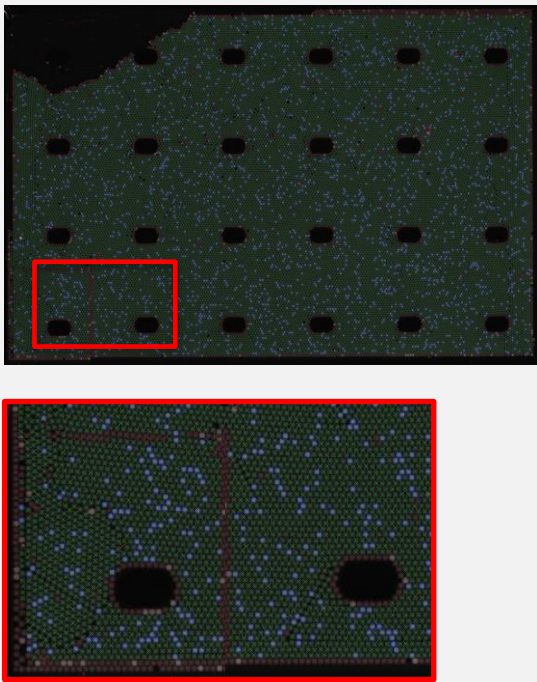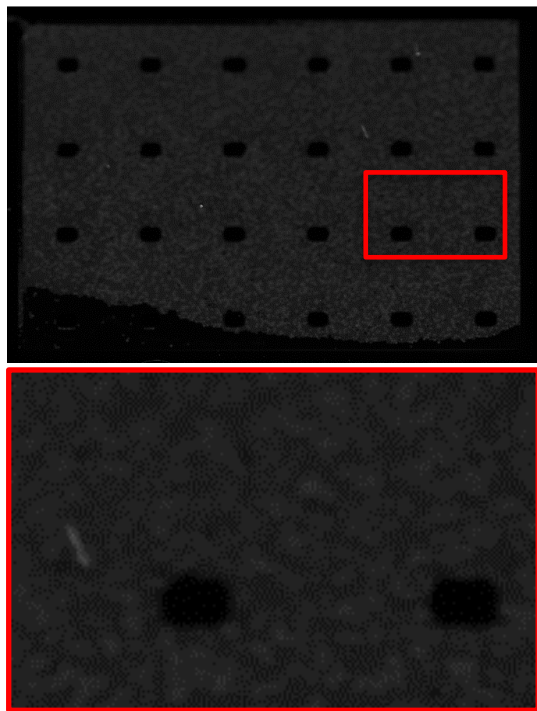
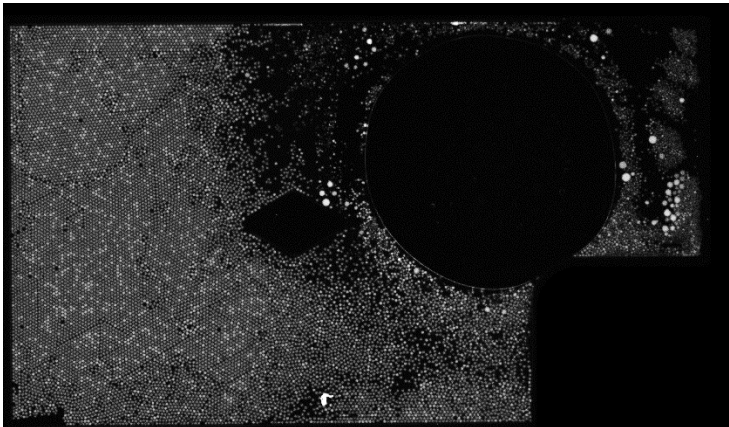
<p>Image stitching event</p> 	<p>The Crystal Reader software stitches several images of the chamber taken subsequently into one final image. Droplets along the stitching lines are occasionally excluded to avoid over/under quantification of the final droplet numbers.</p>	<p>No action is required. Occasional image stitching events are expected with Crystal Reader software imaging. Image stitching events can occur with both chip types- Sapphire Chip and Ruby Chip.</p>
<p>The whole chamber is out of focus</p> 	<p>The chip and/or the holder is not well positioned</p>	<p>Verify that the chip and/or chip holder is positioned flat.</p> <p>A foreign particle might be present between the chip and the holder, forcing the chip to be tilted. Please clean the foil of the chip and the holder with a Precision Wipe and re-scan the concerned chamber.</p>
	<p>The focus value is not well calibrated</p>	<p>Please contact the Technical Support team for assistance.</p>




Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip
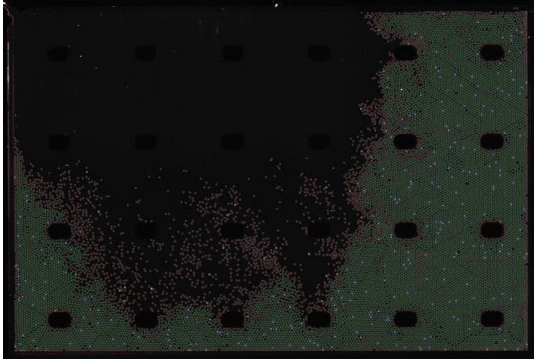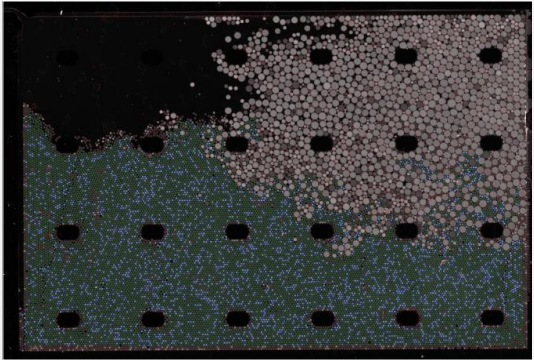
<p>The chamber is not optimally filled</p> 	<p>The sample volume loaded in the chamber was low.</p>	<p>The chamber is still analyzable as the partitioning step was successful. Depending on the assay, it is at the user's discretion whether the number of analyzable droplets is sufficient.</p>
<p>Different sizes of droplets present in the chamber (polydispersity)</p> 	<p>The procedure using the antistatic treatment was omitted or not properly performed.</p>	<p>The chamber is still analyzable as the partitioning step was successful. Depending on the assay, it is at the user's discretion whether the number of analyzable droplets is sufficient.</p> <p>Please refer to the IFU, for the Sapphire Chip or the Ruby Chip, for this step.</p>



Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip



6.3. Ruby Chip

OBSERVATION	POSSIBLE CAUSE	RECOMMENDATION
<p>Empty Chamber</p> 	<p>The sample was not pipetted in the correct well.</p> <p>The sample was ejected from a tip that was introduced too deep (and could touch the bottom of the well) or insufficiently deep (i.e. it could have been at the air-oil interface).</p> <p>When pipetting with a multichannel pipet, the 8 tips were not perfectly horizontal, hence some of them were either above the oil or touching the bottom of the well.</p>	<p>Take stock of the numbering on the wells to ensure that samples are always pipetted in the correct well.</p> <p>When pipetting your sample, stabilize the pipet tip about 1mm above the bottom of the well just by lifting it slightly from the bottom of the well (i.e. neither in contact with the bottom of the well, nor outwith the oil) before ejecting your sample from the tip</p> <p>For more information about how to properly pipet your sample in the Ruby Chip, please refer to the IFU for Ruby Chip.</p>
<p>Bubble</p> 	<p>Small bubbles can sometimes occur during the PCR run, due to heat and pressure cycles.</p> <p>If you observe big and frequent bubbles, the inflated bag containing the Ruby Chip might have been damaged during transport.</p>	<p>It is recommended to check that the inflated pouch is still airtight upon reception.</p>

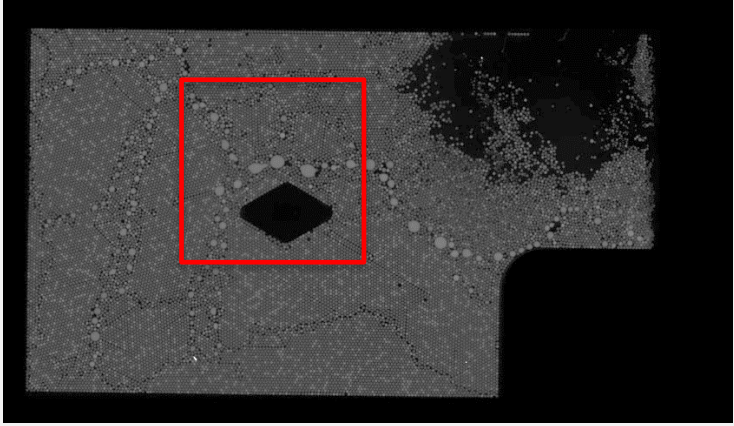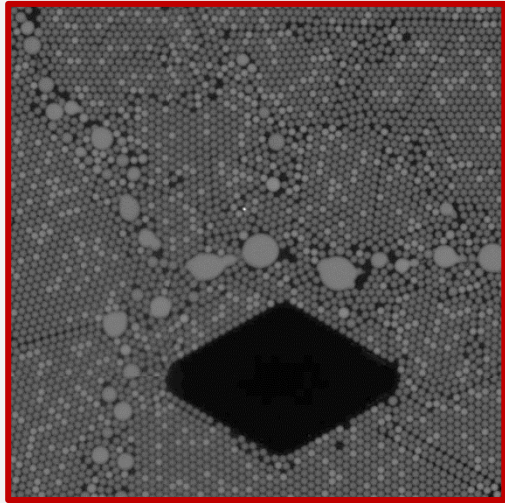
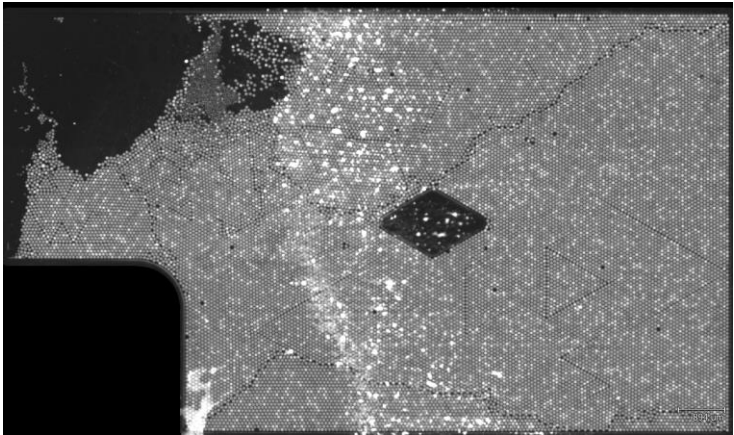
<p>Electrocoalescence</p>  	<p>The procedure using the antistatic wipes was omitted or not properly performed.</p>	<p>Apply the antistatic wipes on all the chips as described in the IFU for the chips. In case of further questions or support, please contact Technical Support</p>
<p>Traces of antistatic product</p> 	<p>The presence of residual antistatic product under the chip generates artifacts in the blue channel.</p>	<p>After applying the antistatic treatment, wipe off the excess of antistatic product with a piece of Precision Wipe according the IFU instruction.</p>

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip
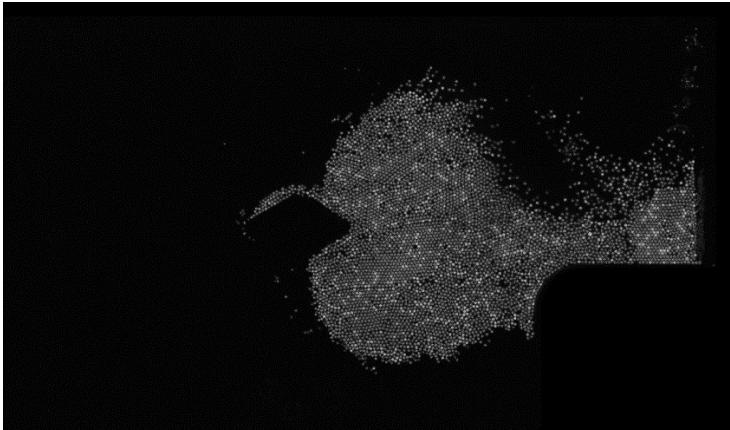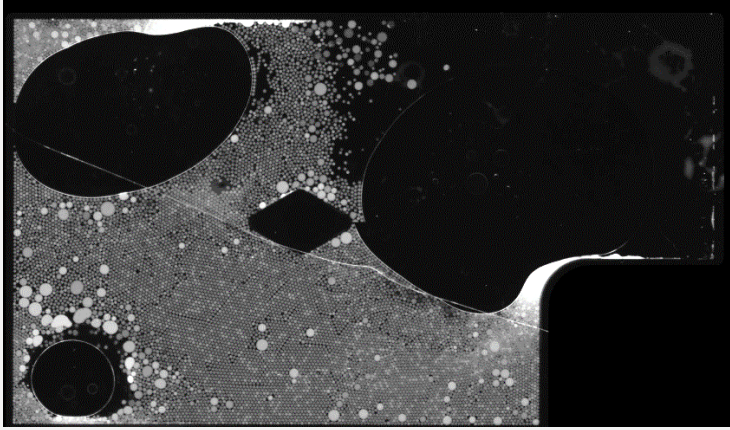
<p>Low number of droplets</p> 	<p>The sample was ejected from a tip that was introduced too deep (and could touch the bottom of the well) or insufficiently deep (i.e. it could have been at the air-oil interface).</p> <p>When pipetting with a multichannel pipet, the 8 tips were not perfectly horizontal, hence some of them were either above the oil or touching the bottom of the well.</p>	<p>When pipetting your sample, stabilize the pipet tip about 1mm above the bottom of the well just by lifting it slightly from the bottom of the well (i.e. neither in contact with the bottom of the well, nor out with the oil) before ejecting your sample from the tip</p> <p>For more information about how to properly pipet your sample in the Ruby Chip, please refer to the IFU for Ruby Chip.</p>
<p>Tension on transparent foil leading to oil traces</p> 	<p>Incorrect pipette tips have been used to pierce the Ruby Chip</p> <p>Not all the 16 wells were pierced before launching the PCR program.</p> <p>The heating plate was not clean, and a dust particle generated tension on the transparent foil.</p>	<p>Only use 10µL and 20µL universal pipette tips as specified in the IFU.</p> <p>Always pierce all 16 wells of each chip, even when some wells will be left empty, before launching the PCR program.</p> <p>Keep the heating plate clean from oil and dust. Clean traces of oil on any part of the Geode.</p> <p>For more information about how to properly clean the Geode, please refer to the cleaning and decontamination standard procedure.</p>
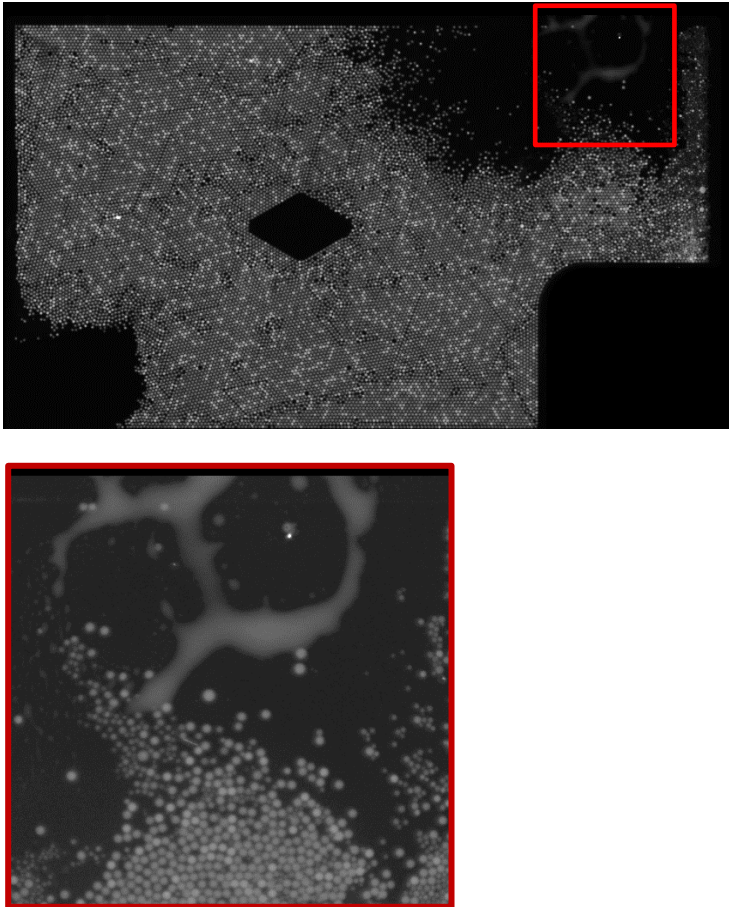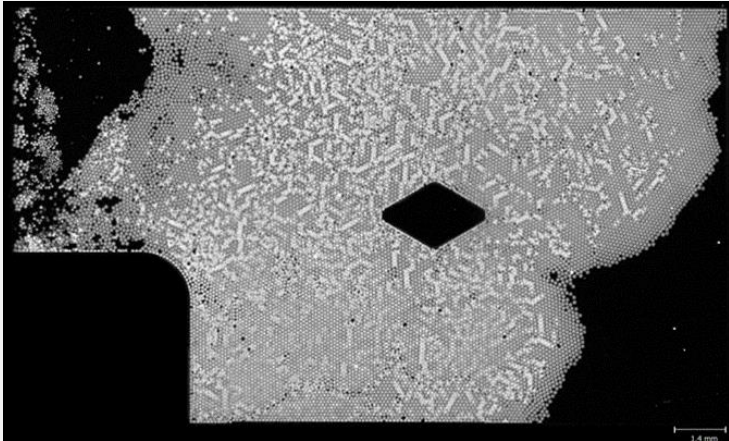
<p>Oil Traces</p> 	<p>The chip contained oil that could condense outside of the wells in normal conditions of use.</p>	<p>Keep the heating plate clean from oil and dust. Clean traces of oil from any part of the Geode.</p> <p>For more information about how to properly clean the Geode, please refer to the cleaning and decontamination standard procedure.</p>
<p>Wetting Droplets</p> 	<p>A higher concentration of qScript XLT 1-Step RT-qPCR ToughMix may cause droplet to wet the surface of the chamber rather than staying spherical during RT-PCR experiments.</p>	<p>These droplets are excluded during the analysis. Use the recommended XLT mix concentration (1 x). Avoid leaving the RT-PCR mix too long in the open air (e.g. in a microplate) before injection into the chip.</p>



Bigger Droplets

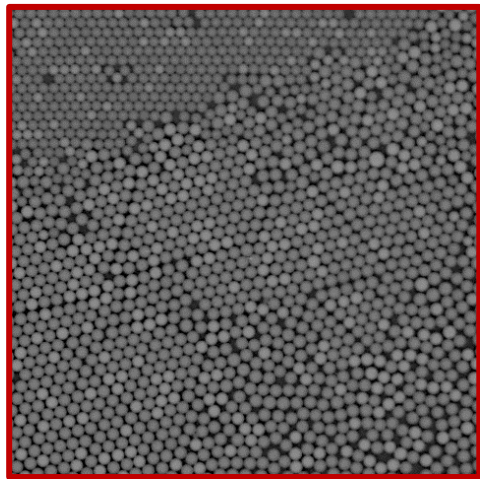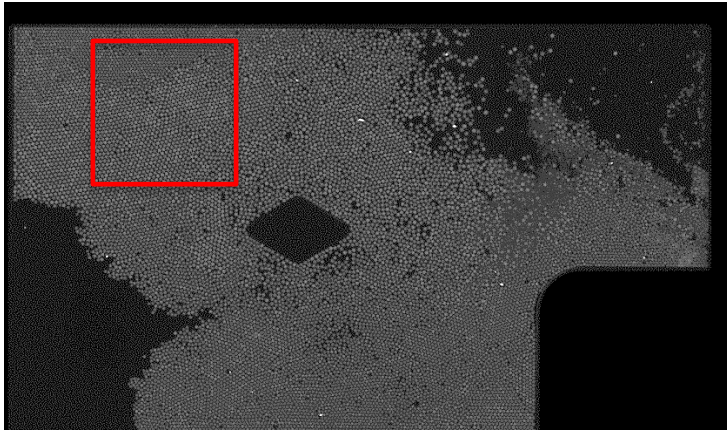

The cycling of temperature and pressure could cause, in some rare cases, the presence of these artifacts.

This population of droplets is excluded during the analysis. The software recognizes the different sizes of the droplets and considers them as artifacts. This phenomenon does not normally affect the output of the assay.

Small Droplets

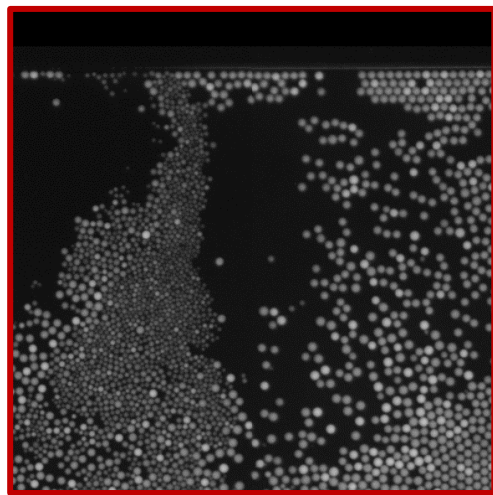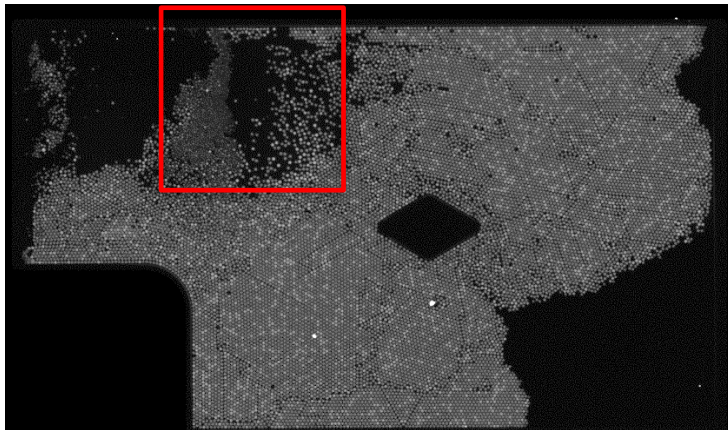


The cycling of temperature and pressure could cause, in some rare cases, the presence of these artifacts.

This population of droplets is excluded during the analysis. The software recognizes the different sizes of the droplets and considers them as artifacts. This phenomenon does not normally affect the output of the assay.

Crystal Reader software v4.0.5.3 / Sapphire Chip & Ruby Chip

Non homogenous distribution of positives

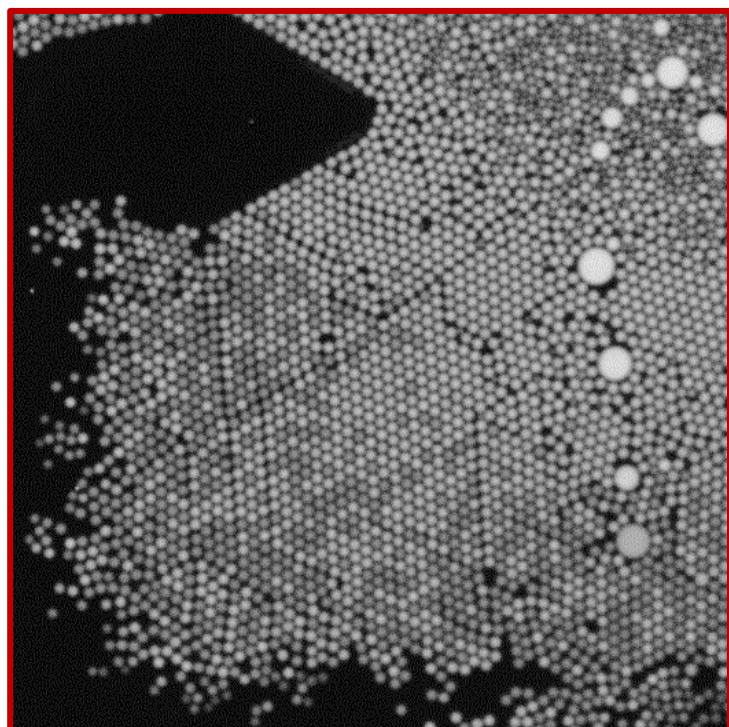

A part of the sample was injected in the chamber after the PCR cycles had begun. Hence, they were amplified in bulk before being partitioned and they present a higher ratio of positive droplets.

The possible reasons for this late injection are:

The sample was ejected from a tip that was introduced too deep (and could touch the bottom of the well) or insufficiently deep (i.e. it could have been at the air-oil interface).

When pipetting with a multichannel pipet, the 8 tips were not perfectly horizontal, hence some of them were either above the oil or touching the bottom of the well.

When pipetting your sample, stabilize the pipet tip about 1mm above the bottom of the well, by just lifting it slightly from the bottom of the well (i.e. neither in contact with the bottom of the well, nor outwith the oil) before ejecting your sample from the tip

For more information about how to properly pipet your sample in the Ruby Chip, please refer to the IFU for Ruby Chip.

7. Software license Information

Crystal Reader software license


©2015-2022 Stilla Technologies. All rights reserved.

For Research Use Only.

Not for use in diagnostic procedures unless specified otherwise by domestic registration.

7.1. How to view the software version

- From the HOME page of the Crystal Reader software, click on the “ABOUT” button. The software version as well as the third-party license information will be displayed:

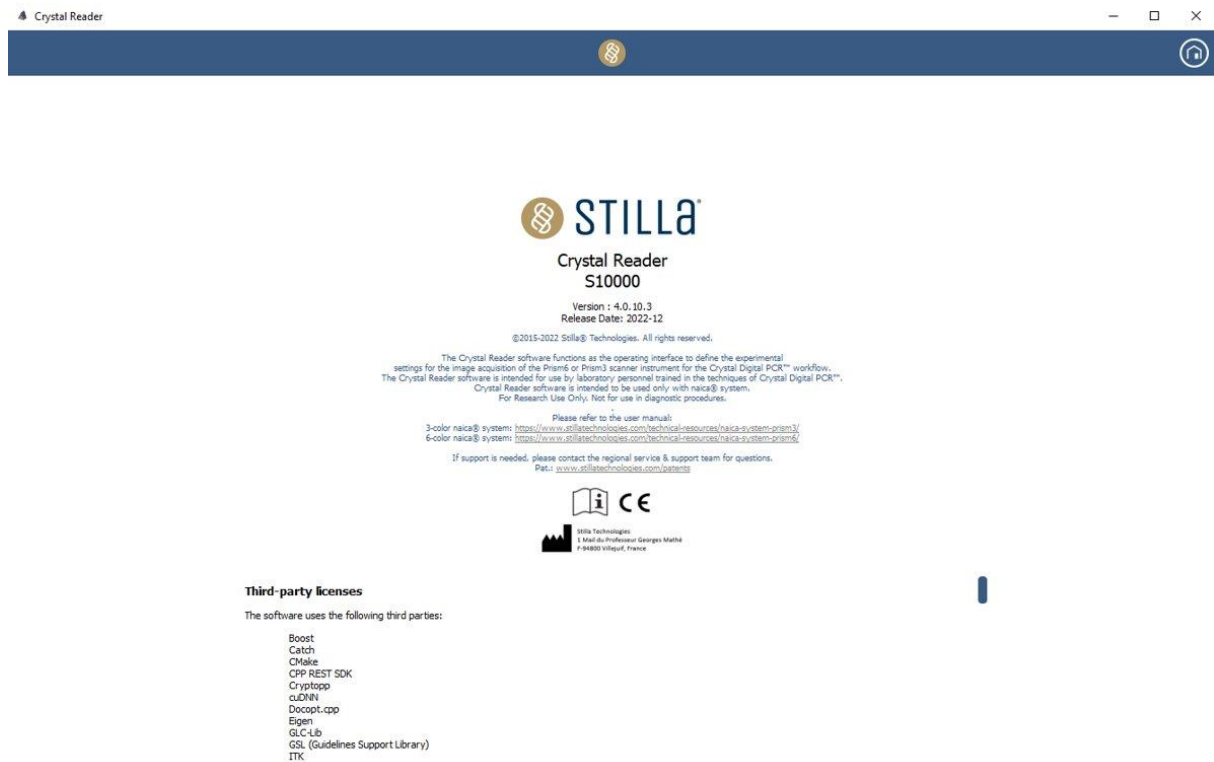

The “About” page of the Crystal Reader software.

7.2. Third-party licenses

The Crystal Reader software uses the following third-party software components:

- Boost
- Catch
- CMake
- cpprestsdk
- cryptopp
- cuDNN
- Docopt.cpp
- Eigen
- GLC-Lib
- GSL (Guidelines Support Library)
- ITK
- JSON for Modern C++
- jwt-cpp
- onnxruntime
- OpenCV
- Python
- Qt
- Quazip
- Qwt
- stlab
- xInt
- Yaml-cpp
- ZLib

To view the license information, please visit the third-party websites or check the Crystal Reader software and Crystal Miner software installation directory (by default in “C:\Program Files\Stilla\CrystalMiner\licenses”).

Note: The license information of all the third-party software components is also accessible in the Crystal Reader software, by clicking on the “About” button on the Home page.

Prism6 and Crystal Reader software for the 6-color naica® system

Headquarters

Stilla Technologies

Villejuif Biopark

1, Mail du Professeur Georges Mathé

94800 Villejuif, FRANCE

Phone: (+33) 09 82 29 50 50

North American Office

Stilla Technologies Inc.

20 Blanchard Road, Suite 9,

Burlington, MA 01803, USA

Tel : 1-833-888-0150

www.stillatechnologies.com

Pat.: www.stillatechnologies.com/patents



MKT-00142 Rev.B