

ddPCR™ Multiplex Supermix

Catalog #	Description
12005909	ddPCR Multiplex Supermix , 1.2 ml (2 x 0.6 ml vials), 200 x 20 µl reactions
12005910	ddPCR Multiplex Supermix , 3 ml (5 x 0.6 ml vials), 500 x 20 µl reactions
12005911	ddPCR Multiplex Supermix , 12.5 ml (5 x 2.5 ml vials), 2,500 x 20 µl reactions

For research purposes only.

Description

ddPCR Multiplex Supermix is a 4x concentrated, ready-to-use reaction mix containing all components — except primers, probe(s), and template — required for probe-based Droplet Digital™ PCR (ddPCR). The hot-start feature of the polymerase in the supermix enables partitioning of sample into droplets while keeping the enzyme inactive at ambient temperature. The supermix has been optimized to provide higher capacity and empower higher-order multiplexing when you use probe-based assays.

Storage and Stability

ddPCR Multiplex Supermix is stable at –20°C through the expiration date printed on the label. Once thawed, it can be stored at 4°C for up to 2 weeks. Repeated freezing and thawing of the supermix is not recommended.

Bio-Rad sells Dithiothreitol (DTT) (catalog #12012171) separately at a 300 mM concentration. DTT tubes should be thawed only once and can be stored at –20°C for 1 year. After thawing the tube, it can be stored at 4°C for up to 2 weeks.

Quality Control

ddPCR Multiplex Supermix is free of contaminating DNase and RNase. Stringent specifications are maintained to ensure lot-to-lot consistency.

Recommendations for Optimal Results

DNA digestion with a restriction endonuclease that does not cut target or reference amplicon and is insensitive to methylation is recommended for optimal performance and precision of ddPCR Multiplex Supermix (see guidelines in DNA Digestion section).

Required Equipment

The QX200™ Droplet Digital PCR System (#1864001), QX200 AutoDG™ Droplet Digital PCR System (#1864100), QX600™ Droplet Digital PCR System (#17007769), QX600 AutoDG Droplet Digital PCR System (#17008371), or QX ONE™ Droplet Digital PCR System (#12006536) is required.

Reaction Setup

- Bring the supermix and samples to room temperature. Mix vigorously by vortexing the tube(s) at **maximum speed (3,200–3,500 rpm) for 15 sec**. It is extremely important to mix the 4x concentrated supermix. This process will ensure its homogeneity and help avoid the formation of a concentration gradient during –20°C storage.

- Centrifuge briefly to collect contents at the bottom of the tube(s).
- Prepare DNA samples at the desired concentration before setting up the reaction mix.
- Prepare reaction mix for the number of reactions needed according to the guidelines in Table 1.

Table 1. Example of multiplex copy number variation assay.

Component	QX200	QX ONE	QX600	Final Concentration
	ddPCR System	ddPCR System	ddPCR System	
	Volume per Reaction, µl	Volume per Reaction, µl	Volume per Reaction, µl	
ddPCR Multiplex Supermix	5	5	5	1x
20x target 1 primers/probe (FAM)	1	1	1	900 nM/250 nM
20x target 2 primers/probe (HEX)	1	1	1	900 nM/250 nM
20x target 3 primers/probe (Cy5)	N/A	1	1	900 nM/250 nM or N/A
20x reference 4 primers/probe (Cy5.5)	N/A	1	1	900 nM/250 nM or N/A
20x target 5 primers/probe (ROX)	N/A	N/A	1	900 nM/250 nM or N/A
20x target 6 primers/probe (ATTO 590)	N/A	N/A	1	900 nM/250 nM or N/A
Diluted restriction enzyme (see DNA Digestion section)	1	1	1	2–5 units
300 mM DTT	0.27 (optional)	0.27 (required)	0.27 (highly recommended)	4 mM
Sample	Variable	Variable	Variable	Up to 330 ng
RNase-/DNase-free water	Variable	Variable	Variable	—
Total volume*	20	20	20	—

* For the Automated Droplet Generator, prepare 22 µl per reaction.

Note: ddPCR Multiplex Supermix is recommended for any assays that use Cy5, Cy5.5, ROX, or ATTO 590 channels.

- To complete the reaction setup after adding DNA sample, it is extremely important to mix vigorously by vortexing each reaction tube at **maximum speed (3,200–3,500 rpm) for 10 sec**.
- Centrifuge briefly to ensure that all components are at the bottom of the reaction tubes.
- Allow reaction tubes to equilibrate at room temperature for about 3 minutes.

8. Transfer the reaction mix from the reaction tubes to the appropriate ddPCR Cartridge as follows:

- For the QX600 or QX200 Droplet Digital PCR System, load 20 µl of each reaction mix into a sample well of a DG8 Cartridge. Follow subsequent instructions as specified in the QX200 Droplet Generator Instruction Manual (10031907)
- For the QX600 or QX200 AutoDG Droplet Digital PCR System, follow instructions in the Automated Droplet Generator Instruction Manual (10043138)
- For the QX ONE Droplet Digital PCR System, load 20 µl of each reaction mix into a sample well of a GCR96 Cartridge. Follow subsequent instructions as specified in the QX ONE Droplet Digital PCR System and QX ONE Software User Guide (10000116512)

DNA Digestion (recommended)

DNA digestion prior to droplet generation enables optimal accuracy by separating tandem gene copies, reducing sample viscosity, and improving template accessibility for input samples >66 ng per well. Two strategies may be used to perform restriction digestion of DNA samples: digestion directly in the ddPCR reaction during setup or conventional digestion prior to Droplet Digital PCR.

Digestion in ddPCR Reaction

- Dilute the restriction enzyme using the recommended diluent buffer according to the manufacturer's instructions and add 1 µl to the ddPCR reaction as outlined in Table 1
- Approximately 2–5 units of restriction enzyme per 20 µl ddPCR reaction are recommended
- The addition of restriction enzyme buffers with high salt can inhibit Droplet Digital PCR and should be avoided
- HaeIII, MseI, AluI, HindIII, and CviQI have been observed to work well for digestions in ddPCR reactions

Digestion Prior to Droplet Digital PCR

- Use 5–10 enzyme units per microgram DNA and 10–20 enzyme units per microgram genomic DNA
- Incubate the reaction for 1 hr at the temperature recommended for the restriction enzyme
- Heat inactivation is not required, but can be considered if long-term storage is required; do not heat inactivate above 65°C
- DNA purification is not necessary after restriction digestion
- Use a minimum 10-fold dilution of the digest to reduce the salt content of the sample in the ddPCR reaction
- Store digested DNA at –20°C or below
- Prepare the ddPCR reaction mix, as outlined in Table 1, without the diluted restriction enzyme

Thermal Cycling Conditions

Follow instructions based on the system in use:

- For the QX600 or QX200 Droplet Digital PCR System, carefully transfer droplets into a clean 96-well plate. Seal the plate using the PX1 PCR Plate Sealer (#1814000) at 180°C for 5 sec. Proceed to thermal cycling (see Table 2)
- For the QX600 or QX200 AutoDG Droplet Digital PCR System, seal the 96-well plate using the PX1 PCR Plate Sealer at 180°C for 5 sec. Proceed to thermal cycling (see Table 2)
- For the QX ONE Droplet Digital PCR System, use appropriate thermal cycling conditions as specified in Table 2 during plate setup. Refer to the QX ONE Droplet Digital PCR System and QX ONE Software User Guide (10000116512) for plate setup instructions

Table 2. Thermal cycling conditions.*

Cycling Step		Temperature, °C	Time	Number of Cycles
Hold (QX ONE ddPCR System only)		25	3 min	1
Enzyme activation		95	10 min	1
Denaturation		94	30 sec	40
Annealing/extension		50–65**	1 min***	40
Enzyme deactivation		98	10 min	1
Hold	QX600 or QX200 ddPCR System	4	30 min	1
	QX ONE ddPCR System	25	1 min	1

* For the PTC Tempo Deepwell Thermal Cycler or C1000 Touch Thermal Cycler with 96-Deep Well Reaction Module, use a heated lid set to 105°C and set the sample volume to 40 µl.

** Annealing/extension temperature depends on application type.

*** Check/adjust ramp rate settings to ~2°C/sec.

Data Acquisition and Analysis

Follow instructions based on the system in use:

- For the QX600 Droplet Digital PCR System and the QX600 AutoDG Droplet Digital PCR System, refer to the QX600 Droplet Reader and QX Manager Software Standard Edition User Guide (10000153877)
- For the QX200 Droplet Digital PCR System and the QX200 AutoDG Droplet Digital PCR System, refer to the QX200 Droplet Reader and QX Manager Software Standard Edition User Guide (10000107223)
- For the QX ONE Droplet Digital PCR System, refer to the QX ONE Droplet Digital PCR System and QX ONE Software User Guide (10000116512) and the QX ONE Software User Guide for Standard Edition (10000116655) or Regulatory Edition (10000116656)

Visit [bio-rad.com/ddPCRMultiplexsmx](https://www.bio-rad.com/ddPCRMultiplexsmx) for more information.

BIO-RAD, AUTODG, DDPCR, DROPLET DIGITAL, QX200, QX600, and QX ONE are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions. All trademarks used herein are the property of their respective owner. © 2023 Bio-Rad Laboratories, Inc.

Purchase of Digital PCR and/or Single-Cell NGS Sample Preparation products (the "Products") from Bio-Rad Laboratories is subject to Bio-Rad Laboratories, Inc. Standard Terms and Conditions of Sale, which can be accessed at <https://www.bio-rad.com/en-us/terms-conditions>. Unless we expressly state otherwise in additional Terms and Conditions, no rights are granted for you to distribute or resell the Products. Unless we expressly state otherwise in additional Terms and Conditions, no rights are granted for the development or commercialization of diagnostic assays for use with the Products without a license from Bio-Rad. It is the user's obligation to obtain a commercial license from Bio-Rad for (i) all commercial uses (not just diagnostic uses) and (ii) sale of assays for use on Bio-Rad's dPCR and ddSEQ instruments. The Products and/or their use are covered by U.S. and foreign patents and/or pending patent applications owned by or under license to Bio-Rad Laboratories, Inc. See <https://www.bio-rad.com/en-us/trademarks>.