


# CFX Duet Real-Time PCR System Quick Start Guide

# **Unpacking and Installing the CFX Duet System**

The CFX Duet System is packed in a single box with the accessories included in the box. The accessories kit includes the power cord, USB cable, shipping screw plug, and an Ethernet cable. Set the box on the benchtop and follow the instructions per the labels affixed to the shipping box. For more detailed instructions, refer to the instrument guide. The CFX Duet System **requires** a connected PC running the most current version of CFX Maestro Software to operate.

## Unpacking the CFX Duet System

1. Remove the straps from the shipping box. DO NOT cut the top of the box open. Go to step 2 to continue unpacking the instrument.

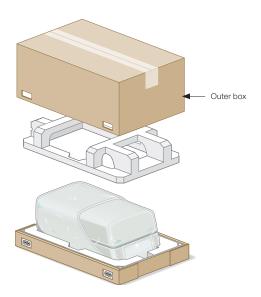

2. Pinch and pull out the clips. After pulling out the four clips, lift up the outer box to expose the instrument.



3. Pull down for tray access.




4. Slide tray to pull out the instrument.





### Removal of the Shipping Screw

- 1. Use the supplied power cord to plug the CFX Duet System into an appropriate electrical outlet.
- 2. Turn on the CFX Duet System using the power switch located on the right-hand side on the back of the instrument.
- 3. Ensure the CFX Duet System is connected via the USB cable to a PC running CFX Maestro Software.
- 4. In CFX Maestro Software, confirm that the CFX Duet System icon appears in the Detected Instruments pane.
- 5. In the Detected Instruments pane of CFX Maestro Software, right click the CFX Duet System icon and select Properties.
- Select the Shipping Screw tab and click Remove Shipping Screw.
- 7. Follow the instructions displayed in the CFX Maestro Software window to remove the shipping screw, which is located on the right-hand side of the instrument.

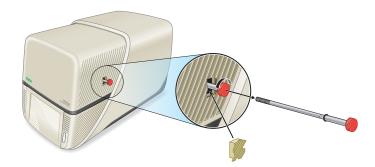

# **Operating the CFX Duet System**

Operating the CFX Duet System requires connection to a PC running CFX Maestro Software. For details, see the user guide for CFX Maestro Software. For details about the CFX Duet System, refer to the instrument guide (10000148048).

Visit bio-rad.com/CFXDuet for more information.

BIO-RAD is a trademark of Bio-Rad Laboratories, Inc.
All trademarks used herein are the property of their respective owner.



Bio-Rad Laboratories, Inc.

Life Science Group Website bio-rad.com USA 1 800 424 6723 Australia 61 2 9914 2800 Austria 00 800 00 24 67 23 Belgium 00 800 00 24 67 23 Brazil 4003 0399 Canada 1 905 364 3435 China 86 21 6169 8500 Czech Republic 00 800 00 24 67 23 Denmark 00 800 00 24 67 23 Finland 00 800 00 24 67 23 France 00 800 00 24 67 23 Gramany 00 800 00 24 67 23 Gramany 00 800 00 24 67 23 Gramany 00 800 00 24 67 23 India 91 124 4029300 Israel 0 3 9636050 Italy 00 800 00 24 67 23 Japan 81 3 6361 7000 Korea 82 2 3473 4460 Luxembourg 00 800 00 24 67 23 Mexico 52 555 488 7670

The Netherlands 00 800 00 24 67 23 New Zealand 64 9 415 2280 Norway 00 800 00 24 67 23 Poland 00 800 00 24 67 23 Portugal 00 800 00 24 67 23 Russian Federation 00 800 00 24 67 23 Singapore 65 6415 3188 South Africa 00 800 00 24 67 23 Spain 00 800 00 24 67 23 Sweden 00 800 00 24 67 23 Switzerland 00 800 00 24 67 23 Taiwan 886 2 2578 7189 Thailand 66 2 651 8311 United Arab Emirates 36 1 459 6150 United Kingdom 00 800 00 24 67 23

10000147782 Ver A (12017192) US/EG 21-0752 1021 Sig 0121

