
QX600™ Droplet Reader and QX Manager Software Standard Edition

User Guide



BIO-RAD

QX600 Droplet Reader and QX Manager Software Standard Edition

User Guide

Version 2.2

For Research Use Only

Bio-Rad Technical Support Department

The Bio-Rad Technical Support department in the U.S. is open Monday through Friday, 5:00 AM to 5:00 PM, Pacific time.

Phone: 1-800-424-6723, option 2

Email: Support@bio-rad.com (U.S./Canada Only)

For technical assistance outside the U.S. and Canada, contact your local technical support office or click the Contact Us link at bio-rad.com.

Legal Notices

No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopy, recording, or any information storage or retrieval system, without permission in writing from Bio-Rad Laboratories, Inc.

Bio-Rad reserves the right to modify its products and services at any time. This guide is subject to change without notice. Although prepared to ensure accuracy, Bio-Rad assumes no liability for errors or omissions, or for any damage resulting from the application or use of this information.

AUTODG, BIO-RAD, DDPCR, DROPLET DIGITAL, DROPLET DIGITAL PCR, QUANTASOFT, QX200, and QX600 are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions.

EvaGreen is a trademark of Biotium, Inc. Bio-Rad Laboratories, Inc. is licensed by Biotium, Inc. to sell reagents containing EvaGreen Dye for use in real-time PCR, for research purposes only.

Purchase of Digital PCR and/or Single-Cell NGS Sample Preparation products (the "Products") from Bio-Rad Laboratories is subject to Bio-Rad Laboratories, Inc. Standard Terms and Conditions of Sale, which can be accessed at <https://bio-rad.com/en-us/terms-conditions>. Unless we expressly state otherwise in additional Terms and Conditions, no rights are granted for you to distribute or resell the Products. Unless we expressly state otherwise in additional Terms and Conditions, no rights are granted for the development or commercialization of diagnostic assays for use with the Products without a license from Bio-Rad. It is the user's obligation to obtain a commercial license from Bio-Rad for (i) all commercial uses (not just diagnostic uses) and (ii) sale of assays for use on Bio-Rad's ddPCR and ddSEQ instruments. The Products and/or their use are covered by U.S. and foreign patents and/or pending patent applications owned by or under license to Bio-Rad Laboratories, Inc. See <https://bio-rad.com/en-us/trademarks> for trademark information.

All trademarks used herein are the property of their respective owner.

Copyright © 2023 by Bio-Rad Laboratories, Inc. All rights reserved.

Revision History

Document	Date	Description of Change
QX600 Droplet Reader and QX Manager Software Standard Edition User Guide Software Version 2.2 DIR No. 10000153877 Ver D	December 2023	Update content for Version 2.2
Software Version 2.1	September 2023	Add EcoTank functionality to Instrument Maintenance section
QX600 Droplet Reader and QX Manager Software Standard Edition User Guide Software version 2.1 DIR No. 10000153877 Ver C	June 2023	Update DIR version
QX600 Droplet Reader and QX Manager Software Standard Edition User Guide Software version 2.1 DIR No. 10000153877 Ver B	May 2023	Update the user guide with new and improved features for Version 2.1; add new legal statement and Managing Waste section
QX600 Droplet Reader and QX Manager Software Standard Edition User Guide Software version 2.0 DIR No. 10000153877 Ver A	August 2022	Create new document for QX Manager with QX600 instrument functionality

This page intentionally left blank.

Table of Contents

Front Cover	a
Revision History	iii
Safety and Regulatory Compliance	11
Regulatory Compliance	11
Safety Warning Labels	13
Safe Use Specifications	14
Personal Protective Equipment	15
Hazards	16
Biohazards	16
Chemical Hazards	17
Explosive or Flammability Hazards	17
Electrical Hazards	17
Surface Decontamination	18
Decommissioning and Disposal	18
Transport	19
Warranty	19
Chapter 1 Introduction to Droplet Digital PCR	21
ddPCR Workflow	22
Finding Out More	22
Chapter 2 About the QX600 Droplet Reader	23
Instrument Specifications	25
Environmental Requirements	26
Included Components	27
Additional Components	28
Compatible ddPCR Process Instruments	30
Connecting and Starting the Instrument	31
Status Indicator Lights	32

Chapter 3 About QX Manager Software	33
Enabling User Management	35
Chapter 4 Getting Started	37
Signing In, Viewing Privileges, and Changing Preferences	37
Signing into the Software	38
Selecting the Instrument	40
Viewing User Privileges	41
Managing Your Preferences	43
Partial Preferences	44
Signing Out or Changing Users	45
Instrument Status Bar	46
Tabs to Functional Windows	47
Compatible File Types	48
Chapter 5 Running Experiments	49
Assembling a Plate	49
Inserting an Assembled Plate into the Instrument	50
Adding the Plate in QX Manager Software	51
Configuring the Plate	52
Plate Information Tab	52
Well Selection Tab	54
Well Information Tab	55
Starting the Run	57
Run Prerequisites Checklist	58
Running the Experiment	61
Tracking the Run Status	62
Using Live Analysis	64
Canceling a Run	65
Chapter 6 Creating or Editing Templates	67
Creating or Modifying Plate Templates	68
Template Plate Editor Window	70
Creating and Storing Plate Templates from CSV Files	79
Creating or Modifying Report Templates	81

Chapter 7 Data Analysis Module Overview	85
Opening a Data File for Analysis	87
Saving a Data File	89
Analysis Display and Output Options	90
Universal Options	91
Expanding or Reducing the Chart	94
Selecting Chart Display Preferences	95
Selecting Chart Scale and Label Options	96
Using Chart Menu Options	98
Using Table Menu Options	100
Analysis Dashboard	103
Modifying the Dashboard Display	104
Dashboard Options	105
Plate View Windows	107
Plate Editor — Analysis View	107
Plate View	108
2D Plate View	109
Data Windows	110
Data Table Tab	110
Plate Editor Well Data Table	117
Analysis Windows	118
Viewing Run Information	119
Chapter 8 Data Analysis Methodology	121
Amplitude Plot Analysis Options	122
1D Amplitude	127
2D Amplitude	132
Data Visualization Chart Analysis Options	135
Concentration	138
Copy Number	141
Ratio	144
Fractional Abundance	146
Event Counts	148
Linkage Analysis	149

Chapter 9 Analysis Module Reporting	151
Creating Analysis Reports	151
Chapter 10 Gene Study Module	153
Gene Study Options	155
Setting up the Gene Study	156
Adding Data Files	158
Removing Data Files	158
Selecting Data Files for Analysis	158
Showing the Plate View	159
Using the Study Analysis Features	160
Gene Study Reports	162
Creating a Gene Study Report	163
Chapter 11 System Utilities	165
System Settings Tab	166
Preferred Location	167
Shared Settings	168
Instrument Settings	169
Software Settings	170
System Log Files	171
Event Log	171
Maintenance Log	173
Maintenance Reports	174
System Data	175
Archiving Raw Data	175
Troubleshooting Data	177
Recovering Data	179
Reprocessing Data	181
System Calibration and Tools	182
Chapter 12 Managing Users	183
Adding Users	184
Adding or Removing User Privileges	186
Modifying User Preferences	188
Removing Users	188

Appendix A Experiment Examples	189
Direct Quantification Multiplexed	190
Setting Up Experiment Parameters	190
Viewing and Adjusting Results	191
Copy Number Variation Triplexed	193
Setting Up Experiment Parameters	193
Viewing and Adjusting Results	195
Drop-Off Assays for Genome Edit Detection	196
Setting Up Experiment Parameters	197
Viewing and Adjusting Results	198
Direct Quantification with the Advanced Classification Method	199
Setting Up Experiment Parameters	199
Viewing and Adjusting Results	201
Appendix B Instrument Maintenance	203
Fluid Maintenance and Waste Management	204
Changing the Droplet Reader Oil Bottle	205
Emptying the Waste Container	206
Using an EcoTank Oil and Waste Container	207
Priming and Flushing the Instrument	213
Managing Waste	214
Recovery From Idle Mode	215
Cleaning or Decontaminating the Instrument	215
Instrument Maintenance Log	216
Appendix C Troubleshooting	217
Event Log	217
Instrument Errors	217
Appendix D Ordering Information	219
Back Cover	227

Table of Contents

This page intentionally left blank.

Safety and Regulatory Compliance

This section cites regulatory requirements for laboratory and electrical equipment, as well as requirements for working with chemicals and hazardous substances, and also explains safety precautions and recommendations.

Important: Only trained personnel should use this instrument.




Regulatory Compliance

The instrument has been tested and found to be in compliance with all applicable requirements of the following safety and electromagnetic standards:

- IEC 61010-1:2010 (3rd:2016), EN61010-1:2010 (3rd ed.). Electrical equipment for measurement, control, and laboratory use - Part 1: General requirements
- IEC 61326-1:202012 (Class A), EN 61326-1:2013 (Class A). Electrical equipment for measurement, control, and laboratory use. EMC requirements, Part 1: General requirements
- IEC 61010-2-081:20195, 3.0 edition, UL 61010-2-081:2015, CAN/CSA C22.2 No. 61010-2-081:19. Safety requirements for electrical equipment for measurement, control, and laboratory use. Part 2-081: Particular requirements for automatic and semi-automatic laboratory equipment for analysis and other purposes (includes Amendment 1)
- CAN/CSA 22.2 No 61010-1-04, Safety requirements for electrical, equipment for measurement, control, and laboratory use, Part 1: General requirements
- CAN/CSA C22.2 No. 61010-2-101:2019 IEC 61010-2-101: 2018 (3rd ed.)
- EN 61326-1:2006 20201(Class A). Electrical equipment for measurement, control, and laboratory use. EMC requirements, Part 1: General requirements
- UL 61010-1:2004, Laboratory equipment, Test & Measurement Equipment and Industrial Process Controls
- UL 61010-2-101:2019. Safety requirements for electrical equipment for measurement, control, and laboratory use. Particular requirements for in vitro diagnostic (IVD) medical equipment
- Restriction of hazardous substances (ROHS) directive (European Union)

- Registration, evaluation, authorization and restriction of chemicals (REACH). European Chemicals Agency (ECHA) June 1, 2007
- Waste electrical and electronic equipment (WEEE) directive 2012/19/EU




Table 1. Regulatory symbols

Symbol	Definition
	The CE symbol indicates that the manufacturer ensures the product conforms with the essential requirements of the applicable EN directives.
	The CSA symbol indicates that a project has been tested to Canadian and U.S. standards, and it meets the requirements of those applicable standards. This equipment has been tested and found to comply with the limits for a Class A digital device pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment.
 2012/19/EU www.bio-rad.com	The Waste Electrical and Electronic Equipment (WEEE) Directive symbol indicates that when the end-user wishes to discard this product, it must be sent to separate collection facilities for recovery and recycling.

Safety Warning Labels

Warning labels alert you about sources of injury or harm. [Table 2](#) defines each safety warning label.

Table 2. Meaning of safety warning labels

Icon	Meaning
	<p>Warning about risk of harm to body or equipment</p> <p>Operating the system before reading this manual can constitute a personal injury hazard. For safe use, do not operate this instrument in any manner unspecified in this manual. Only qualified laboratory personnel trained in the safe use of electrical equipment should operate this instrument. Always handle all components of the system with care and with clean, dry hands.</p>
	<p>Warning about handling biohazardous materials</p> <p>When handling biohazardous samples, adhere to the recommended precautions and guidelines and comply with any local guidelines specific to your laboratory and location.</p>
	<p>Warning about risk of electric shock</p> <p>In order to prevent electric shock, use caution when plugging and unplugging the instrument. Always turn off and unplug the instrument when performing maintenance procedures.</p>

Safe Use Specifications

For safe operation of the instrument, Bio-Rad™ strongly recommends that you comply with instructions listed in this section and in [Appendix B, Instrument Maintenance](#).

Important: This instrument is intended for laboratory use only. Bio-Rad is not responsible for any injury or damage caused by use of this instrument for purposes other than those for which it is intended, or by modifications to the instrument not performed by Bio-Rad or an authorized agent.

- This instrument is for use only by trained personnel.
- Use only the power and connection accessories supplied with the instrument, and the plug adapter corresponding to the electrical outlets in your region.
- For adequate ventilation and easy access to the back of the instrument, place the instrument on a solid, stable surface and provide adequate room at the back and on each side. For information, see [Environmental Requirements on page 26](#).
- This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the provided instructional documentation, may cause harmful interference to radio communications. Operation of the systems in a residential area is likely to cause harmful interference, in which case users will be required to correct the interference at their own expense.

Note: Bio-Rad recommends maintaining a backup power source in case of power outages. A universal power supply (UPS) can protect from brown outs and power surges, while a regular backup generator does not.

For information on environmental requirements, see [Environmental Requirements on page 26](#).

Personal Protective Equipment

Proper use of gloves is recommended with use of oils and sample plates. OSHA requirements for personal protective equipment (PPE) are set forth in the Code of Federal Regulations (CFR) at 29 CFR 1910.132 (General requirements); 29 CFR 1910.138 (Hand protection); 29 CFR 1926.95 (Criteria for standard personal protective equipment). Any gloves with impaired protective ability should be discarded and replaced. Consider the toxicity of the chemicals and factors such as duration of exposure, storage, and temperature when deciding to reuse chemically exposed gloves.

Listed below are features to aid glove selection for handling of machines, assays, oils, and cleaning solvents:

- Butyl gloves are made of a synthetic rubber and protect against peroxide, hydrofluoric acid, strong bases, alcohols, aldehydes, and ketones
- Natural (latex) rubber gloves are comfortable to wear and feature outstanding tensile strength, elasticity, and temperature resistance
- Neoprene gloves are made of synthetic rubber and offer good pliability, finger dexterity, high density, and tear resistance; they protect against alcohols, organic acids, and alkalis
- Nitrile gloves are made of copolymer and provide protection from chlorinated solvents such as trichloroethylene and tetrachloroethene; they offer protection when working with oils, greases, acids, and caustic substances

Hazards

The instrument is designed to operate safely when used in the manner prescribed by the manufacturer. If the instrument or any of its associated components is used in a manner not specified by the manufacturer, the inherent protection provided by the instrument may be impaired.

Bio-Rad Laboratories, Inc. is not liable for any injury or damage caused by the use of this equipment in any unspecified manner, or by modifications to the instrument not performed by Bio-Rad or an authorized agent. Only trained Bio-Rad personnel should perform service on the system.

Biohazards

The instrument is a laboratory product. However, if biohazardous samples are present, adhere to the following guidelines and comply with any local guidelines specific to your laboratory and location.

Note: No biohazardous substances are expended during normal operations of this instrument.

General Precautions

- Always wear laboratory coat, laboratory gloves, and safety glasses with side shields or goggles.
- Keep your hands away from your mouth, nose, and eyes.
- Completely protect any cut or abrasion before working with potentially infectious materials.
- Wash your hands thoroughly with soap and water after working with any potentially infectious material before leaving the laboratory.
- Store all infectious or potentially infectious material in unbreakable leak-proof containers.
- Before leaving the laboratory, remove protective clothing.
- Change gloves frequently. Do not use a gloved hand to write, answer the telephone, turn on a light switch, or touch anything that other people may touch without gloves. Remove gloves immediately when they are visibly contaminated.
- Do not expose materials that cannot be properly decontaminated to potentially infectious material.
- Upon completion of an operation involving biohazardous material, decontaminate the work area with an appropriate disinfectant (for example, a 1:10 dilution of household bleach).

Disposal of Biohazardous Material

Dispose of the following potentially contaminated materials in accordance with laboratory local, regional, and national regulations:

- Clinical samples
- Reagents
- Used reaction vessels or other consumables that may be contaminated

Chemical Hazards

The instrument contains no potentially hazardous chemical materials.

Explosive or Flammability Hazards

The instrument poses no uncommon hazard related to flammability or explosion when used in a proper manner as specified by Bio-Rad.

Electrical Hazards

The instrument poses no uncommon electrical hazard to operators if installed and operated properly without physical modification and connected to a power source of proper specification.

Surface Decontamination



WARNING! To prevent electrical shock, always turn off and unplug the instrument before performing decontamination procedures.

Important: Do not use abrasive or corrosive detergents or strong alkaline solutions. These agents can scratch surfaces and damage the system.

The following areas can be cleaned with 10% bleach solution:

- Outer area and touch screen
- Inner plate holders
- Droplet reading surfaces

To prepare and apply the disinfectant, refer to the instructions provided by the product manufacturer. For questions regarding the use of other cleaning agents, contact Bio-Rad Technical Support.

For more information on surface cleaning, see [Appendix B, Instrument Maintenance](#).

Decommissioning and Disposal

The instrument contains electrical materials that should be disposed of as unsorted waste and must be collected separately, according to European Union Directive 2012/19/EU on waste electrical and electronic equipment — WEEE Directive. The purpose of decommissioning is to make sure that the equipment is electrically and environmentally safe for disposal. Before disposal, contact your local Bio-Rad representative for country-specific instructions.

Transport

You must perform the specified decontamination procedures before moving or shipping the system. Always move or ship the instrument with the supplied packaging materials, which will protect the instrument from damage. If appropriate containers cannot be found, contact your local Bio-Rad office.

Warranty

Important: The instrument is for research use only, and not for use in diagnostic procedures.

The instrument and its associated accessories are covered by a standard Bio-Rad warranty. Contact your local Bio-Rad office for the details of the warranty. Follow the safety specifications listed in this chapter and throughout this guide.

The instrument is intended for laboratory use only. Bio-Rad Laboratories, Inc. is not responsible for any injury or damage caused by use of this instrument for purposes other than those for which it is intended, or by modifications of the instrument not performed by Bio-Rad Laboratories or an authorized agent. Alteration of this instrument voids the warranty and safety certification, as it creates a potential safety hazard.

Use of unapproved supermixes or additives may harm the instrument and voids the warranty.

Use only the power cord supplied with the instrument, using only the plug adaptor that corresponds to the electrical outlets in your region.

This page intentionally left blank.

Chapter 1 Introduction to Droplet Digital PCR

Droplet Digital™ PCR (ddPCR™) is a digital polymerase chain reaction method based on water-oil emulsion droplet technology, which uses a combination of microfluidics and proprietary surfactant chemistries to divide each sample into water-in-oil droplets.

The technology uses reagents and workflows similar to those used for most standard TaqMan probe-based assays, and provides absolute quantification of nucleic acid target sequences by counting nucleic acid molecules encapsulated in discrete, volumetrically defined water-in-oil droplet partitions.

ddPCR is highly effective in the following areas:

- **Absolute quantification** — ddPCR provides a concentration of target DNA copies per input sample without the need for running standard curves, making this technique ideal for target DNA measurements, viral load analysis, and microbial quantification.
- **Genomic alterations such as gene copy number variation (CNV)** — CNVs result in too few or too many dosage-sensitive genes responsible for phenotypic variability, complex behavioral traits, and disease. ddPCR enables measurement of 1.2x differences in gene copy number.
- **Detection of rare sequences** — researchers must amplify single genes in a complex sample, such as a few tumor cells in a wild-type background. ddPCR is sensitive enough to detect rare mutations or sequences.
- **Gene expression and microRNA analysis** — ddPCR provides stand-alone absolute quantification of expression levels, especially low-abundance microRNAs, with sensitivity and precision.
- **Next-generation sequencing (NGS)** — ddPCR quantifies NGS sample library preparations to increase sequencing accuracy and reduce run repeats. Validate sequencing results such as single nucleotide polymorphisms or copy number variations with absolute quantification.
- **Single cell analysis** — the high degree (10-fold to 100-fold) of cell-cell variation in gene expression and genomic content among homogeneous post-mitotic, progenitor, and stem cell populations drives a need for analysis from single cells. ddPCR enables low copy number quantification and gene expression of individual cells.
- **Genome edit detection** — ddPCR enables fast, precise, and cost-effective assessment of HDR (Homology directed repair) and NHEJ (non-homologous end joining) generated by CRISPR-Cas9 or other genome editing tools.

ddPCR has the following benefits for nucleic acid quantification:

- **Unparalleled precision** — The massive sample partitioning afforded by ddPCR enables small fold differences in target DNA sequence between samples to be reliably measured.
- **Increased signal-to-noise ratio** — High-copy templates and background are diluted, effectively enriching template concentration in target-positive partitions. This allows for the sensitive detection of rare targets and enables a $\pm 10\%$ precision in quantification.
- **Removal of PCR efficiency bias** — Error rates are reduced by removing the amplification efficiency reliance of qPCR, enabling accurate quantification of targets to near zero.
- **Simplified quantification** — There is no requirement for a standard curve for absolute quantification.


ddPCR Workflow

The ddPCR process adheres to the following workflow:

- Prepare samples for PCR by combining DNA or RNA with primers, probes dye, and Bio-Rad ddPCR supermix.
- A droplet generator fractionates a sample into thousands of nanoliter-sized droplets with target and background DNA distributed randomly into the droplets during the partitioning process.
- Following droplet generation, the droplets are run through a thermal cycler, which performs PCR amplification of the nucleic acid target in each individual droplet.
- A droplet reader reads each droplet to determine the fraction of positive droplets in the original sample and uses Poisson statistical formulas to determine the absolute quantity.

Note: Positive droplets containing at least one copy of the target DNA molecule exhibit increased fluorescence compared to negative droplets.

Finding Out More

Click the Help tab  on the left, and then click the Bio-Rad Website link to access links to technical notes, manuals, videos, product information, and technical support. The website also provides many technical resources on a wide variety of methods and applications related to PCR, Droplet Digital PCR, and gene expression.

Chapter 2 About the QX600 Droplet Reader

Using the QX600™ Droplet Reader, you can detect rare DNA target copies, determine copy number variation, and precisely measure gene expression levels. Applications and uses include:

- Copy number variation
- Rare sequence detection
- Next generation sequencing (NGS) library quantification
- Viral load determination, including SARS-CoV-2 assays
- mRNA analysis
- Gene expression analysis and single cell gene expression analysis
- Absolute quantification
- Rare mutant detection
- NGS sample preparation
- GMO detection
- Contaminant testing
- High-order multiplexing

Following the droplet generation and thermal cycling processes, the QX600 Droplet Reader collects binary data for analysis as it singulates the nanoliter-sized droplets in each sample, and then streams them in single-file past a six-color detector to determine which droplets contain target molecules (positive reading) and which do not (negative reading).

The QX600 Droplet Reader detects multiple targets in six fluorescent dyes using three interrogation points across defined spectrums. The first interrogation point detects targets in FAM or EvaGreen® and HEX or VIC, the second in Cy5 and Cy5.5, and the third in ROX and ATTO 590.

Important: Use only the dyes listed above (FAM or EvaGreen®, HEX/VIC, Cy5, Cy5.5, ROX, and ATTO 590).

Important: Your QX600 Droplet Reader performs the post-PCR droplet reading function only. A droplet generator and a thermal cycler are required to perform a full ddPCR cycle. You can purchase a droplet generator and thermal cycler, as well as associated accessories and consumables, from Bio-Rad. For information, see [Appendix D, Ordering Information](#).

The following graphic and legend illustrate the functional areas on the instrument exterior:



LEGEND

- 1** Button to open and close lid

- 2** Status indicator lights (see [Status Indicator Lights on page 32](#))

- 3** On the back of the instrument, power and USB outlets

- 4** Oil and waste compartment (see [Instrument Maintenance on page 203](#))

Instrument Specifications

This section contains information on the instrument size and weight of the instrument. Safety information is cited in [Safe Use Specifications on page 14](#) and the proper environmental maintenance is cited in [Environmental Requirements on page 26](#).

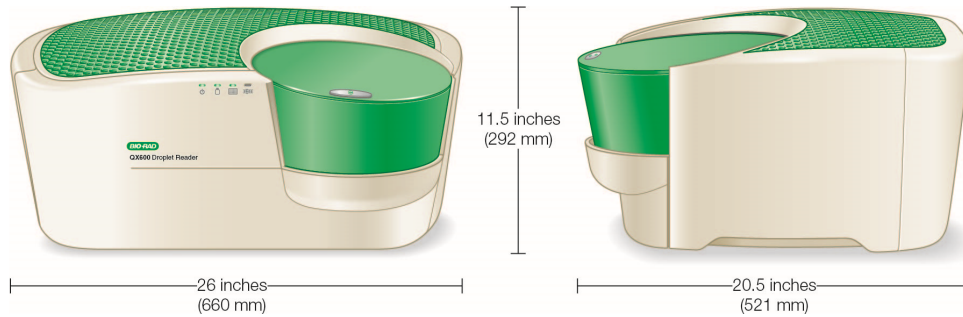


Table 3. Instrument specifications

Element	Specifications
Weight	56.6 lbs (26 kg)
Size (W x D x H)	26 x 20.5 x 11.5" (660 x 521 x 292 cm)

Environmental Requirements

Table 4 lists the environmental requirements for the QX600 Droplet Reader.

Important: You must use the supplied shielded cables with your instrument to ensure compliance with the Class A FCC limits.

Table 4. Conditions for safe use

Usage aspect	Requirements
Rated input power	100–240 V, 50–60 Hz, 2A (plugs into standard AC receptacle)
Fuses	10A 250 V SLOW BLOW
Voltage fluctuations	± 10% for the included external power supply Note: Use only the power cord supplied with the equipment.
Pollution degree	2
Usage temperature	18–30° C
Relative humidity	85% max non-condensing
Altitude	0 to 8,200 ft (0 to 2,500 m)
Installation category	II (external power supply plugs into a standard AC receptacle) Indoor use only
Ventilation and access clear space requirements	The following distances should be unobstructed for proper ventilation and service access <ul style="list-style-type: none"> ■ 5 in (12 cm) on the left and right sides of the instrument ■ 10 in (25 cm) behind the instrument

Included Components

This section identifies the functional items included with your droplet reader purchase, as well as additional components and consumables. For information on items you may need for a full ddPCR cycle, see [Appendix D, Ordering Information](#).

[Table 5](#) explains the components that ship with the QX600 Droplet Reader.

Table 5. QX600 Droplet Reader and included components

Component	Description	Catalog Number
QX600 Droplet Reader	Droplet reading instrument	12013328
USB cable and power cord	USB cable that connects the computer to the instrument and power cord that connects the instrument to the power source Note: To obtain cables or power cords, contact Bio-Rad technical support.	Included
Droplet reader plate holders (2)	Positions and secures the 96-well plate in the droplet reader plate compartment	12006834

Additional Components

The following tables identify additional components you can purchase for use with your system or instrument.

Table 6. Accessories and consumables

Component	Description	Catalog Number
Computer	Computer that connects to the QX600 Droplet Reader for data collection and analysis	12017458
Software purchase	Bio-RadQX Manager Software, Premium Edition Standard Edition is included with your droplet reader purchase.	12012172
ddPCR 96-Well Plates	Clear well/clear shell semi-skirted plates	12001925 (pkg of 25)
Droplet Reader Oil	ddPCR droplet reader oil for QX600 and QX200 instruments and systems	1863004
Rainin Pipets	20 µl pipet for sample loading 50 µl pipet for droplet transfer 8-channel 200 µl pipet for oil	Rainin L-20 Rainin L-50, L8-50 Rainin L8-200
Rainin Pipet Tips	Filtered	Rainin GP-L10F Rainin GP-L200F

Table 7. Thermal cyclers and plate sealer

Product	Description	Catalog Number
C1000 Touch Thermal Cycler with 96-Deepwell Reaction Module	Includes the C1000 Touch Thermal Cycler chassis, 96-deepwell reaction module, USB flash drive	1851197
PTC Tempo Deepwell Thermal Cycler	Includes the thermal cycler, power cord, USB cable, and Ethernet cable	12015392
PX1 PCR Plate Sealer	PCR plate sealer, includes heat sealing instrument, plate support block that holds 96-well and 384-well plates, sealing frame, power cord	1814000
Pierceable Foil Heat Seal	Heat seal for the PX1 PCR Plate Sealer	1814040

Table 8. ddPCR supermixes

Supermix	Description	Catalog number	
ddPCR EvaGreen® Supermix	2x supermix, for use in nucleic acid sample preparation with the QX600 and QX200 ddPCR systems	1864033	2ml (2 x 1ml)
		1864034	5ml (5 x 1ml)
		1864035	25ml (25 x 1ml)
		1864036	50ml (50 x 1ml)
ddPCR Supermix for Probes (no dUTP)	2x supermix, for use in nucleic acid sample preparation with the QX600 and QX200 ddPCR systems	1863023	2ml (2 x 1ml)
		1863024	5ml (5 x 1ml)
		1863025	25ml (25 x 1)
ddPCR Supermix for Probes	2x supermix, for use in sample preparation with the QX600 and QX200 ddPCR systems	1863026	2ml (2 x 1ml)
		1863010	5ml (5 x 1ml)
		1863027	25ml (25 x 1ml)
		1863028	50ml (50 x 1ml)
ddPCR Multiplex Supermix	4x supermix especially suited for probe-based detection of multiple targets in DNA samples using the QX600 and QX200 ddPCR systems	12005909	1.2ml (2 x 0.6ml)
		12005910	3 ml (5 x 0.6 ml)
		12005911	12.5 ml (5 x 2.5 ml),

Table 8. ddPCR supermixes, continued

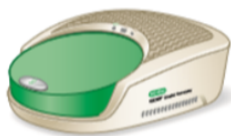
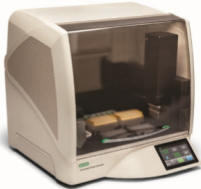
Supermix	Description	Catalog number	
ddPCR Supermix for Residual DNA Quantification	2x supermix, for use in residual DNA detection with the QX600 and QX200 ddPCR systems	1864037	2ml (2 x 1ml)
		1864038	5ml (5 x 1ml)
		1864039	25ml (25 x 1ml)
		1864040	50ml (50 x 1ml)
One-step RT-ddPCR Advanced Kit for Probes	200 or 500x 20 µl reaction kit, for absolute quantification of target RNA in a one-step format with the QX600 and QX200 ddPCR systems	1864021	2ml (2 x 1ml)
		1864022	5ml (5 x 1ml)

Table 9. Buffer controls

Buffer	Catalog Number
ddPCR Buffer Control Kit for Probes	1863052
QX200 Buffer Control Kit for EvaGreen®	1864052

Compatible ddPCR Process Instruments

Along with the QX600 Droplet Reader, you can purchase the droplet generators shown below from Bio-Rad.

QX200 Droplet Generator**Automated Droplet Generator (left)****QX600 Droplet Reader**

Important: A thermal cycler, which can be purchased separately, is required to perform a full ddPCR cycle. For information on additional instruments, accessories, and consumables, see [Appendix D, Ordering Information](#).

Connecting and Starting the Instrument

To execute runs on the QX600 Droplet Reader, you must connect the instrument to a computer on which the QX Manager Software is installed and configured. Bio-Rad recommends purchasing at least one configured computer that can remain connected to the QX600 Droplet Reader and be used for adding and running plates in QX Manager Software. For information on additional items you can purchase for your instrument, see [Additional Components on page 28](#).

To connect and start the instrument

1. Connect the USB cable provided by Bio-Rad to the USB port on the back of the QX600 Droplet Reader, and then to the computer.
2. Press the Power button on the instrument.

The Power indicator light turns solid green to indicate that power is on.

3. Turn on the connected computer and log in with a Windows account. For information, see your system administrator.
4. Open and log into QX Manager Software.

Important: The first time you log into the software, the Instrument window appears and you must select the instrument you are using. For information on selecting the correct instrument in QX Manager Software, see [Selecting the Instrument on page 40](#).













Status Indicator Lights

This section describes the indicator lights on the front of the QX600 Droplet Reader.



Before you start a run, check the indicator lights to ensure the instrument is ready. For information, see [Table 10](#).

Table 10. Droplet Reader Status Indicator Lights

Indicator	 (Power)	 (Fluids)	 (Plate)	 (Run Status)
Solid Green 	Power is on	Fluid levels OK	Plate loaded properly	Complete
Flashing Green 	N/A	<ul style="list-style-type: none">  Less than 40% oil is available  More than 70% waste in container 	N/A	In progress
Flashing Amber 	N/A	<ul style="list-style-type: none">  Less than 10% oil available  More than 90% waste in container 	N/A	Error during run
No color 	Power is off	N/A	No plate loaded	Instrument is idle

If the bottle indicator light is solid green or flashing green, there is enough oil in the bottle to complete a run on 96 wells. You can configure the full plate in QX Manager Software for your run. Once the low level sensor is triggered, the Add Plate button is disabled and you cannot continue to run plates (less than 10%). You must replace the oil and waste bottles with new ones before the instrument can start the run.

Chapter 3 About QX Manager Software


QX Manager Software, when connected to your QX600 Droplet Reader, provides the necessary functionality to create, run, and analyze droplet reading experiments on your samples.

During the installation of Standard Edition, you can enable the User Management module for all user management functions and preferences. If you do not enable User Management, you can use the software with a single generic user.

Using the instrument and software, you can

- Set up customized, real-time ddPCR experiments for plate runs
- Read standard droplets
- Choose between two droplet volumes
- See the available disk space and number of plates you can process
- Use the live analysis function during the droplet reading process
- Analyze your data files in a variety of charts and tables in the Analysis module
- Reprocess data with updated color calibration values
- Create and store plate and report templates, and produce reports on your data

The following image highlights functional areas.

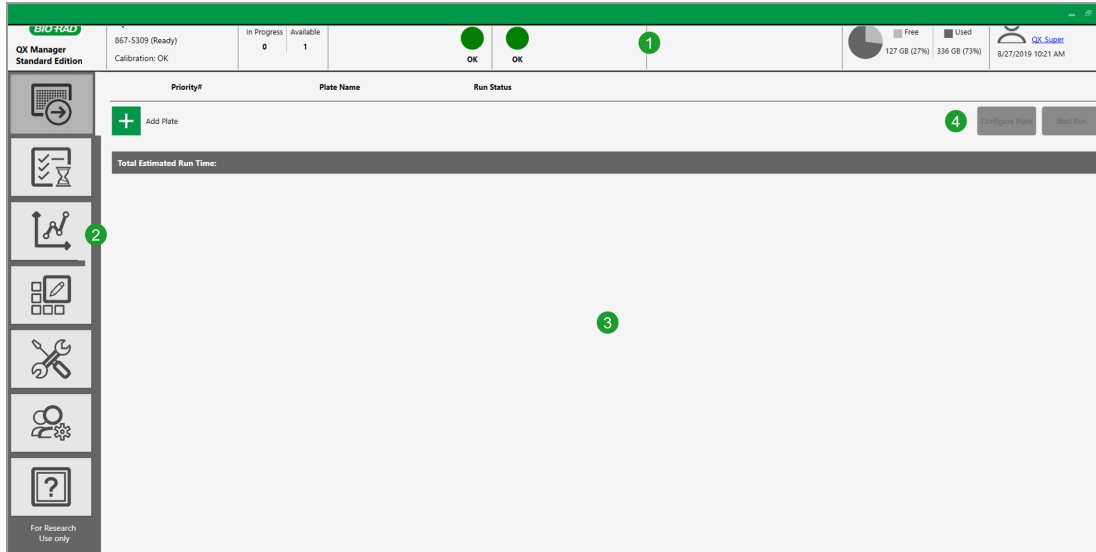

LEGEND

- 1 The status bar displays information about the instrument and user. For information on specific functionality, see [Instrument Status Bar on page 46](#).
- 2 Tabs provide access to the main functional windows.
- 3 The main pane displays the details of the selected tab.
- 4 The buttons allow you to configure the plate and start the run.

Enabling User Management

During the installation of QX Manager Software, Standard Edition, you can choose to enable or bypass the User Management module:

- If the Enable User Management checkbox is selected, all user functions are included in the installation.
- If the checkbox is not selected, User Management is not enabled and the corresponding tab is not included in the side pane. A single generic user, with a small set of user preferences, is always logged in. For information, see [Partial Preferences on page 44](#).

Important: Before you can set up users in the software, all users must be set up on the computer as local Windows users or in Active Directory on your network.

To bypass the User Management module

- ▶ Ensure the checkbox is cleared in the User Management dialog box, and then click Continue to complete the installation.

To enable the User Management module

1. Ensure the Enable User Management checkbox is selected.
2. Enter the user name and click Check Name to validate the user.
3. Click Continue to add the user and complete the installation.

This page intentionally left blank.

Chapter 4 Getting Started

Use the information in this chapter to

- Sign into the software or change users
- Select the instrument
- Become familiar with the instrument status bar and main functional windows
- Understand the compatible file types for your software edition

Tip: If QX Manager Software goes into standby mode after being idle, click in the lockout screen to log into the software again.

Signing In, Viewing Privileges, and Changing Preferences

Important: If User Management was not enabled during the software installation, QX Manager Software creates a single generic user, which is already logged in and is available to all users for setting up and running plates, and analyzing the results. The user can access a set of partial preferences using the generic user link. For more information, see [Partial Preferences on page 44](#).

Typically, a system or software administrator sets up users and privileges for QX Manager Software, and communicates the information to you.

Note the following:


- When you sign in the first time after the software is installed, you must select the instrument to which the computer is connected. See [Selecting the Instrument on page 40](#) for information.
- You cannot change your user privileges unless you are assigned the Add/Manage Users privilege. For information, see [Viewing User Privileges on page 41](#).
- You can change your user preferences, which are set to defaults. For information, see [Managing Your Preferences on page 43](#).

Signing into the Software

Important: If User Management was not enabled during the installation, the generic user is automatically signed in when you open the software.


To sign in

1. Right-click the software icon on the desktop and select Open to display the Sign in dialog box.
2. Do one of the following to enter your user name:
 - If the Sign in to: label shows the correct domain name, enter only the user name and password.

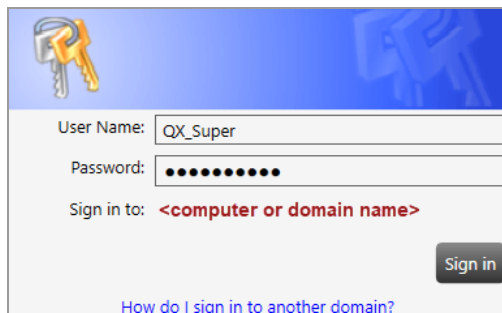


The screenshot shows a sign-in dialog box with a blue header and a key icon. It contains three input fields: 'User Name' with 'QX_Super', 'Password' with masked characters, and 'Sign in to:' with '<computer or domain name>'. A 'Sign in' button is at the bottom right, and a link 'How do I sign in to another domain?' is at the bottom.

- If the Sign in to: label is blank or shows a different domain, then enter it as **<computer or domain name>****user name**, followed by the password.

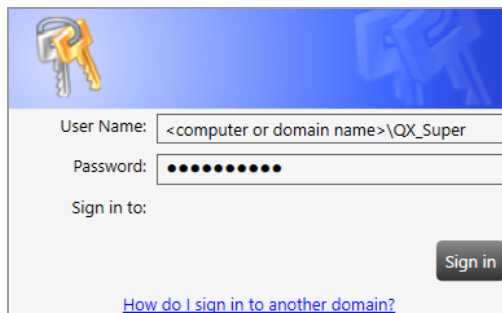

The screenshot shows the same sign-in dialog box. The 'User Name' field now contains '<computer or domain name>\QX_Super'. The 'Sign in to:' field is empty. The 'Sign in' button and the link 'How do I sign in to another domain?' are also present.

For more information, click the [How do I sign into another domain?](#) link.

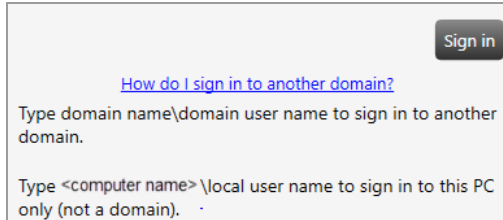

Note: If you are logging into a network domain, but don't know the name, contact your system administrator.

3. Ensure the Do not show EULA on sign in checkbox is selected and tap I Agree.



Each user must agree to the end user license agreement on initial login in order to use the software. If you clear the checkbox, the End User License Agreement appears the next time you log into the software, and you must agree again to the terms before the software opens. If you do not agree to the EULA, the application closes immediately.

When you agree to the EULA, the dialog box closes and the application opens. Your user name appears in the Instrument Status bar.

If you are connected to the QX600 Droplet Reader, the Add Plate window opens.

If you are opening the software on a computer not connected to an instrument, the Data Analysis window opens.


Selecting the Instrument

When you open the software and sign in for the first time, the instrument connection window appears and you must select the connected instrument.

This procedure assumes the instrument is connected to the computer and QX Manager Software is installed. If you are using the QX200 Droplet Reader and the QX600 Droplet Reader in your lab, you can change instruments and then select the connected instrument in System Settings.

To select the instrument

1. Select QX600, and then click Save.

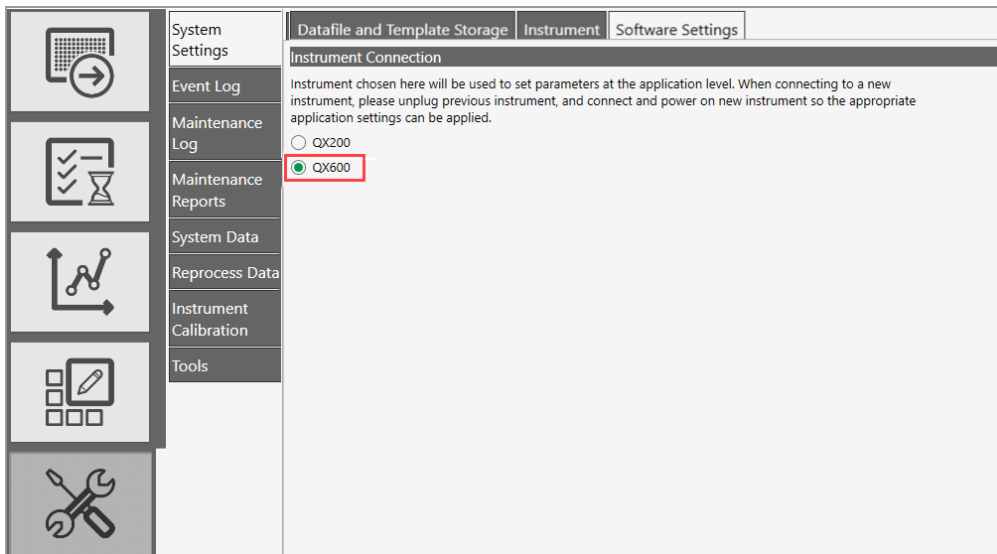

Viewing User Privileges

Important: If you did not enable the User Management module during the installation, the related functions are not available.

In the User Management window, your privileges are identified by the selected checkboxes.

To view your assigned privileges

- ▶ Select the User Management tab.



Unless you are assigned the Add/Manage Users privilege, checkboxes for your assigned privileges are selected but not enabled, as shown in the following graphic. To request changes to your privileges, contact your software administrator.

User Name:	<input type="text" value="(computer or domain name\QX_User1)"/>	<input type="button" value="Check Name"/>
Full Name:	QX_User1 (full name)	
<u>Privileges</u>		
<input type="checkbox"/>	Add/Manage Users	
<input checked="" type="checkbox"/>	Create New Templates	
<input checked="" type="checkbox"/>	View Datafiles (created by other users)	
<input type="checkbox"/>	Overwrite Existing Datafile name	
<input checked="" type="checkbox"/>	System Settings	
<input checked="" type="checkbox"/>	Maintenance	
<input type="checkbox"/>	System Data	
<input type="checkbox"/>	Manage Plate	

[Table 11](#) lists the available privileges, which can be assigned in any combination.

Table 11. User privileges

Privilege	Purpose
Add and manage users	Add or remove users, set privileges, and change preferences. Important: Users who are assigned the Add/Manage Users privilege can create users and update privileges and preferences for any user. Only the superuser can remove this privilege from user accounts.
Create new templates	Save a design for a plate or analysis report as a template.
View data files (created by other users)	See run files that were created from runs executed by other users.
Overwrite existing data file name	Save the file with a new file name. Files are always saved with the extension for the application you are using. For information, see Compatible File Types on page 48 . Note: Users without this permission can open files and perform analysis but cannot create a new file.
System settings	View logs, reprocess data with new calibration values, and change file locations. All users can view the file locations. Important: If your administrator selected the Preferred Locations checkbox in System Settings, individual user location preferences are overridden.
Maintenance	View the event log and the system data settings. Note: All users can reprocess data, view system folder settings, as well as the maintenance log and maintenance reports.
System Data	Move raw data from the computer to free up disk space. Access troubleshooting data.
Manage Plate	Cancel a run while it is in process on the QX600 Droplet Reader.

Managing Your Preferences

Important: If you did not enable User Management during the software installation, the User Management module is not included and a truncated set of preferences is available. See [Partial Preferences on page 44](#) for information.

Use the User Management window to modify your personal user preferences. Default folder locations appear in your preferences, but you can change them. You can also choose to keep all data files and templates private. The software prompts you to choose either your personal folder or the shared folder each time you save a file or template.

To change your preferences

- Change any of the following preferences:
 - Enter a different file path for your data files and templates.
 - Select or clear the check boxes to change your data file and template privacy settings.

Important: The paths and checkboxes are disabled if your administrator has set Preferred Locations in System Settings.
 - Enter a different system timeout period.
 - Enter a different total of completed plates to show in the Run Status window, up to a maximum of 100.
 - Select the Acquire all Wells By Default checkbox. If you do not select the checkbox, all wells are excluded by default and you must manually include the wells for each droplet reading run.
- Do one of the following:
 - If you are setting preferences for a new user, click Add.
 - If you are updating preferences, click Save.
- When the confirmation message appears, click Yes, and then click OK.

Partial Preferences


If User Management was not enabled during the software installation, you can still access partial functionality.

To open the menu

1. Click the user name link in the upper-right corner to display the pop-up.

Important: The Service Sign in option is used by Bio-Rad service engineers only.

2. Click User Preferences to display the dialog box.

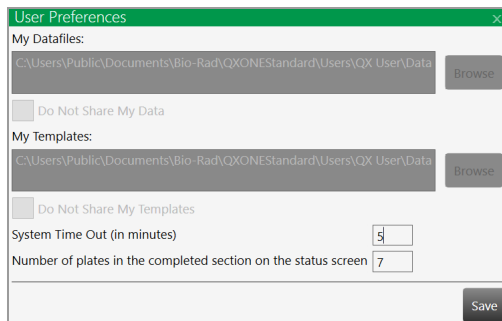

The screenshot shows a 'User Preferences' dialog box with the following fields and options:

- My Datafiles:** A text field containing the path 'C:\Users\Public\Documents\Bio-Rad\QXONEStandard\Users\QX User\Data' and a 'Browse' button.
- Do Not Share My Data:** An unchecked checkbox.
- My Templates:** A text field containing the path 'C:\Users\Public\Documents\Bio-Rad\QXONEStandard\Users\QX User\Data' and a 'Browse' button.
- Do Not Share My Templates:** An unchecked checkbox.
- System Time Out (in minutes):** A text input field with the value '5'.
- Number of plates in the completed section on the status screen:** A text input field with the value '7'.
- Save:** A button at the bottom right.

3. You can change the following:

- System time out in minutes (default is 5 minutes)
- Number of plates in the completed section of the status screen (default is 7)

4. After you have made your changes, click Save.

Signing Out or Changing Users

While the instrument is idle, one user can log out and another user can log in.

Important: If User Management was not enabled during the software installation, the sign in, sign out, and change users functions are not available.

Note the following:

- Do not change users while a run is in progress or data might be affected.
- If the screen is locked and the initial user left unsaved changes, an advisory prompt appears. Click Save before proceeding or the previous user's work is lost.

To sign out

1. Click the user name link in the upper-right corner and select Sign Out.
2. Click Yes to confirm.

To change users

1. Click anywhere in the lockout screen to display the Sign in window.
2. Enter a user name.
 - The domain name appears below the user login fields. If you are on the same domain as the previous user, you can enter your user name only.

- If the user is on a different domain, enter the domain name followed by a backslash and the user name.

<domain name>\<user name> (for example, global\john_smith)


3. Enter the password and click Sign in.

Instrument Status Bar

An instrument status bar appears above all windows except the analysis displays, which open in a separate module.

Important: If User Management was not enabled during the installation, the generic user name appears on the far right.

When a computer is connected to the instrument, the status bar appears as shown below. If a computer is not connected to the instrument, the status bar shows only the software, disk space, and user information.

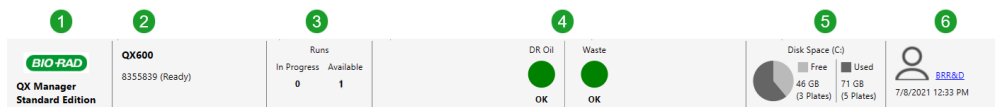

If you have installed an EcoTank, the status bar appears as shown below:



LEGEND

- 1 Software name and edition

- 2 Information and status of the connected instrument

- 3 Instrument availability

- 4 Oil and waste levels

- 5 Used and free disk space
Available storage is displayed in GB, and also in number of plates that you can run before running out of space on the computer.
Note: When QX Manager Software detects low disk space, an advisory prompt to archive data appears.








- 6 Name of the current user, along with the current date and time.

- 7 The status bar with an EcoTank loaded in lieu of individual oil and waste bottles.
If an EcoTank is loaded into the droplet reader, oil and waste levels are displayed as a single circle with one status for both. For information on EcoTanks, see [Using an EcoTank Oil and Waste Container on page 207](#).

Tabs to Functional Windows

Table 12 describes the functional areas and primary windows in QX Manager Software.

Table 12. Window Tabs

Tab	Name	Purpose
	Add Plate	<ul style="list-style-type: none"> ■ Add and configure a plate for droplet reading
	Run Status	<ul style="list-style-type: none"> ■ In the top pane, view the run in progress ■ In the bottom pane, view up to 100 completed runs
	Data Analysis	<ul style="list-style-type: none"> ■ Access the Data Analysis and Gene Study modules
	Templates	<ul style="list-style-type: none"> ■ Set up templates for plates and reports, and search for existing template files
	System	<ul style="list-style-type: none"> ■ View storage locations, logs, and maintenance reports ■ Archive data, reprocess data with new calibration values, access troubleshooting information <p>Important: Available functionality depends on your assigned user privileges. Instrument calibration is available only to the Bio-Rad service engineer.</p>
	Users	<ul style="list-style-type: none"> ■ View the user privileges you have been assigned, and view and modify your personal preferences ■ Create, edit, or remove QX Manager Software users (if you are assigned the Add/Manage Users privilege) <p>Important: In Standard Edition, user functionality is available only if User Management was enabled during the installation.</p>
	Help	<ul style="list-style-type: none"> ■ Access software version information, current calibration values, End User License Agreement company website, and open source software license information <p>Tip: You can also link the user and instrument guides to the labeled buttons.</p>

Compatible File Types

This section describes the file types you can open in QX Manager Software.

The following applies if your organization is using more than one edition of the software:

- You can open files created using Standard Edition only in Standard Edition.
- If you open a .qlp file in QX Manager, Standard Edition, the file opens and is saved as a .ddpcr file.

Table 13. Compatible file types

File Type	Extension	Details
Plate template	.ddplt	Plate template file containing setup details to perform experiments; this file type opens in the Plate Editor.
Plate template	.csv	Importable plate template file containing setup details in CSV format; this file type opens in the Plate Editor.
Data	.ddpcr	Contains the results of an experiment run performed on the droplet reading instrument and opened in QX Manager Software, Standard Edition.
Data	.qlp	File type from earlier QuantaSoft™ versions, for two-channel experiment runs and data analysis; this file type opens in the Analysis module.


Chapter 5 Running Experiments

When the Start Run button is enabled, you can start the run. You can also cancel a run after it has started if you are assigned the Manage Plate user privilege. For information, see [Canceling a Run on page 65](#).

Assembling a Plate


The QX600 Droplet Reader accepts a solid 96-well plate that fits into a secure holder. You must assemble the plate components before you can insert the plate into the QX600 Droplet Reader. Three separate components comprise the plate:

- A base section to hold the plate
- The plate itself
- A top section to secure the plate

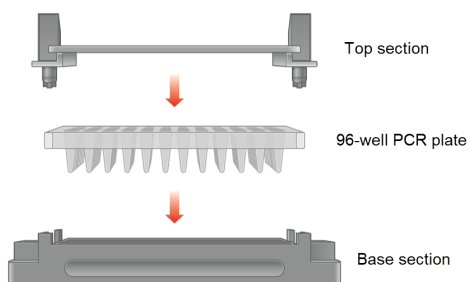

To assemble the plate

1. Place the plate into the base section, and then raise the release tabs on the top section.
2. Place the top section over the plate, and then lower the release tabs to secure the plate in the holder.

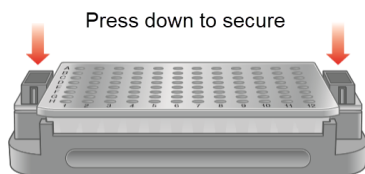


Inserting an Assembled Plate into the Instrument

This section explains how to insert an assembled plate into the QX600 Droplet Reader.


Note: You can insert the plate into the droplet reader before or after you add the plate in QX Manager Software, but you cannot configure the plate until the software recognizes the plate in its compartment.

To insert a plate into the QX600 Droplet Reader


1. Press the button to open the compartment.

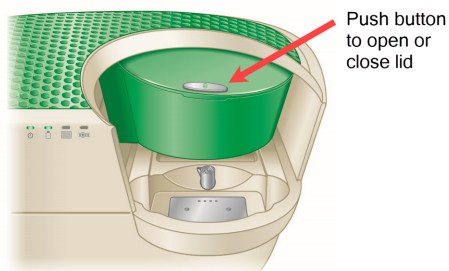

2. Fit the assembled plate into the compartment. Well A1 must be in the top-left position.

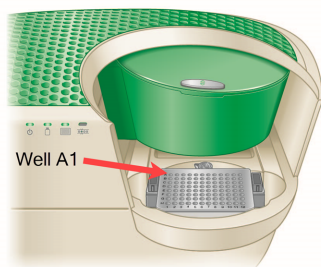

Important: Make sure the tabs on the plate holder are securely clamped and the plate lays flat. Otherwise, the sampling needle can strike the lid and damage the instrument.

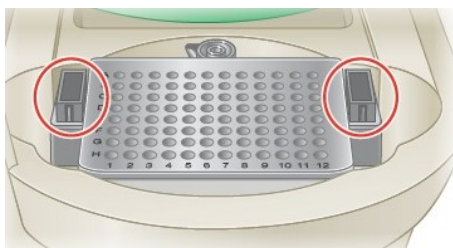

3. Press the button again to close the lid.

Adding the Plate in QX Manager Software

This section explains how to add the plate in QX Manager Software, Standard Edition when using the QX600 Droplet Reader for your droplet reading run.


To add a plate in QX Manager Software

1. Click the Add Plate tab.




If the Runs Available number in the Instrument Status Bar is 1, the Add Plate button is enabled.



2. Click  next to Add Plate.



When QX Manager Software recognizes the plate holder in the compartment, the software enables the Configure Plate button.

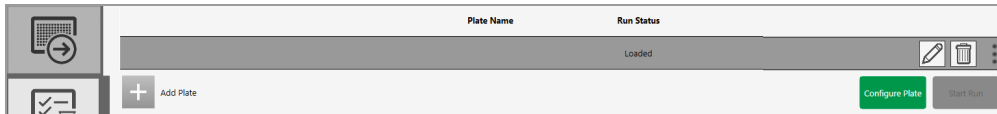

Note: You can insert the plate into the droplet reader before or after you add the plate in QX Manager Software, but you cannot configure the plate until the software recognizes the plate in its compartment. If the software does not recognize the plate holder, an advisory message appears. Remove and reinsert the plate. If necessary, contact Bio-Rad Technical Support.

3. Click Configure Plate.
4. When the Plate Configuration window appears, continue to [Configuring the Plate on page 52](#).

Configuring the Plate

The Plate Configuration window features the Plate Information, Well Selection, and Well Information tabs.

All users can start runs using the initial plate configuration, an altered configuration, or just the plate and data file names and the supermix. Users with the Create New Template user privilege can save the configuration as a template before the droplet reading run.

Plate Information Tab

The Plate Information tab is selected by default when the Plate Configuration window opens.

In the Plate Information window, you can

- Select an existing template from the dropdown list
 - Create a new template
 - Browse for an existing template
- Note:** When using the Browse button to select a template, the template is applied to the run, but is not imported or saved in QX Manager Software.
- Name your plate and select a supermix
 - QX Manager Software automatically enters the plate name into the data file name field, but you can change it.
 - You can start the run with the plate name and supermix only, and then you can identify the remainder of your well information parameters in the Analysis module Plate Editor after the run has concluded.
 - Select to acquire wells by columns or rows

Identifying Plate Information

To identify the plate

1. Do one of the following:

- To choose an existing template, select a template file from the Choose Existing dropdown list. QX Manager Software automatically populates the fields under Create New Plate.
- To create a new template, click Create New, and see [Creating or Editing Templates on page 67](#).
- To browse for a template, click Browse and navigate to the templates folder (if necessary), and then select a template file. QX Manager Software automatically populates the fields under Create New Plate.


Note: When using the Browse button to select a template, the template is applied to the run, but is not imported or saved in QX Manager Software.

- To identify the plate, enter a plate name and select a supermix under Create New Plate. When you enter the plate name, the Save data file as field is automatically populated with the same name, but you can change it.

2. (Optional) Under Acquire Wells By, select the alternative well acquisition method. The default is Column.

Well Selection Tab

When you select the Well Selection tab, the Plate Editor grid appears. The wells in the plate layout are automatically excluded and you must manually include the wells to be processed. If you typically run full plates, you can change the setting in User Preferences to acquire all wells by default. For information, see [Managing Your Preferences on page 43](#).

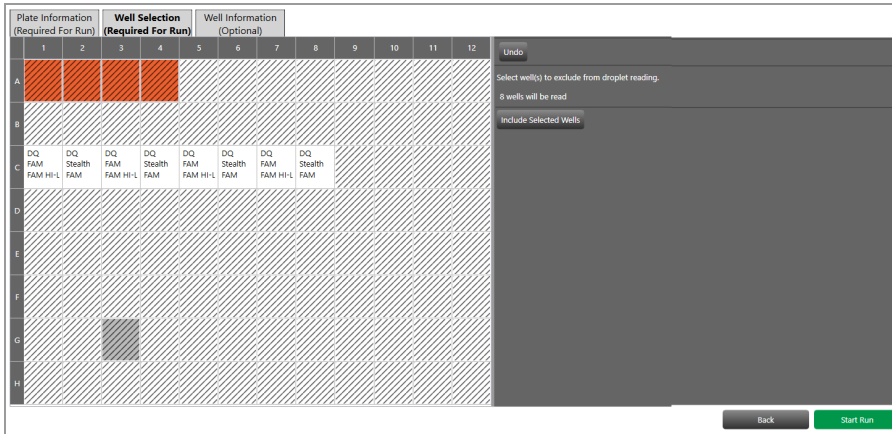


Including or Excluding Wells in the Plate Layout

All wells are excluded by default when the plate editor appears.

To include wells in the run

1. Select the Well Selection tab.
2. In the Plate Editor layout, select the wells to include in the run (those containing sample), and then click the Include Selected Wells toggle button.
3. On the right, QX Manager Software specifies the number of wells that will be included in the droplet reading run and the button text changes to Exclude Selected Wells.

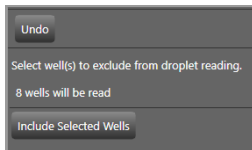


Well Information Tab

When you select the Well Information tab, the Plate Editor opens. You can configure your experiment parameters for selected wells before or after the run.

Note: QX Manager Software can calculate automatic thresholds for positive control wells, but you must select Pos Ctrl from the Sample Type dropdown list. You can make the selection before or after the run. For information, see [Sample Types on page 74](#).

Defining or Editing Well Information

You can assign different experiment parameters to each well or groups of wells.

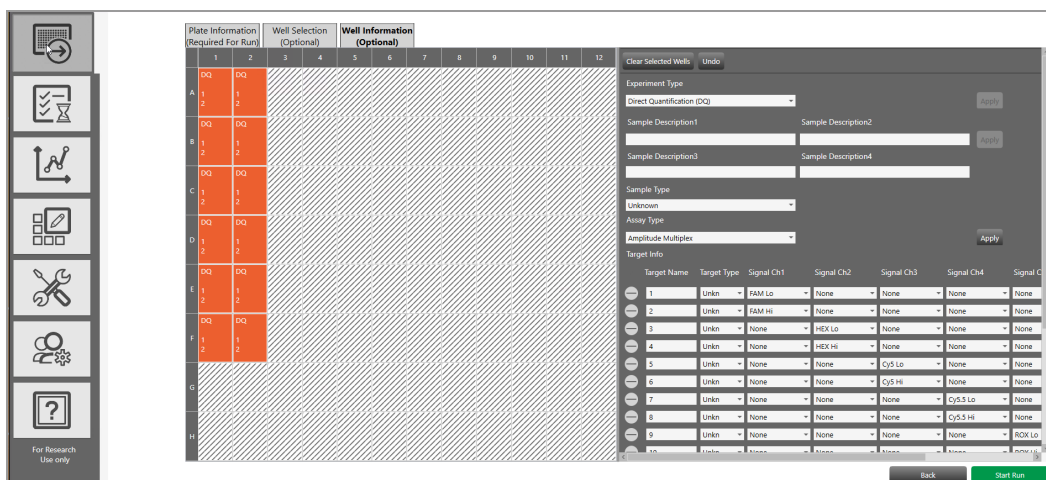

You can apply your entries at any time. You must click Apply for the software to recognize your new or edited information. If you enter information in a field and need to apply earlier changes first, QX Manager Software displays an advisory message.

To define the configuration for the wells


1. Select a well or group of wells.
2. Select an experiment type; see [Experiment Types on page 73](#).
3. In the Sample Description fields, enter up to four words or phrases that describe the sample; see [Sample Descriptions on page 74](#).
4. Select a sample type; see [Sample Types on page 74](#).

Note: For automatic thresholding of positive control wells, you must select Pos Ctrl.

5. Select a supermix; see [Supermixes on page 75](#).

Chapter 5 Running Experiments

6. Select an assay type.
7. Available assay types vary, depending on the selected experiment type; see the fluorophore options in [Assay Types and Fluorophores on page 76](#).
8. (Optional) Under Target Info, change the fluorophore assigned to a channel. The fluorophore information populates automatically for each target, with the maximum number of rows allowed for the assay method. However, you can modify the default fluorophore assignments, if applicable.
 - a. Click the dropdown arrow and select an alternative.
 - b. To delete a row, click the minus (-) icon next to the target.
 - c. To add a row, click the plus (+) icon.
9. (Optional) To clear one or more wells at any time, select the wells and click Clear Selected Wells.
10. Click Apply. After you have applied your entries and selections, you can pause the cursor on a well to see the well information.

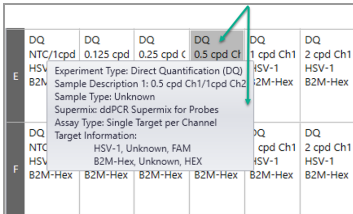

E	DQ	DQ	DQ	DQ	DQ	DQ
	NTC/1cpd	0.125 cpd	0.25 cpd	0.5 cpd	1 cpd Ch1	2 cpd Ch1
	Experiment Type: Direct Quantification (DQ)				HSV-1	HSV-1
	Sample Description: 1: 0.5 cpd Ch1/1cpd Ch2				B2M-Hex	B2M-Hex
	Sample Type: Unknown					
	Supermix: ddPCR Supermix for Probes					
	Assay Type: Single Target per Channel					
F	DQ	Target Information:			XQ	DQ
	NTC	HSV-1, Unknown, FAM			1 cpd Ch1	2 cpd Ch1
	HSV	B2M-Hex, Unknown, HEX			HSV-1	HSV-1
	B2M-Hex	B2M-Hex	B2M-Hex	B2M-Hex	B2M-Hex	B2M-Hex

11. When you finish configuring the plate, click Save.

Starting the Run

When all required elements to set up a plate for a run are satisfied, QX Manager Software enables the Start Run button. You can configure the elements of the plate before you start the run, or you can run the plate with just the plate name and supermix, and define the configuration from the Analysis module.

To start the droplet reading run


1. Tap Start Run.



The software performs a mandatory prerequisites check to ensure that the instrument is ready.


One of the following occurs:

- If the instrument passes the check, the run starts normally. Continue to [Running the Experiment on page 61](#).
- If the check finds issues with the instrument, the software displays a dialog box in which the issues are identified. The instrument cannot start the run until all issues are resolved. For information, see [Run Prerequisites Checklist on page 58](#).

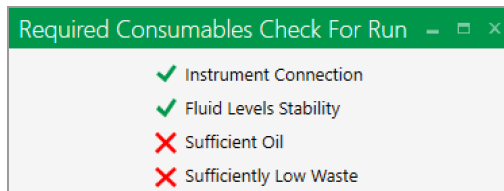

Run Prerequisites Checklist

If the prerequisites check on the instrument status finds one or more problems with instrument readiness, QX Manager Software identifies them with an **✗** in the Check for Run dialog box. You must resolve all issues before the instrument can start the run.

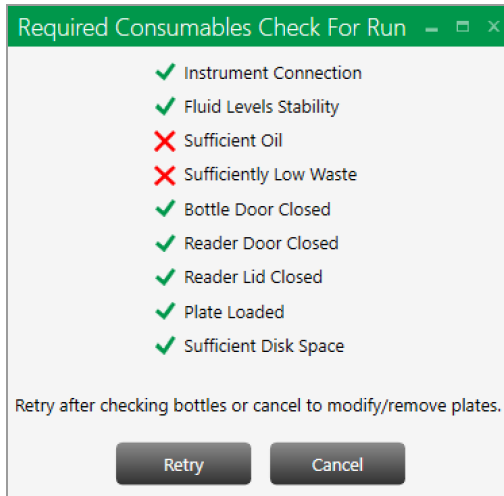

Note: If you encounter a problem with the instrument or software that is not described here, see [Appendix C, Troubleshooting](#).

For information on each issue that can prevent the instrument from starting the run, see [Table 14 on page 60](#).

Table 14. Pre-Run Error Indicators

Status Check Item	Issue	Suggested Resolution
Instrument Connection	Power connection not recognized.	Ensure the power outlet is working. Disengage the power cord from the outlet, and then reinsert it into the outlet.
	USB cable connection not recognized.	Disconnect, and then reconnect, the USB cable between the instrument and the computer.
	Instrument connection not recognized.	Ensure the correct instrument is selected in the System Settings tab > Instrument tab.
Fluid Levels Stability	The oil or waste is moving within the bottle so the level cannot be accurately determined.	Wait until the liquid is settled, then try again. Ensure the instrument rests on a solid, stable surface.
Sufficient Oil	Oil level is below the acceptable level for a run.	Remove the oil bottle and replace it with a full one.
Sufficiently Low Waste	Waste level is above the acceptable level for a run.	<ol style="list-style-type: none"> 1. Empty any remaining oil from the bottle and then place a Waste label (provided by Bio-Rad) over the Oil label. 2. Insert the bottle into the Waste slot.
Bottle Door Closed	Bottle door is not closed properly.	Open the bottle door and ensure there are no blockages, and then close the door again.
Reader Door Closed	Reader door is not closed properly.	Open the reader door and ensure there are no blockages, and then close the door again.
Reader Lid Closed	Reader lid is not closed properly.	Open the reader lid and ensure there are no blockages, and then close the lid again.

Table 14. Pre-Run Error Indicators, continued

Status Check Item	Issue	Suggested Resolution
Plate Loaded	Plate is not loaded properly.	<ol style="list-style-type: none">1. Open the reader door and remove the plate.2. Ensure the tabs of the plate cover are securely clamped.3. Insert the plate again and ensure it lays flat.
Sufficient Disk Space	There is not enough storage space available for the run data.	Archive older data files to clear space.

Running the Experiment

The instrument allows you to do the following:

- During the run, perform real-time analysis on the droplets in each well after the well has been read
- When the run concludes, QX Manager Software moves the run to the Completed list in the Run Status window, and saves the analysis data file. Data files are always saved to your personal file location or a preferred location, and can also be saved to a shared location. For information, see [Managing Your Preferences on page 43](#).

Important: if your system administrator has designated Preferred Locations in System Settings, they override any file paths indicated in your preferences.

- During and after the run, QX Manager Software displays run information in the Run Status window.
- Cancel a run (with the Manage Plate permission).

Note: Canceling a run also shuts down the instrument.

Plate Name	Run Status	Location	Time Remaining	Time Elapsed
Test		Droplet Reader	01:58:37	00:00:12

Total Time Remaining: 1:58:37

Plate Name	Run Status	Run Completed	File Name
Test		07/21/2021 10:38:46 AM	Test_20210721_103740_583.ddpcrs
.TestFile		07/19/2021 04:16:15 PM	.TestFile_20210719_161528_357.ddpcrs

Tracking the Run Status

You can track each run from the Run Status window.

Plate Name	Run Status	Location	Time Remaining	Time Elapsed
Test		Droplet Reader	01:58:37	00:00:12

Total Time Remaining: 1:58:37

Plate Name	Run Status	Run Completed	File Name
.Test		07/21/2021 10:38:46 AM	.Test_20210721_103740_583.ddpcrs
.TestFile		07/19/2021 04:16:15 PM	.TestFile_20210719_161528_357.ddpcrs

The top pane shows the run in progress. The run remains in the top pane for the duration of the processing.

Icons displayed in the Run Status column indicate each process point, as shown in [Table 15 on page 62](#).

Table 15. Run status indicators

Status Icon	Description
	Ready to begin
	Droplet reading run
	Run completed
	Run completed, plate removed
	Run stopped or canceled by the system, or canceled by the user before it began

When the run is finished, the software moves the run to the bottom pane, where up to 100 completed runs are visible.

Tip: You set the number of completed runs to display in your user preferences. See [Managing Your Preferences on page 43](#).

The content of each column is explained in [Table 16](#).

Table 16. Run Information Columns

Column Name	Location	Description
Plate name	Top and bottom panes	Name entered under Create New Plate Note: If you selected an existing template, the template name appears.
Run status	Top and bottom panes	Icons for each process step (see Table 15 on page 62)
Time remaining	Top pane only	Hours and minutes until the run will be finished
Time elapsed	Top pane only	Time, in hours and minutes the run has been in progress
Total time remaining	Below the top pane	Total hours and minutes remaining for all in-progress runs
Run completed	Bottom pane only	Date and time stamp showing when the instrument completed the run
File name	Bottom pane only	Plate name and system-generated identifier


After run completion, the software saves the analysis data file in a location designated by you or a system administrator.

Using Live Analysis


When the QX600 Droplet Reader initiates droplet reading, QX Manager Software provides an option for real-time analysis of the wells that have been read, while the process as a whole is still active.

QX Manager Software initially shows each well in the Plate Editor as disabled. As the software confirms that the instrument has acquired a well and read the droplets, the software enables the well for live analysis in read-only mode. The process continues until all wells are enabled and the process is finished.

To use live analysis

1. Tap the droplet reading icon () for the run in progress.

The Droplet Reading dialog box opens.

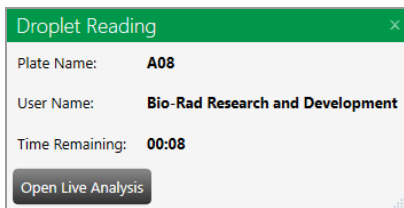

2. Click Open Live Analysis.
3. Select an analysis window.
4. In the well selector, select one or more wells that are enabled.


Important: QX Manager Software displays read-only data in the selected analysis window. You can view the data but to make changes, you must wait for the run to complete, and then change the data in the Analysis module.

Canceling a Run

You can cancel a run if you are assigned the Manage Plate user privilege.


To cancel the run

1. In the Run Status window, click the X on the right to cancel the run.

Location	Time Remaining	Time Elapsed	
Droplet Reader	00:03	00:00	


Note: Canceling a run also shuts down the instrument.

2. When the Are you sure message appears, click Yes.

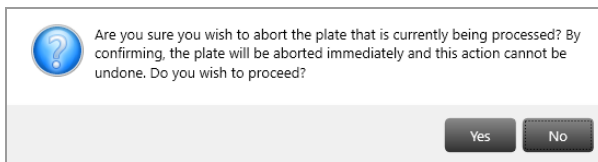

The instrument finishes reading the current well, stops the process and reverts to its pre-run state.

Note: If you click No, or if you do not respond to the confirmation message within 30 sec, the process continues without interruption.

The  icon appears in the Run Status column.

QX Manager Software creates the data file for the wells that were processed on the canceled plate and the status in the Event Log is specified as “run aborted by user.”

This page intentionally left blank.

Chapter 6 Creating or Editing Templates

If you have been assigned the Create New Templates user privilege, you can create, edit, and save plate and report templates for your common experiments. Templates that are available to you appear in the dropdown list in the Plate Information window, and also in the default view of the Template Setup tab. If the list is long, you can use the Search field to narrow the results.



All users can configure a plate template in a CSV file, and then import the template into QX Manager Software. For information, see [Creating and Storing Plate Templates from CSV Files on page 79](#).

Note: You can select a CSV template from the dropdown list in the Plate Information window only. They do not appear in the My Templates list.

For information on report templates, see [Creating or Modifying Report Templates on page 81](#).

Templates you created are saved to one of the following folders:

- The path specified in your user preferences; these appear under My Templates in the Templates window.
- The paths specified by your administrator in the shared or preferred location fields in System Settings.

Important: Preferred locations override personal folder settings and templates from all users are saved there.

You can also browse for templates when you are setting up your plate for a run.

Creating or Modifying Plate Templates

Important: You must be assigned the Create New Template user privilege to create or edit, and then save, a template.

To create or edit a template

1. Do one of the following:
 - Navigate to the Plate Information window (click Plate Setup > Add Plate > Configure Plate) and continue to Step 2
 - Open the Templates window and then go to Step 3.
2. From the Plate Information window, under Plate Template, do one of the following:
 - Select an existing template from the Choose Existing dropdown list.
 - Click Create New. The plate layout opens in Edit mode.
 - Click Browse and navigate to an existing template.

Continue to Step 4.

3. From the Templates window, do one of the following:
 - Select the Plate Setup tab and click Add New Plate > Edit Plate Layout.
 - Double-click an existing template from the list.

The Plate Editor opens in Edit mode. The Instrument Type field is grayed out.

Note: CSV templates do not appear in the Templates window, but are available from the dropdown list in the Plate Information tab.

Continue to Step 4.

4. Click the Exclude button, and select the wells to include in the plate layout, and then click Include selected wells.

By default, all wells in the Plate Editor are automatically excluded. You must select and include new wells in the template. If you typically run full plates, you can change the setting to Acquire all wells by default in your user preferences. For information, see [Managing Your Preferences on page 43](#).

5. Click the Edit button to display the layout for experiment and sample identification.

The fields are enabled when you include wells.

6. For one or more wells at a time, complete any or all of the following substeps to include the selections and entries that should be static. Leave blank any field that should be completed at the time of the experiment run.

- a. Select an experiment and sample type, supermix, assay type, and enter sample descriptions.

- [Experiment Types on page 73](#)
- [Sample Types on page 74](#)
- [Sample Descriptions on page 74](#)
- [Supermixes on page 75](#)
- [Assay Types and Fluorophores on page 76](#)

Note: For automatic thresholding of positive control wells, you must select the Pos Ctrl sample type.

- b. If you selected Residual DNA Quantification as the experiment type, enter a conversion factor.
- c. Enter up to four sample descriptions, and any relevant well or plate notes.
- d. Edit the Target Name and Target Type fields.

7. Click Apply.

8. Click Save and select one of the following:

- Shared Templates. The template is saved in the folder specified in the System Settings.
- My Templates (saves the file to the path in your user preferences) or System Templates (saves the file to the preferred location specified by your system administrator).
- If your system administrator has specified preferred locations for all users, templates are automatically saved to that location.

Template Plate Editor Window

Use the Template Plate Editor to

- Create new or edit existing plate template files, and save new or modified templates for reuse
- Add up to four sample descriptions per well
- Exclude wells from droplet reading


Important: When you create an experiment from a template with wells excluded in the plate layout, the instrument pierces the foil for all wells in the physical plate, but reads droplets for active wells only.

To open the template plate editor


1. Click the Templates tab on the left.



2. In the Templates window, click Edit Plate Layout

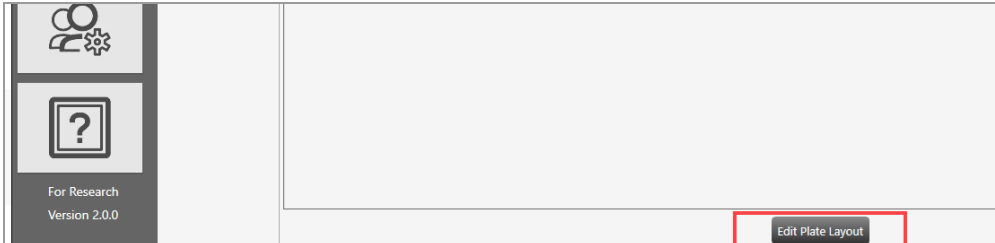

You can also open the template plate editor from the Add Plate window.

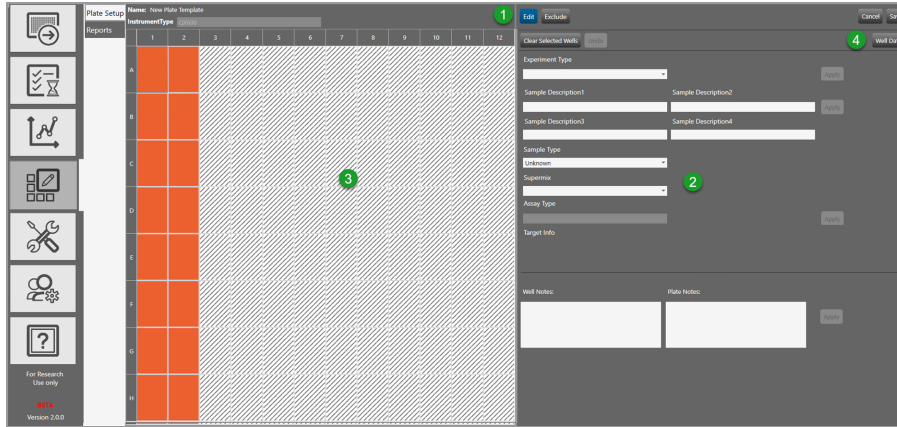

LEGEND

- 1 The Edit and Exclude buttons are accessible from the top of the screen.

- 2 The right pane contains the interface for defining your experiment parameters.

- 3 The left pane displays the plate grid and configuration information in each well.

- 4 The Well Data button toggles to the Well Data table, where you can see your plate setup in a tabular format, and then back to the Plate Editor. When the Well Data table appears, the button name changes to Plate Layout.

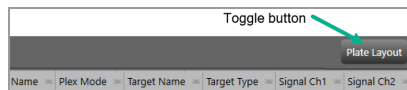


Plate Editor Tools

Use the buttons described in [Table 17](#) when you are creating a template or defining an experiment in the Plate Template Editor.

Table 17. Plate Editor Buttons

Button	Description
Edit	<p>Allows you to</p> <ul style="list-style-type: none"> ■ Set the experiment type, sample name and type, supermix, and assay type ■ Set the target name, type, and fluorophore for each channel ■ Add well or plate notes
Exclude	<p>Toggle button that allows you to include or exclude wells from droplet reading</p> <p>Note: Wells in the plate editor are initially excluded by default.</p>
Cancel	Allows you to cancel and return to the Template Setup window
Save	Opens a dialog box where you can enter a file name and storage location
Clear Selected Wells	Allows you to delete information from selected wells
Undo	<p>Allows you to reverse your last action</p> <p>Note: This button is enabled only after you have applied at least one experiment parameter (for example, the selection for Experiment Type).</p>
Well Data or Plate Layout	Toggle button that allows you to access the Well Data table and return to Plate Layout view
Apply	Applies your entries and choices for selected wells

Experiment Types

QX Manager Software offers the experiment types described in [Table 18](#), each enabling different assay options and analysis tools. You can assign multiple experiment types across the plate, but you can assign only one experiment type to each well.

Table 18. Experiment types

Type	Description
Direct quantification (DQ)	DQ uses absolute quantification to determine the concentration (copies/ μ l) of target DNA copies in a sample. DQ measures the number of positive and negative droplets for each target in a well based on their fluorescence amplitude, and uses a Poisson algorithm to calculate the starting concentration of each target DNA molecule. Note: The ABS experiment type created in QuantaSoft Software versions 1.4 through 1.7 is automatically mapped into the software as a DQ experiment type.
Copy number variation (CNV)	CNV determines concentration and calculates the copy number of an unknown (CNV) target relative to a known reference or references within the same well.
Mutation detection (MUT)	MUT determines concentration and calculates the fractional abundance of an unknown mutant present at a low frequency in a wild-type background.
Rare event detection (RED)	RED determines the concentration of a known mutant or a rare target species relative to a given reference species within a large pool of background DNA.
Drop-off (DOF)	DOF determines absolute quantification of targets for assays designed to detect non-wild type sequences, such as indels and genome edits; DOF is designed to support an assay strategy where one probe counts all alleles and one “drop-off” probe sits on top of a predicted cut site or mutation site.
Gene expression (GEX)	GEX determines concentration (as with DQ) and calculates relative expression levels of an unknown target relative to a known reference or references within the same well.
Residual DNA quantification (RDQ)	RDQ provides a precise method for residual host cell DNA quantification. Note: Users can enter a conversion factor (1/C-value) for a given genome (human, mouse, etc.) to calculate the mass for a well based on the number of copies measured, up to three decimals. The software displays the result in pg/ μ l of the ddPCR well in the Molecular Weight column of the Data Table. Starting sample concentration can then be back-calculated based on dilution factor.

Sample Descriptions

QX Manager Software allows you to enter descriptive words or phrases for your sample in up to four fields per well. Descriptors can include information such as research type, dilution factors, and so forth.

Tip: Only the information from the first field appears in the Plate Editor layout. However, you can see entries in the second, third, and fourth fields in the Well Data table, or you can pause on a well to display a tool tip, as shown in the following graphic.

4	5	6	7	8
th	DQ S.a. 1cpd Sa822	DQ Stealth FAM	DQ S.a. 1cpd Sa822	DQ Stealth FAM
ti	Experiment Type: Direct Quantification (DQ) Sample Description 1: Stealth Sample Type: Unknown Supermix: ddPCR Supermix for Probes Assay Type: Single Target per Channel Target Information: FAM, Unknown, FAM HEX, Unknown, HEX			X

Sample Types

You can assign multiple sample types within the plate, but you can assign only one sample type to each well.

QX Manager Software provides four sample types for selection:

- Unknown
- NTC (no template control)
- Pos Ctrl (positive control) where one or more targets are expected
- Neg Ctrl (negative control), where no response is expected

Important: Beginning in v1.2, QX Manager Software can automatically calculate thresholds for positive control wells, but you must select Pos Ctrl from the Sample Type dropdown list. You can select Pos Ctrl during plate setup or in the Analysis module after the run is finished.

Supermixes

QX Manager Software offers five PCR supermixes. Each supermix is optimized to deliver maximum PCR efficiency and sensitivity for the amplification and detection of DNA and RNA targets.

Important: Use of unapproved supermixes can harm the instrument and void the warranty.

ddPCR Supermix	Description
Multiplex Supermix	For use in nucleic acid sample preparations to amplify and detect DNA targets in probe-based experiments involving multiple targets. Note: Bio-Rad recommends using ddPCR Multiplex Supermix for Cy5 and Cy5.5 probes assays.
Supermix for Probes (No dUTP)	For use in nucleic acid sample preparations, with sensitivity for the amplification and detection of DNA targets using hydrolysis probe-based assays.
Supermix for Probes	For use in sample preparation with uracil N-glycosylase (UNG) decontamination protocols to prevent the reamplification of carryover PCR products between experiments.
Supermix for Residual DNA Quantification	For use in residual DNA detection.
One-Step RT Advanced Kit for Probes	For use in absolute quantification of target RNA molecules.
EvaGreen® Supermix	For use in gene expression.

Assay Types and Fluorophores

QX Manager Software offers the assay methods explained in [Table 19](#).

Table 19. Assay types

Assay type	Description
Single target per channel	<p>Method assuming up to six probe colors (FAM/EvaGreen®, HEX/VIC, Cy5, Cy5.5, ROX, and ATTO 590) and up to six targets per channel.</p> <p>This is the default assay type, with the maximum six rows displayed. You can delete up to three rows.</p>
Amplitude multiplex	<p>Method to increase multiplexing up to twelve targets per well, with one or two targets detected per channel. The maximum of eight rows are displayed. You can delete up to seven rows.</p>
Probe mix triplex	<p>Triplex mode that allows three groups of three targets each.</p> <p>Group 1 has one target in FAM, one target in HEX/VIC, and one target in both channels. Group 2 has one target in Cy5, one target in Cy5.5, and one target in both. Group 3 has 1 target in ROX, 1 target in ATTO 590, and one target in both groups. You can delete a group, but not the individual targets within the group.</p> <p>Note: The channels within a group are not interchangeable; therefore FAM is always paired with HEX/VIC and so forth.</p>
Advanced classification	<p>Method to increase multiplexing rate in any combination of targets detected per channel. This method requires manual classification of droplets.</p>
Basic drop-off	<p>This mode allows two groups of two targets each. The first group appears by default, with one target in FAM <u>and</u> HEX/VIC and one target in FAM <u>or</u> HEX/VIC.</p> <p>This assay type is available only for the DOF experiment type.</p> <p>You can add one more group of two targets, with one target in Cy5 <u>and</u> Cy5.5, one target in Cy5 <u>or</u> Cy5.5. You can delete a group, but not the individual targets in the group.</p> <p>Note: The channels within a group are not interchangeable; therefore FAM is always paired with HEX/VIC and so forth.</p>

Table 19. Assay types, continued

Assay type	Description
Double drop-off	<p>This assay type is available only for the DOF experiment type, and allows three groups of three targets each.</p> <p>The first group appears by default, and looks for one target in FAM <u>and</u> in HEX/VIC (target type reference) one target in FAM (target type unknown), and one target in HEX/VIC (target type unknown).</p> <p>You can add the second group to look for one target in Cy5 <u>and</u> Cy5.5 (target type reference, with two reference copies), one target in Cy5 (target type unknown), and one target in Cy5.5 (target type unknown).</p> <p>You can also add a third group to look for one target in ROX <u>and</u> ATTO 590 (target type reference), one target in ROX (target type unknown), and one target in ATTO 590 (target type unknown).</p> <p>By default, all reference copies are set to 2.</p> <p>The channels are static for each group:</p> <ul style="list-style-type: none"> ■ Channel 1 FAM is always paired with Channel 2 HEX/VIC). ■ Channel 3 Cy5 is always paired with Channel 4 Cy5.5. ■ .Channel 5 ROX is always paired with Channel 6 ATTO 590

The experiment type you select determines the available assay methods. [Table 20](#) displays the sample type associated with each assay method.

Table 20. Assay types by experiment type

Assay type	Experiment type						
	DQ	CNV	MUT	DOF	GEX	RDQ	RED
Single target per channel	✓	✓	✓		✓	✓	✓
Amplitude multiplex	✓	✓	✓		✓	✓	✓
Probe mix triplex	✓	✓	✓		✓	✓	✓
Advanced classification	✓						✓
Basic drop-off				✓			
Double drop-off				✓			

QX Manager Software automatically populates the fluorophore in each channel for each target but you can change the default selections.

Use [Table 21](#) to select from the fluorophore options for each assay method listed.

Table 21. Fluorophore options

Assay method	Signal Ch1	Signal Ch2	Signal Ch3	Signal Ch4	Signal Ch5	Signal Ch6
Single target per channel, for 1 to 6 targets	FAM	HEX	Cy5	Cy5.5	ROX	ATTIO590
	EvaGreen®	VIC	None	None	None	None
	None	None				
Amplitude multiplex, for 1 to 12 targets	FAM Hi	HEX Hi	Cy5 Hi	Cy5.5 Hi	ROX Hi	ATTO 590 Hi
	FAM Lo	HEX Lo	Cy 5 Lo	Cy 5.5 Lo	ROX Lo	ATTO 590 Lo
	EvaGreen® Hi	VIC Hi	None	None	None	None
	EvaGreen® Lo	VIC Lo				
	None	None				
Probe mix triplex, for 9 targets	FAM	HEX	Cy5	Cy5.5	ROX	ATTO 590
	None	VIC	None	None	None	None
		None				
Advanced classification	FAM	HEX	Cy5	Cy5.5	ROX	ATTO 590
	1–10	1–10	1–10	1–10	1–10	1–10
	EvaGreen®	VIC	None	None	None	None
	1–10	1–10				
	None	None				

Creating and Storing Plate Templates from CSV Files

You can create a plate template in a spreadsheet application, save it in .csv format (CSV), and then store it for use in your droplet reading runs.

Note the following:

- CSV templates do not appear in the Templates window; however, when you set up a run, you can access them from the Plate Information window using the following methods:
 - Click the Choose Existing dropdown arrow and select a CSV template file from the list
 - Click Browse and navigate to the applicable storage folder, and then select the CSV template file
- You can store CSV template files in any location and navigate to the file using the Browse function, but the Choose Existing dropdown list automatically populates with templates that are stored in the following locations:
 - The personal templates storage folder defined in your user preferences (see [Managing Your Preferences on page 43](#))
 - The shared or preferred location specified by your system administrator (see [Shared Settings on page 168](#))

As shown in [Fig. 1](#) and [Fig. 2 on page 80](#), when you create a CSV template, the headings for the columns containing the experiment information must match the fields and order in the Plate Editor. If they do not match, an error message appears when you select the template.

Important: After you specify the information in row 1, do not change the field entries until you install a new software version or edition.

To create and save a CSV template

1. Open a spreadsheet program and enter the run settings into the template in the format shown below.

Fig. 1: Template CSV Layout Example, Columns A through H

	A	B	C	D	E	F	G	H
1	ddplate - DO NOT MODIFY		ApplicationName=QX Manager	ApplicationVersion=	ApplicationEdition=	User=QLDR-XXXX\	CreatedDate=7/24/2022	
2	THIS LINE	Version=1	Standard Edition	2.0.0.11098	ResearchEmbedded	QXManager_user	1:54:15 PM	
3	PlateSize=GCR96							
4	PlateNotes=Plate note text							
5	Well	Perform Droplet Reading	ExperimentType	Sample description 1	Sample description 2	Sample description 3	Sample description 4	SampleType
6	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
7	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
8	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
9	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
10	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
11	A01	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
12	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
13	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
14	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
15	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
16	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown
17	A02	Yes	Direct Quantification (DQ)	Description text 1	Description text 2	Description text 3	Description text 4	Unknown

Fig. 2: Template CSV Layout Example, Columns I through R

	I	J	K	L	M	N	O	P	Q	R	S	T	U	V
	SupermixName	AssayType	TargetName	TargetType	Signal Ch1	Signal Ch2	Signal Ch3	Signal Ch4	Signal Ch5	Signal Ch6	Reference Copies	Well Notes	Plot?	RdgConversion Factor
	ddPCR Multiplex Supermix	Single Target per Channel	1	Unknown	FAM	None	None	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	2	Unknown	None	HEX	None	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	3	Unknown	None	None	Cy5	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	4	Unknown	None	None	None	Cy5.5	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	5	Unknown	None	None	None	None	ROX	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	6	Unknown	None	None	None	None	None	ATTO590		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	1	Unknown	FAM	None	None	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	2	Unknown	None	HEX	None	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	3	Unknown	None	None	Cy5	None	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	4	Unknown	None	None	None	Cy5.5	None	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	5	Unknown	None	None	None	None	ROX	None		Well note text	FALSE	
	ddPCR Multiplex Supermix	Single Target per Channel	6	Unknown	None	None	None	None	None	ATTO590		Well note text	FALSE	

2. When you finish populating the fields, navigate to your templates storage location and save the template.

If you save the CSV file in

- The personal or shared templates storage location defined in your user preferences or shared settings, the CSV template is available from the Choose Existing dropdown list.
- Another storage location folder, you can navigate to the template file using the Browse feature in the Plate Setup Plate Information tab.

Creating or Modifying Report Templates

If you are assigned the Create New Templates user privilege, you can create and save a report template by editing the default report structure or by opening and editing an existing report.

To create or edit a report template

1. Select the Template Setup tab.




2. Click Reports and do one of the following:

- To create a new template, click Add New Template. and click Edit Report Template.
- To open an existing template, select the template and click Edit Report Template.



3. Click the arrow to the left of each category to expand the list of reportable items.

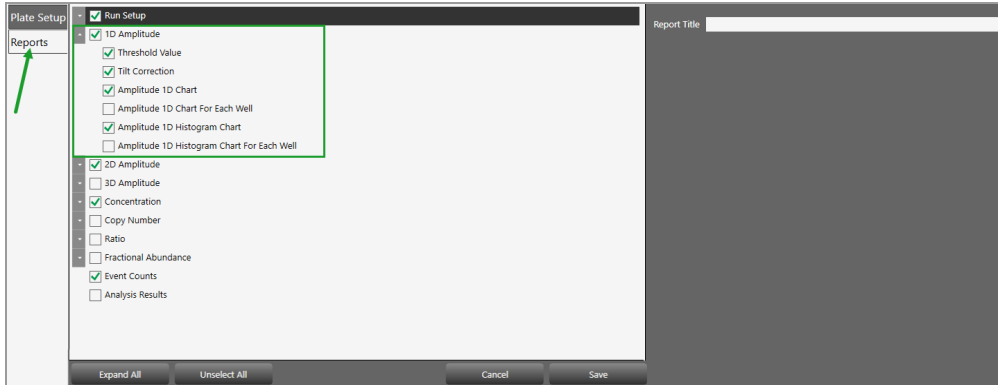

4. Select and clear checkboxes to create the report structure.


Note: Use the buttons at the bottom of the screen to expand or collapse all categories at once, or to select or clear all checkboxes at once. You can also drag and drop the section to rearrange the order of appearance in the report, and drag and drop items in the section to reorder them.



5. Enter a name for your report in the Report Title field and click Save.



6. When the Save dialog box appears, enter a file name.

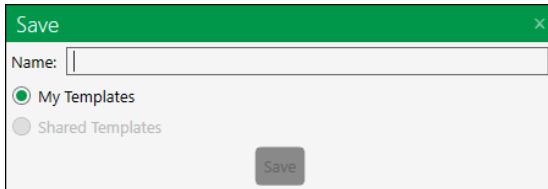

7. Select a storage location. Choose from the following:
 - System Templates (preferred location, if specified by the administrator)
 - My Templates (personal storage location specified in user preferences)
 - Shared Templates (files saved to this location are available to all users)
8. Click Save again to save the report format as a template.


This page intentionally left blank.

Chapter 7 Data Analysis Module Overview

As the instrument reads droplets in each well, the QX600 Droplet Reader collects raw data. At the end of each run, QX Manager Software uses the raw data to produce a data file, which contains analysis information based on the number of positive and negative droplets identified for each target in each sample. The fraction of positive droplets is fit to a Poisson algorithm to determine the starting concentration of the target DNA molecule in units of copies/ μ l input.

This section summarizes the analysis functionality, and also describes the general options governing your chart displays. More detailed information, including specific information and uses for each analysis window, as well as information on data calculations, is contained in [Chapter 8, Data Analysis Methodology](#).

When you select the Data Analysis tab, you can access the Analysis module and the Gene Study module.

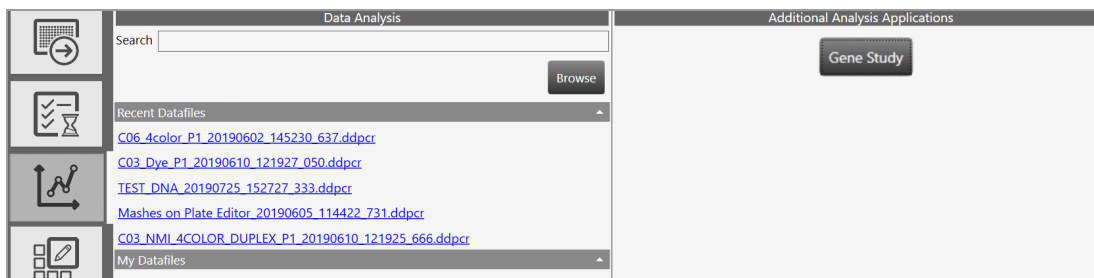

When you open the data file in the Analysis module, you can view the data with the automatic analysis thresholds and calculations, or you can choose different options to recalculate, replot, or recolor the results.

The Data Analysis module functionality allows you to

- View your droplet data in 1D, 2D, and 3D plots
- Analyze multiple targets within a well using amplitude multiplexing and probe mixing strategies
- Detect genomic editing events and non-wild type events with a “drop-off assay” analysis option
- Visualize and export data with improved flexibility
- Generate reports for different analysis types
- The Gene Study module allows you to
- Collect data by adding one or more data files
- View gene expression calculations across multiple experiments

Important: The Gene Study module is accessible from a standalone computer only. The maximum number of samples you can analyze in a gene study is limited by the size of the RAM and virtual memory of the standalone computer.

Opening a Data File for Analysis

You can open the following file types in the Data Analysis module:

- .qlp
- .ddpcr

To open data files, you must be on a computer on which QX Manager Software is installed and have access to the storage locations.

To open a data file for analysis

1. In QX Manager Software, select the Data Analysis tab.



Available data files are displayed under the following headings.

- **Recent Datafiles**

Displays a list of files you have recently opened.

- **System Datafiles or My Datafiles**

Files are always saved to one of the following locations:


- If preferred locations are enabled, then files for all users are saved there and they appear under System Datafiles.
- If preferred locations are not enabled for all users, your files are saved to your personal folder and appear under My Datafiles.

- **Shared Datafiles**

You can also save your data files to a shared location; these appear under Shared Datafiles.

- Click a file to open it in the Analysis module.

Tabs for each analysis and data window appear on the left of the Analysis window.

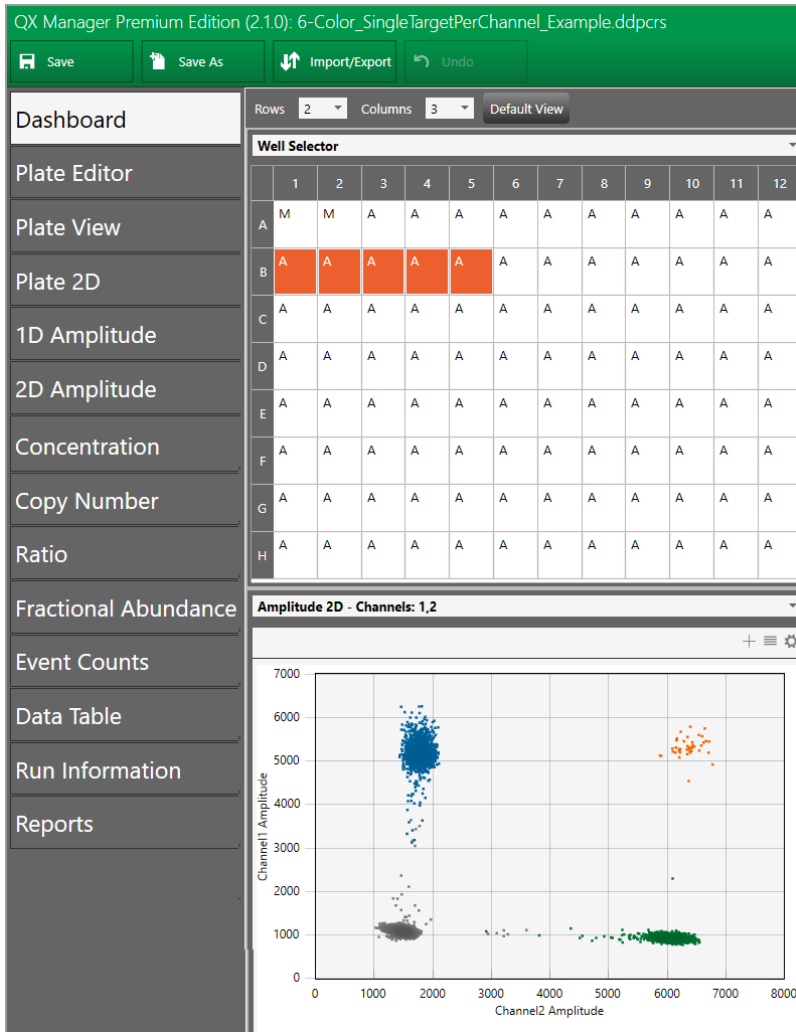

Tip: From the computer attached to the QX600 Droplet Reader, you can also open the data file from the Run Status window. In the completed files section, click the completed run icon (✔) for the file, and then click Launch Analysis.

Saving a Data File

You can save data files with or without modifications, with the same or different file names, and in the same or different locations.



Important: When you save a .qlp.qlps file type in QX Manager Software, the software automatically saves the file as a .ddpcr file. This creates a new file and does not overwrite the original.

To save a data file

1. Do one of the following:
 - To save the file under the same file name and in the same location, click Save.
 - To save the file under a different file name or in a different location, click Save as, and then do one of the following:
 - If you are saving the file under the same name and in the same location, QX Manager Software saves the file immediately.
 - If you are saving the file under a different file name or in a different location, navigate to the new location (if applicable) and enter the new file name. You must be assigned the Overwrite existing file name user privilege to change the file name.
2. Click Save.

Analysis Display and Output Options

QX Manager Software provides the following option that you can choose or set:

- **Universal Options** — available from the primary toolbar to save, import, or export data files
- **Chart Display Preferences** — to change how your chart appears in the analysis window
- **Chart Scale Options** — to change the scale represented on an axis
- **Chart Menu Options** — available in chart views (from the  icon) to copy, save, and print chart images. In Amplitude charts showing data points, you can show or hide
- **Table Menu Options** — available in table views (from the  icon) to copy data, export data to Excel or CSV files, and show or hide columns

You can also expand the chart display to full-screen size and then return to the original display.

Universal Options

Universal options appear on the top menu bar, and are available from all windows, although functionality within the main option might differ according to the analysis window that appears.



Universal options are explained in [Table 22](#), and are available in all Analysis module windows.

Table 22. Universal options

Option	Description
Save	Saves changes to the analysis file under its current file name or file extension
Save as	Saves the file with a different name If you open a file in the latest application that was created in an earlier application, the file assumes the file extension of the new application. See Saving a Data File on page 89 for information.
Import/export	<ul style="list-style-type: none"> ■ Imports or exports plate setup data to a CSV file ■ Imports or exports plate setup data to a plate template file ■ Exports visible data to a CSV file ■ Exports amplitude, cluster, and linkage data ■ Exports and prints the visible charts for the 2D plate layout
Undo	Enabled in the Plate Editor only, after you have made a change
User Settings	Allows you to restore your user preferences to the default settings and to show or hide color matrix change messages

To import or export setup or analysis data

- ▶ Click Import/Export.

[Table 23](#) describes the available import/export options.

Note: If there is unsaved data, the software prompts you to save the file before its contents can be imported or exported.

Table 23. Import/export options

Import/export option	Description
Import plate setup CSV	Import plate layout data from a CSV file to the Plate Editor Note: This option is only available in the Plate Editor. The new plate layout overwrites the existing layout.
Export plate setup CSV	Export a configured plate layout to a CSV file
Import plate setup	Import a plate layout from a plate template (.ddplt) file Note: This option is only available in the Plate Editor.
Export plate setup	Export the plate layout to a plate template file Note: This option is only available in the Plate Editor.
Export visible data to CSV	Export only the data from the wells you selected in the Well Selector to a CSV file
Export amplitude data	Export the data to an Excel spreadsheet, and indicates whether or not the target was found (1 = yes, 0 = no)
Export cluster data	Export the data to an Excel spreadsheet Note: Includes (for each well) whether or not the target was found (1 = yes, 0 = no), the droplet count, and the mean and standard deviation for each dye, and the cluster ID.
Export linkage data	Exports to a CSV file the number of copies of a target that are present in overabundance, versus. the expected value in copies per μL
Export visible charts	Exports all charts that are displayed
Print visible charts	Prints all charts that are displayed

Table 24 shows where the options are available.

Table 24. Option availability

Import/export option	Plate editor	Plate view	Plate 2D	Ampl plots	Data visual charts	Event counts	Data table
Import plate setup from CSV	✓						
Export plate setup to CSV	✓						
Import plate setup	✓						
Export plate setup	✓						
Export visible data to CSV				✓	✓	✓	✓
Export amplitude data	✓	✓	✓	✓	✓	✓	✓
Export cluster data	✓	✓	✓	✓	✓	✓	✓
Export linkage data	✓	✓	✓	✓	✓	✓	✓
Export visible charts			✓	✓	✓	✓	
Print visible charts			✓	✓	✓	✓	

Note: Amplitude plots are available in 1D, 2D and 3D, and Data Visualization charts are Concentration, Copy Number, Ratio, and Fractional Abundance. In the Dashboard view corresponding options are available for each plot or chart displayed.


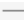
Expanding or Reducing the Chart

You can expand or reduce the chart size in the following windows:

- Dashboard
- 1D Amplitude
- 2D Amplitude
- 3D Amplitude
- Concentration
- Copy Number
- Ratio
- Fractional Abundance
- Event Counts

If you expand an amplitude or data visualization chart to full size, the Well Selector and the Well Data table are hidden from view until you collapse the chart again.

To expand or reduce the chart size

1. To increase the chart size to the full window, click the  icon in the upper-right corner of the chart.
2. To reduce the window to its original size, click the  icon.

Selecting Chart Display Preferences

The Chart Display preferences primarily control font sizes and positioning. [Table 25](#) explains each preference, as well as the windows in which it is available.

Table 25. Chart display preferences

Preference	Window	Description
Channel or Dye	Plate 2D 1D Amplitude 2D Amplitude 3D Amplitude	Select Channel or Dye to identify your chart accordingly <ul style="list-style-type: none"> ■ If you select Channel, the channel number appears as the chart title ■ If you select Dye, FAM, HEX, Cy5, C5.5, ROX, or ATTO 590 appears as the chart title
Axis Data Label Size	1D Amplitude 2D Amplitude 3D Amplitude Concentration Copy Number Ratio Fractional Abundance Event Counts	Select a different font size
Chart Data Label Size	Concentration Copy Number Ratio Fractional Abundance Event Counts	Select a different font size
Chart Marker Size	Concentration Copy Number Ratio Fractional Abundance	Select a different marker size
Chart Marker Label Position	Concentration Copy Number Ratio Fractional Abundance	Reposition the marker label

1. Click the Chart Display button.

To select chart display preferences


1. Click the Chart Display button.
2. Select your options.

The display changes immediately to reflect your choices.

3. Do one of the following:
 - To save your choices as your default settings, click Save as my default. When the success message appears, click OK to close the dialog.
 - To close the menu without saving the settings, click anywhere in the chart display.

Selecting Chart Scale and Label Options

Chart scale options primarily control chart scale and label displays.

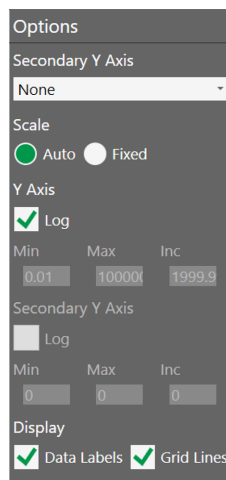


Each preference, as well as the windows in which it is available, is explained in [Table 26 on page 97](#).

Table 26. Chart scale options

Option	Description
Scale	<p>Appears in options for all charts</p> <p>The Scale option determines the numerical range for your chart data:</p> <p>Auto — Selected by default; QX Manager Software automatically displays a scale to match the data distribution.</p> <p>Fixed — Selecting Fixed enables the Axis fields, and allows you to manually adjust the range and increment.</p>
Axis	<p>Appears in options for all charts</p> <p>When the Axis fields are enabled, you can manually adjust the minimum (Min) and maximum (Max) points for the data display on each axis, as well as the increments (Incr).</p> <p>X — Horizontal axis Y — Vertical axis</p> <p>Note: You can modify the x axis scale in the 2D Amplitude chart only.</p>
Secondary Y Axis	<p>Appears in options for Concentration only</p> <p>From the dropdown list, you can select from Channel 2 concentration, ratio, copy number, or fractional abundance to display the corresponding data in a secondary row on the y axis.</p>
Log	<p>Appears in options for all Data Visualization charts (Concentration, Copy Number, Ratio, and Fractional Abundance)</p> <p>You can select the Log checkbox to display the y axis in base 10 logarithmic format.</p> <p>Note: In the Concentration chart, you can also select Log for the secondary y axis.</p>
Inverse	<p>Appears in options for Ratio and Fractional Abundance only</p> <p>The Inverse option equals the opposing number of the Ratio data point to a total of 2, and the opposing number of the Fractional Abundance data point to a total of 100.</p>
Data labels	<p>Appears in options for all Data Visualization charts</p> <p>A data label appears by default for each point in the chart, but you can clear the checkbox to remove the labels.</p>
Grid Lines	<p>Appears in options for all charts</p> <p>Gridlines appear by default. You can clear the check box to remove them.</p>

To select chart scale and label options

1. Click the  icon in the upper-left corner of the chart to display the corresponding chart options.
2. To manually define the chart scale, select Fixed.
Note: Auto is the default setting.
3. In the Min, Max, and Inc fields, enter your lower and upper limits, and the increment, for the y axis. For 2D Amplitude, you can also change the scale on the x axis.
4. For Concentration, Copy Number, Ratio, Fractional Abundance, and Event Counts, select an option for a secondary y axis display. To change to a base 10 logarithmic scale on the y axis, select the Log checkbox.
5. For Ratio and Fractional Abundance, to view an inverse calculation, select the Inverse checkbox.
6. To remove the data labels or grid lines from the display, clear the Data Labels or Grid Lines checkbox.

Using Chart Menu Options

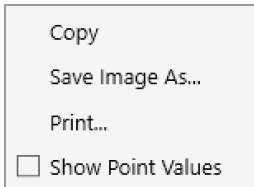


Chart menu options are explained in [Table 27](#).

Table 27. Chart menu options

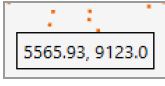
Option	Description
Copy	Copies the chart to the clipboard, which you can paste into a separate document, such as an MS Word file
Save Image as	Saves the chart as an image file
Print	Prints the image
Show Point Values	Shows corresponding point values when you pause on a data point  Tip: Select the Show Point Values checkbox to see the data. Data is available for 1D and 2D amplitude only.

Chart menu options are available in the following windows:

- 1D Amplitude
- 2D Amplitude
- 3D Amplitude
- Concentration
- Copy Number
- Ratio
- Fractional Abundance
- Event Counts


To use the chart menu options

1. Select one or more wells in the Well Selector.

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E	A	A	A	A	A	A						
F	A	A	A	A	A	A						
G	A	A	A	A	A	A						
H	A	A	A	A	A	A						

Corresponding data appears in the Well Data table and the chart.

Well	Target	Conc(copies/μL)	Status	Experiment	SampleType
F01	1	534	OK	DQ	Unknown
F01	2	531	OK	DQ	Unknown
F01	3	539	OK	DQ	Unknown
F01	4	551	OK	DQ	Unknown
F02	1	551	OK	DQ	Unknown
F02	2	551	OK	DQ	Unknown
F02	3	550	OK	DQ	Unknown

2. In the upper-right corner of the chart, click the menu icon () and select from the dropdown list.
 - Select Copy to copy the chart to the Clipboard. To paste the data, navigate to the location (such as a Word file), right-click and select Paste.
 - Select Save Image As and follow the prompts to save the file.
 - Select Print and follow the prompts to print the image.
3. Click anywhere in the display to close the dialog box.

Using Table Menu Options

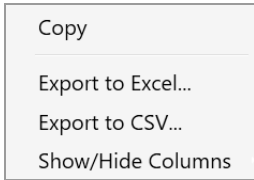


Table menu options are explained in [Table 28](#).

Table 28. Table options

Option	Description
Copy	Copies the information from the selected wells to the clipboard
Export to Excel	Exports the data to an .xlsx file
Export to CSV	Exports the data to a .csv file
Show/Hide Columns	<p>Add or remove columns from a table display</p> <p>When you select Show/Hide Columns, the following dialog box opens. Select or clear check boxes to show or hide columns in the data table.</p>

Table menu options are available in the following windows:

- 1D Amplitude
- 2D Amplitude
- 3D Amplitude
- Concentration
- Copy Number
- Ratio
- Fractional Abundance
- Event Counts

Tip: You can access the menu icon from the individual windows, or from the windows in the Dashboard display.

To use the table menu options

If you are using these options in the Data Table, skip to Step 3.

1. Select wells in the Well Selector.

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E	A	A	A	A	A	A						
F	A	A	A	A	A	A						
G	A	A	A	A	A	A						
H	A	A	A	A	A	A						

Corresponding data appears in the Well Data table.

Well Data							+	☰
	Well	Target	Conc(copies/μL)	Status	Experiment	SampleType		
▾	F01	1	534	OK	DQ	Unknown		▲
	F01	2	531	OK	DQ	Unknown		
	F01	3	539	OK	DQ	Unknown		
	F01	4	551	OK	DQ	Unknown		
	F02	1	551	OK	DQ	Unknown		
	F02	2	551	OK	DQ	Unknown		
	F02	3	550	OK	DQ	Unknown		▼


2. In the upper-right corner of the table, click the menu icon (☰) and select from the dropdown list.
 - To copy the data from the selected wells to the Clipboard, select Copy.

Note: To paste the data, navigate to the location (for example, an Excel file) and then right-click and select Paste.
 - To export your data, select Export to Excel or Export to CSV, and then follow the prompts to save the file.

Note: Exporting to Excel exports the data in spreadsheet format, while Exporting to CSV exports the data in comma-delimited format.
 - To show or hide the corresponding column in the table, select Show/Hide Columns, and then select or clear checkboxes.
3. Click anywhere in the display to close the dialog box.

Analysis Dashboard

When you open a data file, the Analysis module opens to a Dashboard window containing default analysis and data panes in a summary view. You can customize the view to change panes, or increase or decrease the number of panes.

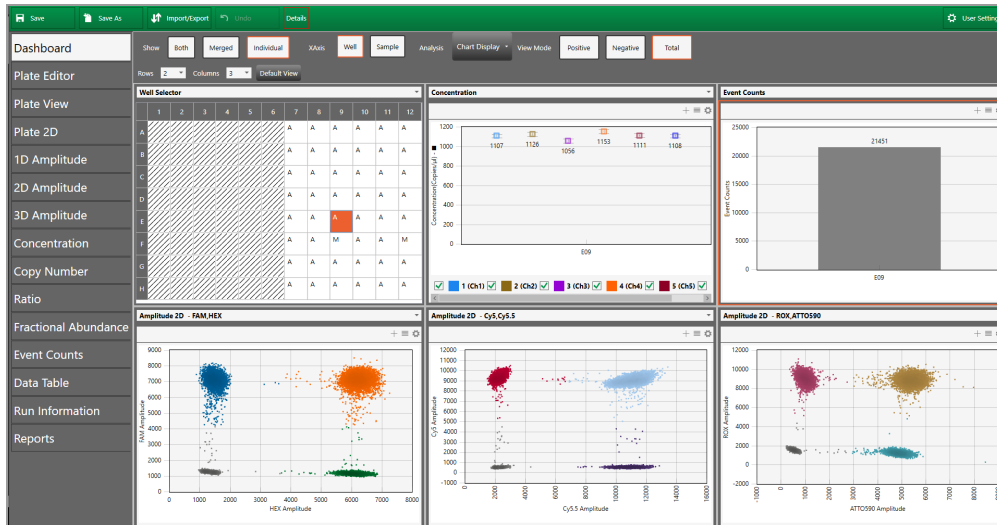

Note: If you open a .qfp file, the Plate Editor opens as the default view and you must select the Dashboard tab to open it.


From the tabs on the left, you can access plate views, amplitude charts, data visualization distribution charts, droplet counts, and a complete data table. You can also view data specific to the run in the Run Information window, and run reports on any of the analysis charts from the Reports window.

Modifying the Dashboard Display


You can modify the column number and row number to display any or all of the analysis windows simultaneously.

To change the number of windows in the display

1. Click the dropdown arrow for Rows and then select a different number.

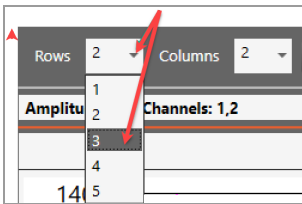


2. Click the dropdown arrow for Columns, and then select a different number.

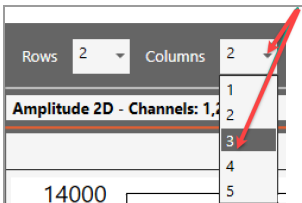

You can also use the dropdown menus to change the windows that appear in the Dashboard.

To display a different window in a particular Dashboard slot


1. Click the dropdown arrow in the upper-right corner to display the list of available windows, and then select a window. The dropdown list closes and the window changes immediately.

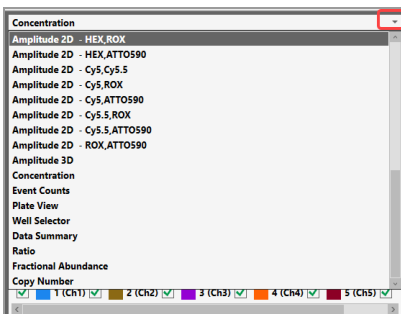


Dashboard Options

From the Dashboard window, depending on the individual dashboard window you have selected, you can use a range of options as described below:

- When you select a window, the options that appear in the menu bar correspond to the options you see if you open the same window from the side menu. For example, if you select the Concentration chart in the Dashboard, the Concentration menu options appear.

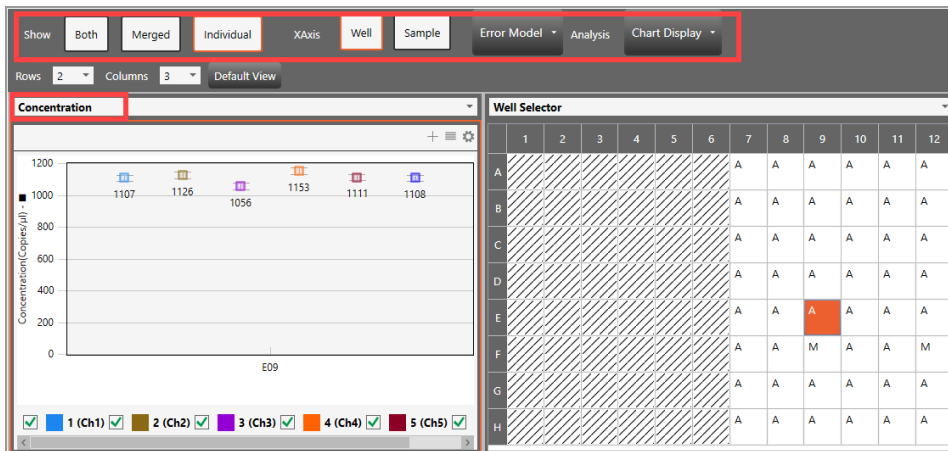

- You can reset the entire dashboard display by adding or removing rows or columns to display more or fewer individual windows. For information, see [Modifying the Dashboard Display on page 104](#).

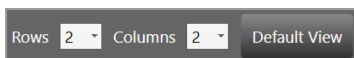

Click Default View to return to the default display of two rows and two columns.

- From any window you can
 - Expand or collapse the window
 - Copy, save, or print the image
 - Select display, scale, axis, and logarithmic options
 - Use the dropdown list to replace the current display with a different display
 - Use the menu bar options for the particular chart type

Table 29 describes the option sets identified in the following graphic.

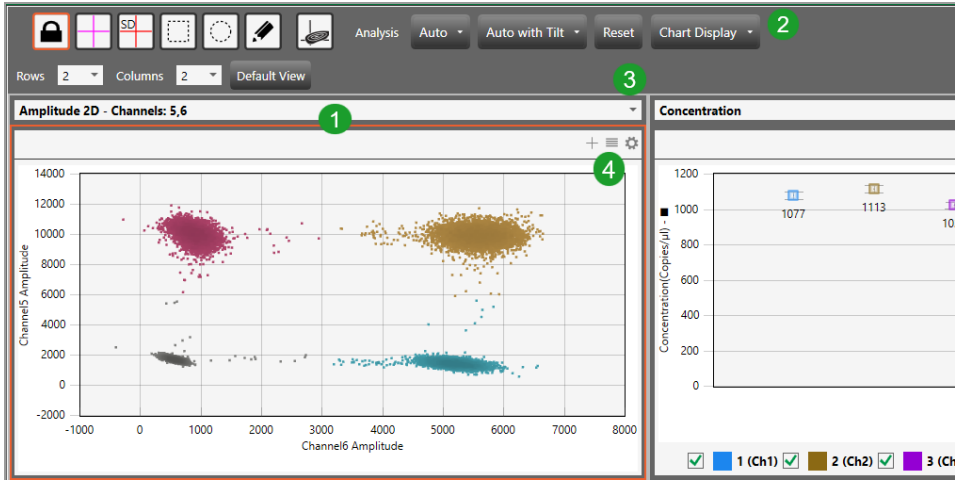


Table 29. Dashboard options

Item	Description
1	Border to identify selected window
2	Toolbar corresponding to the selected window type in the Dashboard
3	Dropdown arrow to replace the current display with a new display
4	Chart options

Plate View Windows

In the QX Manager Software Analysis module you can open the following windows that contain data on all wells in the run:

- **Plate Editor (analysis view)** — Displays the experiment type, sample names, sample types, supermixes, and assay type for each well in the run (similar to the Template Setup Plate Editor)
- **Plate View** — Displays concentration results in a text format for each processed well
- **2D Plate View** — Displays a two-dimensional amplitude chart for each processed well

Plate Editor — Analysis View

You can use the Plate Editor tab in the Analysis module to select wells, and then view or change the experiment parameters. Similar to other Plate Editor displays, the Analysis Plate Editor displays the experiment information for each well in the run.

Wells you select in the Analysis Plate Editor determine the data initially displayed in other analysis windows.

The screenshot displays the 'Plate Editor' window in the QX Manager Software Analysis module. The interface is divided into a main grid and a right-hand configuration panel.

Main Grid: A 96-well plate layout with columns numbered 1-12 and rows labeled A-H. Each well contains a 'DQ 6 Color Cl' label. Well C9 is highlighted in orange.

Right Panel Configuration:

- Clear Selected Wells:** Button
- Experiment Type:** Direct Quantification (DQ) [Apply]
- Sample Descriptions:** Four input fields for Sample Description1-4. Sample Description2 is set to '(Standard)'. [Apply]
- Sample Type:** Unknown [Apply]
- Assay Type:** Single Target per Channel [Apply]
- Target Info Table:**

Target Name	Target Type	Signal CH1	Signal CH2	Signal CH3	Signal CH4	Signal CH5
1	Unkn	FAM	None	None	None	N
2	Unkn	None	HEX	None	None	N
3	Unkn	None	None	Cy5	None	N
4	Unkn	None	None	None	Cy5.5	N
5	Unkn	None	None	None	None	ROX
6	Unkn	None	None	None	None	N
- Well Notes:** Input field
- Plate Notes:** Input field [Apply]

Plate View

The Plate View tab provides, within the plate layout grid, a text view of the calculated concentration for each target in each well processed in the run. For information on the concentration results in chart format, see [Concentration on page 138](#).


	1	2	3	4	5	6	7	8	9	10	11	12
A	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	971.01	959.87	959.87	959.87	959.87	959.87	959.87	959.87	959.87	959.87	959.87	959.87
	1015.5	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7	1007.7
	986.50	989.9	989.9	989.9	989.9	989.9	989.9	989.9	989.9	989.9	989.9	989.9
	1047.7	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4	1038.4
B	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9	1029.9
	1004.5	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3	1006.3
	1045.9	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2	1054.2
	1020.4	991.46	991.46	991.46	991.46	991.46	991.46	991.46	991.46	991.46	991.46	991.46
C	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	1021.5	984.14	984.14	984.14	984.14	984.14	984.14	984.14	984.14	984.14	984.14	984.14
	1020.9	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3	1019.3
	1007.8	964.53	964.53	964.53	964.53	964.53	964.53	964.53	964.53	964.53	964.53	964.53
	1067.8	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5	1033.5
D	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	991.27	1018	1018	1018	1018	1018	1018	1018	1018	1018	1018	1018
	1041.4	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9	1033.9
	978.97	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8	1023.8
	1033.3	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2	1052.2
E	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	941.47	959.99	959.99	959.99	959.99	959.99	959.99	959.99	959.99	959.99	959.99	959.99
	914.91	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9	1015.9
	945.91	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7	1006.7
	988.11	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1	1044.1
F	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ	DQ
	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check	Color Check
	931.9	993.8	993.8	993.8	993.8	993.8	993.8	993.8	993.8	993.8	993.8	993.8
	967.49	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2	1024.2
	957.87	993.69	993.69	993.69	993.69	993.69	993.69	993.69	993.69	993.69	993.69	993.69
	994.92	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9	1069.9

2D Plate View

To facilitate comparisons of amplitude in the processed wells, the Plate 2D tab provides a two-dimensional amplitude view in a separate chart for each well in a single display. To see a larger representation of a well, select the well in the 2D Amplitude view.

To view different color arrays

- ▶ Select the channel combination from the dropdown on the left, and then click Update Grid.

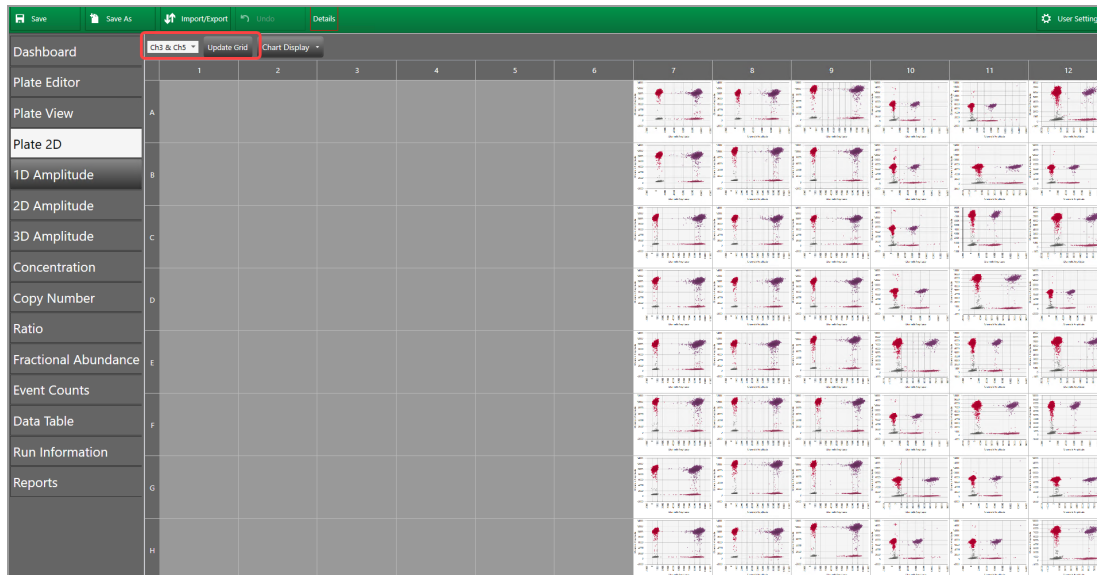

Data Windows

QX Manager Software provides the following table views for your well and run data:

- **Data Table** — available from the main tabs on the left, shows your configured experiment parameters, plus all calculations from the run
 - Note:** You can display the data in individual well format, merged format, or both.
- **Plate Well Data Table** — available in plate setup windows (including the Analysis module Plate Editor), shows only your configuration information for each individual well
- **Analysis Well Data Table** — available in the Analysis module, shows a subset of configuration and run data for the wells selected in the Well Selector

Data Table Tab

The Data Table window contains columns of data generated for each target in each well, as droplet clusters are identified through thresholding or manual clustering tools.

Well	SampleDescr	Target	Conc(copies/μL)	Status	Status Reason	Experiment	SampleType	TargetType	Supermix	DyeName(s)	Copies/20μLWell	TotalC
E01	NTC/1cpd Ch2	HSV-1	No Call	CHECK	Automatic Analysis Unsuccessful	DQ	Unknown	Unknown	ddPCR S...	FAM		
E01	NTC/1cpd Ch2	B2M-...	1050	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	21005	
E02	0.125 cpd Ch...	HSV-1	121	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	2420	
E02	0.125 cpd Ch...	B2M-...	1008	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20158	
E03	0.25 cpd Ch1...	HSV-1	238	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	4768	
E03	0.25 cpd Ch1...	B2M-...	1046	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20920	
E04	0.5 cpd Ch1/...	HSV-1	482	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	9636	
E04	0.5 cpd Ch1/...	B2M-...	1018	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20357	
E05	1 cpd Ch1/1c...	HSV-1	956	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	19127	
E05	1 cpd Ch1/1c...	B2M-...	1020	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20400	
E06	2 cpd Ch1/1c...	HSV-1	1939	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	38786	
E06	2 cpd Ch1/1c...	B2M-...	997	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	19938	
F01	NTC/1cpd Ch2	HSV-1	No Call	CHECK	Automatic Analysis Unsuccessful	DQ	Unknown	Unknown	ddPCR S...	FAM		
F01	NTC/1cpd Ch2	B2M-...	1022	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20435	
F02	0.125 cpd Ch...	HSV-1	126	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	2529	
F02	0.125 cpd Ch...	B2M-...	995	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	19890	
F03	0.25 cpd Ch1...	HSV-1	250	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	4991	
F03	0.25 cpd Ch1...	B2M-...	1008	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20156	
F04	0.5 cpd Ch1/...	HSV-1	475	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	9492	
F04	0.5 cpd Ch1/...	B2M-...	1031	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20621	
F05	1 cpd Ch1/1c...	HSV-1	943	OK		DQ	Unknown	Unknown	ddPCR S...	FAM	18865	
F05	1 cpd Ch1/1c...	B2M-...	1048	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20962	
F06	2 cpd Ch1/1c...	HSV-1	No Call	CHECK	Droplet Count < 10000	DQ	Unknown	Unknown	ddPCR S...	FAM		
F06	2 cpd Ch1/1c...	B2M-...	No Call	CHECK	Droplet Count < 10000	DQ	Unknown	Unknown	ddPCR S...	HEX		
G01	NTC/1cpd Ch2	HSV-1	No Call	CHECK	Automatic Analysis Unsuccessful	DQ	Unknown	Unknown	ddPCR S...	FAM		
G01	NTC/1cpd Ch2	B2M-...	1012	OK		DQ	Unknown	Unknown	ddPCR S...	HEX	20245	

Data Table Column Definitions

Table 30 defines each field in the Data Table.

Table 30. Data Table columns

Column name	Content description
Well	The well location in the plate Note: The instrument accepts 96-well plates only.
Sample description 1 Sample description 2 Sample description 3 Sample description 4	Words or phrases to describe the sample in each well Note: Only the sample description 1 data appears in plate layout views but you can pause on a well in the Plate Editor to display a tool tip containing the entries in all completed Sample description fields.
Target	Target name from Plate Editor tab
Conc(copies/ μ L)	Concentration of the target molecules recorded as copies per microliter
pg/ μ L	Mass concentration of DNA in the well in pg/ μ L
Status	Contains either Check, OK, Multi, or Manual <ul style="list-style-type: none"> ■ If Check, either the automatic analysis for the well failed or fewer than 10,000 droplets were counted in the well ■ If OK, enough droplets were counted to trigger automatic analysis ■ If Multi, the data was automatically analyzed as part of a multi-well selection ■ If Manual, the droplets were analyzed manually
Status Reason	Identifies the reason for the status (unsuccessful automatic analysis, not enough droplets counted, or droplets analyzed manually)
Experiment	Experiment type selected for the well in the Plate Editor
SampleType	Sample type selected for the well in the Plate Editor
TargetType	Target type selected for each dye (FAM or EvaGreen®, HEX or VIC, Cy5, Cy5.5, ROX, and ATTO 590) in the Plate Editor

Table 30. Data Table columns, continued

Column name	Content description
Supermix	Supermix selected for the well in Plate Editor
DyeName(s)	Dye assigned to channels 1, 2, 3, 4, 5, and 6, respectively, for the well in the Plate Editor
Copies/20 μ LWell	Concentration of the target normalized to a volume of 20 μ L
TotalConfMax	For merged wells the high error bar for the target concentration of the combined wells at a 95% confidence interval
TotalConfMin	For merged wells the low error bar for the target concentration of the combined wells at a 95% confidence interval
PoissonConfMax	Maximum target concentration normalized for the high error bar of the droplet Poisson distribution for the 95% confidence interval
PoissonConfMin	Minimum target concentration normalized for the low error bar of the droplet Poisson distribution for the 95% confidence interval
Accepted Droplets	Total number of droplets accepted by quality algorithm
Positives	Number of droplets that contain the target
Negatives	Number of droplets that do not contain the target
Copies/ μ l linked molecules	Shows number of copies of a target that are present in overabundance versus the expected value in copies per μ L
CNV	Copy number calculated for the target relative to the reference
TotalCNVMax	For merged wells, the high error bar for the copy number of the combined wells at a 95% confidence interval
TotalCNVMin	For merged wells, the low error bar for the copy number of the combined wells at a 95% confidence interval
PoissonCNVMax	Maximum copy number normalized for the high error bar of the droplet Poisson distribution for the 95% confidence interval
PoissonCNVMin	Minimum copy number normalized for the low error bar of the droplet Poisson distribution for the 95% confidence interval

Table 30. Data Table columns, continued

Column name	Content description
ReferenceCopies	Copy number identified for the reference target in the Plate Editor Default is 2, indicating 2 copies per diploid genome.
UnknownCopies	This field is currently unused in QX Manager Software
Threshold1	Threshold value of first threshold line from left to right and bottom to top
Threshold2	Threshold value of second threshold line from left to right and bottom to top
Threshold3	Threshold value of third threshold line from left to right and bottom to top
ThresholdSigmaAbove	Distance of the threshold from the mean of the negative cluster, as a multiple of the negative cluster standard deviation
ThresholdSigmaBelow	Distance of the threshold from the mean of the positive cluster, as a multiple of the positive cluster standard deviation
ReferenceUsed	Identifies which target was used as a reference
Ratio	The ratio of the target against the identified reference
TotalRatioMax	For merged wells the high error bar for the ratio of the unknown against the reference of the combined wells at a 95% confidence interval
TotalRatioMin	For merged wells the low error bar for the ratio of the unknown against the reference of the combined wells at a 95% confidence interval
PoissonRatioMax	Maximum ratio of the unknown against the reference normalized for the high error bar of the droplet Poisson distribution for the 95% confidence interval
PoissonRatioMin	Minimum ratio of the unknown against the reference normalized for the low error bar of the droplet Poisson distribution for the 95% confidence interval

Table 30. Data Table columns, continued

Column name	Content description
FractionalAbundance	Calculation of fractional abundance of this unknown target vs. the reference target
TotalFractionalAbundanceMax	For merged wells the high error bar for the fractional abundance of the combined wells at a 95% confidence interval
TotalFractionalAbundanceMin	For merged wells the low error bar for the fractional abundance of the combined wells at a 95% confidence interval
PoissonFractionalAbundanceMax	Maximum fractional abundance normalized for the high error bar of the droplet Poisson distribution for the 95% confidence interval
PoissonFractionalAbundanceMin	Minimum fractional abundance normalized for the low error bar of the droplet Poisson distribution for the 95% confidence interval
MeanAmplitudeOfPositives	Mean amplitude value of all droplets that contain the target
MeanAmplitudeOfNegatives	Mean amplitude value of all droplets that contain no target
MeanAmplitudeTotal	Mean amplitude value of all droplets
ExperimentComments	Data populates this field only from files created in QuantaSoft versions 1.4 – 1.7, if comments were added.
MergedWells	Identifies which wells were merged together
TotalConfidenceMax68	For merged wells the high error bar for the target concentration of the combined wells at a 68% confidence interval
TotalConfidenceMin68	For merged wells the low error bar for the target concentration of the combined wells at a 68% confidence interval
PoissonConfidenceMax68	Maximum target concentration normalized for the high error bar of the droplet Poisson distribution for the 68% confidence interval
PoissonConfidenceMin68	Minimum target concentration normalized for the low error bar of the droplet Poisson distribution for the 68% confidence interval

Table 30. Data Table columns, continued

Column name	Content description
TotalCNVMax68	For merged wells the high error bar for the copy number of the combined wells at a 68% confidence interval
TotalCNVMin68	For merged wells the low error bar for the copy number of the combined wells at a 68% confidence interval
PoissonCNVMax68	Maximum target concentration normalized for the high error bar of the droplet Poisson distribution for the 68% confidence interval
PoissonCNVMin68	Minimum target concentration normalized for the low error bar of the droplet Poisson distribution for the 68% confidence interval
TotalRatioMax68	For merged wells the high error bar for the ratio of the unknown against the reference of the combined wells at a 68% confidence interval
TotalRatioMin68	For merged wells the low error bar for the ratio of the unknown against the reference of the combined wells at a 68% confidence interval
PoissonRatioMax68	Maximum ratio of the unknown against the reference normalized for the high error bar of the droplet Poisson distribution for the 68% confidence interval
PoissonRatioMin68	Minimum ratio of the unknown against the reference normalized for the low error bar of the droplet Poisson distribution for the 68% confidence interval
TotalFractionalAbundanceMax68	For merged wells the high error bar for the fractional abundance of the combined wells at a 68% confidence interval
TotalFractionalAbundanceMin68	For merged wells the low error bar for the fractional abundance of the combined wells at a 68% confidence interval
PoissonFractionalAbundanceMax68	Maximum fractional abundance normalized for the high error bar of the droplet Poisson distribution for the 68% confidence interval
PoissonFractionalAbundanceMin68	Minimum fractional abundance normalized for the low error bar of the droplet Poisson distribution for the 68% confidence interval plate list after correcting the issue

Table 30. Data Table columns, continued

Column name	Content description
TiltCorrected	Indicates if tilt correction was applied to the well Note: Tilt correction rotates the angles from the double-negative cluster centroid to each droplet, such that the two single positive cluster centroids are orthogonal with the double-negative cluster centroid.
Ch1+Ch2+	Number of droplets that contain both channel 1 and channel 2 targets
Ch1+Ch2-	Number of droplets that contain just the channel 1 target
Ch1-Ch2+	Number of droplets that contain just the channel 2 target
Ch1-Ch2-	Number of droplets that contain neither channel 1 nor channel 2 targets
Ch3+Ch4+	Number of droplets that contain both channel 3 and channel 4 targets
Ch3+Ch4-	Number of droplets that contain just the channel 3 target
Ch3-Ch4+	Number of droplets that contain just the channel 4 target
Ch3-Ch4-	Number of droplets that contain neither channel 3 nor channel 4 targets
Ch5+Ch6-	Number of droplets that contain just the channel 5 target
Ch5-Ch6+	Number of droplets that contain just the channel 6 target
Ch5-Ch6-	Number of droplets that contain neither channel 5 nor channel 6 targets
Ch5+Ch6+	Number of droplets that contain both channel 5 and channel 6 targets

Plate Editor Well Data Table

A well data table containing both configuration and run data appears in the analysis windows, next to the Well Selector. This table displays a subset of data columns for only the wells selected in the Plate Editor or Well Selector. [Table 31](#) defines each column in the consolidated table.

Table 31. Fields in the Analysis Well Data table

Column name	Content description
Well	The well location in the plate the sample is from
Perform Droplet Reading	Yes or No
Experiment Type	Experiment type from the Plate Editor
Sample description 1 – 4	The sample name used to identify the well in the Plate Editor tab
SampleType	Sample type from the Plate Editor
Supermix Name	Supermix from the Plate Editor
Assay type	Assay type selected in the Plate Editor
Target Name	Target Name from Plate Editor tab
TargetType	Target type assigned to the target in the Plate Editor
Signal Ch1	Selected fluorophore in Channel 1
Signal Ch2	Selected fluorophore in Channel 2
Signal Ch3	Selected fluorophore in Channel 3
Signal Ch4	Selected fluorophore in Channel 4
Signal Ch5	Selected fluorophore in Channel 5
Signal Ch6	Selected fluorophore in Channel 6
Reference Copies	Copy number identified for the reference target in the Plate Editor tab
Well Notes	Entries by the user
Plot?	If selected, target will be used to calculate the copy number for the well
RDQ Conversion Factor	The Residual DNA Quantification conversion factor for the well Note: RDQ is only available in Residual DNA Quantification experiments.

Analysis Windows

In its Analysis module, QX Manager Software provides amplitude displays and data visualization charts in the Analysis module. For more information on each chart, see [Chapter 8, Data Analysis Methodology](#).

Amplitude Charts

Amplitude charts illustrate the droplets by degree of dye fluorescence. QX Manager Software displays fluorescence for positive and negative targets identified with FAM/EvaGreen®, HEX/VIC, Cy5, Cy5.5, ROX, and ATTO 590 fluorophores in the following analysis windows:

- 1D Amplitude
- 2D Amplitude
- 3D Amplitude

Data Visualization Charts

Data visualization charts show calculations based on positive droplet concentrations. QX Manager Software counts the number of positive and negative droplets for each fluorophore in each sample, and then fits the fraction of positive droplets to a Poisson algorithm to determine the starting concentration of the target DNA molecule in units of copies/μl input in the following analysis windows:

- Concentration
- Copy Number
- Ratio
- Fractional Abundance

When you open the analysis data file in QX Manager Software and select one or more wells in the grid, the software plots each calculated Poisson distribution average as a data point in the grid. You can pause on the data point to see the high, median, and low numbers.

Note: You can display the data using a confidence level of either 68% or 95% (default setting).

Event Counts

Event counts record the number of droplets in which the target was found (positive) and was not found (negative). The Event Counts tab shows a bar chart view of the number of positive, negative, and total droplets for each of the targets in each of the selected wells.

Viewing Run Information

Use the Run Information window to view the information about the plate and run. You can also view or add post-run notes regarding the plate. .

Run Details	
Instrument Type:	CX300
Instrument Serial Number:	867-5309
Plate Name:	DQ_Test
Date of Run:	9/2/2021 10:46:18 AM
Run Started by:	USPLC01K1YFZ/QX_Super
Supermix Type:	ddPCR Supermix for Probes
Droplet Size:	Standard
Number Of Wells Read:	8
Oil Name:	Probes
DR Elapsed Time:	0h 0m 31s
Run Status:	Success
Software Version:	2.0.0
MCU Version:	8
DLL Version:	2.0.0.462
FPGA Version:	16
Fluidics Version:	3
Motion Version:	10
System Version:	7
Calibration State:	OK

Post Run Plate Notes:
N/A

This page intentionally left blank.

Chapter 8 Data Analysis Methodology

As summarized in [Chapter 7, Data Analysis Module Overview](#), the software presents analysis data in several different formats. When you open the data file in the Analysis module, you can view the data with the automatic analysis thresholds and calculations or you can choose different options to recalculate, replot, or recolor the results.

Data is shown in the following formats:

- Amplitude, which shows each droplet in each well in a scatter plot, in views
 - Data visualizations, which show estimated target molecule concentration, ratio, and fractional abundance of one target against another target based on selected error model and confidence interval
- Note:** Copy number variation is also calculated if the user ran a CNV experiment.
- Event counts, which are precise counts of positive and negative droplets for each target in each well

Amplitude Plot Analysis Options

QX Manager Software provides three amplitude plot displays in which you can view the data with automatic thresholding where applicable, or choose different options to replot or recolor the results.

In the 1D amplitude and 2D amplitude plots, you can

- Show well calculations

Note: The 1D amplitude plot shows the data for each selected well in its own section, while the 2D amplitude plot shows all selected wells together.

- Modify automatic thresholds calculated by QX Manager Software
- View data in a heat map
- View histograms (1D amplitude tab only)
- Re-apply automatic analysis
- Apply tilt correction to automatic analysis

In the 3D amplitude plot, you can

- Rotate the plot to gain a better view of the droplets in 3D orientation
- View thresholds previously set in 1D or 2D amplitude display
- View 2D data behind 3D droplets in clusters or heat maps

Table 32 explains the analysis options that are available in both 1D amplitude and 2D amplitude plots.

Table 32. 1D and 2D amplitude tool options




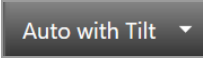

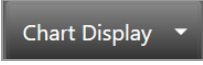
Icon	Mode	Description
	View mode	Data appears with the thresholds or clusters assigned by QX Manager Software. When this icon is selected, the view is locked and you cannot modify the thresholds.
	Heat map	Shows the distribution of droplets, with higher density of droplets shown as hot.
	Automatic analysis	Disables tilt correction and automatically reanalyzes the data.
	Automatic analysis with tilt correction	Enables tilt correction and automatically reanalyzes the data. Note: Tilt correction rotates the angles from the double-negative cluster centroid to each droplet, such that the two single positive cluster centroids are orthogonal with the double-negative cluster centroid.
	Reset	Resets the data to the original automatically analyzed data.
	Chart display options	Choose whether to display axis with channel number or fluorophore, and change the size of axis markers, data labels, and axis labels.

Table 33 explains the analysis options available only for 1D amplitude plots.

Table 33. 1D amplitude tool options











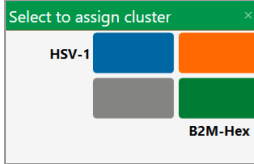
Icon	Mode	Description
	Automatic thresholding	<p>Individual Wells: Clusters are generated for each selected well separately.</p> <p>Combined Wells: Droplets from all selected wells are combined, and clusters are generated on that combined set.</p> <p>Positive Control Wells: Thresholds are generated for each channel based on any positive controls currently selected, and applies those thresholds to all currently selected wells.</p> <p>Note: If you change the thresholds manually, you can click Reset to return to the automatic thresholds.</p>
	Threshold, single well	<p>For a single well or multiple wells, the user can manually input thresholding values in the 1D amplitude plot.</p> <p>You can click anywhere in the plot to apply threshold lines to the data.</p> <p>To adjust the threshold values</p> <ul style="list-style-type: none"> ■ Drag the lines horizontally in the plot area ■ Enter a new threshold value in the boxes at the end of each threshold line ■ Click the up and down arrows to incrementally adjust the numbers
	Threshold, multiple wells	
	Threshold, single well SD	<p>For a single well or multiple wells, QX Manager Software automatically calculates and populates thresholding values with standard deviations (SD).</p> <p>Important: SD thresholds are not recommended for RMD, since these clusters are not typically located at right angles to each other.</p> <p>Note: The SD thresholds are set at a limit that is less than the mean number of positive targets, and more than the mean number of negative targets. Starting from where the user clicks in the plot, SD locates the threshold between the positive and negative clusters such that the classification error is minimized.</p>
	Threshold, multiple wells SD	
	Show histogram	Hides or displays histogram plots for 1D amplitude

Table 34 explains the analysis options available only for 2D amplitude plots.

Table 34. 2D amplitude tool options


Icon	Mode	Description
	Automatic thresholding	<p>Individual Wells: Clusters are generated for each selected well separately.</p> <p>Combined Wells: Droplets from all selected wells are combined, and clusters are generated on that combined set.</p> <p>Positive Control Wells: The software generates thresholds for each channel based on any positive controls currently selected, and applies those thresholds to all currently selected wells.</p> <p>Note: If you change the thresholds manually, you can click Reset to return to the automatic thresholds.</p>
	Threshold line	<p>Use the Threshold line icons to establish thresholds across a pair of channels, dividing droplets into distinct clusters separated by different colors.</p> <p>You can click anywhere in the plot to apply threshold lines to the data.</p> <p>To adjust the threshold values, you can</p> <ul style="list-style-type: none"> ■ Drag the lines horizontally or vertically in the plot area ■ Enter a new threshold value in the boxes at the end of each threshold line ■ Click the up and down arrows to incrementally adjust the numbers
	Threshold line SD	<p>QX Manager Software automatically calculates and populates threshold line values with standard deviations (SD).</p> <p>Important: SD thresholds are not recommended for RMD, since these clusters are not typically located at right angles to each other.</p> <p>Note: The SD thresholds are set at a limit that is less than the mean number of positive targets, and more than the mean number of negative targets. Starting from where the user clicks in the plot, SD locates the threshold between the positive and negative clusters such that the classification error is minimized.</p>

Icon	Mode	Description
	Threshold cluster square	Use the Threshold Cluster icons to establish thresholds across a pair of channels, dividing droplets into distinct clusters separated by different colors.
	Threshold cluster circle	To classify your droplets, you can
	Threshold cluster freeform	<ul style="list-style-type: none"> ■ Click any of the three Threshold Cluster Mode buttons (square, circle, or free form) ■ Draw the chosen shape around the cluster of droplets <p>Note: When using the cluster mode tools to draw a shape around a cluster of droplets, a dialog box with color selections by target appears in the window to help you select the correct target combination.</p>  <p>Target names are shown along the two axes based on the signal values assigned during plate setup. Click the colored box that best represents the location of the cluster of droplets selected to automatically classify the droplets as positive or negative for each given target.</p> <p>Note: The Cluster dialog box display varies according to the plex mode.</p>

1D Amplitude


Use the 1D Amplitude analysis window to view droplet amplitudes for each channel, in selection order, for each well selected in the Well Selector.

The following graphic shows six colors in standard-size droplets.

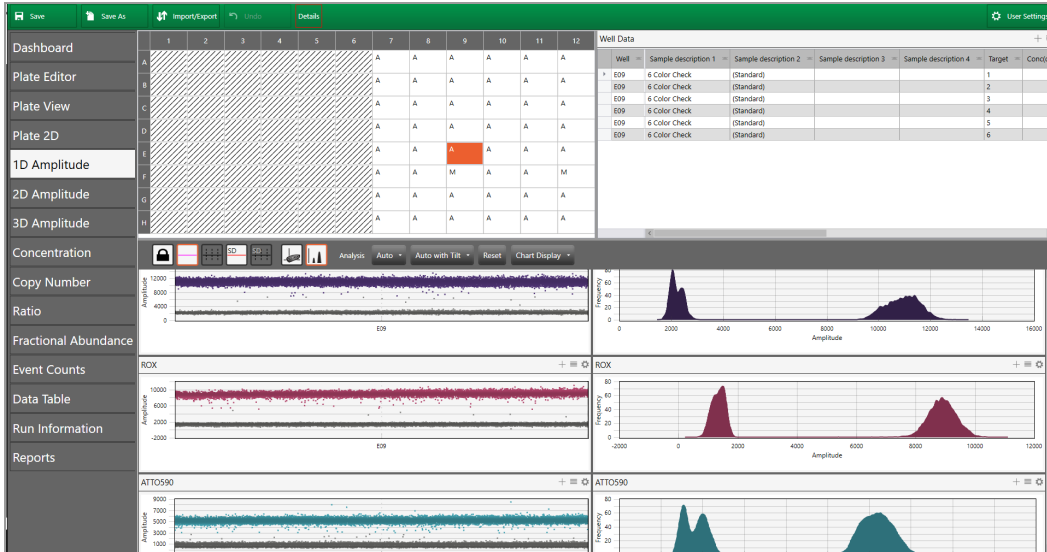


The 1D Amplitude plots illustrate

- All accepted droplets, with a visual representation of the positive and negative droplets in each well
 - Positive droplets appear in color, while negative droplets appear in gray.
 - The Well Selector displays an A in each well where QX Manager Software automatically thresholded the droplet data. An M appears in any well in which you manually changed the threshold.
- Amplitude in a histogram plot

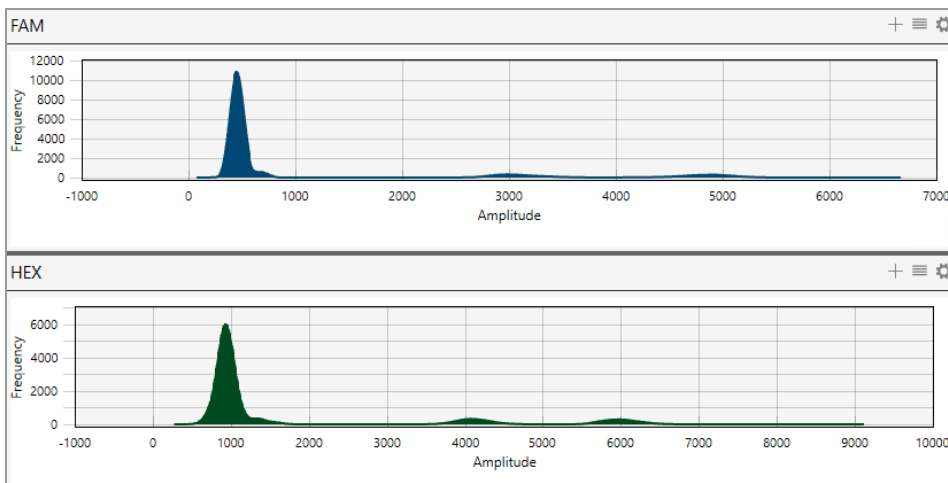


Note: The Histogram icon is enabled by default so a Histogram plot appears automatically for each fluorophore (channel).

- Amplitude in a Heat Map view

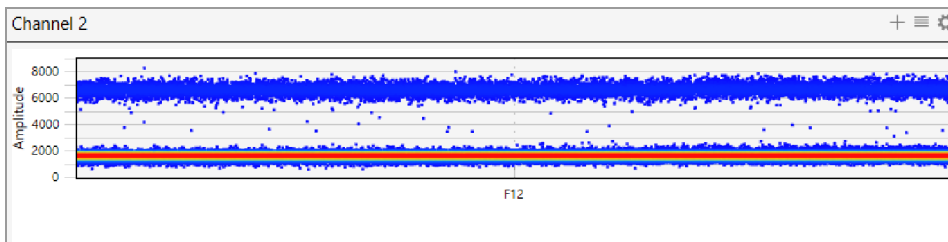

When you select the 1D Amplitude tab, the toolbar options that allow you to change thresholds and show or hide the heat map and histogram appear above the plots.

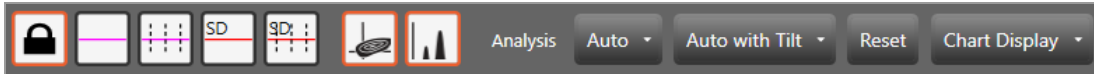


Note: For detailed information on each toolbar option, see [Amplitude Plot Analysis Options on page 122](#).

Default color assignments for 1D amplitude plots are

Positive FAM	=	Blue
Positive HEX	=	Green
Positive Cy5	=	Red
Positive Cy5.5	=	Violet
Positive ROX	=	Maroon
Positive ATTO 590	=	Aqua
Negative	=	Gray


The images in this section display data from a direct quantification experiment involving six targets:


Negative droplets appear in grayscale in each plot.

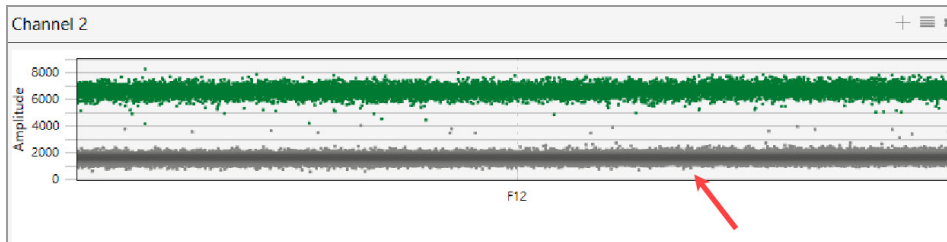


For experiments where enough droplets are recognized, QX Manager Software automatically configures an amplitude threshold, which distinguishes between positive and negative droplets.


Note: Well data must meet certain quality metrics (minimum of 10,000 droplets) before the software will automatically calculate a threshold above which droplets are considered positive. You can use the manual thresholding tools to ensure correct quantification of positive and negative droplets.

By default, the Histogram icon () is enabled so the Histogram plot automatically appears with the amplitude plot.

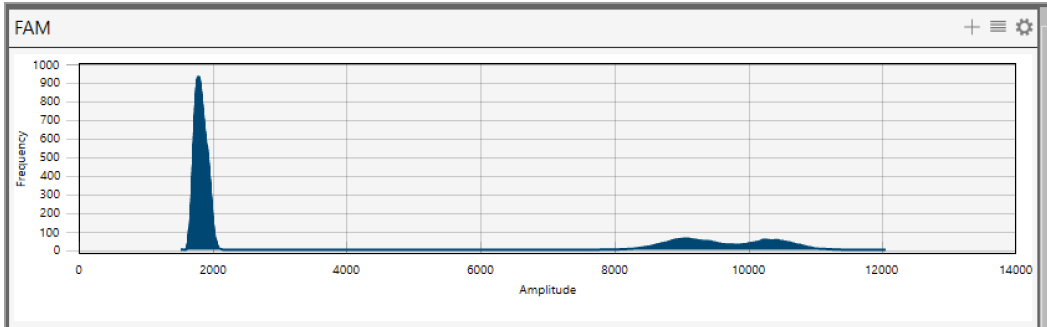

To manually adjust the thresholds

1. Select one or more wells.
2. Click any of the threshold icons () to set manual thresholds:
 - Regular threshold for a single well; when multiple wells are selected, you can set a different threshold for each well
 - Regular threshold for multiple wells, where you can set one threshold for all selected wells
 - Standard deviation threshold for a single well; when multiple wells are selected, you can set a different threshold for each well
 - Standard deviation for multiple wells, where you can set one threshold for all selected wells
3. Click on the location in the plot where you want to draw the new threshold.
4. When the line appears, use the up or down arrows, or click and drag the threshold, to adjust the location.

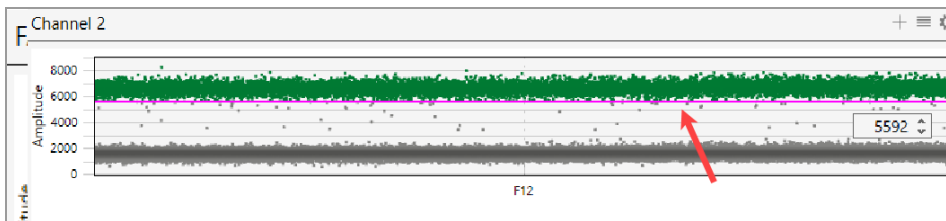

QX Manager Software displays the manual threshold and changes the color of the droplets that are now above or below the threshold.

Note: You can also manually enter a threshold number.


When you manually modify a threshold, the well selector displays an M for manual instead of the A for Automatic.

/ / / / / / / /	A	A	A	A	A	A
/ / / / / / / /	A	A	A	A	A	A
/ / / / / / / /	A	A	A	A	A	M
/ / / / / / / /	A	A	A	A	A	A

5. (Optional) To reset to the automatic threshold, click Auto.
6. (Optional) To adjust the automatic threshold to consider tilt correction, click Auto with Tilt.
7. (Optional) To reset everything to the original display, click Reset.

To view a heat map representation of the plot data


- ▶ Click the heat map () icon.

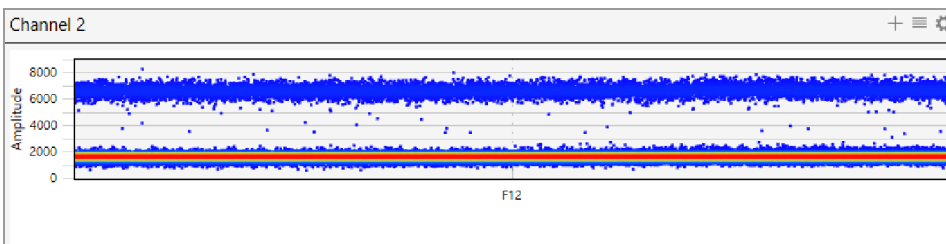

2D Amplitude

For the wells you select in the analysis view of the Plate Editor, the 2D Amplitude window illustrates fluorescence for six targets. The images in this section display data from an experiment involving six targets. Each plot shows individual cluster representations for first and second target positive droplets, droplets containing both targets, and droplets containing neither target.

The following graphic shows six colors in standard-size droplets.

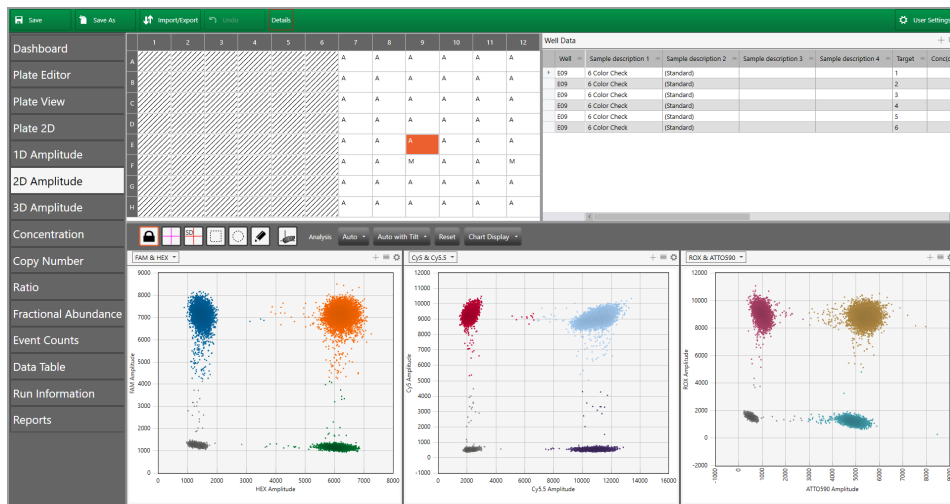

When you select the 2D Amplitude tab, the toolbar options allow you to change thresholds, manually cluster wells, and show or hide the heat map above the plots.



For detailed information on each toolbar option, see [Amplitude Plot Analysis Options on page 122](#).

Default color assignments for 2D amplitude plots are:


Positive FAM	=	Blue
Positive HEX/VIC	=	Green
Positive FAM and HEX/VIC	=	Orange (may appear brown on some monitors)
Positive Cy5	=	Red
Positive Cy5.5	=	Violet
Positive Cy5 and Cy5.5	=	Light Blue
Positive ROX	=	Maroon
Positive ATTO 590	=	Gold
Positive ROX and ATTO 590	=	Aqua
Double negative	=	Gray


Note: Well data must meet certain quality metrics (minimum of 10,000 droplets) before QX Manager Software will automatically calculate a threshold above which droplets are considered positive. You can use the manual thresholding tools to ensure correct quantification of positive and negative droplets.




To use the 2D amplitude tools


1. Open your data file in the Analysis module and select the 2D Amplitude tab.
2. Choose a threshold line or cluster option.

Note: Use the cluster option when the data is not orthogonal and the threshold lines do not adequately divide the populations of droplets.

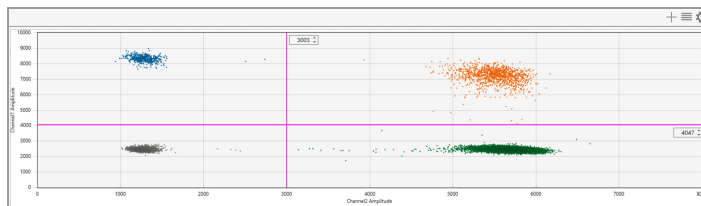- To adjust threshold lines, select a threshold line type () and then click in the plot to apply crosshair lines. Use any of the following methods:
 - Drag the lines horizontally or vertically in the plot area.
 - Enter a new threshold value in the boxes at the end of each threshold line.
 - Tap the up and down arrows to incrementally adjust the numbers.



- To adjust thresholds using the cluster option, select a cluster type (  ) and draw the chosen shape around the cluster of droplets to classify.

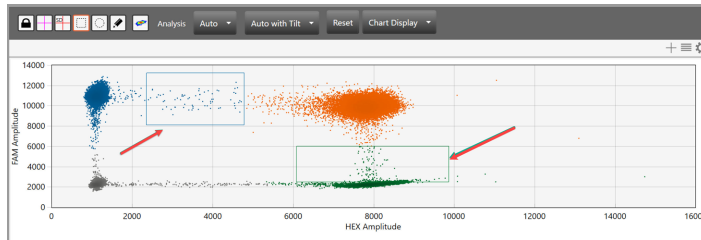

3. (Optional) To reset to the automatic threshold, click Auto.
4. (Optional) To adjust the automatic threshold to consider tilt correction, click Auto with Tilt.
5. (Optional) To reset everything to the original display,click Reset.

Data Visualization Chart Analysis Options

QX Manager Software also provides chart displays where you can view the data from the original calculations, or choose different options to recalculate the results.





In charts reflecting Data Visualization distributions, you can

- Show individual or merged well data
- Change the error model and confidence level

Note: For event counts, you can also alter the view mode to see positive, negative, or all droplet counts.

[Table 35](#) explains the available options.

Table 35. Chart tool options for Data Visualization distribution charts

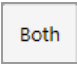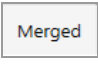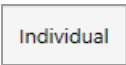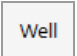Icon	Mode	Description
	Show both	Shows the calculated total or percentage as a merged calculation for all selected wells and as an individual calculation for each selected well
	Show merged	Shows the calculated total or percentage as a combined number for all selected wells with the same parameters Note: Well setup must match for each well before the software can merge them.
	Show individual	Shows the calculated total or percentage as an individual calculation for each selected well
	X-axis well	Shows the well number on the x axis





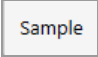

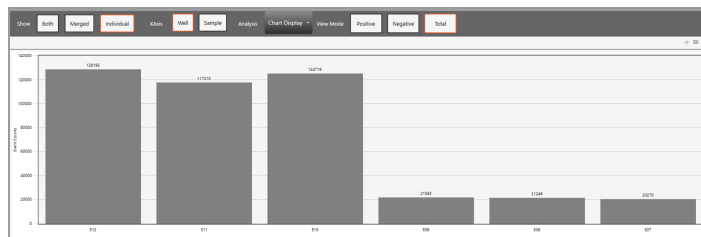
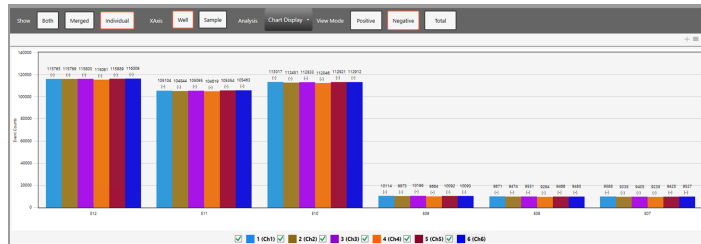
Icon	Mode	Description
	X-axis sample	Shows a partial sample name on the x axis Note: You can select Well and Sample simultaneously.
	Error model and confidence interval	Select an error model and confidence level Note: The default choices are Poisson and 95%, to yield the highest confidence in the results.  Important: The software uses the following multipliers in its calculations: <ul style="list-style-type: none"> ■ For 95% confidence interval, the sigma multiplier is 1.96 ■ For 68% confidence interval, the sigma multiplier is 1.0

Table 36 identifies the additional options available in the Event Count window. These options are available for event counts only.

Table 36. Additional options for event counts


Icon	Mode	Description
	View mode positive	Displays in a bar chart the number of positive droplets for each target in each well
	View mode negative	Displays in a bar chart the number of negative droplets for each target in each well
	View mode total	Displays in a bar chart the number of total droplets in each well

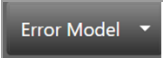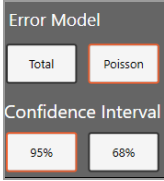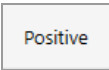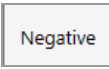


Concentration

Concentration measures the number of copies of the target molecules in each μl of sample based on the default settings. Use the toolbar to merge or separate the calculated results for each well, and to recalculate your results based on a different error model and confidence interval.


For information on toolbar options, see [Data Visualization Chart Analysis Options on page 135](#).

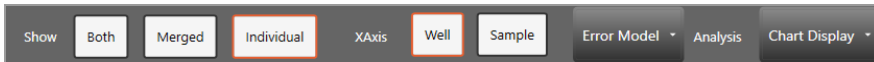

The following graphic shows the concentration for six colors in standard-size droplets.

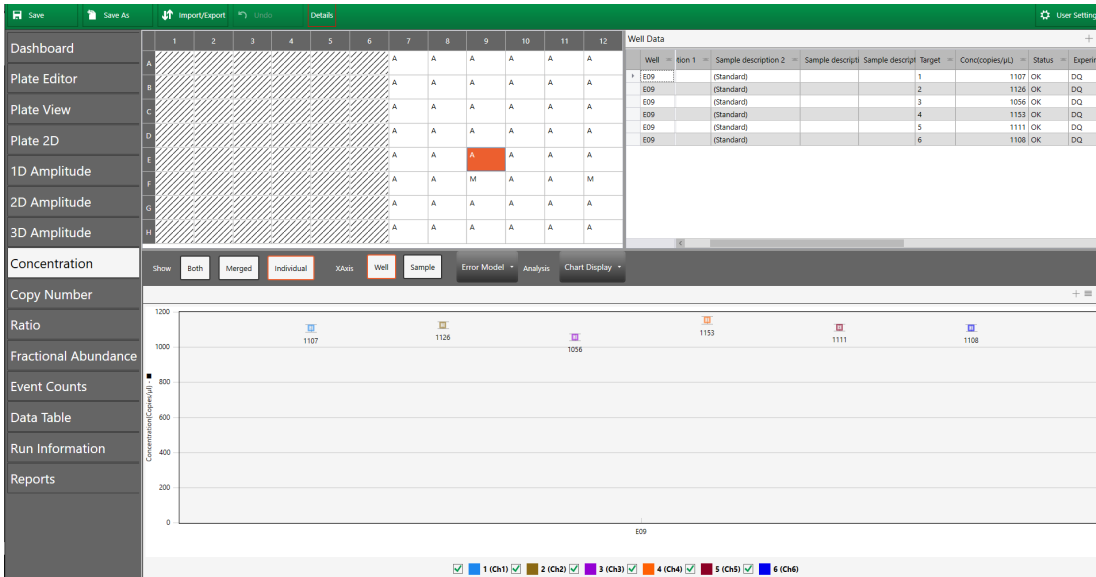

QX Manager Software plots the averaged concentration for each target as a data point. Pause on a data point to see the high, median, and low numbers.

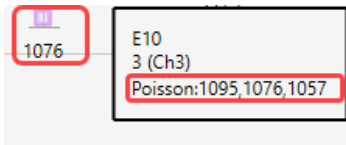

Concentration Calculation

The software uses the formula and variables shown below.

$$c = -\ln\left(\frac{N_{neg}}{N}\right) / V_{droplet}$$

Where:

-ln	Negative natural logarithm
V_{droplet}	Volume of droplet
N_{neg}	Number of negative droplets
N	Total number of droplets

Confidence Interval Calculation

Note: If you select a 68% confidence interval, the multiplier is changed from 1.96 to 1.00.

$$p = \frac{N_{neg}}{N}$$

$$\partial_{N_{neg}} = \sqrt{Np(1-p)}$$

$$N_{neg.min} = N_{neg} - 1.96 * \partial_{N_{neg}}$$

$$N_{neg.max} = N_{neg} + 1.96 * \partial_{N_{neg}}$$

Where:

$V_{droplet}$	Volume of droplet
N_{neg}	Number of negative droplets
N	Total number of droplets
p	p-value (probability of finding the target in the sample)
∂	Delta


Copy Number

Important: To see copy number variation data, you must run a copy number variation (CNV) ddPCR experiment.


Copy number measures the count of the target gene, relative to the reference gene, to calculate the number gene targets per genome. Use the toolbar to merge or separate the calculated results for each well, and to recalculate your results based on a different error model and confidence interval. For information on toolbar options, see [Data Visualization Chart Analysis Options on page 135](#).



The following graphic shows copy number data for six colors in standard-size droplets.

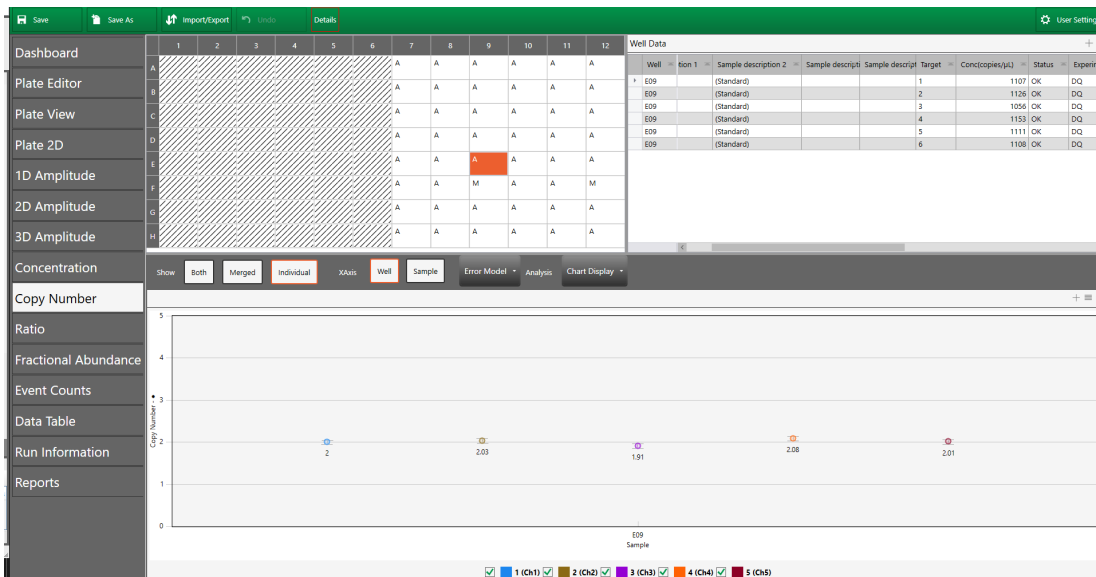

QX Manager Software plots each copy number variation as a data point. Pause on a data point to see the high, median, and low numbers.

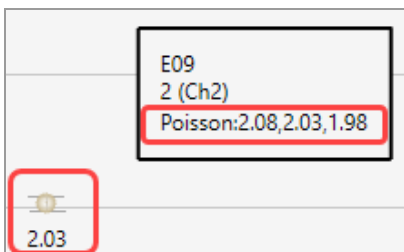

Copy Number Calculation

The software uses the formula and variables shown below.

$$CNV = \frac{A}{B} N_B$$

Where:

A	Target concentration
B	Reference concentration
N_B	Number of reference copies in genome (usually 2)

Confidence Interval Calculations

Note: If you specify the lesser confidence interval when you define your error model, the software uses 68% in its calculations instead of 95%.

$$CI_{CNV} = N_B \cdot \frac{A}{B} \sqrt{\frac{CI_A^2}{A^2} + \frac{CI_B^2}{B^2}}$$

$$(CNV_{low_95\%}, CNV_{hi_95\%}) = (CNV - (1/2)CI_{CNV}, CNV + (1/2)CI_{CNV})$$

Where:

A	Target concentration
B	Reference concentration
N_B	Number of reference copies in genome (usually 2)
CI_A	Poisson confidence interval of target (95% max - 95% min)
CI_B	Poisson confidence interval of reference (95% max - 95% min)

Notes:

Copy Number analysis involves determining the number of copies of a given target DNA sequence as compared to an invariant reference DNA sequence. Use the CNV analysis to determine structural variations, such as ploidy changes (addition or deletion of chromosomes), deletions (missing partial chromosomes) or duplications (repeated target sequences) in the genome.

The ability to discriminate between consecutive CN states has been a technical challenge in copy number assessment. Fundamentally, as CN state increases, the percentage difference in target genomic material between states decreases. For example, for a given target locus, a CN of 3 is 50% more abundant in concentration per genome than a CN of 2, while a CN of 5 is only 25% more abundant per genome than a CN of 4, and so on.


Due to the massive partitioning of the reaction across up to 20,000 droplets per well, and the absolute nature and precision of the ddPCR concentration measurement, ddPCR CNV analysis enables the quantitative discrimination required to resolve consecutive copy number states beyond CN 3. For example, a 20% difference in target concentration between a CN 5 versus CN 6 sample can be reproducibly resolved using ddPCR. This discrimination is both a function of concentration differences between consecutive CN states, as well as the amount of DNA assayed per well.

Ratio


Ratio determines the quotient of two target molecules in each μl of sample based on the default settings.

T1:T2


Use the toolbar to merge or separate the calculated results for each well, and to recalculate your results based on a different error model and confidence interval. For information on toolbar options, see [Data Visualization Chart Analysis Options on page 135](#).

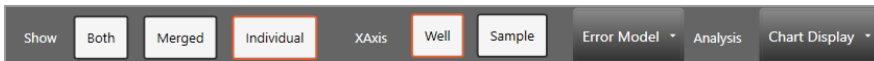

The following graphic shows the ratio for standard-size droplets.

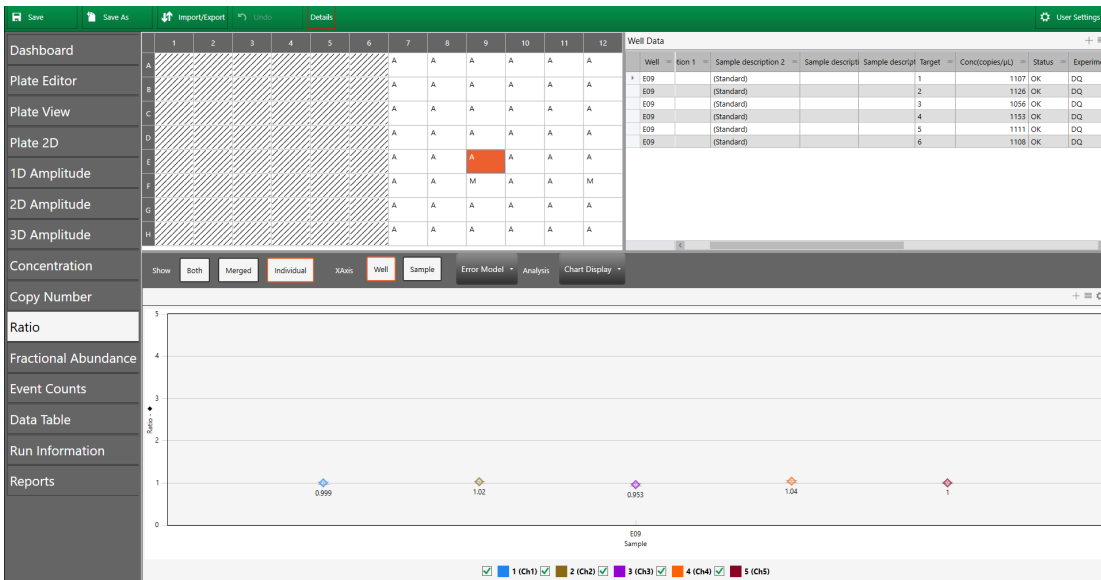

QX Manager Software plots each ratio as a data point. To see the confidence level numbers, pause on the data point.

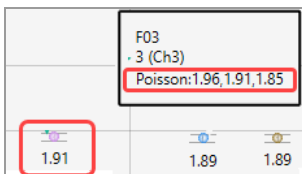

Ratio Calculation

The software uses the formula and variables shown below.

$$r = \frac{A}{B}$$

Where:

A	Target concentration
B	Reference concentration

Note: If no reference is selected, the target with the highest dye concentration in the highest channel will automatically be considered as the reference. If multiple references are selected, the geometric mean is used.

Confidence Interval Calculations

Note: If you specify the lesser confidence interval when you define your error model, the software uses 68% in its calculations instead of 95%.

$$CI_r = \frac{A}{B} \sqrt{\frac{CI_A^2}{A^2} + \frac{CI_B^2}{B^2}}$$


$$(r_{low_95\%}, r_{hi_95\%}) = (r - (1/2)CI_r, r + (1/2)CI_r)$$

Where:


A	Channel 1 target concentration
B	Channel 2 target concentration
CI_A	Confidence interval of Channel 1 target (95% max - 95% min)
CI_B	Confidence interval of Channel 2 target (95% max - 95% min)

Fractional Abundance

Fractional Abundance determines the relative abundance of your targets in each μl of sample based on the default settings. Use the toolbar to merge or separate the calculated results for each well, and to recalculate your results based on a different error model and confidence interval. For information on toolbar options, see [Data Visualization Chart Analysis Options on page 135](#).

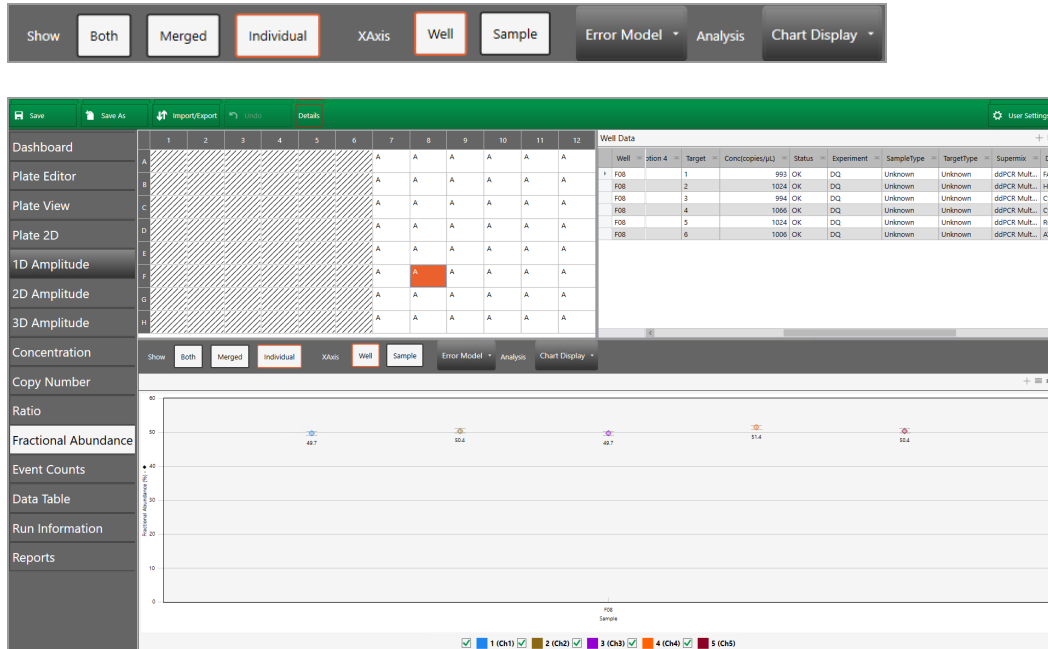

QX Manager Software plots the fractional abundance as a data point. To see the high and low confidence interval numbers, pause on the data point.

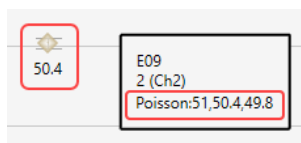

Fractional Abundance Calculation

The software uses the formula and variables shown below.

$$f = \frac{A}{A+B}$$

Where:

A	Target concentration
B	Reference concentration

Note: If no reference is selected, the target with the highest dye concentration in the highest channel will automatically be considered as the reference. If multiple references are selected, the geometric mean is used.

Confidence Interval Calculations

Note: If you specify the lesser confidence interval when you define your error model, the software uses 68% in its calculations instead of 95%.

$$CI_f = \frac{1}{(A+B)^2} \sqrt{B^2 CI_A^2 + A^2 CI_B^2}$$


$$(f_{low_95\%}, f_{hi_95\%}) = (f - (1/2)CI_f, f + (1/2)CI_f)$$

Where:


A	Channel 1 target concentration
B	Channel 2 target concentration
CI_A	Confidence interval of Channel 1 target (95% max - 95% min)
CI_B	Confidence interval of Channel 2 target (95% max - 95% min)

Event Counts

The Event Counts tab shows a bar chart view of the number of positive and negative droplets for each of the targets in each of the selected wells.

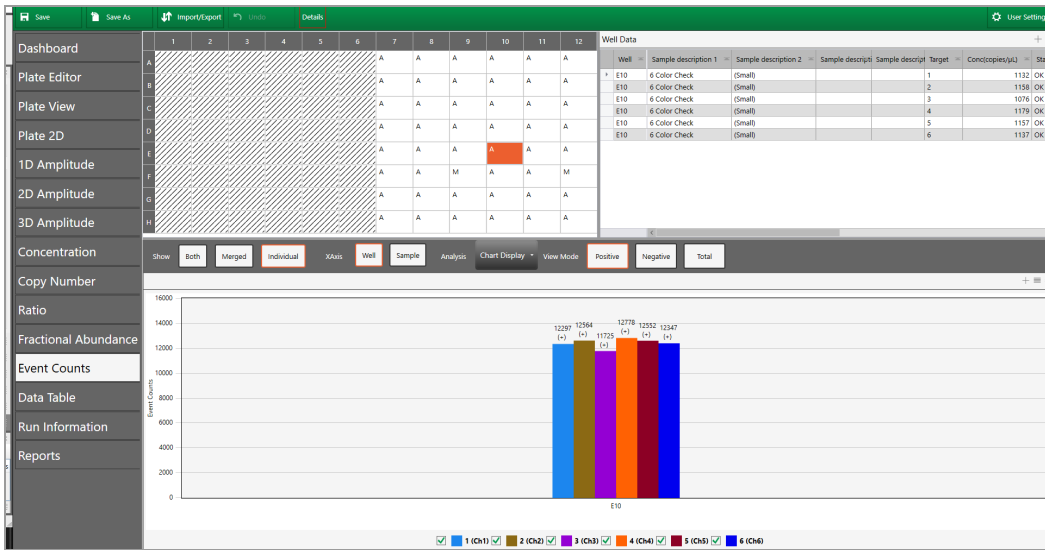

Use the toolbar to change the display by combining or separating data for each well. You can also show data by well name or by sample name, and you can show any combination of positive, negative, or combined droplet counts.

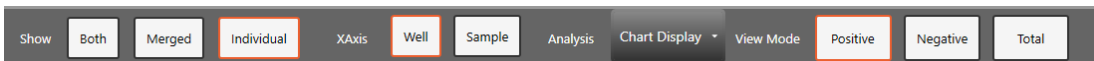

Linkage Analysis

In ddPCR, physical linkage of two targets can occur in one of the following patterns:

- Tandem repeats of the same target sequence
- Proximity of two target sequences physically linked on the same piece of DNA

QX Manager Software uses the formulas and variables shown below.

- The following formula presents the linkage in copies per μL :

$$\text{linkage} = \ln(N) - \ln\left(n_{00} + n_{01} + n_{10} + \frac{n_{01}n_{10}}{n_{00}}\right) \left(\frac{1000}{\text{droplet volume}}\right)$$

- The following formula presents the linkage in copies per droplet.

$$\text{linkage} = \ln(N) - \ln\left(n_{00} + n_{01} + n_{10} + \frac{n_{01}n_{10}}{n_{00}}\right)$$

Where:

N	Total number of droplets
ln	Natural logarithm
-ln	Negative natural logarithm
n_{00}	Number of negative droplets for both channels
n_{01}	Number of negative droplets in channel 1 and positive droplets in channel 2
n_{10}	Number of positive droplets in channel 1 and negative droplets in channel 2

Where:

Confidence Interval Calculation

$$CI_{\text{linkage}} = (\text{linkage} - (1.96)\sigma_{\text{linkage}}, \text{linkage} + (1.96)\sigma_{\text{linkage}})$$

Where:

σ	Standard deviation
----------	--------------------

This page intentionally left blank.

Chapter 9 Analysis Module Reporting

Use the Reporting function to create a variety of reports in the QX Manager Software Analysis module.

You can create reports on the fly from any analysis file or, if you need to generate specific reports on a regular basis, you can set up report templates for reuse.


This section contains information on how to create analysis reports.

Creating Analysis Reports

You can create reports with or without a template. For information on creating templates for reports you generate often, see [Creating or Modifying Report Templates on page 81](#).

To create a report

1. Select the Reports tab to display the Reports window.

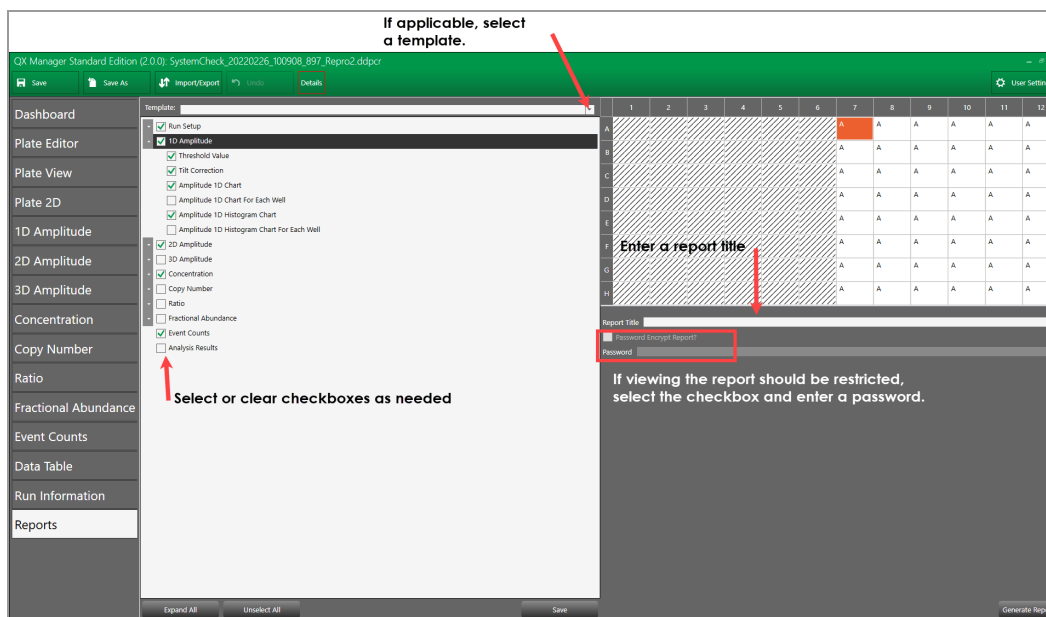

2. Do one of the following:

- To create a report from a template, click the Template dropdown arrow and select a template from the list
- To create a custom report, click the arrows to expand reporting categories, and then select the checkbox for each report element to include, and clear checkboxes for elements to eliminate.


Note: Use the buttons at the bottom of the screen to expand or collapse all categories at once, or to select or clear all checkboxes at once. You can also drag and drop the section to rearrange the order of appearance in the report, and drag and drop items in the section to reorder them.



3. Enter a report title.
4. (Optional) To restrict viewing, select the Password Encrypt Report checkbox and then enter a password.
5. Click Generate Report to generate a PDF report.
6. To save the report, click Save.

Chapter 10 Gene Study Module

You access the Gene Study module from the Data Analysis window.

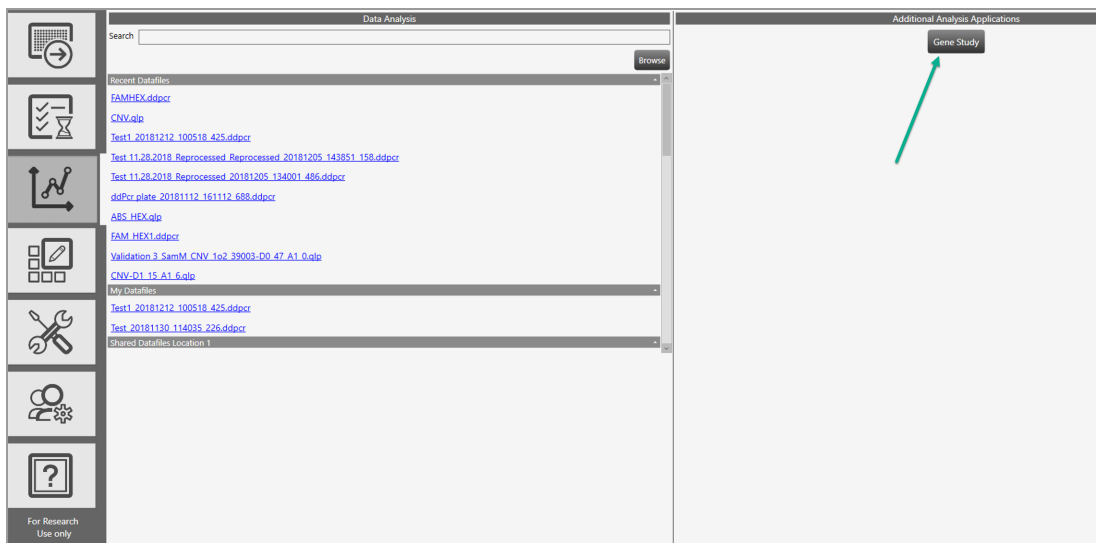


To create a gene study, you add one or more data files and then compare gene expression data from one or more ddPCR experiments.

Note: The maximum number of samples you can analyze in a gene study is limited by the size of the computer's RAM and virtual memory.

To open the Gene Study module

- ▶ Click Gene Study.

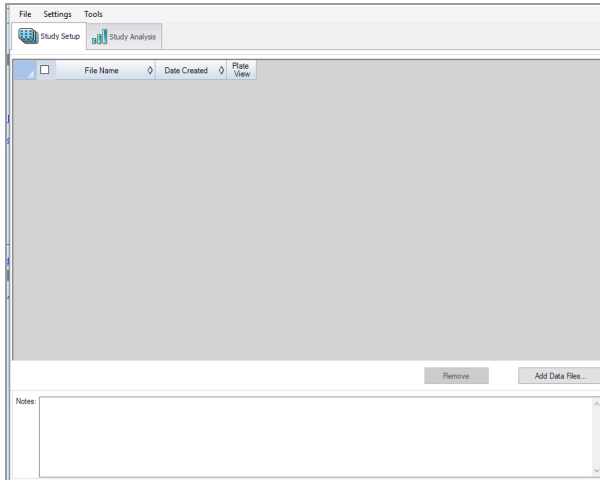

The Gene Study dialog box includes two tabs:

- **Study Setup** — manages the run files in the gene study
 - Important:** Adding or removing data files in a gene study does not change the data in the original file.
- **Study Analysis** — displays the gene expression data for the combined runs.

From these tabs you can add files to set up a gene study, analyze your data in charts and plots, and run reports on the gene study data.

Gene Study Options

The Gene Study module provides the menus and corresponding options described below.

- **File** — Open, save, and close the gene study, and add files
- **Settings** — Restore the default layout
- **Tools** — Create a report

To open a gene study file

1. Click File → Open.
2. In the Open dialog box, navigate to and select the data file, and then click Open.

To save a gene study file

1. Click File and select Save or Save As.
2. If prompted, enter a file name and click Save.
3. If prompted to overwrite an existing file, click Yes or No.

To close a gene study file

- ▶ Click File → Close.

Note: If you have not saved the file, an advisory prompt appears.

To add data files

1. Click File → Add Data Files.
2. In the Open dialog box, navigate to the data files to be added to the study.
3. Select one or more files and then click Open to add the files to the Gene Study window.

To restore default settings

- ▶ Click Settings → Restore default settings.

Note: At least one modified data file must be open.

To generate a report

- ▶ See [Gene Study Reports on page 162](#).

Setting up the Gene Study

Use the Study Setup window to

- Add data files to include in the study
- Remove data files from the study
- Select data files for viewing and analysis
- Show the plate view for a data file


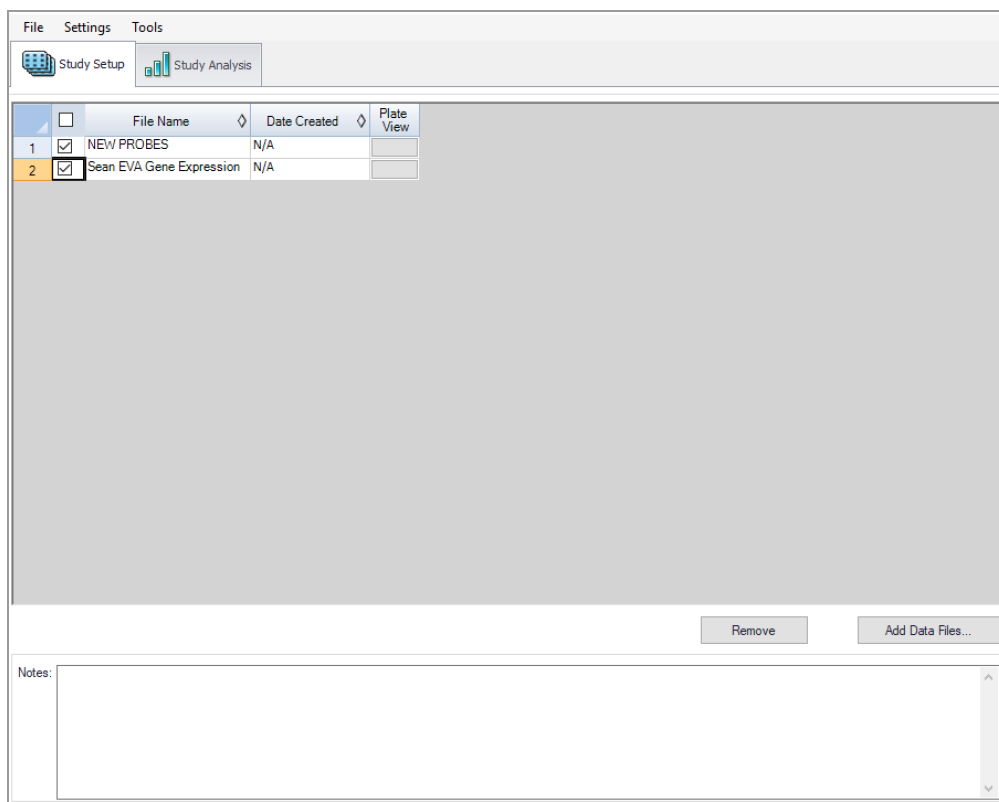


Table 37 defines each column in Study Setup.

Table 37. Gene Study Setup Fields

Column Title	Description
File Name	Name of the data file
Date Created	Date the file was imported into the Gene Study module
Plate View	Opens a map of the plate with the data for each dye in each well included in the Gene Study

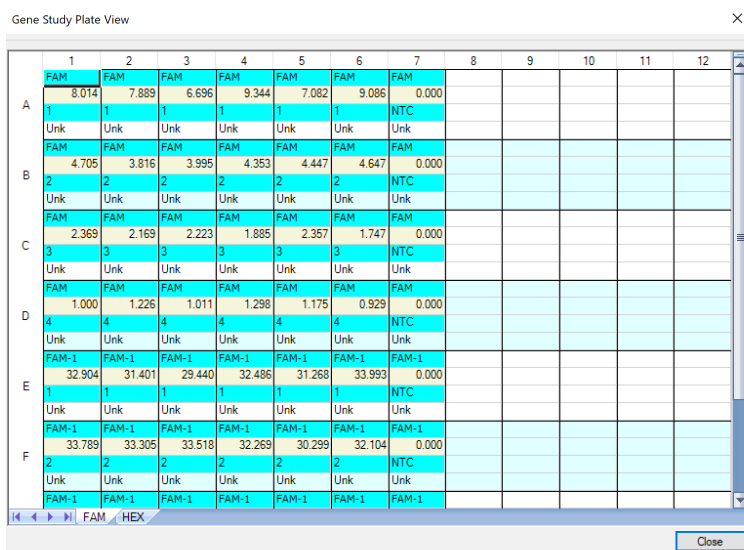


Adding Data Files

To add data files

1. On the Study Setup tab, click Add Data Files.
2. In the Open dialog box, navigate to the data files to be added to the study.
3. Select one or more files and then click Open.

The files are added to the Study Setup window.

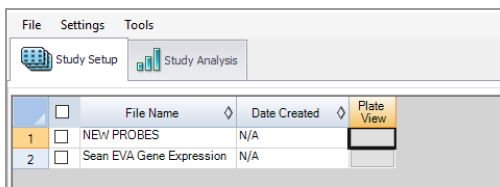

Removing Data Files

To remove data files

- ▶ Select the checkbox next to the file and then click Remove.


The file is removed from the Study Setup window.

Note: The Remove button is disabled until you select a file.

Selecting Data Files for Analysis

To select data files for analysis

1. Select the checkboxes next to the files to be included in the analysis.

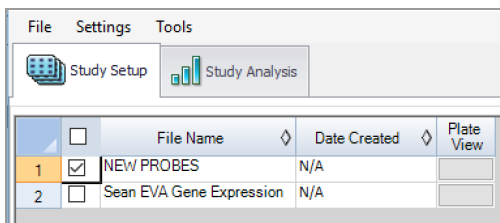


2. Select the Study Analysis tab to see the analysis results.

Showing the Plate View

To show the plate view

1. In the Plate View column, click the button for the plate view to be displayed.

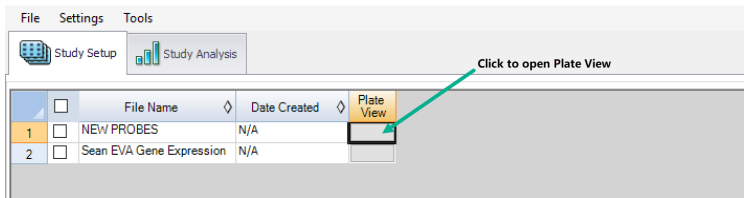

The Plate view appears.

The screenshot shows a window titled 'Gene Study Plate View' containing a grid of data. The grid has 12 columns (wells) and 6 rows (A-F). Each cell contains a value or label. The values are: Row A: FAM (8.074, 7.889, 6.696, 9.344, 7.082, 9.086, 0.000), Unk (1, 1, 1, 1, 1, 1, NTC); Row B: FAM (4.705, 3.816, 3.995, 4.353, 4.447, 4.647, 0.000), Unk (2, 2, 2, 2, 2, 2, NTC); Row C: FAM (2.369, 2.169, 2.223, 1.885, 2.357, 1.747, 0.000), Unk (3, 3, 3, 3, 3, 3, NTC); Row D: FAM (1.000, 1.226, 1.011, 1.298, 1.175, 0.929, 0.000), Unk (4, 4, 4, 4, 4, 4, NTC); Row E: FAM-1 (32.904, 31.401, 29.440, 32.486, 31.268, 33.993, 0.000), Unk (1, 1, 1, 1, 1, 1, NTC); Row F: FAM-1 (33.789, 33.305, 33.518, 32.269, 30.299, 32.104, 0.000), Unk (2, 2, 2, 2, 2, 2, NTC). The status bar at the bottom shows 'FAM / HEX' and a 'Close' button.

	1	2	3	4	5	6	7	8	9	10	11	12
A	FAM 8.074	FAM 7.889	FAM 6.696	FAM 9.344	FAM 7.082	FAM 9.086	FAM 0.000					
	Unk 1	Unk 1	Unk 1	Unk 1	Unk 1	Unk 1	NTC					
B	FAM 4.705	FAM 3.816	FAM 3.995	FAM 4.353	FAM 4.447	FAM 4.647	FAM 0.000					
	Unk 2	Unk 2	Unk 2	Unk 2	Unk 2	Unk 2	NTC					
C	FAM 2.369	FAM 2.169	FAM 2.223	FAM 1.885	FAM 2.357	FAM 1.747	FAM 0.000					
	Unk 3	Unk 3	Unk 3	Unk 3	Unk 3	Unk 3	NTC					
D	FAM 1.000	FAM 1.226	FAM 1.011	FAM 1.298	FAM 1.175	FAM 0.929	FAM 0.000					
	Unk 4	Unk 4	Unk 4	Unk 4	Unk 4	Unk 4	NTC					
E	FAM-1 32.904	FAM-1 31.401	FAM-1 29.440	FAM-1 32.486	FAM-1 31.268	FAM-1 33.993	FAM-1 0.000					
	Unk 1	Unk 1	Unk 1	Unk 1	Unk 1	Unk 1	NTC					
F	FAM-1 33.789	FAM-1 33.305	FAM-1 33.518	FAM-1 32.269	FAM-1 30.299	FAM-1 32.104	FAM-1 0.000					
	Unk 2	Unk 2	Unk 2	Unk 2	Unk 2	Unk 2	NTC					
	FAM-1	FAM-1	FAM-1	FAM-1	FAM-1	FAM-1	FAM-1					

2. Click Close to close the view.

Using the Study Analysis Features

The Study Analysis tab displays the data from all runs in the gene study in the views described in [Table 38](#).

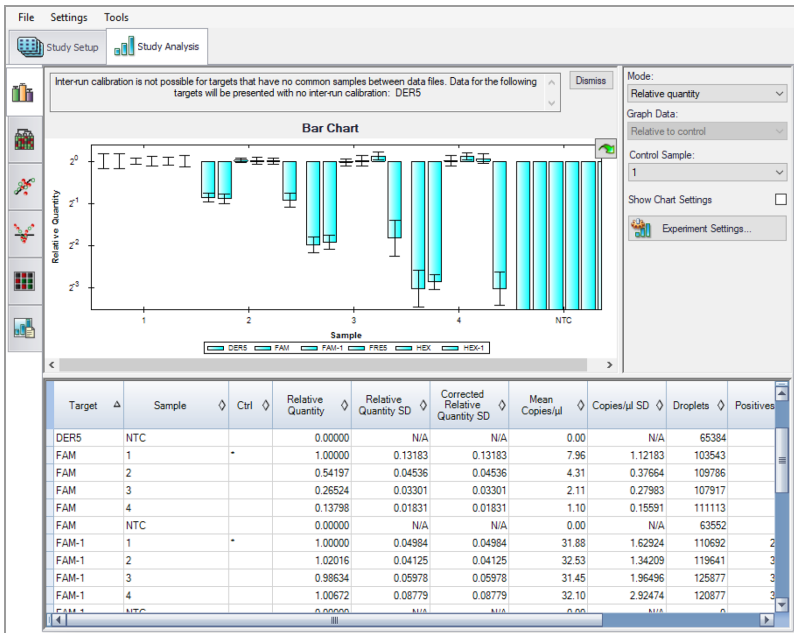


Table 38. Analysis Toolbar







Button	Description
	Size of bars correspond to normalized expression or relative quantity
	Data is arranged in a hierarchy based on the degree of similarity of expression for different targets and samples
	Dot plot of normalized expression of targets for a control versus an experimental sample

Table 38. Analysis Toolbar, continued

Button	Description
Volcano Plot 	Dot plot of change in expression (regulation) of a target for an experimental sample compared to a control
Heat Map 	Plate layout grid showing target regulation for an experimental compared to a control sample based on relative normalized expression Note: If the same targets are in the same location on multiple plates but with different samples, use the dropdown menu to select a particular plate for analysis.
Results Table 	Provides results in a tabular format

Gene Study Reports

Use the Gene Study Report dialog box to arrange the gene study data into a report. Table 34 lists all the options available for a gene study report.

Table 39. Gene Study Report Categories

Category	Option	Description
Header		
	Report Information	Date, user name, data file name, data file path, and the selected well group
	Gene Study File List	List of all the data files in the gene study
	Notes	Notes about the data report
Study Analysis Bar Chart		
	Analysis Settings	List of the selected parameters
	Chart	Gene expression bar chart showing the data
	Target Names	List of targets in the gene study
	Sample Names	List of samples in the gene study
	Data	Spreadsheet that shows the data
	Target Stability	Target stability data
	Inter-run Calibration	Inter-run calibration data
Study Analysis Clustergram, Scatter Plot, Volcano Plot, and Heat Map		
	Analysis Settings	Settings for each chart type
	Chart	Gene expression chart showing the data
	Data	Spreadsheet listing the data in each target

Creating a Gene Study Report

Before creating a report, make any needed adjustments to the gene study data and charts.

To create a gene study report

1. Select Tools → Reports in the Gene Study menu to open the Report dialog box.
2. Select the options you want to include in the report.

The report opens with default options selected. Select or clear the checkboxes to change whole categories or individual options within a category.

For information on the available options to display, see [Gene Study Reports on page 162](#).

3. Change the order of categories and items in a report.

Drag the options to the required position. Note that Items can be reordered only within the categories to which they belong.

4. Click Update Report to update the Report Preview window with any changes.
5. Print or save the report:
 - a. Click the Print Report button on the toolbar to print the current report.
 - b. Select File → Save to save the report in PDF, MHT or MHTML file format and select a location to save the file.

Note: MHT and MHTML are Microsoft formats.
 - c. Select File → Save As to save the report with a new name or in a new location.
6. (Optional) To save the current report settings as a template, select Template → Save or Save As.

This page intentionally left blank.

Chapter 11 System Utilities

From the System tab, you can manage tasks related to the instrument, software, data, and storage.

- **System Settings** — View shared and assigned storage locations, select an alternate graphics rendering method, and change the instrument.
- **Event Log** — View logged data about all software activities.
- **Maintenance Log** — Record maintenance activities.
- **Maintenance Reports** — PDF maintenance records from Bio-Rad maintenance.
- **System Data** — Archive raw data to clear up space in your primary storage areas, access data for troubleshooting, and use the data recovery function.
- **Reprocess Data** — Reprocess a .data file using new calibration values.
- **Tools** — Flush or prime the instrument.

System Settings Tab

When you select the System tab, the System Settings window opens by default.

The screenshot shows the 'System Settings' window with the 'Datafile and Template Storage' tab selected. The window is divided into three main sections: 'Preferred Location', 'Shared Settings', and 'Save'. The 'Preferred Location' section includes a checkbox for 'Use for all users' and two rows for storage locations: 'Datafile Storage Location' and 'Template Storage Location', each with a 'Browse' button. The 'Shared Settings' section includes three rows for storage locations: 'Datafile Storage Location', 'Datafile Storage Location 2', and 'Template Storage Location', each with a 'Browse' button. A 'Save' button is located at the bottom of the 'Shared Settings' section.

- Use the Datafile and Template Storage tab to view storage locations:
 - **Preferred Locations** — if the checkbox selected, storage for all users is limited to a single folder for data files and a single folder for templates.
 - **Shared Settings** — if the Preferred Locations checkbox remains cleared, then shared locations can be defined, but remain optional. Users can choose to save their files to their personal folders specified in User Preferences or to the shared location.

Note: The system administrator specifies shared and preferred storage locations, and can also define a backup location for data file storage if the priority folder is full.

- Use the Instrument tab to select the instrument model that is connected to QX Manager Software.
- Use the Software Settings tab to change the graphics rendering method.

Preferred Location

If the administrator sets Preferred Location file paths, then data files and templates for all users are automatically saved to the specified folders. Preferred Locations override selections in individual user preferences.

Important: Only users assigned the System Settings user privilege can set preferred locations.

System Settings	Datafile and Template Storage	Instrument
Event Log	Preferred Location ⓘ	
Maintenance Log	<input checked="" type="checkbox"/> Use for all users	
Maintenance Reports	Datafile Storage Location:	
System Data	<input type="text" value="C:\Users\Public\Documents\Bio-Rad\QXMgrStandard\AppData\Data"/>	
Reprocess Data	<input type="button" value="Browse"/>	
	Template Storage Location:	
	<input type="text" value="C:\Users\Public\Documents\Bio-Rad\QXMgrStandard\AppData\Templates"/>	
	<input type="button" value="Browse"/>	

To specify preferred locations

1. Select the Use for all users checkbox.
2. (Optional) Keep the default paths or modify them to alternative secure storage locations.
3. Click Save.
4. When the confirmation message appears, click OK.

Shared Settings

Under Shared Settings, you can choose to save your files to locations that are available to all users.

The screenshot shows the 'System Settings' window with a sidebar on the left containing categories like System Settings, Event Log, Maintenance Log, Maintenance Reports, System Data, Reprocess Data, Instrument Calibration, Instrument Check, and Tools. The main area is titled 'Datafile and Template Storage' and 'Instrument'. The 'Shared Settings' section is highlighted with a red border and contains the following fields:

- Preferred Location** (with a help icon)
- Use for all users
- Datafile Storage Location:** C:\Users\Public\Documents\Bio-Rad\QX\Mgr\Standard\AppData\Data (with a 'Browse' button)
- Template Storage Location:** C:\Users\Public\Documents\Bio-Rad\QX\Mgr\Standard\AppData\Templates (with a 'Browse' button)
- Shared Settings** (with a help icon)
- Datafile Storage Location:** <location available to all users> (with a 'Browse' button)
- Datafile Storage Location 2:** <backup location available to all users> (with a 'Browse' button)
- Template Storage Location:** (with a 'Browse' button)
- Save** button

The software allows the software administrator to set a primary storage folder and a backup storage folder. Bio-Rad recommends routinely checking available storage space and if necessary, using the Archive function.

Important: Only users assigned the System Settings user privilege (typically your software administrator) can change storage locations in the System Settings window.

To specify shared folder locations

1. In each of the storage path fields, enter a file path or click Browse to search for the folder.
2. Click Save.
3. When the confirmation message appears, click OK.

Instrument Settings

When you open QX Manager Software for the first time after installation, the Instrument connection window appears and you must select the instrument you are using.

Changing the Instrument


You can choose between the QX200 Droplet Reader and the QX600 Droplet Reader.

To change the instrument

1. Select the System Settings tab.



2. In the System Settings window (default), select the Instrument tab.

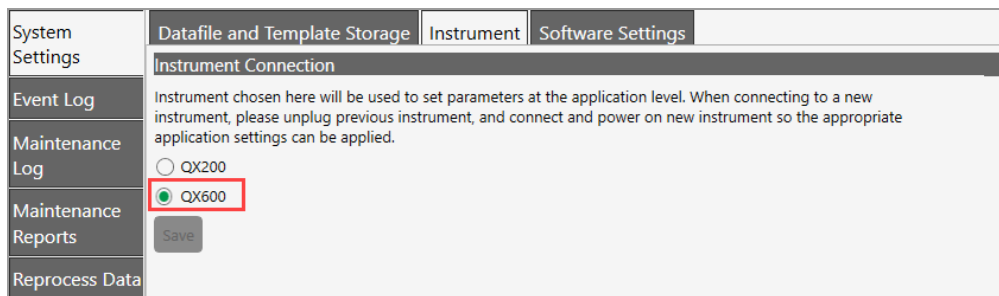


3. Select the instrument and then click Save.

Features associated with the selected instrument become available when the instrument is connected.

Software Settings

To improve chart appearance, the Software Settings tab provides an alternate option for chart rendering.

Important: Poor rendering typically occurs with some Intel graphics cards only. Contact Bio-Rad Technical Support for information and assistance. Do not change the setting unless instructed to do so by Bio-Rad Technical Support.

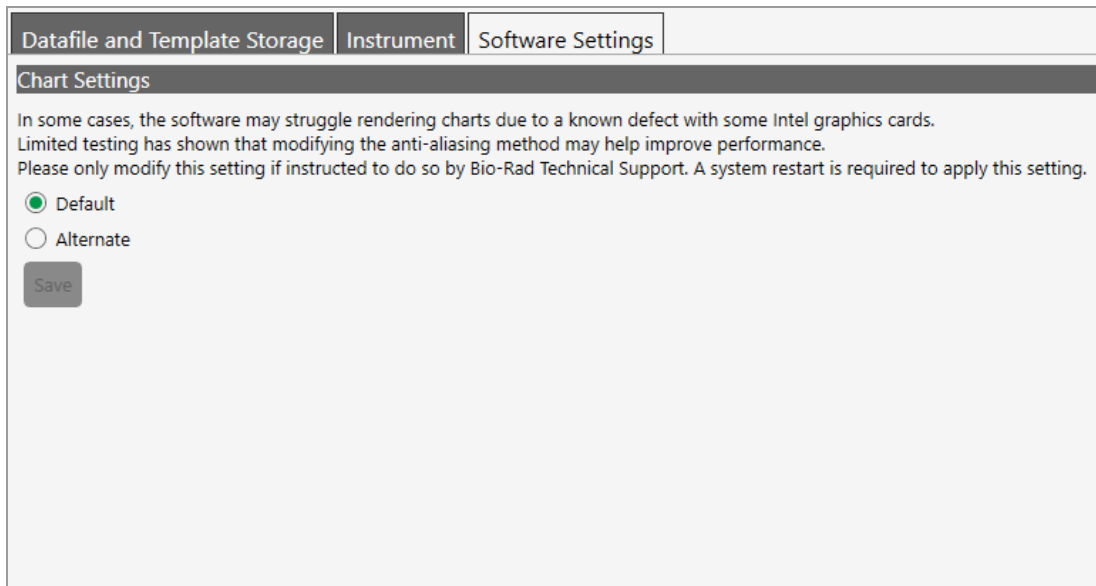

The screenshot shows a software interface with three tabs: 'Datafile and Template Storage', 'Instrument', and 'Software Settings'. The 'Software Settings' tab is selected. Below the tabs is a section titled 'Chart Settings'. The text in this section reads: 'In some cases, the software may struggle rendering charts due to a known defect with some Intel graphics cards. Limited testing has shown that modifying the anti-aliasing method may help improve performance. Please only modify this setting if instructed to do so by Bio-Rad Technical Support. A system restart is required to apply this setting.' Below the text are two radio button options: 'Default' (which is selected) and 'Alternate'. A 'Save' button is located below the radio buttons.

Complete the following steps if directed to do so by Bio-Rad Technical Support.

To change the graphics rendering method

1. Under Chart Settings, click Alternate.
2. Click Save.
3. Close QX Manager Software.
4. To enable the change, restart the connected computer, and then reopen the software.

System Log Files

From the System tab, you can access log files for system and software events (event log), and maintenance activities (maintenance log).

Event Log

Users who are assigned the Maintenance user privilege can view the system event log, which contains timestamped information on all software activities.

System Settings	Log File: QXMgr< edition> History.log			Open Log	PDF Password	PDF Report
Event Log	Date/Time	Message	Severity	User Name	Search	
Maintenance Log	8/2/2021 11:54:59 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	USPLECOTKYFZ\QX_!		
Maintenance Reports	8/2/2021 11:20:32 AM	User Privileges Changed For User: QX_Super (USPLECOTKYFZ). New Privileges are: Add/Manage Users,Create New Templates,V...	Info	QX_Super		
Reprocess Data	8/2/2021 11:20:13 AM	User Privileges Changed For User: QX_Super (USPLECOTKYFZ). New Privileges are: Add/Manage Users,Create New Templates,V...	Info	QX_Super		
Instrument Calibration	8/2/2021 11:19:52 AM	User Privileges Changed For User: QX_Super (USPLECOTKYFZ). New Privileges are: Add/Manage Users,Create New Templates,V...	Info	QX_Super		
Tools	8/2/2021 11:16:10 AM	User Preferences Changed For User: QX_Super. Timeout Value Changed. Old Value: 5 New Value: 30	Info	QX_Super		
	8/2/2021 11:16:09 AM	User Privileges Changed For User: QX_Super (USPLECOTKYFZ). New Privileges are: Add/Manage Users,Create New Templates,V...	Info	QX_Super		
	8/2/2021 11:15:46 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	8/2/2021 10:58:16 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	8/2/2021 10:52:15 AM	Gene Study Application Opened. Version: 2.0.0.462	Info	QX_Super		
	8/2/2021 10:52:08 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/30/2021 12:18:32 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/30/2021 9:59:50 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/30/2021 9:59:44 AM	Failed to Login User QX_Super. Error code: InvalidPassword	Info	QX_Super		
	7/30/2021 9:59:37 AM	Failed to Login User QX_Super. Error code: InvalidPassword	Info	QX_Super		
	7/30/2021 9:49:25 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/30/2021 8:45:16 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 4:16:09 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 3:49:37 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 3:43:09 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 2:54:49 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 2:47:10 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 2:25:21 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 1:00:31 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 12:44:20 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 12:27:28 PM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 11:09:43 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		
	7/29/2021 10:56:22 AM	User Logged in. User Name: QX_Super, User Full Name:	Info	QX_Super		

To view the Event Log

1. Select the System Utilities tab.



2. Select Event Log.
3. Click Open Log to choose either the debug log or the history log.
4. Review the log entries.
5. (Optional) To automatically generate and display a PDF document, click PDF Report.

Tip: To restrict access to the file, you can assign a password. Select the PDF Password check box and then enter a password in the field. The file is saved in the Log Reports folder and can be opened only by entering the password.

Maintenance Log

All users can view the maintenance log, which contains a list of maintenance records. Users who are assigned the Maintenance user privilege can set up maintenance records, specify associated activities, and generate PDF reports.

Tip: This guide also contains a form on which you can document completion of maintenance activities. For information, see [Instrument Maintenance Log on page 216](#).

The screenshot shows a software interface for the Maintenance Log. On the left, there is a sidebar with tabs: System Settings, Event Log, Maintenance Log, Archive Data, and Reprocess Data. The 'Maintenance Log' tab is active, displaying a table with the following data:

Date Time	User	Product Version	Firmware Version	Activity
12/10/2018	lbr&D	1.0.66		
12/10/2018	lbr&D	1.0.66		

Below the table is a form for adding a new activity. It includes fields for Date Time (12/10/2018 10:50:16 AM), User (lbr&D), Product Version (1.0.66), and Firmware Version. There is also a large text area for the Activity description and a PDF Report button at the bottom right.

To add or view an activity

1. Select the System Utilities tab.



2. Select Maintenance Log. Note that activities that are already scheduled appear in the grid, sorted by newest activity first.
3. Do one of the following:
 - Scroll to view a particular activity.
 - **Important:** You can view or add static activities; you cannot modify them.
 - Add an activity; click the + icon next to New Activity.

A line item appears in the bottom section, with all fields except Activity completed automatically.


4. In the Activity field in the bottom section, type the activity into the Activity field, and click Save.
5. Repeat to add another activity.
6. (Optional) Click PDF Report to create a report of the added activities.

Maintenance Reports

Reports generated from the Bio-Rad service engineer when performing maintenance on your instrument are stored on the Maintenance Reports tab.

To open a report

- ▶ Click the list item.

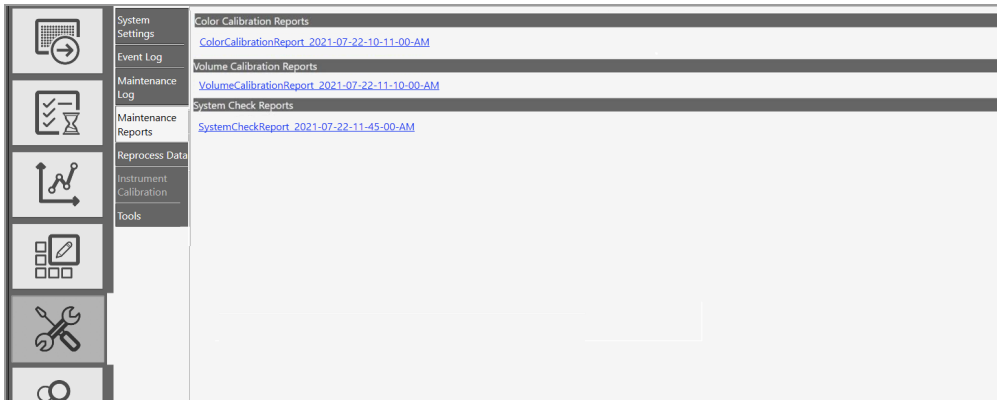

Following is an example of a Color Calibration report:

Color Calibration and Carryover Report		
<p>BIO-RAD</p> <p>Color Calibration and Carryover Report</p> <p>Username of the Person who did the Run: BBRAD</p> <p>Report Generation Date and Time: 7/22/2019 09:03:29 PM</p> <p>Instrument Serial Number: CHARLE-1</p> <p>Firmware Version: 1.2790</p> <p>Software Version: 1.0.774.0628</p> <p>Software: QX ONE Software Standard Edition</p> <p>Lot Information</p> <p>GC836 Cartridge: 18230</p> <p>Droplet Generation: 08.BKDR101.BKDR101</p> <p>Droplet Reader: 08.BKDR101.BKDR101</p> <p>Calibration Effective Date: 27/2019 03:20:11 PM</p>		
Color Calibration Result	Carryover Test Result	# of carryover events
Channel 1 (FAM)	Pass	0
Channel 2 (HEX/NEC)	Pass	0
Channel 3 (GFP)	Pass	0
Channel 4 (R-5-F)	Pass	0
Low Mix	Pass	0

Copyright © 2019 Bio-Rad Laboratories, Inc. All rights reserved. Page 51

System Data

In the System Data window, you can

- Archive raw data
- Access troubleshooting data
- Recover data

Archiving Raw Data

When the storage space on the computer connected to the QX600 Droplet Reader approaches the limit displayed under Disk Space, you can free up space by moving raw data to an archive storage folder in a different location, such as a network or USB drive.

Important: You must be assigned the System Data user privilege to archive data.


To archive raw data

1. Select the System Utilities tab.



2. Select System Data.

The Archive Raw Data tab is selected by default.

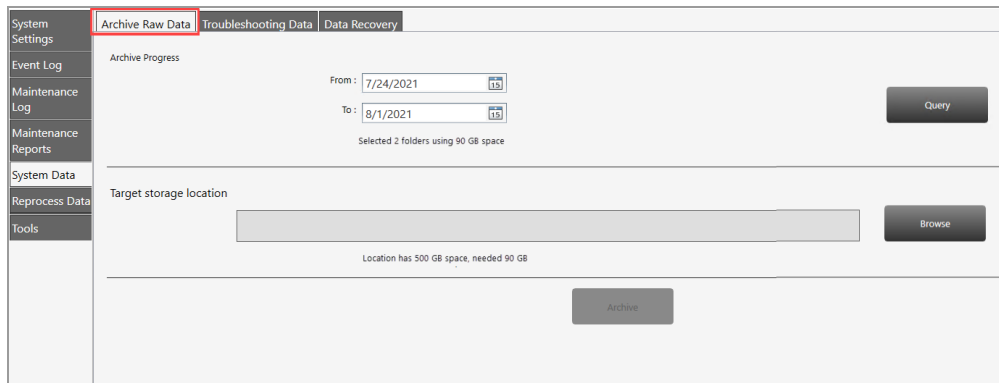

3. Under Archive Progress, click the calendar icons and select a From date and a To date to define a date span.
4. Click Query to search for files dated within the date span.
5. Click Browse to search for and select the target storage location.
6. Click Archive.

Troubleshooting Data

If issues occur with the instrument or software, you can access data to assist in troubleshooting the problem.

Note: You must be assigned the System Data user privilege to access troubleshooting data.

Important: To minimize the size of the zip file and ensure a smoother transfer of troubleshooting data, delete the contents in C:\ProgramData\Bio-Rad\QXMgrStandard\Logs before you perform the following procedure.

To collect data for troubleshooting


1. Select the System Settings tab.



2. Select System Data.

The window opens to the Archive Raw Data screen by default.

3. Select the Troubleshooting Tab.

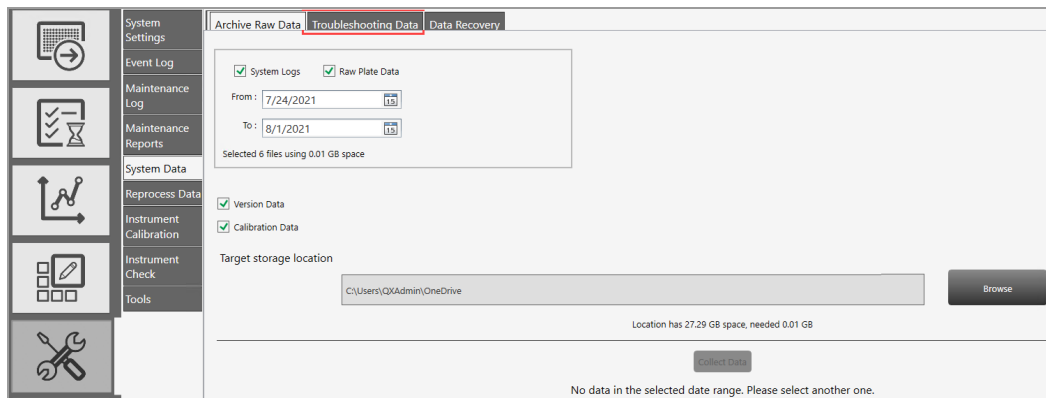

4. Click the From and To calendar icons to select a date range.

5. Select any of the following checkboxes:


- System Logs
- Raw Plate Data
- Version Data
- Calibration Data

6. Click Browse to navigate to a folder where the data will be stored.

QX Manager Software advises how much space you need and how much is available, and also advises if there is no data from your specified date range.

7. Click Collect Data.

A progress bar appears.

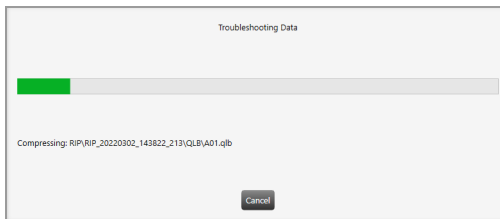

QX Manager Software creates a subfolder containing the data you specified and places it in the target location.

Recovering Data

If a run fails on the QX600 Droplet Reader, you can use the Data Recovery tab in System Data to recover data that has been saved to a temporary file.


Important: You must be assigned the System Data user privilege to access the Data Recovery tab.

To create a.ddpccr file from the data in the temporary file

1. Click the System Utilities tab.



2. Click System Data, and then click Data Recovery.

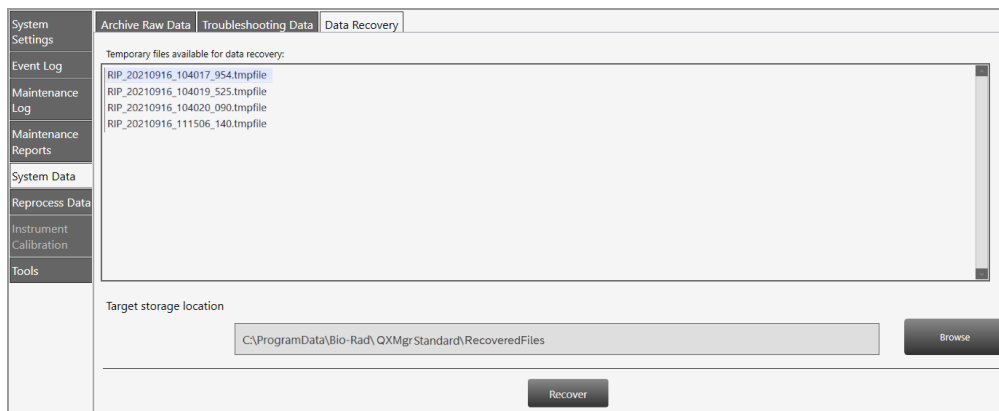


Temporary files that are available for conversion are listed in the top pane, which is scrollable.

3. Click a file, and then click Browse


4. Navigate to the target storage location on your computer or network, and then click Recover.

The following advisory message appears.

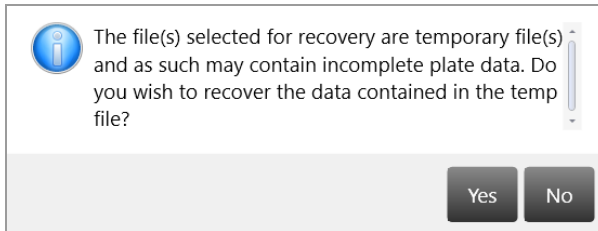

5. Click Yes.

A dialog box appears with a progress bar, and automatically closes when the conversion is complete.

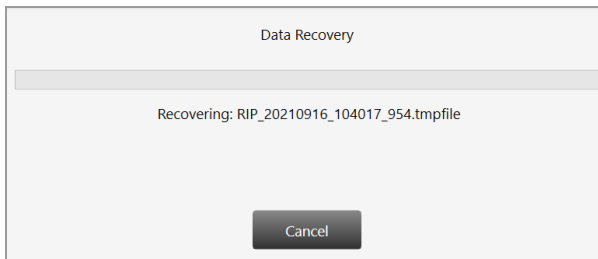

The .ddpcr file is stored in the folder you specified.

Reprocessing Data

After your instrument has been recalibrated, you can reprocess a data file from a run using different color calibration values.

Note: The instrument must be in an idle state before you can reprocess the data in QX Manager Software. If the file was created in an older software version, the software cannot reprocess the file and displays an advisory message.

To reprocess a data file

1. Select the System Utilities tab.



2. Click Reprocess Data.
3. Under Datafile, click Browse and navigate to the analysis data file.
4. In the Select Datafile dialog box, select the file and click Open.

The calibration values appear in the first Color Calibration table.

Dye	Channel 1	Channel 2	Channel 3	Channel 4	Channel 5	Channel 6
FAM	10000/10000	0/0	0/0	0/0	0/0	0/0
EVAGREEN	10000/10000	0/0	0/0	0/0	0/0	0/0
HEX	0/0	10000/10000	0/0	0/0	0/0	0/0
TRIC	0/0	10000/10000	0/0	0/0	0/0	0/0
CY5	0/0	0/0	10000/10000	0/0	0/0	0/0
CY5.5	0/0	0/0	0/0	10000/10000	0/0	0/0
ROX	0/0	0/0	0/0	0/0	10000/10000	0/0
APTES990	0/0	0/0	0/0	0/0	0/0	10000/10000

5. In the table, enter new values for the fluorophores.
6. Click Reprocess Data.

System Calibration and Tools

You cannot execute runs on the QX600 Droplet Reader unless the instrument is properly calibrated. Only a Bio-Rad service engineer can calibrate your instrument and use the Instrument Calibration options in QX Manager Software.


If your instrument requires service, contact Bio-Rad Technical Support.

The Tools functions are currently applicable only to instrument maintenance. For information, see [Appendix B, Instrument Maintenance](#).

Chapter 12 Managing Users

Important: This chapter applies only if you enabled the User Management module during the installation of QX Manager Software Standard Edition. If you did not enable the module, QX Manager Software creates a generic user with the ability to set up, run, and analyze experiments.

Use the User Management window to add, change, or remove QX Manager Software users, and set or change user privileges and preferences.

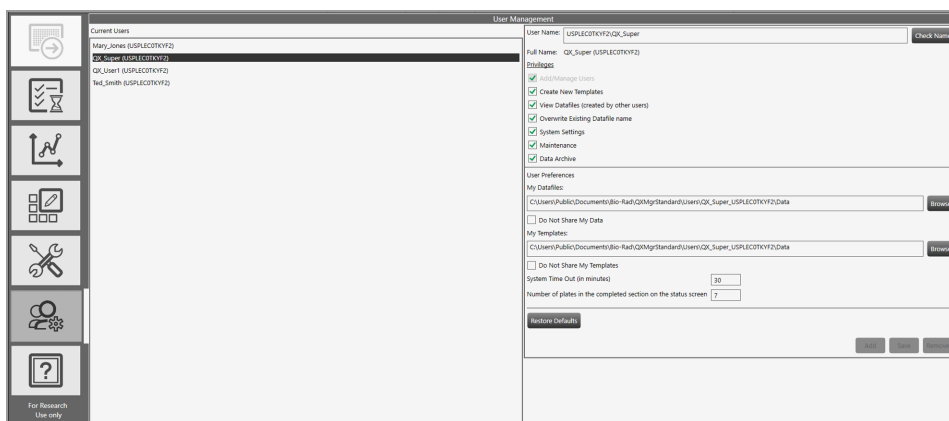

An administrative user account, with all privileges, should be set up first. You will use this user account to create additional users and assign user privileges. If you need assistance, contact Bio-Rad Technical Support. Users who are not assigned any of the specified privileges can still execute runs and use the Analysis module.

You can set up local or domain accounts. If the computer connected to the QX600 Droplet Reader

- Cannot access your company's network, your IT department or system administrator can set up user names as local Windows accounts.
- Can access your company's network, QX Manager Software can recognize the Active Directory user names.

You can create user accounts with any user name convention and password character structure, as long as it complies with your organizational requirements. The software is designed to validate each user against the name it finds in either the Active Directory or local domain, so users must log in with the proper domain and user name.

Adding Users

This section applies only if the User Management module was enabled during the installation of QX Manager Software Standard Edition.

Before you can add users to QX Manager Software, you must be logged in with the initial superuser account, or another account that is assigned the Add/Manage Users privilege.

Important: Before you begin, ensure that all users either have local Windows user accounts on the local computer or user accounts in Active Directory (network domain).

To add a user

1. Open QX Manager Software and sign into the application.
2. Click the User Management tab on the left



3. In the User Name field on the right, enter the user name as follows:
 - If the computer is connected to a network and you are using Active Directory accounts, enter <domain name>/<user name>. For example, **global/johnsmith**.

Important: Active Directory users must log in every time with the domain name preceding the user name.

 - If the computer is not connected to a network, enter <local computer name>/<user name>. For example, **analysiscomputer/johnsmith**.

Important: To locate the computer name, right-click on the yellow folder in the lower menu bar and select File Explorer. When the dialog box appears, the computer name is located at the top of the left panel. Local users are not required to reenter the computer name after the initial login.

4. Click Check Name.

One of the following occurs:

- If the software recognizes the network domain or computer name and the user name, the software displays a validation message.
- If the user is not recognized, an error message appears. Do one of the following:
 - For network users, ensure there is a working connection to the network and then verify your entry is accurate.
 - For local users, verify the entry is accurate. If the user is not set up locally or on the network, contact your system administrator for assistance.

5. When the This user is valid message appears, click OK, and then click Add at the bottom of the pane.

An Are you sure you want to add this user? message appears.

6. Click Yes.

A confirmation message appears.

7. Click OK.

8. Continue until all users are added.

Adding or Removing User Privileges

This section applies only if the User Management module was enabled during the installation of QX Manager Software Standard Edition.

Only a user assigned the Add/Manage users privilege can add or remove user privileges. By default, individual users without privileges can set up and execute a run, open files, view and analyze results.

For descriptions of all user privileges, see [Viewing User Privileges on page 41](#).

User Name: (computer or domain name)\QX_User1 Check Name

Full Name: QX_User1 (full name)

Privileges

- Add/Manage Users
- Create New Templates
- View Datafiles (created by other users)
- Overwrite Existing Datafile name
- System Settings
- Maintenance
- System Data
- Manage Plate

[Table 40](#) contains a sample set of privileges assigned by user roles common to a laboratory. User roles can be defined and used in any scenario, but might be required in regulatory environments.

Table 40. Example User Roles

Privilege	Superuser	Lab Manager	Group Lead	Technician/ Student Intern
Add/manage users	✓	✓		
Create new templates	✓	✓	✓	
View data files created by other users	✓	✓	✓	✓
Override existing ddPCR file name	✓	✓	✓	
System settings	✓			
Maintenance	✓			

Table 40. Example User Roles, continued

Privilege	Superuser	Lab Manager	Group Lead	Technician/ Student Intern
Archive data files	✓	✓	✓	✓
Access the module recovery tool	✓	✓	✓	✓
Manage Plates	✓	✓	✓	

You can assign the Create new templates privilege to non-administrative roles if anyone can create and save a template. If template formats are restricted, then you should assign this permission to administrative or supervisory roles only.

Note: Bio-Rad recommends defining at least one additional superuser as a backup, as well as other administrative and standard roles based on your needs. Guest users can open shared templates, execute runs, and perform analysis but are not typically assigned other privileges.

To add or modify privileges

1. Click the Add/Manage Users tab to open the User Management window.
2. Select a user from the Current Users list. The name appears in the User Name field. A network or computer name and backslash might appear before the user name.
3. Select or clear the checkbox for each privilege in accordance with the user's role in using the software. You can assign user privileges in any combination.
4. When the confirmation message appears, click Yes to save the changes, and then click OK to close the pop-up.

Modifying User Preferences

This section applies only if the User Management module was enabled during the installation of QX Manager Software Standard Edition. If User Management was not enabled, only the options for system timeout duration and the number of completed plates displayed in the Runs window are available. For information, see [Partial Preferences on page 44](#).

Users assigned the Add/Manage Users privilege can change the preferences for other users. Individual users can also change their own preferences.

To modify a user's preferences

1. Click the Add/Manage Users tab, and then enter the user name.
2. Change any of the following preferences for the user:
 - Enter a different file path for user data files and templates, and choose to keep them private.
Important: You can also specify Preferred Locations in System Settings, to override all file paths specified in individual preferences. For information, see [System Settings Tab on page 166](#).
 - Enter a different default system timeout period.
 - Enter a different number of completed plates to show in the Status section, up to a maximum of 100.
3. Click Save. QX Manager Software displays a confirmation message.
4. Click Yes to save the changes, and then click OK to close the dialog box.

Removing Users

This section applies only if the User Management module was enabled during the installation of QX Manager Software Standard Edition.

Only a user assigned the Add/Manage Users privilege can remove users. Removing a user from the software does not remove the user from other systems, such as Windows, or databases.

To remove a user

1. Click the Add/Manage Users tab to open the User Management window.
2. In the Current Users pane, select the user and click Remove. QX Manager Software displays a confirmation message.
3. Click Yes to remove the user, and then click OK to close the dialog box.

Appendix A Experiment Examples

This section contains setup and analysis information for the following sample experiments:

- Direct Quantification, using the Amplitude Multiplex assay type
- Copy Number Variation, using the Probe Mix Triplex assay type
- Drop-Off, using the Basic Drop-off Assay type for genome edit detection
- Direct Quantification, using the Advanced Classification Method assay type

Direct Quantification Multiplexed

You can increase the number of targets to be evaluated by using the Amplitude Multiplex option. Optionally, to expedite plate setup you can copy and paste data from one well to another, and you can click Undo to undo an applied well setting. You cannot change the supermix that was used during data collection.

Setting Up Experiment Parameters

To set experiment parameters

1. Open the analysis data file and select the Plate Editor tab.
2. Select one or more wells to be assigned well analysis attributes.
3. Change the experiment type to Direct Quantification (DQ), and then click Apply.
4. (Optional) Change the sample description identifiers.
5. (Optional) Change the sample type. The default type is Unknown.

Tip: For automatic thresholding of positive control wells, you must select Pos Ctrl.

6. Under Assay Type, select Amplitude Multiplex and click Apply. Twelve target rows appear.
7. Change the target name and type (unknown or reference) for each target in the well.

The software defaults to FAM Lo, FAM Hi, HEX Lo, HEX Hi, Cy5 Lo, Cy5 Hi, Cy5.5 Lo, Cy5.5 Hi, ROX Lo, ROX Hi, ATTO 590 Lo, and ATTO 590 Hi, respectively, for each target.

8. (Optional) Change the high and low target identifiers.

Amplitude Multiplex allows for up to two assays in each channel, indicated by Lo and Hi labels. In this multiplex mode, each target may have a signal for only one channel. You can

- Select a different dye, if the channel supports two dyes (for example change HEX to VIC or VIC to HEX)
- Select a different channel for the target (for example, change FAM to HEX)
- Change Hi to Lo or Lo to Hi

9. Click Apply.


Viewing and Adjusting Results

Automatic thresholding is not available for Amplitude Multiplex experiments.

For clusters that are easily identifiable across the chart, use Threshold Lines mode to set manual thresholds. If the data is not orthogonal and the threshold lines do not adequately divide the populations of droplets, use Threshold Clusters mode to set manual thresholds.


To set thresholds in the 2D Amplitude window

1. Select the 2D Amplitude tab.

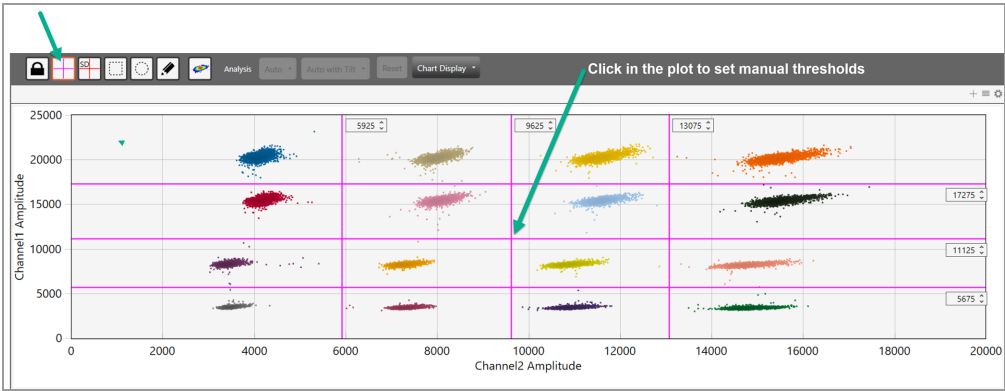

2. On the toolbar, click the Threshold Line Mode button.
3. Click anywhere on the plot to apply “best fit” threshold lines to the data.
4. To adjust the thresholds, drag the lines horizontally or vertically in the plot area.


Tip: You can also enter a specific threshold value in the boxes at the end of each threshold line.

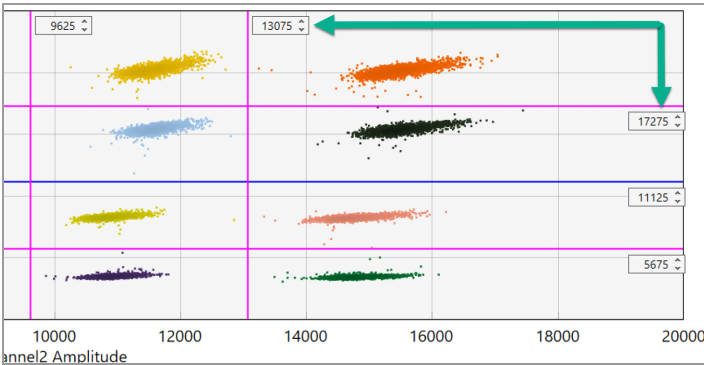

To apply and adjust manual clusters

1. Click any of the three cluster mode buttons (square, circle, or freehand).
2. Press the left mouse button and draw the chosen shape around the cluster of droplets to be classified first.

Tip: When you release the mouse button, a pop-up tool appears to help you select the correct target combination. Target names are shown along the two axis based on the signal values assigned during plate setup.

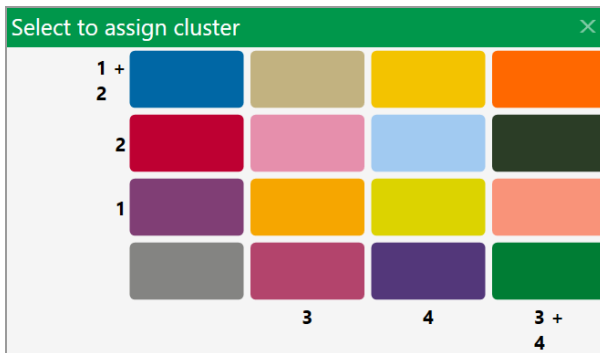

3. To classify droplets as positive or negative for each given target, select the colored box that best represents the location of the cluster of droplets.

Copy Number Variation Triplexed

A CNV experiment with probe mix triplexing uses two targets and a reference in three groups of three to produce analysis data. Optionally, to expedite plate setup you can copy and paste data from one well to another, and you can click Undo to undo an applied well setting.

Setting Up Experiment Parameters

To set experiment parameters

1. Open the analysis data file and select the Plate Editor tab.
2. Select one or more wells to be assigned well analysis attributes.
3. Set the experiment type to Copy Number Variation (CNV), and click Apply.
4. (Optional) Change the sample description identifiers.
5. (Optional) Change the sample type. The default type is Unknown.

Tip: For automatic thresholding of positive control wells, you must select Pos Ctrl.

Note: You cannot change the supermix that was used during data collection.

6. Click Apply.
7. From the Assay Type dropdown list, select Probe Mix Triplex and click Apply.

QX Manager Software displays three targets.

Important: You cannot add or delete triplex targets.

8. Enter target names and select the target type (unknown or reference) for each target in the well.

This creates three groups of three targets each.

9. Assign the targets as follows:

Group 1

Target	Target Type	Signal Ch1	Signal Ch2	Ref Copies	Plot?
Target 1	Unknown	FAM	None	N/A	N/A
Target 2	Unknown	None	HEX/VIC	N/A	N/A
Target 3	Reference	FAM	HEX/VIC	2	Yes

Group 2

Target	Target Type	Signal Ch3	Signal Ch4	Ref Copies	Plot?
Target 4	Unknown	Cy5	None	N/A	N/A
Target 5	Unknown	None	Cy5.5	N/A	N/A
Target 6	Reference	Cy5	Cy5.5	2	Yes

Group 3

Target	Target Type	Signal Ch3	Signal Ch4	Ref Copies	Plot?
Target 7	Unknown	ROX	None	N/A	N/A
Target 8	Unknown	None	ATTO 590	N/A	N/A
Target 9	Reference	ROX	ATTO 590	2	Yes

10. Click Apply.


Viewing and Adjusting Results

Note: Linear thresholding is not available for Probe Mix Triplex experiments.

To see the results

1. After all wells are configured, select the 2D Amplitude tab to evaluate the data and manually set clusters.
2. Click any of the three cluster mode buttons (square, circle, or freehand).
3. Press the left mouse button and draw the chosen shape around the cluster of droplets to be classified first.

Tip: When you release the mouse button, a pop-up tool appears to help you select the correct target combination. Target names are shown along the two axis based on the signal values assigned during plate setup.

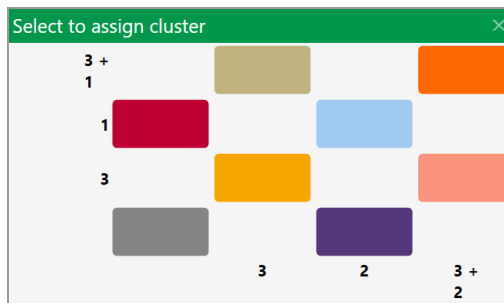

4. To classify droplets as positive or negative for each given target, select the colored box that best represents the location of the cluster of droplets.


Drop-Off Assays for Genome Edit Detection

Optionally, to expedite plate setup you can copy and paste data from one well to another, and you can click Undo to undo an applied well setting.

Two configurations of drop-off assays are possible and are shown below. Assign a signal (expected cluster position) to wild-type and NHEJ indel alleles. The signal produced by a wild-type allele will always be positive for both FAM and HEX, while the signal produced by an NHEJ indel allele depends on the assay configuration used (as shown below left and right).


Drop-Off Probe in HEX

NHEJ Probe = HEX and Reference Probe =FAM

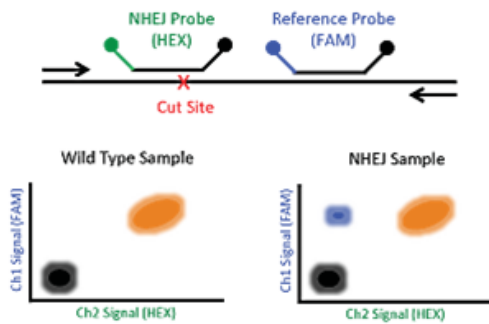

Drop-Off Probe in FAM

NHEJ Probe = FAM and Reference Probe =HEX

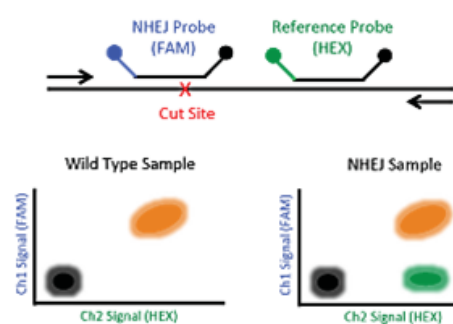

This type allows three groups of two targets per well; Group 1, shown by default, with 1 target in FAM and HEX (reference), and 1 target in FAM or HEX (unknown). You can add a second group with 1 target in Cy5 and Cy5.5 (reference) and 1 target in Cy5 or Cy5.5 (unknown) and a third group with 1 target in ROX and ATTO 590 (reference) and 1 target in ROX and ATTO 590 (unknown).

Note: The channels within each group are not interchangeable therefore FAM is always paired with HEX/VIC. Cy5 with Cy5.5. and ROX with ATTO 590.

Setting Up Experiment Parameters

To set experiment parameters

1. Open the analysis data file and select the Plate Editor tab.
2. Select one or more wells to be assigned well analysis attributes.
3. Set the experiment type to Drop-Off (DOF) and click Apply.

The screenshot shows the 'Experiment Type' dialog box with 'Drop-Off (DOF)' selected. Below this are four 'Sample Description' fields. The 'Sample Type' is set to 'Unknown'. The 'Assay Type' is set to 'Basic Drop-Off'. The 'Target Info' table is as follows:

Target Name	Target Type	Signal Ch1	Signal Ch2	Signal Ch3	Signal Ch4	Signal Ch5	Signal Ch6	Ref Copies
Target-1	Ref	FAM	HEX	None	None	None	None	2
Target-2	Unkn	FAM	None	None	None	None	None	
3	Ref	None	None	Cy5	Cy5.5	None	None	2
4	Unkn	None	None	Cy5	None	None	None	
5	Ref	None	None	None	None	ROX	ATTO590	2
6	Unkn	None	None	None	None	ROX	None	

4. (Optional) Change the sample description identifiers.
5. (Optional) Change the sample type. The default type is Unknown.

Tip: For automatic thresholding of positive control wells, you must select Pos Ctrl.

Note: You cannot change the supermix that was used during data collection.

6. Click Apply.

QX Manager Software automatically assigns the assay method as Basic Drop-Off.

7. Enter the target names, and then designate one target as Reference and the other as Unknown.

The Reference target enables ratio calculation of results.

8. Click Apply.


Viewing and Adjusting Results

Thresholding is not available for Drop Off experiments.

To apply and adjust clusters

1. After all wells are configured, select the 2D Amplitude tab to evaluate the data and manually set clusters.
2. Click any of the three cluster mode buttons (square, circle, or freehand).
3. Press the left mouse button and draw the chosen shape around the cluster of droplets to be classified first.

Tip: When you release the mouse button, a pop-up tool appears to help you select the available cluster assignments. Target names are shown along the two axis based on the signal values assigned during plate setup.

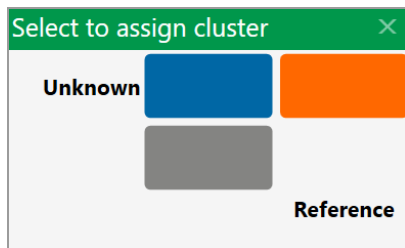

4. To classify droplets as positive or negative for each given target, select the colored box that best represents the location of the cluster of droplets.

Direct Quantification with the Advanced Classification Method


The Advanced Classification Method analysis mode allows the user complete control of droplet identification, and does not force or preassign any clusters. This allows a higher degree of multiplexing through radial dye mixing. Optionally, to expedite plate setup you can copy and paste data from one well to another, and you can click Undo to undo an applied well setting.

Important: This mode is intended for experienced users of Droplet Digital PCR.

Setting Up Experiment Parameters

To use the advanced classification method

1. Open the analysis file and select the Plate Editor tab.

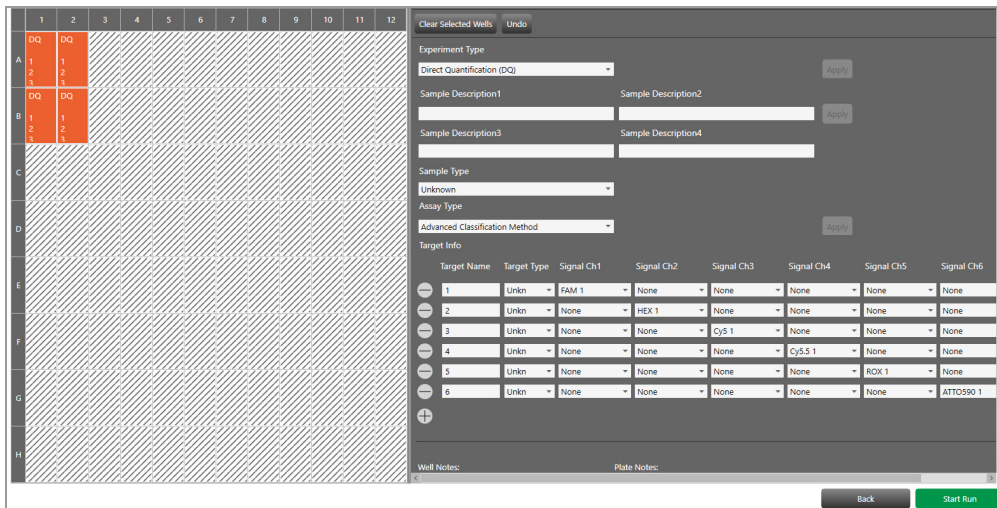

2. Select one or more wells to be assigned well analysis attributes.
3. Set the experiment type to Direct Quantification (DQ) or RED and click Apply.
4. (Optional) Change the sample description identifiers and click Apply.
5. (Optional) Change the sample type and click Apply. The default type is Unknown.

Important: For automatic thresholding of positive control wells, you must select Pos Ctrl. You cannot change the supermix that was used during data collection.

6. Under Assay Type, select Advanced Classification Method and use the + icon to add up to ten targets.

7. Enter the target name and type (unknown or reference) for each target in the well, and then click Apply.

If you select Reference, you can change the number of reference copies. The default is 2.

- If no reference is selected, the target with the highest dye concentration in the highest channel is used as the reference. For example, if there are targets in all six channels, the target in Channel 6 is used; however, if there are targets only in the first five channels, then Channel 5 is the reference target.
- If more than one reference is selected in the well, checkboxes appear to the right for using the reference in the plot.

Note: Each Reference target enables ratio calculations in the results.

8. Assign the first target an expected signal in a channel.
9. Assign your last target a signal in a channel that was not selected for your first target.
Each target in between will be assigned a signal depending on mixture ratio of the target.
10. Click Apply to confirm all well settings.

Viewing and Adjusting Results


To view and adjust results

1. Click the 2D Amplitude tab.

All droplets are unclassified by default and are colored red. Only the Threshold Cluster mode is available for setting thresholds in Advanced Classification experiments in the 2D plot.

2. Use one of the Threshold Cluster mode tools (box, circle or freehand) to assign droplet clusters to targets.

When you select a droplet cluster, a cluster assignment box appears.

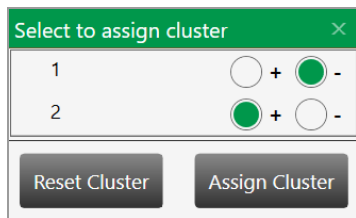

3. Select a radio button to identify the composition of the cluster.

For example, if the cluster is composed of positive droplets for Target-2, select the positive radio button for Target-2 and the negative radio buttons for the remainder.

Important: For calculations to occur, you must have at least one positive cluster for each target, as well as the assigned negative cluster.

Note: The software sets aside unassigned for calculation purposes. If you are unsure of droplet/cluster composition, leave it unassigned (red).

You can view the concentration results in the data table and in the Concentration tab.

If one of your samples was identified as a Reference under AssayType, you can view a ratio of concentration for unknown versus reference in the Ratio tab.

This page intentionally left blank.

Appendix B Instrument Maintenance

Bio-Rad recommends regular maintenance of your equipment, which includes

- Keeping the surface areas clean, post-run or as necessary

Note: Wipe down external panels and the plate holder with 70% ethanol. See [Cleaning or Decontaminating the Instrument on page 215](#) for information.

- Regularly checking the fluid levels
- Inspecting the equipment for damaged parts or wiring

Bio-Rad also recommends that a Bio-Rad service engineer perform annual preventive maintenance on your instrument. Contact Bio-Rad if your instrument requires service or recalibration.

Fluid Maintenance and Waste Management

You must ensure there is enough oil for at least one full run, and you must empty the waste container when it approaches a full state.

You can verify your fluid levels by checking the following:

- Instrument status bar
- Instrument indicator lights
- The oil and waste bottles or EcoTank

[Table 41](#) explains fluid level maintenance for your instrument.

Table 41. Fluid Levels

Fluid Level	Status
Oil greater than 40% or waste less than 70%	Bottle levels OK. You can start your run.
Oil less than 40% or waste greater than 70%	<p>You may be able to start a run with fewer than 96 wells.</p> <p>If there is not enough oil in the dispensers to perform a run on the number of wells you set up in the Plate Editor, the software prevents the run until you fill the oil dispenser and empty the waste container.</p> <p>Pay attention to the fluid level by checking instrument status bar in the QX Manager Software or by checking the oil bottles.</p>
Oil less than 10% or waste greater than 90%	The software prevents the run until you fill the oil dispenser and empty the waste container.

Changing the Droplet Reader Oil Bottle

To replace the oil bottle

1. Slide the container carriage out using the handle on the side of the instrument.
2. Remove the oil bottle and set it aside, to use when changing the waste bottle.

Once opened, a pop-up appears on your screen asking you to indicate which type of bottle you are installing on the system.

Bottle door has opened

Please select the action taking place:

New EcoTank is being loaded into instrument. (Note: The EcoTank can be used for 6 months from load date).

Batch number:

Expiration date: YYYY-MM-DD (YYYY-MM-DD)

New Oil and/or Waste bottle is being loaded into instrument.

Nothing is changing.

Please close the bottle door to continue.

OK

Note: If using an EcoTank, see [Using an EcoTank Oil and Waste Container on page 207](#) for information.

3. Insert a new bottle of oil and tighten the cap into place.
4. Prime and flush the instrument at least once after the new oil bottle is installed. See [Priming and Flushing the Instrument on page 213](#) for information.

Note: If you are replacing reagents in the middle of a run, the run resumes when the bottle door is closed and you have clicked OK in the dialog box. You can replace with the same type of bottle only. For example, if the run started with an EcoTank loaded, you must replace the spent EcoTank with another EcoTank.

After an EcoTank is loaded into the system, the software tracks the number of plates that have been run with the EcoTank. Once you have run 30 plates, the software begins to warn you to replace the EcoTank. After you have run 40 plates, you are required to replace the EcoTank before proceeding with a new run.

Emptying the Waste Container

Important: Use each empty oil bottle as a new waste bottle once and then discard it. Before discarding, add 50 ml of 10% bleach solution to the bottle to prevent microbial growth, and then label the bottle as waste. If you are using an EcoTank, see [Using an EcoTank Oil and Waste Container on page 207](#) for information.

To replace the waste bottle

1. Slide the container carriage out using the handle on the side of the instrument and remove the waste bottle.
2. Place the new Waste label over the Oil label on the oil bottle.
3. Insert the last empty oil bottle into the waste container slot.
4. Flush the instrument after the new waste bottle is installed. See [Priming and Flushing the Instrument on page 213](#) for information.

For information on disposing of waste, see [Managing Waste on page 214](#).

Using an EcoTank Oil and Waste Container

EcoTanks are high-capacity reservoirs that you can use instead of individual oil and waste bottles in your QX600 Droplet Reader.

The ddPCR Droplet Reader Oil EcoTank for Probes distributes and recirculates droplet reader oil, and also collects waste generated by the instrument into a single partitioned container. This ensures that floating waste droplets remain on the waste side of the bottle, while the droplet reader oil can be redistributed through the instrument for another droplet reading cycle.

Usage Advisories

Note the following when using the Droplet Reader Oil EcoTank for Probes:

- The EcoTank can read up to 40 plates.
- Do not add bleach to the EcoTank before or while it is installed in the instrument. You can add bleach after you remove the EcoTank from the instrument and before storage or disposal.
- Do not reuse a spent EcoTank.
- An unused EcoTank is stable for up to 12 months when stored at room temperature (25°C).
- You must replace the EcoTank before or at 6 months after date you install it in the instrument. QX Manager Software tracks the EcoTank and locks out the user after the EcoTank passes the 6-month limit.
- The EcoTank for Probes is not compatible with QX200 ddPCR EvaGreen® Supermix (Catalog No. 1864033, 1864034, 1864035, 1864036).

Installing an EcoTank Bottle in the Droplet Reader

To install an EcoTank bottle

1. Open QX Manager Software.
2. Open the reagent door on the left side of the QX200 Droplet Digital PCR System and place an EcoTank in front of the currently installed oil bottles.

Note: The EcoTank is symmetrical and new bottles can be loaded in either orientation. Bio-Rad does not recommend changing the EcoTank orientation after it has been installed in the droplet reader.

3. Remove the caps from the EcoTank.
4. Remove from the currently installed bottles the caps that are attached to the instrument, and attach them to the EcoTank.
5. Cap and remove the currently installed bottles.
6. Place the EcoTank in the bottle tray. Maintain the orientation to prevent twisting of the oil lines.

After the bottle door has opened, QX Manager Software prompts you to select the action taking place. Continue to the next section.

Selecting the EcoTank in QX Manager Software

To specify an EcoTank is being used

1. Select *New EcoTank is being loaded into instrument*.
2. Enter the batch number and expiration date, and then click OK.

Important: You must close the bottle door to enable the OK button.

Bottle door has opened

Please select the action taking place:

New EcoTank is being loaded into instrument. (Note: The EcoTank can be used for 6 months from load date).

Batch number:

Expiration date: (YYYY-MM-DD)


New Oil and/or Waste bottle is being loaded into instrument.

Nothing is changing.

Please close the bottle door to continue.

OK

In the Instrument Status bar, QX Manager displays the EcoTank for the droplet reader and waste oils, as well as the number of plates remaining.

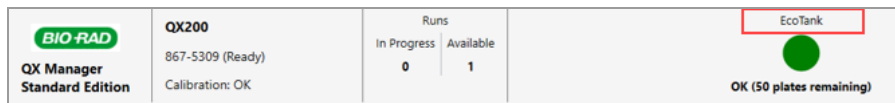

Important: If the instrument has been idle for more than a week, you must prime the instrument before running a plate. From the Setup menu, click Tools > Prime to fill the lines with oil before starting a run.

Reinstalling the Standard Bottles

You can return the QX600 Droplet Reader to the initial two bottle configuration at any time, except when a run is in progress.

Complete the steps in the following sections to change the bottle configuration.

To exchange the EcoTank for the two-bottle configuration

1. Open QX Manager Software.
2. Complete the following substeps to prepare a new 1 L waste bottle for the QX600 Droplet Reader:
 - a. Add 50 ml 10% bleach to the Waste bottle to prevent microbial growth.
 - b. Place an identifying label on the Waste bottle.
3. Open the reagent door on the left, and then place a droplet reader oil bottle and the new waste bottle into the space in front of the EcoTank bottle.
4. Remove the caps from the EcoTank and bottles.
5. Replace the oil and waste caps on the used EcoTank bottle, and then use the handle on the EcoTank bottle to move it out of position and remove it from the instrument.
6. Place the oil and waste bottles into the corresponding slots and close the reagent door.

After the bottle door has opened, QX Manager Software prompts you to select the action taking place. Continue to the next section.

To specify that separate oil and waste bottles are being used in QX Manager Software

- ▶ Select *New Oil and/or Waste bottle is being loaded into the instrument* and click OK.

Bottle door has opened

Please select the action taking place:

New EcoTank is being loaded into instrument. (Note: The EcoTank can be used for 6 months from load date).

Batch number:

Expiration date: (YYYY-MM-DD)

New Oil and/or Waste bottle is being loaded into instrument.

Nothing is changing.

Please close the bottle door to continue.

OK

QX Manager displays the DR Oil and Waste indicators.

BIO-RAD QX Manager Standard Edition	QX600 8355839 (Ready)	Runs In Progress: 0 Available: 0	DR Oil OK	Waste OK	Disk Space (C:) Free: 46 GB (3 Plates) Used: 71 GB (5 Plates)	 BESAC 7/8/2021 12:33 PM
--	---------------------------------	--	------------------	-----------------	---	--------------------------------

Important: If the instrument has been idle for more than a week, you must prime the instrument before running a plate. From the Setup menu, click Tools > Prime to fill the lines with oil before starting a run.

Disposing of Waste

Droplet generation and reader fluids from Bio-Rad are based on fluorinated hydrocarbon chemistry and should be disposed of in accordance with institutional, state, and local regulations. For information, see [Managing Waste on page 214](#).

Disposing of the EcoTank and EcoTank Waste

When QX Manager indicates 40 plates have been run, you must replace the used EcoTank bottle. After removing the EcoTank from the instrument, you can install a new EcoTank or reinstall the standard two-bottle configuration. If the EcoTank bottle is removed before 40 plates are run, discard the EcoTank bottle and oil in accordance with institutional, state, and local regulations.

To remove and dispose of the EcoTank and EcoTank Waste

1. Uncap one side of the EcoTank and take it to the liquid waste receptacle.
2. Hold the uncapped opening slightly lower than the capped opening and slowly pour out as much of the contents from the uncapped side as possible.
3. Return the bottle to its upright position to allow the fluid trapped in the closed cap side of the bottle to move to the open cap side, and repeat the process.
4. Pour entire contents of spent bottle into the waste receptacle.
5. Place the cap on the empty EcoTank bottle and then discard per waste disposal regulations.

Cleaning the EcoTank Waste Cap

Repeat the following process once per month at a minimum.

To clean the waste cap

1. Dampen a lint-free wipe with fresh 10% bleach solution.
2. Remove the cap on the waste side of the EcoTank.
3. Wipe down anywhere microbial growth might occur, including along the threading and inside of the cap .
4. If the cleaning leaves excess liquid on the cap, dry the cap using a dry lint-free wipe.
5. After the cleaning is done and the cap is dry, place and tighten the cap on the EcoTank.

Priming and Flushing the Instrument

You should prime the instrument after you replace the oil bottle and you should prime and flush the instrument when you change the waste bottle.

Important: Before you initiate a prime or flush, *the lid must be closed*. Do NOT press the button to open the lid while a prime or flush is in progress.

To prime the instrument

1. Select the System Utilities tab.
2. Tap Prime.

A message appears, asking if you want to prime the instrument.

3. Tap Yes to begin.
4. A “Performing Prime” message appears.

When the priming process is finished, an operation complete message appears.

5. Tap OK.

To flush the instrument

1. Select the System Utilities tab.
2. Tap Flush.

A message appears, asking if you want to flush the instrument.

3. Tap Yes to begin.
4. A “Performing Flush” message appears.

When the flushing process is finished, an operation complete message appears.

5. Tap OK.

Managing Waste

You must apply standard Material Safety Data Sheet (MSDS) and OSHA practices when handling and disposing of generated waste. A typical waste profile should contain the following:

Table 42. Managing Waste

Chemical name	Weight %
Fluorinated fluids	>50%
Water	<50%
Bleach	<10%
Other materials (proteins, nucleic acids, fluorescent dye)	<10%

Note: The chemical identities and exact percentages in the above table are withheld as trade secrets.

Bio-Rad droplet generation and reader fluids are based on fluorinated hydrocarbon chemistry, and should be disposed of in accordance with institutional, state and local regulations. Avoid release to the environment and prevent entry into sewer systems or bodies of water. As a disposal alternative, incinerate in a permitted high-temperature waste incineration facility.

These nonflammable fluids are inert and have low environmental impact and low toxicity.

Toxicological Information

Chemical name: (mixture)

Oral LD50: >5000 mg/kg (Rat)

Dermal LC50: >5000 mg/kg (Rabbit)

Inhalation LD50: >8838 ppm (Rat/8-hr)

Ecological Information

Acute aquatic hazard: LC50 > 500 mg/l (96-hr./static; Pimephales promelas/Fathead minnow)

Chronic aquatic hazard: No information available (GHS Chronic 4)

Recovery From Idle Mode

Complete the steps below if the instrument has been idle:

1. Check the expiration date on the oil bottle. If expired, replace the oil bottle.
2. Ensure the waste bottle is not full.
3. If needed, add 50 ml of 10% bleach to the waste bottle.
4. If you have changed the bottles, or if you have not used the system for a week or more, perform the prime and flush processes as specified in [Priming and Flushing the Instrument on page 213](#).

Cleaning or Decontaminating the Instrument

Use the protocols specified below for the following:

- Regular post-run instrument cleanup
- Instrument decontamination, as needed

If you have questions regarding the use of cleaning agents other than those specified, contact Bio-Rad Technical Support.

To clean the instrument

1. Remove the plate after a run is complete and ensure it is properly discarded.
2. Using 70% ethanol, wipe down the external panels, as well as the plate holder and its compartment.

To decontaminate the instrument

1. Remove the plate after a run is complete and ensure it is properly discarded.
2. Wipe down the plate holder and its compartment with 10% bleach.

Important: Do not use bleach inside the system.

3. Wipe down all external surfaces with 10% bleach.
4. Identify the equipment as *Decontaminated*.

Instrument Maintenance Log

Serial Number	Last Annual Maint Date																																														
		QX600 Droplet Reader System Maintenance Log		Month	Year																																										
As necessary																																															
Restart the droplet reader																																															
Restart the connected computer and software																																															
Replace droplet reader oil																																															
Wipe droplet reader plate holder (if wet)																																															
Discard waste bottle after one-time use																																															
Use empty oil bottle as new waste bottle																																															
Add 50 mL of 10% bleach to new waste bottle																																															
Initials																																															
Daily																																															
Clean work area (before and after use)																																															
Discard used consumables																																															
Initials																																															
Weekly (at start of week)																																															
Check droplet reader oil expiration date																																															
Prime and flush the system (if the system has been idle)																																															
Initials																																															
Monthly																																															
Clean metal surface of instrument plate holder with 70% ethanol																																															
Initials																																															
Quarterly																																															
Perform a database backup																																															
Initials																																															
Annually																																															
Call Bio-Rad																																															
Initials																																															

Appendix C Troubleshooting

This section contains information on troubleshooting and correcting minor issues you may experience with your instrument or software.

Event Log

The Event Log contains records for all system and user activities, including system errors. If you do not understand, or cannot find an immediate remedy to, the error event, contact Bio-Rad Technical Support.

Instrument Errors

If an error occurs during a run, the software cancels the run and displays an X in the Status column, along with a Run Aborted by System message as the Run End status. [Table 43](#) contains error grouped by type.

1. Click the X to display the corresponding error code and description.
2. Perform the steps in the solution column before starting a new run.
3. If the suggested solution does not resolve the problem, contact Bio-Rad Technical Support immediately.

Note: The error code and description are also displayed in the Run Information window .

Communication errors can be caused by instrument, connectivity, or power failures.

Table 43. Communication errors

Error Code	Description	Solution
102, 103, 104*, 110*, 111	Communication loss between the instrument and software, detector board error, loss during a run, or unable to reestablish a connection Note: For codes marked with an asterisk (*), some data may be usable.	Disconnect and reconnect the USB cable, and then power cycle the instrument.
302, 308	Communication error involving pump Note: Some data might be usable.	Power cycle the instrument.

Table 43. Communication errors, continued

Error Code	Description	Solution
502	Invalid checksum value	Retry.
900	Invalid command	

An electrical error is typically triggered by an invalid setting.

Table 44. Electrical error codes

Error Code	Description	Solution
406	Electrical Error; invalid power value setting. Note: Some data might be usable.	Power cycle the instrument.

Mechanical errors can be caused by fluid issues (low oil, high waste), instrument operation errors, or binary data collection errors.

Table 45. Mechanical error codes

Error Code	Description	Solution
105, 106	Fluidics or motion response was invalid. Note: Some data may be usable.	Power cycle the instrument.
201	Motors disabled Note: Some data may be usable.	<ol style="list-style-type: none"> 1. Remove and reinsert the plate and plate lid. 2. Check for incorrect or multiple heat seal foil layers. 3. Check for strain on all motor axes. 4. Power cycle the instrument.
205, 207, 210, 211, 212, 218	X, Y, or Z-axis, unable to reach target location or home in allotted time. Note: Some data might be usable.	Check for motor obstructions, and then power cycle the instrument.
301, 303, 304, 305, 307, 312, 316, 317	Pump errors Note: Some data might be usable.	Power cycle the instrument.

Appendix D Ordering Information

This appendix contains descriptions and catalog numbers for new or replacement instruments, accessories, and consumables for Bio-Rad ddPCR products, including your QX600 Droplet Reader.

ddPCR System and Instrument Packages

Table 46. ddPCR system and instrument packages

Product	Description	Catalog Number
QX200 Droplet Digital PCR System	Includes the QX200 Droplet Generator, QX200 Droplet Reader, computer, software, and consumables	1864001
QX600 Droplet Digital PCR System	Includes the QX200 Droplet Generator, QX600 Droplet Reader, computer, software, and consumables	17007769
QX200 AutoDG™ Droplet Digital PCR System	Includes the AutoDG, QX200 Droplet Reader, computer, and QX Manager Software Standard Edition	1864100
QX600 AutoDG Droplet Digital PCR System	Includes the AutoDG, QX600 Droplet Reader, computer, QX Manager Software Standard Edition, and consumables	17008371
Automated Droplet Generator	Includes the AutoDG (1), cooling block accessory (1), oil purge reservoir (1), and power cord (1)	1864101
QX200 Droplet Generator	Includes the droplet generator (1), cartridges and gaskets (24 each), cartridge holders (2), and power cord (1)	1864002
QX200 Droplet Reader	Includes the droplet reader, two plate holders, USB cable, and power cord	1864003
QX600 Droplet Reader	Includes the droplet reader, two plate holders, USB cable, and power cord	12013328

Thermal Cyclers and Plate Sealer

Table 47. Thermal cyclers and plate sealer

Product	Description	Catalog Number
C1000 Touch Thermal Cycler with 96-Deepwell Reaction Module	Includes the C1000 Touch Thermal Cycler chassis, 96-deepwell reaction module, USB flash drive	1851197
PTC Tempo Deepwell Thermal Cycler	Includes the thermal cycler, power cord, USB cable, and Ethernet cable	12015392
PX1 PCR Plate Sealer	PCR plate sealer, includes heat sealing instrument, plate support block that holds 96-well and 384-well plates, sealing frame, power cord	1814000
Pierceable Foil Heat Seal	Heat seal for the PX1 PCR Plate Sealer	1814040

AutoDG Accessories

Table 48. Automated Droplet Generator accessories

Product	Description	Catalog Number
Cooling block accessory	Prevents evaporation during droplet generation	12002819
Oil purge reservoir	Collects oil waste from priming and flushing Note: Contact Bio-Rad Technical Support for replacement information.	N/A
Power cord	Connects the AutoDG to the power source Note: Contact Bio-Rad Technical Support for replacement information.	N/A

QX200 Droplet Generator Accessories

Product	Description	Catalog Number
DG8 droplet generator cartridges and gaskets	Microfluidic cartridges used to mix sample and oil for droplet generation, and gaskets to seal the cartridges Note: You must seal the cartridge to prevent evaporation and to apply required pressure for droplet generation.	1864007
Droplet generator cartridge holder	Positions and holds the droplet generator cartridge in the instrument compartment	1863051
Power cord	Connects the QX200 Droplet Generator to the power source Note: Contact Bio-Rad Technical Support for replacement information.	N/A

QX600 Droplet Reader Accessories

Table 49. QX600 Droplet Reader accessories

Product	Description	Catalog Number
Computer	Computer that connects to the QX600 Droplet Reader for data collection and analysis	12017458
USB cable and power cord	Cable connecting the computer to the instrument and power cord connecting the instrument to the power source Note: Contact Bio-Rad Technical Support for replacement information.	Included
Droplet reader plate holders (2)	Used to position the 96-well plate in the droplet reader plate compartment	12006834

QX200 Droplet Reader Accessories

Table 50. QX200 Droplet Reader accessories

Product	Description	Catalog Number
Computer	Computer that connects to the QX200 Droplet Reader for data collection and analysis	12017458
USB cable and power cord	Cable connecting the computer to the instrument and power cord connecting the instrument to the power source Note: Contact Bio-Rad Technical Support for replacement information.	Included
Droplet reader plate holders (2)	Used to position the 96-well plate in the droplet reader plate compartment	12006834

ddPCR Consumables and Other Materials

Table 51. Consumables and other materials

Product	Description	Catalog Number
ddPCR 96-Well Plates	Clear well/clear shell semi-skirted plates	12001925 (pkg of 25)
DG32 Droplet Generation Cartridges	Microfluidic cartridge with gasket for the AutoDG	1864108 (pkg of 30) 1864109 (pkg of 60)
Pipet Tips	Pipet tips for the AutoDG	1864120 (pkg of 20) 1864121 (pkg of 40)
Pipet Tip Bins	Pipet tip waste bins for the AutoDG	1864125
Automated Droplet Generation Oils	Automated droplet generation oils for the AutoDG: ■ Automated Droplet Generation Oil for Probes ■ Automated Droplet Generation Oil for EvaGreen®	1864110 1864112
Droplet Generation Oils	Droplet generation oil for the QX200 Droplet Generator: ■ Droplet Generation Oil for Probes, 10 x 7 ml ■ Droplet Generation Oil for EvaGreen®, 2 x 7 ml ■ Droplet Generation Oil for EvaGreen®, 10 x 7 ml	1863005 1864005 1864006
Rainin Pipets	■ 20 µl for sample loading ■ 50 µl for droplet transfer ■ 8-channel, 200 µl for oil	L-20, L8-20 L-50, L8-50 L8-200
Rainin Pipet Tips	Filtered	GP-L10F GP-L200F
Foil Plate Seals for PX1	Pierceable foil plate seals for PX1 Plate Sealer	1814040
ddPCR Droplet Reader Oil	Droplet reader oil for the QX600 Droplet Reader and QX200 Droplet Reader	1863004
Droplet reader waste bottle	You can use an empty droplet reader oil bottle to collect the waste from droplet reading.	N/A

ddPCR Supermixes

Table 52. ddPCR supermixes

Supermix	Description	Catalog number	
ddPCR EvaGreen® Supermix	2x supermix, for use in nucleic acid sample preparation with the QX600 and QX200 ddPCR systems	1864033	2ml (2 x 1ml)
		1864034	5ml (5 x 1ml)
		1864035	25ml (25 x 1ml)
		1864036	50ml (50 x 1ml)
ddPCR Supermix for Probes (no dUTP)	2x supermix, for use in nucleic acid sample preparation with the QX600 and QX200 ddPCR systems	1863023	2ml (2 x 1ml)
		1863024	5ml (5 x 1ml)
		1863025	25ml (25 x 1)
ddPCR Supermix for Probes	2x supermix, for use in sample preparation with the QX600 and QX200 ddPCR systems	1863026	2ml (2 x 1ml)
		1863010	5ml (5 x 1ml)
		1863027	25ml (25 x 1ml)
		1863028	50ml (50 x 1ml)
ddPCR Multiplex Supermix	4x supermix especially suited for probe-based detection of multiple targets in DNA samples using the QX600 and QX200 ddPCR systems	12005909	1.2ml (2 x 0.6ml)
		12005910	3 ml (5 x 0.6 ml)
		12005911	12.5 ml (5 x 2.5 ml),
ddPCR Supermix for Residual DNA Quantification	2x supermix, for use in residual DNA detection with the QX600 and QX200 ddPCR systems	1864037	2ml (2 x 1ml)
		1864038	5ml (5 x 1ml)
		1864039	25ml (25 x 1ml)
		1864040	50ml (50 x 1ml)
One-step RT- ddPCR Advanced Kit for Probes	200 or 500x 20 µl reaction kit, for absolute quantification of target RNA in a one-step format with the QX600 and QX200 ddPCR systems	1864021	2ml (2 x 1ml)
		1864022	5ml (5 x 1ml)

Buffer Controls

Table 53. Buffer controls

Buffer	Catalog Number
ddPCR Buffer Control Kit for Probes	1863052
QX200 Buffer Control Kit for EvaGreen®	1864052

ddPCR Kits

Table 54. ddPCR Kits

Product	Description	Catalog number
ddPCR Mutation Screening Kits		
ddPCR KRAS G12/G13 Screening Kit		1863506
ddPCR KRAS Q61 Screening Kit		12001626
ddPCR BRAF V600 Screening Kit		12001037
ddPCR NRAS Q61 Screening Kit		12001006
ddPCR NRAS G12 Screening Kit		12001094
ddPCR NRAS G12/G13 Screening Kit		12001627
ddPCR EGFR Exon 19 Deletions Screening Kit		12002392
All kits include 20x multiplex assay and 2x ddPCR supermix for probes (no dUTP).		
ddPCR Residual DNA Quantification Kits		
ddPCR CHO Residual DNA Quantification Kit		17000031
ddPCR E.coli Residual DNA Quantification Kit		17000032
200x 20 µl reactions, includes 20x CHO or E.coli RDQ assay and 2x ddPCR Supermix for Residual DNA Quantification.		
ddPCR Copy Number Determination Kits		
ddPCR SMN1 Copy Number Determination Kit		1863500
ddPCR SMN2 Copy Number Determination Kit		1863503
200x 20 µl reactions, includes assay at 20x concentration, 2x ddPCR Supermix for Probes (no dUTP) and positive controls		
ddPCR Library Quantification Kit		
ddPCR Library Quantification Kit for Illumina TruSeq		1863040
200x 20 µl reactions, includes 1 vial of primers and probes at 20x concentration, 2x ddPCR Supermix for Probes (no dUTP) and positive controls		

ddPCR Assays

Table 55. ddPCR Assays

Product	Catalog number	
ddPCR HDR Gene Edit Assay	12002312	100 rxns
ddPCR HDR Gene Edit Assay	12002313	500 rxns
ddPCR HDR Gene Edit Package	12003796	1,000 rxns
ddPCR HDR Ref Assay, Predesigned	12003805	100 rxns
ddPCR HDR Ref Assay, Predesigned	12003806	500 rxns
ddPCR HDR Ref Package, Predesigned	12003793	1,000 rxns
ddPCR NHEJ Gene Edit Assay	12002314	100 rxns
ddPCR NHEJ Gene Edit Assay	12002315	500 rxns
ddPCR NHEJ Gene Edit Package	12003794	1,000 rxns



**Bio-Rad
Laboratories, Inc.**

Life Science
Group

Website bio-rad.com **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399
Canada 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23
France 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300 **Israel** 0 3 9636050
Italy 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 080 007 7373 **Luxembourg** 00 800 00 24 67 23 **Mexico** 52 555 488 7670
The Netherlands 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23 **Portugal** 00 800 00 24 67 23
Russian Federation 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23 **Spain** 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23
Switzerland 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311 **United Arab Emirates** 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23